Selladora de bolsas SEAL-1000

Bag sealing machine SEAL-1000

SEAL-1000

4101712

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1 Packing list and accessories

Once the unit is unpacked, visually check, the packing list (See 1.1) and that there are no parts damaged during transportation. Save the packing for some days for later sending.

1.1 Package contents

- Sealing machine
- Output tray
- 2x Paper roll stand (black)
- 2x Inox rod Ø 6
- 2x Inox rod Ø 11
- 4x Buffer rules
- 16x Rollers Ø 11
- 16x Rollers Ø 16
- Power cord
- Sealing lever
- 4x Screw M4x7
- 4x Screw M4x25
- 4x Screw M5x10
- 2x Spare cutters

1.2 Accessories

Sterilizing paper roll

From polipropilene and polyester, over paper core.
Sterilization status indication ink.

| Wide | Content | Code |
|-------|---------|----------------|
| 7.5cm | 100m | 4100075 |
| 10cm | 100m | 4100100 |
| 20cm | 100m | 4100200 |

Cutting paper roll stand

Code 4100045

Capacity for 2 rolls of 20cm or 4 rolls of 10cm.
Circular cutter shape.

Size: Height: 130mm Width: 510mm Depth: 320mm

Date and caducity stamp

Code 4100046

Sterilization date and caducity mark.

2 Safety cautions

- The sealing area could reach dangerous hot temperatures. Handle the unit only for the indicated parts, lever and cutter knob.
- Choose a workspace, flat, levelled and let a free space around it.
- Avoid that water splashes drops over the sealing machine.
- Perform some tests to set the sealing temperature. An excess or a defect of temperature could reduce the quality of the sealing.
- This equipment must only be operated by personnel trained and suitable qualified to work in a sterilization facilities and its common hazards.
- Not suitable to use on flammable or explosive ambients.

3 Overview

The sealing machine SEAL-1000 has been designed to seal paper/polipropilene or paper/polyester to wrap goods to be sterilized on an autoclave.

It provides a wide sealed area and temperature can be selected to suit any material.

It has a two-side cutter for a high quality cut.

4 Specifications

| | |
|------------------------|---------------------------------------|
| Power supply: | AC220/110 50Hz (See plate) |
| Power: | 500W |
| Fuse: | T5A |
| Max. paper roll width: | 250mm |
| Sealing wide: | 10mm |
| Ambient conditions: | 5 ... 40°C Humity (Hr) Lower than 85% |
| Dimensions: | 430 x 340 x 250mm |
| Weight: | 5.5Kg |

5 Warranty

The warranty covers the defective components or faulty assembling of the unit for 1 year since the purchasing date. Keep purchasing bill as the only valid document for the purchasing date.

The warranty does not cover damages caused for a unit malfunction. The unit must be suitably qualified before used for valuable samples processing.

Return goods are not accepted without previous J.P. SELECTA, s.a.u. approval.

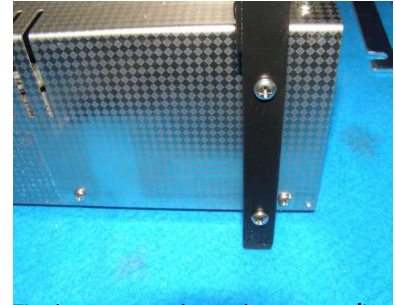
6 Initial assembling

The sealing machine comes disassembly for an easy transportation and packaging.


Locate the two black stands and the 4 screws M4x25 mm.

Fix the two stands on the rear sealing machine body.

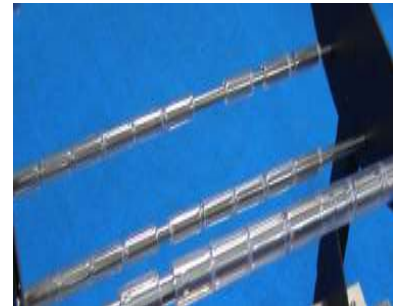Place 16 rollers on each rod.

Fix the two smaller rods on the black stands lower holes with the M4x7 screws.

Fix the two bigger rods on the black stands upper holes with the M5 screws.

Insert the sealing lever on his squared shape fixing hole.

Screw the cutter knob

Fix the output tray into the two front side pivots.

7 Starting-up the sealing machine

Place the unit over a flat surface, levelled, stable and suitable for the unit weight. A mains socket near than 1.5m will be necessary.

To work comfortable, let a free space of about 50cm around the sealing machine.

7.1 Connection to mains

Locate a power supply socket (rated to 16A) near to the unit. Use the power cord supplied or another of similar specifications. For safety reasons, the socket must have ground terminal, since the unit must work grounded.

Before plugging the mains, check than voltage matches the one indicated on the unit plate.

7.2 Sealing machine set up

1. Loading paper roll

Place the paper roll on the stand and the buffer rules to avoid the side moving.

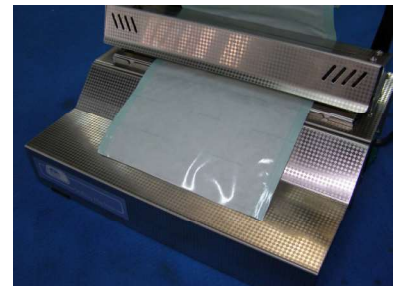
Insert the paper in the sealing head form the rear side.

Place the buffer rules according to roll size.

Inserting the paper to the sealing head.

Pull the paper from the head front side.

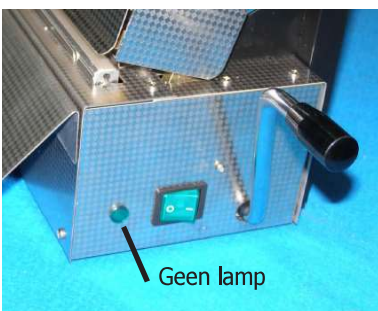
Sealing temperature selection knob.

2. Set the sealing temperature

Use the sealing temperature selection knob to adjust the optimal temperature according paper type. We suggest to run some tests before setting the proper temperature.

As a general rule, set the knob between point number 3 and 4.

It is not recommended to use a temperature knob set over point 4.

Wait for green lamp becomes Off to perform the seal.

3. Wait the indication «Ready to seal»

Once the mains switch is On, the unit starts to heat the sealing area. During this time the green lamp is On.

When the green lamp becomes Off, the unit will be ready to seal.

8 Operation

8.1 Sealing and cutting

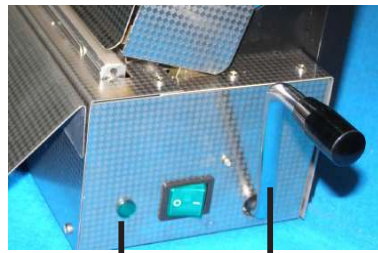
The sealing machine has two main functions:

Sealing:

Once mains switch is ON, wait until the green lamp becomes Off to press the lever. Keep the lever low around 4 second to perform the seal.

Cutting:


To cut the paper, move the cutter knob from one side to the other. It is provided with a double cutting edge.

Green lamp


Sealing lever

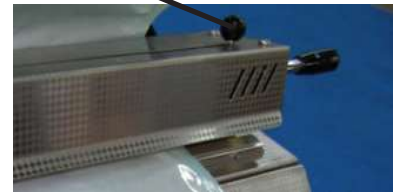Cutter knob.

8.2 Wrapping goods

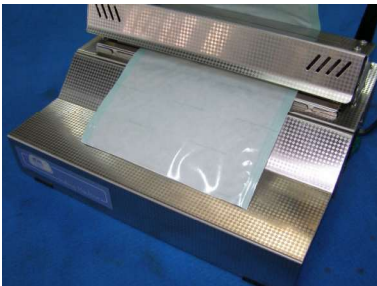
Clean the load properly before wrapping it and proceed as follows:


Pull the paper until it gets the suitable piece of paper to wrap the goods. Seal and cut.

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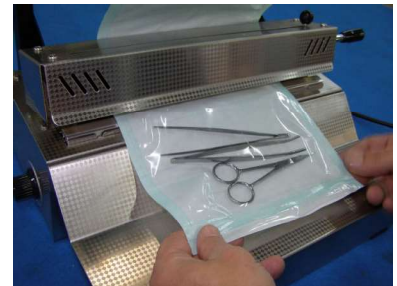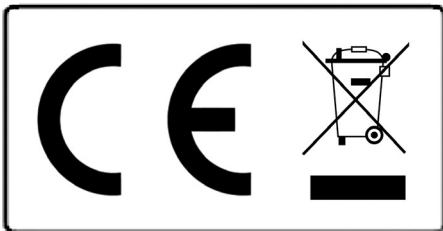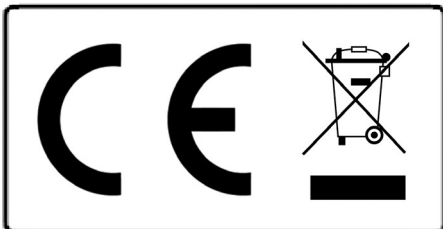
**Notice to customers:**

The product is made up of various components and various materials that must be recycled or, failing that, deposited in the corresponding debris removal sites when the product's life has been completed or when otherwise it is necessary to dispose of it. To do this, the end user who acquires the product must know the current regulations of each municipality and / or locality based on the waste electrical and electronic equipment. The user who acquires this product must be aware of and responsible for the potential effects of the components on the environment and human health as a result of the presence of hazardous substances. Never place the product in a conventional container of citizen scope if a previous dismantling and knowledge of the components that incorporates. If you do not know the procedure to follow, consult with the city council for more information.