**MEDIDOR MULTIPARAMÉTRICO MP-3006**  
***MULTI-PARAMETER MP-3006***

**4120702**

# Contenido

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## 5.4.5 Configuración de parámetros de pH

### 5.4.5.1 Información del electrodo de pH

The screenshot shows the 'Parameter Setting\pH Parameter' screen. The left sidebar contains menu items: pH Electrode, pH Buffer Group, Resolution, pH Alarm Setting, Calibrate Reminder, and pH STD Check. The main area displays the following information:

- Name:** pH Electrode
- Serial No.:** 12345678
- Recognition:** Auto Mode
- CAL Result:** A table showing calibration results for four buffer points (BUF1-BUF2, BUF2-BUF3, BUF3-BUF4, BUF4-BUF5). All show 100.00% slope and 0.00mV offset, with a 'Perfect' comment.
- CAL Time:** 2022/05/08 16:23:43
- User ID:** REX Team

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The screenshot shows a software interface for configuring DO parameters. The title bar reads 'Parameter Setting\DO Parameter'. The user is logged in as 'Admin' at '11:53:19' on '2022/05/20'. The left sidebar has menu items: 'DO Electrode', 'Compensation', 'DO Alarm Setting', 'Calibrate Reminder', 'DO Air-Sat Check', and 'DO Zero Check'. The main area displays the following information:

- Name: DO Electrode
- Serial No.: 12345678
- CAL Result: A table with 4 columns: No., Name, Result, Temp.
  - Row 1: 1, Zero Point, 0.00, 25.0°C
  - Row 2: 2, Air-Sat, 500.0, 25.0°C
- CAL Time: 2022/05/20 11:46:47
- User ID: REX Team

An 'Enter' button is located at the bottom right.

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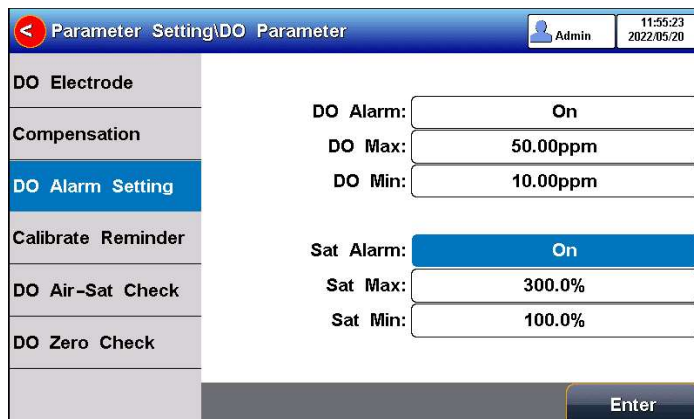
The screenshot shows the same software interface as before, but with the 'DO Alarm Setting' menu item selected in the sidebar. The main area displays the following alarm settings:

- DO Alarm: On
- DO Max: 50.00ppm
- DO Min: 10.00ppm
- Sat Alarm: On
- Sat Max: 300.0%
- Sat Min: 100.0%

An 'Enter' button is located at the bottom right.

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The screenshot shows the 'Parameter Setting\DO Parameter' screen. The left sidebar has a menu with 'Calibrate Reminder' selected. The main area contains the following settings:

DO Electrode	DO CAL Message:	On
Compensation	CAL Interval:	10
DO Alarm Setting	Time Unit:	Hour
Calibrate Reminder	DO Compulsive CAL:	On
DO Air-Sat Check	Validity:	10
DO Zero Check	Time Unit:	Hour

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The screenshot shows the 'Parameter Setting\DO Parameter' screen with 'DO Zero Check' selected in the sidebar. The main area contains the following settings:

DO Electrode	DO Zero Check	On
Compensation	Check Interval:	30
DO Alarm Setting	Time Unit:	Day
Calibrate Reminder	STD Value	0.00ppm
DO Air-Sat Check	Delta Value:	0.50ppm
DO Zero Check	Check Time:	Not checked
	Conclusion:	Not checked

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The screenshot shows a software interface titled "Database" with a user profile "Admin" and a timestamp "14:12:54 2022/05/20". Below the title bar are five tabs: "Filter", "Results", "Report", "Statistic", and "Output". The "Filter" tab is active. The interface contains several input fields and buttons:

- View Type:** Stored Result
- Parameter:** pX
- Filter by:** Location No.
- Start No.:** 0001
- End No.:** 0003
- Stored Num:** 0003/1000
- Display Type:** Display in Table

On the right side, there are three buttons: "Start Search", "Clear", and "Back".

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No.	Time	Sample ID	STD Value	Meas Value	Conclusion
0001	2022/05/20 14:17:46	Sample ID	7.000pH	7.000pH	Pass
0002	2022/05/20 14:17:50	Sample ID	7.000pH	7.000pH	Pass
0003	2022/05/20 14:17:53	Sample ID	7.000pH	7.000pH	Pass

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```

*****
Report Title
-----
Measure Time:2021/01/19 12:27:28
Operator: Operator 1
Model: M500T Multi-parameter Analyzer
Serial Number:
SW Version: Ver 1.00
-----
..... MATCHED INFO
Stored Num:          28
Matched Num:         1
Stored No. :         15
-----
..... CALIB INFO
Calib Operator:      REX Team
Calib Time:          2020/05/13 08:30:00
Calib Num:           3
-----
..... CALIB RESULT
STD 1:  4.01pH      176.8mV   25.0c
STD 2:  7.00pH      0.0mV    25.0c
STD 3: 10.01pH     -178.1mV  25.0c
pH Slope 1:                100.00%
pH EO 1:                    0.0mV
pH Slope 2:                100.00%
pH EO 2:                    0.0mV
-----
..... BRIEF INFO
Reading Mode:           Timed Reading
Stable Type:            Medium
Temp Comp Type:         ATC
-----
..... SAMPLE INFO
Sample ID:              Sample 1
-----
..... RESULT
Result:                 7.000pH
Signal Value:           -0.0mV
Temp Value:             25.0c
-----
Signature:

```

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# 1 Introduction

## 1.1 Introduction

The MP-3006 benchtop multi-parameter analyser is a multifunctional water quality testing device equipped with pH/ISE conductivity, dissolved oxygen, and temperature modules. The meter supports measurement parameters including pH, mV, ORP, pX, ion concentration, conductivity, resistivity, TDS, salinity, dissolved oxygen concentration, dissolved oxygen saturation, and temperature.

Module	Parameters	MP-3006
pH	pH, mV, ORP,	•
ISE	pX, ion concentration	•
Conductivity	Resistivity, conductivity, TDS, salinity, ash	•
Dissolved oxygen	Dissolved oxygen current, dissolved oxygen concentration, saturation	•
Temperature measurement	Temperature	•

### General Features

- Colour High contrast LCD touchscreen, 7.0 inches.
- An intelligent operation system provides features including user management, calibration management, method management, data management, etc.
- The multi-reading feature allows auto-reading, timed-reading, and continuous reading.
- Automatic/Manual temperature compensation ensures accurate results.
- Data Storage 1000 sets for each parameter (GLP-compliant).
- The data analysis feature helps users review, compare and recalculate results.
- Support for USB or RS-232 communication.
- The reset feature automatically resumes all settings back to factory default options.
- IP54 waterproof.

### pH

- 1-6 points calibration with Standard Recognition.
- Selectable pH buffer groups, including DIN, NIST, USA, MERK, JIS, and GB.
- Support for pH standard group customization.
- Automatic electrode diagnosis with pH slope and offset display.

### Ion

- 1-6 points calibration
- Selectable measurement units, including  $\mu\text{g/L}$ ,  $\text{mg/L}$ ,  $\text{g/L}$ ,  $\text{mmol/L}$ , pX, ppm, and ppb.
- Multi-measurement modes are supported, including Direct Reading mode, Standard Addition mode, Sample Addition mode, and GRAN mode.
- Over 10 methods are built-in, including  $\text{F}^-$ ,  $\text{Cl}^-$ ,  $\text{Br}^-$ ,  $\text{I}^-$ ,  $\text{NO}_3^-$ ,  $\text{BF}_4^-$ ,  $\text{NH}_4^+$ ,  $\text{K}^+$ ,  $\text{Na}^+$ ,  $\text{Ca}^{2+}$ ,  $\text{Cu}^{2+}$ ,  $\text{Pb}^{2+}$ ,  $\text{Ag}^+$  and etc., user-defined method is supported.

## Conductivity

- 1-5 points calibration automatically recognizes the standard solutions.
- Settable parameters, including cell constant, reference temperature (5/10/15/18/20/25°C), temperature compensation coefficient, and TDS factor.
- Temperature compensation type (none, linear, pure water, non-linear).

## DO

- Support for air-saturated water or zero oxygen calibration.
- Auto barometric pressure compensation.
- Manual Salinity Factor Correction.
- Selectable pressure unit, including kPa, mbar, Torr, Atm.

## 1.2 Technical Specification

Table 1-1 Instrument Specifications

Model		MP-3006
pH/pX level		0.001 pH/pX
mV	Range	(-2000.00~ 2000.00) mV
	Resolution	0.01mV
	Accuracy	±0.03% or ±0.1mV
	Repeatability	0.1mV
	Input Current	≤1×10 <sup>-12</sup> A
	Input Impedance	≥3×10 <sup>12</sup> Ω
pH	Range	(-2.000~ 20.000) pH
	Resolution	0.001pH
	Accuracy	±0.002pH
	Repeatability	0.001pH
	Measurement Accuracy	±0.01pH
	Measurement Repeatability	0.005pH
pX	Range	(-2.000~ 20.000) pX
	Resolution	0.001pX
	Accuracy	±0.002pX
	Repeatability	0.001pX
	Measurement Accuracy	±0.01pX
	Measurement Repeatability	0.005pX
Ion concentration	Range	( 1.000e <sup>-9</sup> ~9.999e <sup>+9</sup> ), mol/L, mmol/L, g/L, mg/L, µg/L, ppm, ppb
	Resolution	4 significant digits
	Measurement Accuracy	±0.3%

Conductivity	Range	0.000 $\mu$ S/cm $\sim$ 2000mS/cm
	Resolution	0.001 $\mu$ S/cm, automatic switching according to the range
	Accuracy	$\pm$ 0.5% (FS)
	Repeatability	0.17%(FS)
	Measurement Accuracy	$\pm$ 0.80% (FS)
	Measurement Repeatability	0.40% (FS)
Resistivity	Range	5.00 $\Omega$ .cm $\sim$ 100.0M $\Omega$ .cm
	Resolution	0.01 $\Omega$ .cm, automatic switching according to the range
	Accuracy	$\pm$ 0.5% (FS)
TDS	Range	0.000 mg/L $\sim$ 1000g/L
	Resolution	0.001mg/L, automatic switching according to the range
	Accuracy	$\pm$ 0.5% (FS)
Salinity	Range	(0.00 $\sim$ 8.00)%
	Resolution	0.01%
	Accuracy	$\pm$ 0.1%
	Measurement Accuracy	$\pm$ 0.2%
Dissolved Oxygen	Range	(0.00 $\sim$ 99.99)ppm (0.00 $\sim$ 50.00) ppm (measuring range)
	Resolution	0.01ppm
	Accuracy	$\pm$ 0.10ppm
	Repeatability	0.15ppm
	Accuracy at 0% DO	$\leq$ 0.1ppm
	Measurement Accuracy	$\leq$ 20.00ppm: $\pm$ 0.30 ppm, >20.00ppm: $\pm$ 10.0%
	Response time	$\leq$ 45s (90% response at 20°C).
	Salinity Compensation Accuracy	$\pm$ 2%
Saturation	Range	(0.0 $\sim$ 600.0)%
	Resolution	0.1%
	Accuracy	$\pm$ 2.0%
	Measurement Accuracy	$\pm$ 10.0%
Temperature	Range	(-10.0 $\sim$ 135.0) °C/ (14.0 $\sim$ 275.0) °F
	Resolution	0.1 °C/0.1 °F
	Accuracy	$\pm$ 0.1 °C
	Instrument indication error	$\pm$ 0.3°C(0°C $\sim$ 60°C), $\pm$ 1.0 °C(Else)
	Work environment	Ambient temperature: (0 $\sim$ 40) °C Relative humidity: not more than 85%
	Dimensions (LxBxH), weight (kg)	220mm x195mmx68mm, 0.95kg
	Power supply	Power adapter (Input AC 100 $\sim$ 240V. Output DC 24V)

## 1.3 Function Introduction

Table 1-2 Functions Specification

Features		Explanation
Basic Function	Languages	English
	Backlight adjustment	•
	Automatic diagnostics	•
	Factory reset	•
	Default parameter	•
	Prompt Sound	•
	Time settings	•
	Power failure protection	•
	Login password protection	•
	Firmware upgrade	•
	Anti-interference automatic recovery	•
	Automatic shutdown	•
	Protection	IP54
Reading Function	Reading stability settings	•
	Auto-lock reading	•
	End point judgment/reading mode	Continuous, Auto, Timed
	Sample ID	•
	Alarm	•
Data Management	Storage	1000 sets of measurement parameters each 100 sets of calibration data
		1000 sets pH standard solution verification data
	Search - 10 -	Storage number, time, operator, method name, sample ID, and electrode ID.
	View	Sheet and Graph
Data Management	Delete	
	Delete	•
	GLP	•
Communication & external devices	Connect to Printer	RS232
	U Disk	•
	Content and format customization	Standard, GLP, Custom format
	PC Output	•
	Connect to a PC for instrument control	•
	Scanner	•
	Auto sampler	•

pH/mV Measurement	pH electrode status/performance	Slope, Electrode status (Excellent, good, bad)
	Multi-point calibration	6 points
	Automatic standard solutions recognition	6 groups
	Standards customization	•
	Standard groups customization	1 group
	Temperature compensation	Automatic/ Manual
	pH Electrode Diagnostics	•
	pH electrode calibration reminder	•
	pH electrodes mandatory calibration	•
	pH standard solution verification	•
	pH Mandatory Verification	•
	ORP Measurement	•
	ORP electrode calibration reminder	•
	ORP electrodes mandatory calibration	•
pX/ISE Measurement	Built-in ions methods	12 kinds
	Ion's customization	•
	Multi-point calibration	6 points
	Optional units	mol/L, mmol/L, g/L, mg/L, µg/L, ppm, ppb
	Measurement mode	Direct reading method, Standard addition method, Sample addition method, GRAN method
	Ion electrode calibration reminder	•
	Ion electrode mandatory calibration	•
Conductivity Measurement	Conductivity	•
	Resistivity	•
	TDS	•
Conductivity Measurement	Salinity	•
	Conductivity-ash	•
	Reference temperature	6 types, default 25.0°C
	Multipoint calibration	5 points
	Automatic standard solutions	Universal Standard,

Features		Explanation
	recognition	GB Standard
	Cell constant set	•
	Temperature compensation coefficient set	•
	Salinity compensation coefficient set	•
	Compensation mode	Non-compensatory, linear, pure water, no-linear
	Automatic temperature compensation	•
	Manual temperature compensation	•
Dissolved oxygen Measurement	Measurement principle	Polarographic
	Zero calibration	•
	Air calibration	•
	Automatic temperature compensation	(0.0~45.0) °C
	Automatic atmospheric pressure compensation	(600~1100)mbar
	Manual atmospheric pressure compensation	(600~1100)mbar
	Atmospheric pressure units	kPa, mbar, Torr, Atm
	Manual salinity compensation	(0.0~50.0) g/L
Temperature Measurement	Temperature units	°C, °F
	Temperature calibration	•
Management	GMP Mode	•
	System & Running log	•
	User management	Level 3 permissions, 8 users
	Method management	50
	Sample list	•

## 2 Safety Notices

Please read the manual before use. The user **MUST** use the instrument following this manual to avoid damage to the user and equipment.

Before using the meter, **READ** the following notes:

- **DO NOT DISASSEMBLE** the device for inspection or repair.
- To prevent electric shock or damage to the device, **DO NOT** place cables and connectors in any liquid, wet or corrosive environment.
- Please use the defaulted power adapter, **DO NOT** use it if the power cord is damaged (the wire is exposed or broken).
- **DO NOT** use in flammable and explosive environments.
- **DO NOT** use if the user finds any abnormalities such as damage or deformation of the device.

## 3 Terms Explanation

### pH/pX

- **pH/pX Slope:** The amount of potential change generated by each 1 pH/pX change, expressed in mV/pH or by 100% Theoretical Slope (PTS).  $pX = -\log[X]$ , where  $[X]$  means molar concentration (mol/L) of X ions.
- **E0 of pH:** Also known as "zero potential", it usually refers to the potential value at a pH of 7.
- **One-point calibration:** Calibration with a standard solution.
- **Two-point calibration:** Calibration with two standard solutions.
- **Multi-point calibration:** Calibration with more than two standard solutions.

### Redox potential (ORP).

- **Redox potential:** Also known as ORP value, it refers to the potential difference between the potential difference of a measuring battery relative to a standard hydrogen electrode consisting of an indicator electrode, a reference electrode, and a solution under test. The symbol for ORP measurement is "RmV", and the unit is mV.
- **One-point calibration:** Calibration with an ORP standard solution.
- **Offset:**  $= E_{RmV} - E$  ( $E_{RmV}$  potential relative to the hydrogen electrode,  $E$  measured potential). At a certain temperature, when the ORP calibration solution is used for calibration, the potential value "RmV" and the potential difference between the measured system.

### Electrical conductivity

- **Cell Constant:** Also known as the conductivity cell constant. The ratio of the distance to the area of the electrode sheet is expressed in  $\text{cm}^{-1}$ . Usually, there are conductance electrodes with several cell constants such as 0.01, 0.1, 1.0, 10, etc. The conductance electrode with a cell constant of 1.0 is the most used one and has a wide measurement range.
- **Temperature Coefficient:** The change in conductivity caused by a 1°C change in temperature is usually expressed in %/°C, and the default is 0.02, which is 2.00%/°C.
- **TDS Conversion Factor:** The conversion factor between conductivity and TDS, which defaults to 0.5.

### Dissolved oxygen

- **Dissolved oxygen concentration:** The content of oxygen dissolved in water. Expressed in milligrams of oxygen per liter of water, usually denoted as DO.
- **Dissolved Oxygen Saturation:** The ratio of the on-site dissolved oxygen concentration to the saturated dissolved oxygen concentration under the same conditions.
- **Salinity:** Salt content in water, expressed in g/L. When the salinity increases by 1g/L at 15°C, the saturated dissolved oxygen of water decreases by about 0.0559 mg/L.
- **Zero-point calibration:** Electrodes are calibrated in "oxygen-free water" (freshly formulated 5% sodium sulphite solution).
- **Fullness calibration:** The electrodes are calibrated in air or water fully dissolved and saturated with air.
- **Atmospheric pressure compensation:** The atmospheric pressure affects the measurement of dissolved oxygen concentration and dissolved oxygen saturation, and atmospheric pressure compensation is required. Before calibration, the on-site atmospheric pressure needs to be input, expressed in mbar, and the default is 1013 mba.

### General

- **Calibration reminder (Recommended):** "Electrode calibration" prompts on the homepage, user can start measurement without the last calibration.

- **Calibration reminder (Mandatory):** "Electrode calibration" prompts on the homepage, calibration is required for accurate measurement.
- **Upper limit:** The upper limit value is monitored for the measured or calibrated data, and the upper limit value must not be lower than the lower limit value.
- **Lower limit:** Lower limit monitoring of measured or calibrated data.

## 4 Overview and Installation

### 4.1 Overview

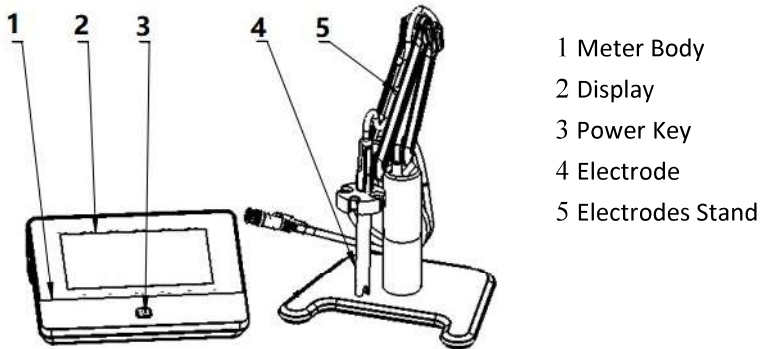
Figure 4-1 Overview-Front View


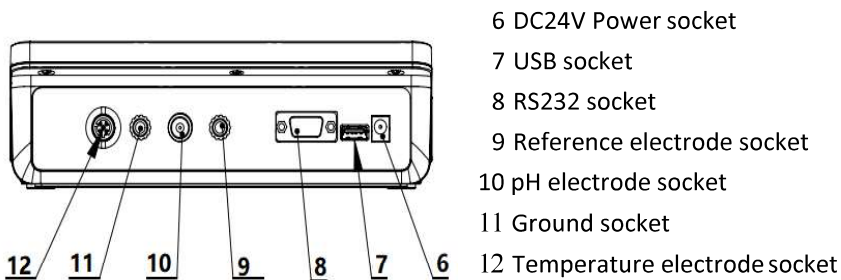
Figure 4-2 Overview- Back View

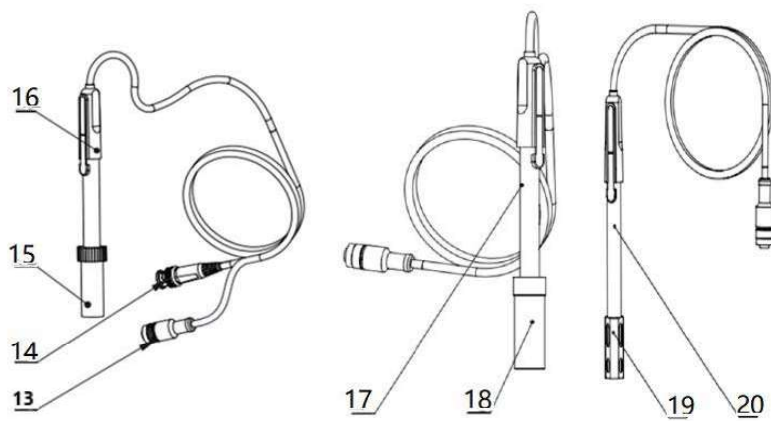- 13 Four-pin aviation connectors
- 14 pH electrode connectors
- 15 Electrode protection cap
- 16 pH electrodes
- 17 Conductivity electrodes
- 18 Conductivity electrode protection cap
- 19 Dissolved Oxygen Electrode protection cap
- 20 Dissolved oxygen electrodes

Figure 4-3 Electrodes and connectors Table 4-1 Connector Specifications

Connector specifications	Electrodes Connection
BNC(Q9)	pH electrode, ORP electrode, ISE probe
Four-pin aviation	DO electrode, ATC probe
Five-pin aviation	Conductivity electrode
Banana	Reference electrode

## 4.2 Instrument Installation

### 4.2.1 Electrodes Stand Installation

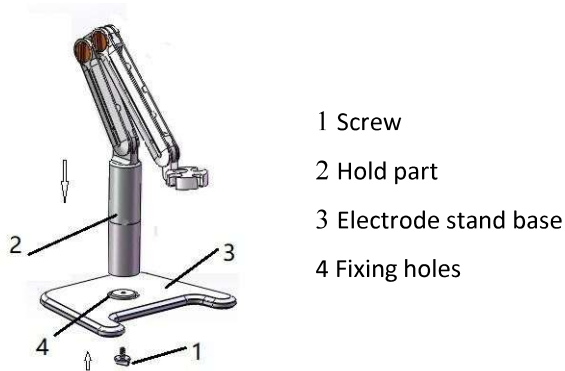
Figure 4-4 Electrode Stand Installation

- 1) Put the electrode base on the table.
- 2) Screw the fixed screw into the fixing hole.
- 3) Place the electrode holder from top to bottom to fix at the vertical shaft of the fixing screw.

### 4.2.2 Electrodes Connection

#### 4.2.2.1 Connection of pH electrodes

Push the pH Electrode into the electrode holder. Remove the socket protector cap of the pH electrode. Connect the pH electrode into the right socket. If the ATC probe is applied, or ATC has been integrated into the pH probe, please connect the ATC probe onto the DO/T electrode socket.

#### 4.2.2.2 Connection of conductivity electrodes

Push the conductivity electrode into the electrode holder. Remove the protector cap of the conductivity electrode. Connect the conductivity electrode into the right socket. Combination conductivity probes integrated with ATC probe. If the separate ATC probe is applied, please connect the ATC probe onto the DO/T electrode socket. At the measurement, please choose the right input of ATC in the meter when you applied a separate ATC probe.

#### 4.2.2.3 Connection of dissolved oxygen electrodes

Push the dissolved oxygen electrode into the electrode holder. Remove the protector cap of the dissolved oxygen electrode socket. Connect the dissolved oxygen electrode into the right socket. All dissolved oxygen electrodes are integrated with the ATC probe.

## 5 Instrument Operation

### 5.1 Switch On/Off

Press and release  to switch on the meter. The start-up screen shows the device model, name, software version, and other related information. After the self-test program, the screen turns to the homepage and the meter is ready to measure.

Also, press and hold  the key for more than 3 seconds and release it to shut down.

### 5.2 Screen Icons

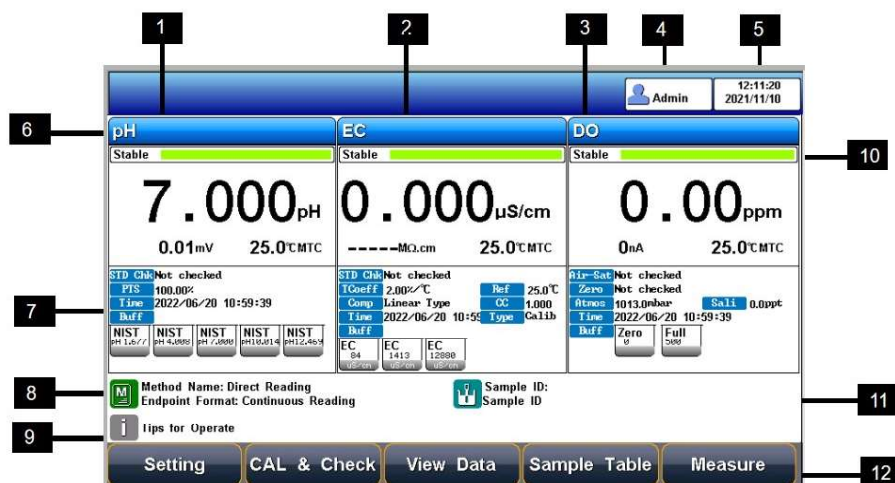
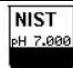
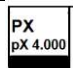
Figure 5-1 Screen icons explanation

1 Measurement box. 2 Measurement boxes. 3 Measurement box. 4 User ID. 5 System time. 6 Measurement parameters. 7 Calibration information. 8 Method management. 9 Operation tips. 10 Reading status. 11 Sample ID. 12 Soft function keys.

The instrument displays symbol identification that has the following functional implications:

Table 5-1 Symbol Explanation

No.	Symbol	Explanation
1		Reading status, display the measurement status of reading, stable, locked, each indicates that the processing, stable, and reading completed.
2	PTS	The percentage slope of the pH electrode calibration data.
3	ATC	Automatic temperature compensation
4	MTC	Manual temperature compensation
5	ORP	Redox potential value, in mV
6	Offset	Offset potential, in mV
7		Measurement method management, display the current method information
8		Standard solution for pH calibration
9		Standard solution for ion calibration
10		User ID
11		Sample ID

## 5.3 Methods Management

The meter provides a library of built-in methods. Select the proper method to measure, the system loads the method's preset information, including the method name, a brief overview, creator, creation date, and measurement parameters.

In the method settings, it supports the navigation setting including the method information, parameters, reading mode, DO parameters setting, temperature setting, and data management setting. For the first use, please follow the guide to settings the measurement parameters.

Figure 5-2 Methods of Information

### 5.3.1 Update Method

When the method is revised by the setting, there is a red mark in the method management logo to remind the user that the current method is different from the method in the library.

When pressing the method in the method library, there is a system tip to give the user a choice to update or create a new method.

### 5.3.2 Create a new Method

There are three modes to create a new method: create the method step by step in the system, save the setting as a new method in the system, and save the method with a new name, description, and location. The location can be the system and U disk.

Figure 5-3 Method setting view

### 5.3.3 Delete a Method

There are 17 methods saved in the meter as the default setting in the library method. The No.001-No.017 method cannot be deleted, and the others can be deleted.

### 5.3.4 Save and Load a Method

It supports storing up to 50 methods in the meter and saves and loads the methods in the U Disk.

## 5.4 Parameter Setting

In the measuring, users can set the instrument parameters by pressing "Parameter Setting" to set measuring parameters including tutorial setting, basic method information, select parameters, endpoint format, pH parameter, pX/ISE parameter, ORP parameter, temperature parameter, data management, output setting, user management, system setting, and GMP mode manage.

Figure 5-4 Parameter settings of the instrument

### 5.4.1 Tutorial setting

For the first use, please follow the guide to settings the measurement parameters. After all the settings, press the "Enter" to return to the previous page.

### 5.4.2 Method basic information

The meter provides a library of built-in methods. Select the proper method to measure, the system loads the method's preset information, including the method name, a brief overview, creator, creation date, and measurement parameters.

### 5.4.3 Select parameters

It could select one measurement parameter from pH, pX, Ion Conc and ORP, resistivity, conductivity, TDS, salinity, ash, DO, and saturation for every test.

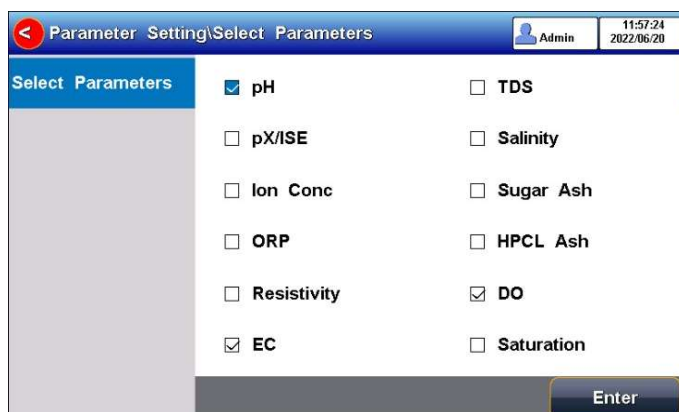
Figure 5-5 Select parameters

The instrument contains up to 3 measurement modules, and each measurement function can select different measurement parameters (Switch the parameters by clicking on the blank area of the measurement box on the screen).

Select the proper measurement parameters to perform a measurement. The screen displays the detail of three measurement parameters when three modules are chosen. Two parameters are shown as two modules are chosen. A single parameter is shown by one measurement method selected.

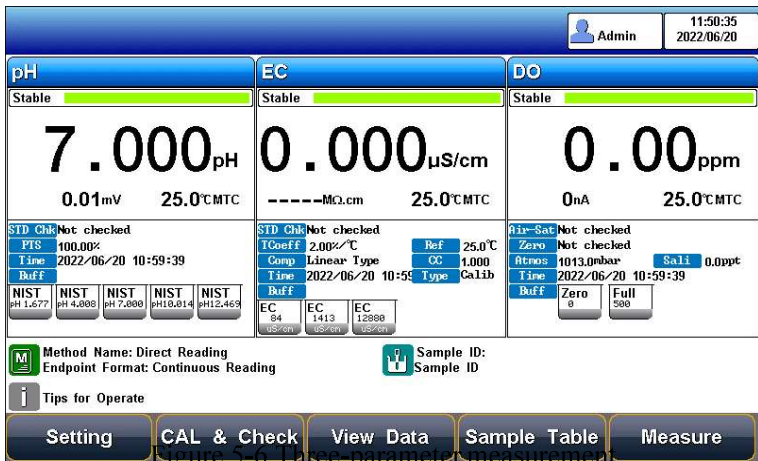
Figure 5-6 Three-parameter measurement

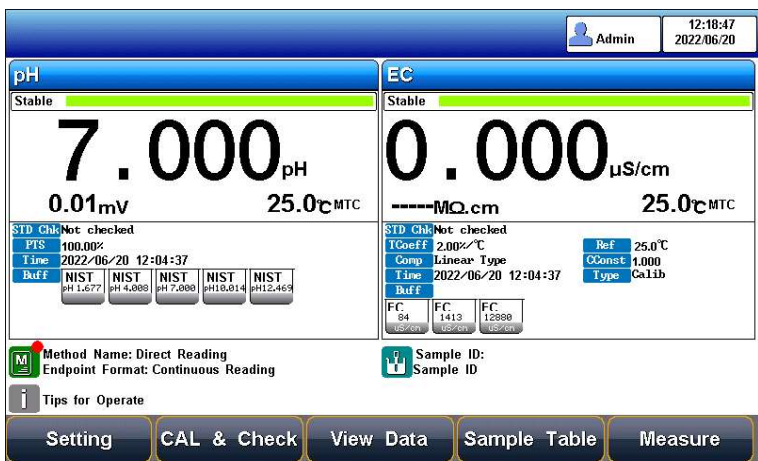
Figure 5-7 Two-parameter measurement

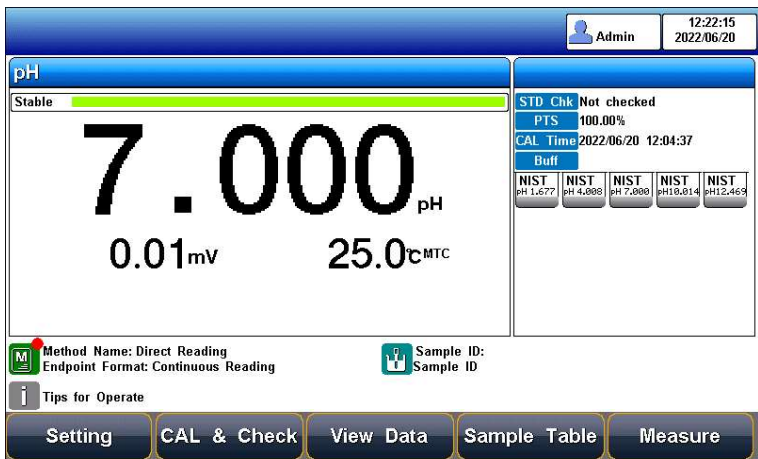
Figure 5-8 Single-parameter measurement

## 5.4.4 Reading Mode Settings

The meter provides three reading modes, including continuous reading, auto reading, and timed reading.

- Continuous reading: The instrument displays real-time measurement results. Users can end the measurement at any time and save the last result.
- Auto-reading: The measurement reached the balance, and the meter locked the reading result. The meter offers "Fast", "Medium", "Strict" and "Custom" four options for endpoint detection conditions.
- Timed reading method: Timed Reading contains two kinds of timed reading methods: "Interval Measurement" and "Timed Measurement". "Interval Measurement" provide measurement results at interval time and "Timed Measurement" provide measurement result after a set time.

Table 5-2 Reading Parameters Settings

Stability Type		pH	pX/ Ion concentratio	Conductivity	DO
Fast	Stable time	4s	4s	5s	5s
	Fluctuation	0.6mV	0.3mV	1.0%	4nA
Medium	Stable time	6s	8s	8s	8s
	Fluctuation	0.1mV	0.08mV	0.4%	3nA
Strict	Stable time	8s	12s	15s	15s
	Fluctuation	0.03mV	0.03mV	0.1%	2nA
Custom (Recommended value)	Stable time	1 to 30s	1 to 30s	1 to 30s	1 to 30s
	Fluctuation	0.03~1mV	0.03~1mV	0.1~2%	2~5nA

## 5.4.5 pH parameter setting

### 5.4.5.1 pH electrode information

The screenshot shows the 'Parameter Setting \ pH Parameter' screen. The interface includes a navigation menu on the left with options like 'pH Electrode', 'pH Buffer Group', 'Resolution', 'pH Alarm Setting', 'Calibrate Reminder', and 'pH STD Check'. The main area displays the following information:

- Name:** pH Electrode
- Serial No.:** 12345678
- Recognition:** Auto Mode
- CAL Result:** A table showing calibration results for four buffer groups (BUF1-BUF2, BUF2-BUF3, BUF3-BUF4, BUF4-BUF5). Each entry shows 100.00% slope, 0.00mV zero-point, and a 'Perfect' comment.
- CAL Time:** 2022/05/08 16:23:43
- User ID:** REX Team

An 'Enter' button is located at the bottom right of the form.

Figure 5-9 pH electrode information

In the meter setting, a set of pH Electrode info defaults. At the measurement, the meter loads the defaulted pH electrode's info including serial number, last calibration result, the number of calibration points, slope value, zero-point potential value and electrode state description, etc., the calibration time and calibrator will also be loaded together.

### 5.4.5.2 Recognition

Auto Mode and Manual Mode.

In some special cases, it is necessary to use some non-standard pH buffer solutions or use two very close pH standard

buffer solutions for electrode calibration. In this case, the manual standard solution identification function can be used. When set to "Manual Mode", the pH value of the current standard solution can be input during and used for electrode calibration.

### 5.4.5.3 pH Buffer groups

The meter provides various Standards Group including GB, DIN, NIST, USA, MERK, and JIS. And allows the user to prepare the customized Standard groups.

Table 5-3 Standard Solution Groups

Groups	Contents
GB	1.680pH, 3.559pH, 4.003pH, 6.864pH, 7.409pH, 9.182pH, 12.460pH
DIN	1.680pH, 2.000pH, 3.557pH, 3.775pH, 4.008pH, 6.865pH, 7.000pH, 7.416pH, 9.184pH, 10.014pH, 12.454pH
NIST	1.677pH, 4.008pH, 6.864pH, 7.000pH, 7.416pH, 10.014pH, 12.469pH
USA	1.680pH, 4.010pH, 7.000pH, 10.010pH
MERK	2.000pH, 4.000pH, 7.000pH, 9.000pH, 12.000pH
JIS	1.680pH, 4.008pH, 6.865pH, 7.413pH, 9.180pH, 10.010pH

Figure 5-10 Selection of standard groups and standard solution

Usually, we use the pH value corresponding to 25.0°C to mark the pH standard buffer solution, such as NIST 7.00pH standard solution, which means the standard pH buffer solution is 7.00pH, and it is 7.00pH at 25.0°C.

After selecting the standard solution group, we need to select the standard buffer solution used for calibration from the standard solution group. The instrument supports up to 6-point calibration, that is, up to 6 standard solutions can be selected. Since the pH values of multiple standard buffer solutions in the standard solution group may be very close, ensuring that the instrument can correctly identify the standard buffer solution, will limit the selection of standard solutions with neighbouring pH values. The check mark indicates the currently used standard solution group and the corresponding standard solution.

### 5.4.5.4 Resolution settings

The pH measurement resolution of the instrument is adjustable. pH Resolution: 0.001pH, 0.01pH and 0.1pH. mV resolution: 0.01mV, 0.1 mV and 1 mV.

Set the result resolution through "pH Parameters" - "Resolution".

### 5.4.5.5 pH Alarm setting

The meter support pH measurement and calibration result monitoring alarm. The setting includes slope limit, potential value limit, and monitoring options. When the measurement results out of the selected range, the meter alarms.

pH Electrode	pH CAL Alarm:	On
pH Buffer Group	Slope Max(%)	120.00%
Resolution	Slope Min(%)	80.00%
<b>pH Alarm Setting</b>	E0 Max:	30.00mV
Calibrate Reminder	E0 Min:	-30.00mV
pH STD Check	pH Alarm:	On
	pH Max:	10.000pH
	pH Min:	4.000pH

Enter

Figure 5-11 pH Alarm setting information

### 5.4.5.6 pH Calibrate reminder settings

The meter provides a calibration prompts function. From the setting, the user can select calibration reminder (recommended) and calibration reminder (mandatory) for future calibration.

Calibration reminder (recommended): When enabling the calibration reminder, the instrument verifies whether the calibration data of the pH electrode is within the validity period. If calibration information expires, a striking calibration reminder will appear at "CAL&Check" on the main interface, but it will not affect the measurement and data saved.

Calibration reminder (mandatory): When enabling the mandatory calibration, the instrument verifies whether the calibration data of the pH electrode is within the validity period. If calibration information expires, a striking calibration reminder will appear at "CAL&Check" on the main interface, the measurement data cannot be saved and output.

pH Electrode	pH CAL Message:	On
pH Buffer Group	CAL Interval:	10
Resolution	Time Unit:	Hour
pH Alarm Setting	pH Compulsive CAL:	On
<b>Calibrate Reminder</b>	Validity:	10
pH STD Check	Time Unit:	Hour

Enter

Figure 5-12 pH calibration reminder setting information

### 5.4.5.7 pH Verification Reminder settings

The meter provides a verification prompts function. The electrode slope and zero potential of pH electrodes drift slightly over time, and this change will have a direct impact on the measurement results. So, the verification of standards is important for accurate results.

Verification reminder (mandatory): When enabling the verification reminder, the instrument verifies whether the verification data of the pH electrode is within the validity period. If verification information expires, a striking verification reminder will appear at "CAL&Check" on the main interface, the measurement data cannot be saved and output.

It allows the user to input the pH standard value and the allowable deviation. If the deviation amount is set to 0, it means that no verification is required.

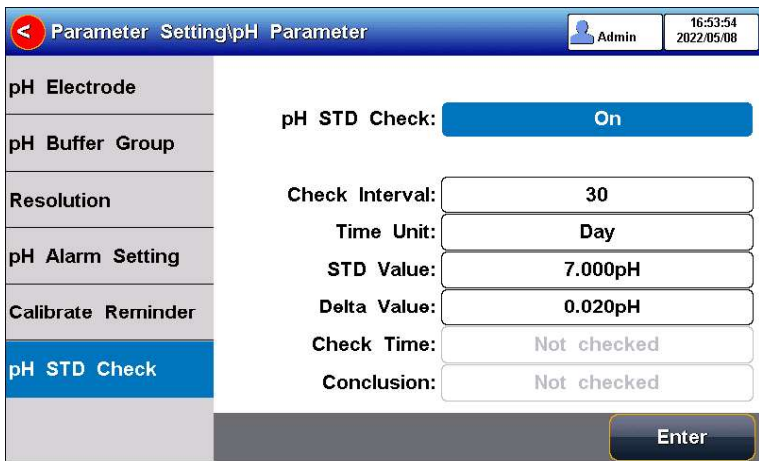
Figure 5-13 pH verification information

### 5.4.6 pX/ISE parameter settings

#### 5.4.6.1 pX electrode information

The instrument has one pX electrode created by default, which is convenient for users. When using the built-in default electrode, the instrument automatically loads the electrode serial number and the last calibration result, including calibration points, slope value, zero-point potential value, etc., and the calibration time and calibrator are also loaded.

#### 5.4.6.2 Ion mode

Ion mode corresponds to pX/ISE, ion concentration measurement. The instrument supports conventional ion modes and user-defined ion modes. The instrument provides a variety of commonly used ion modes which are convenient for use. Users can directly measure the concentration of the corresponding ions as long as they are equipped with the corresponding ion selective electrode.

Custom ion mode: Press the "Create" key, enter the ion name (maximum 8 characters), then enter the molecular weight (molar mass), and select the ionic valency.

Figure 5-14 Ion mode management

#### 5.4.6.3 Select measurement mode

The instrument supports 4 ion measurement modes: direct reading concentration method, standard addition method, sample addition method, and GRAN Measurement method.

Users can choose the following ion measurement methods according to their requirements. In general, when the ion

measurement parameter selects pX or ion concentration measurement, it is recommended to use the direct reading concentration method for measurement, which has a fast measurement speed and is suitable for the measurement of batch samples. Standard addition, sample addition, and GRAN measurements are only valid when the measurement parameter selects ion concentration.

Figure 5-15 Ion Reading mode selection

The direct reading concentration method: also called the standard curve method, is the most commonly used method for measuring ion concentration. According to the Illuminant formula, a linear relationship is established between the ion concentration and the electrode potential, to determine the ion concentration of the sample.

Standard addition method: Also called the known addition method, it is mainly used to measure samples with low content. The standard solution is added to the sample, and the concentration value of the sample is determined by measuring the potential change before and after the addition.

Sample addition method: Similar to the standard addition method, it is mainly used to measure samples with higher content. The sample to be tested is added to the standard solution, and the concentration value of the sample is determined by measuring the potential change before and after the addition.

GRAN measurement method: similar to the standard addition method, it is mainly used to measure samples with low content, and the concentration value of the sample is measured according to the mathematical principle of the GRAN mode.

#### 5.4.6.4 Resolution settings

The pX measurement resolution of the instrument is adjustable. pX resolution: 0.001pX, 0.01pX and 0.1pX.

mV resolution: 0.01mV, 0.1 mV and 1 mV.

Set the result resolution through the "pX/ISE parameter" - "Set Resolution".

#### 5.4.6.5 pX/ISE Alarm setting

The meter supports the pX/ISE measurement alarm. When the measurement results out of the selected range, the meter alarms.

Enter the ion alarm limit setting through "pX/ISE parameter" - "Alarm setting", and you can set the max and min limit values of ion concentration measurement.

Parameter Setting\pX/ISE Parameter		Admin	13:57:20 2022/05/20
pX/ISE Electrode	Ion Type:	F-	
Ion Mode	pX Alarm:	On	
Resolution	pX Max:	10.000pX	
Alarm Setting	pX Min:	4.000pX	
Calibrate Reminder			
		Enter	

Figure 5-16 pX/ISE Alarm setting information

#### 5.4.6.6 pX Calibrate reminder settings

The meter provides a calibration prompts function. From the setting, the user can select calibration reminder (recommended) and calibration reminder (mandatory) for future calibration.

Calibration reminder (recommended): When enabling the calibration reminder, the instrument verifies whether the calibration data of the pH electrode is within the validity period. If calibration information expires, a striking calibration reminder will appear at "CAL&Check" on the main interface, but it will not affect the measurement and data saved.

Calibration reminder (mandatory): When enabling the mandatory calibration, the instrument verifies whether the calibration data of the pH electrode is within the validity period. If calibration information expires, a striking calibration reminder will appear at "CAL&Check" on the main interface, the measurement data cannot be saved and output.

Under "pX/ISE parameter" - "Calibrate Reminder", set calibration reminder and mandatory calibration. Fulfill the electrode calibration reminder function by setting the calibration period and time, and the electrode calibration function by setting the mandatory calibration validity period and time.

Parameter Setting\pX/ISE Parameter		Admin	13:58:00 2022/05/20
pX/ISE Electrode	pX CAL Message:	On	
Ion Mode	CAL Interval:	10	
Resolution	Time Unit:	Hour	
Alarm Setting	pX Compulsive CAL:	On	
Calibrate Reminder	Validity:	10	
	Time Unit:	Hour	
		Enter	

Figure 5-17 pX calibration reminder setting information

### 5.4.7 ORP Parameter Settings

#### 5.4.7.1 ORP electrode information

In the meter setting, a set of ORP Electrode info defaults. At the measurement, the meter loads the defaulted pH electrode's info including serial number, resolution (mV supports 0.01mV, 0.1 mV, and 1mV settings), recognition mode, and last calibration result (nominal and offset), the calibration time and calibrator are also loaded.

### 5.4.7.2 ORP Alarm setting

The meter supports ORP measurement and calibration result monitoring alarm. When the measurement results out of the selected range, the meter alarms.

Enter the ORP alarm limit setting through "ORP parameter" - "Alarm Setting", and you can set the max and min limit values of ORP measurement.

### 5.4.7.3 ORP calibration reminder setting

ORP electrode calibration reminder: If the ORP electrode is not calibrated after the calibration period, the instrument will display a prominent calibration reminder at the "CAL&Check" on the initial interface, but it does not affect the use.

ORP electrode mandatory calibration: After turning on, the instrument verifies whether the ORP electrode calibration data is within the validity period. If the calibration information expires, only the test results can be read before normal calibration, and the data cannot be saved and output.

Under "ORP parameter"- "Calibrate Reminder", set the calibration reminder and mandatory calibration settings. Fulfill the electrode calibration reminder function by setting the calibration period and time, and the electrode calibration function by setting the mandatory calibration validity period and time.

## 5.4.8 Conductivity Parameter Settings

### 5.4.8.1 EC Electrode management

The electrode information includes the name, serial No., user ID, registration time, service life, calibration message, calibration number, and last calibration time. It allows the user to choose one electrode to calibrate and measure samples.

The meter has one created default EC electrode, which is convenient for users. When using the built-in default electrode, the meter automatically loads the electrode serial number and the last calibration result, including calibration points, constant cell, etc., and the calibration time and calibrator will also be loaded.

No.	STD Value	Measured	Cell Cont
STD1	84.0µS/cm	84.0µS/cm	1.000
STD2	141.3µS/cm	141.3µS/cm	1.000
STD3	12.88µS/cm	12.88µS/cm	1.000

Figure 5-18 EC electrode information

The meter allows users to create and save electrode information up to 5. The electrode information contains electrode name, electrode serial number, electrode type, registrant, registration time, shelf life, shelf-life reminder, calibration No., latest calibration time, etc.

### 5.4.8.2 EC calibration type

EC calibration type: Cal by Standards and input manually.

Cal by Standards: Cell constant is calibrated with standard conductivity standard solution.

Input manually: It allows the user to set the cell constant.

Conductivity electrode: Four conductivity cell constant 0.01, 0.1, 1, 10.

The defaulted conductivity cell constant is 1. Users need to enter the cell constant value on the label of the conductivity electrode for accurate measurement.

### 5.4.8.3 Cal Information

The meter provides two EC standards groups including the universal group and GB group. And allows the user to prepare the customized standard groups.

For neighbouring standards, please choose the customization to perform calibration.

Table 5-3 Standard solution groups

Standard group	Standard solution
Universal Group	10 $\mu$ S/cm, 84 $\mu$ S/cm, 500 $\mu$ S/cm, 1413 $\mu$ S/cm and 12880 $\mu$ S/cm.
GB group	146.5 $\mu$ S/cm, 1408 $\mu$ S/cm, 12852 $\mu$ S/cm and 111310 $\mu$ S/cm.

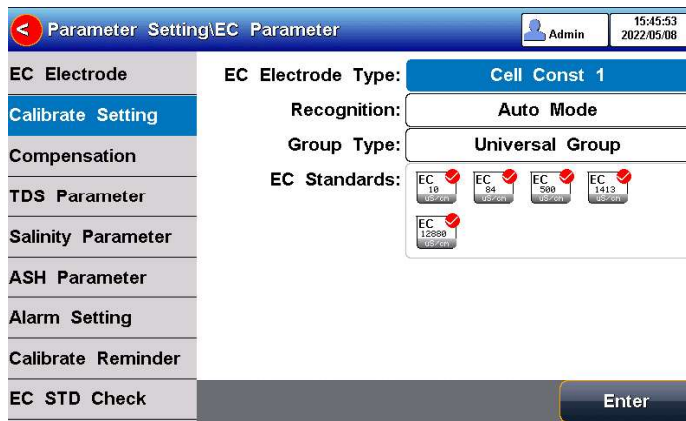
Figure 5-19 Selection of conductivity standards

### 5.4.8.4 EC Reference temperature

Conductivity reference temperature: The conductivity of the solution is greatly affected by temperature, to make the conductivity measurement results at different temperatures comparable, the conductivity and temperature values at the time of measurement are usually recorded and converted into the conductivity value at a certain temperature through temperature compensation, which is the reference temperature.

When the measurement parameter is conductivity, TDS, or resistivity, the meter allows settings of 5.0°C, 10.0°C, 15.0°C, 18.0°C, 20.0°C, 25.0 °C 6 reference temperatures, the default reference temperature is 25 °C.

When the measurement parameter is salinity, the default reference temperature is 18 °C and cannot change.

When the measurement parameter is Conductivity ash, the default reference temperature is 20 °C and cannot change.

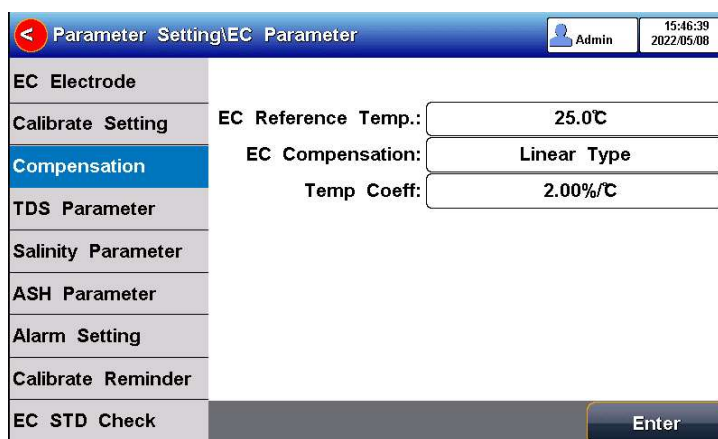
Figure 5-20 Conductivity temperature compensation setting

### 5.4.8.5 Conductivity Compensation

EC Compensation Mode: Three different compensation modes can be used for various applications. The meter supports Linear type, DI water type, and non-comp type.

1) Linear type: Linear compensation is usually used for the measurement of medium and high conductivity solutions. With linear compensation, you can set the temperature compensation coefficient, which defaults to 2.00%/°C (approximately the temperature compensation coefficient of a sodium chloride solution at 25°C). It allows the user to set the temperature coefficient.

2) DI water type: DI water compensation is usually used for the measurement of pure water and ultrapure hydropower conductivity below 5µS/cm. It allows the user to set the temperature coefficient.

3) Non-comp type: Non-compensation is usually used to obtain the true conductivity value at the measured temperature.

4) Non-linear type: Non-linear compensation is usually used for the measurement of surface water.

### 5.4.8.6 TDS Parameter

The TDSF calibration type can be calibrated by standard solution or by setting the TDS Factor. The TDS factor can be adjustable, and the default is 0.500. It allows the user to calibrate the TDS Factor.

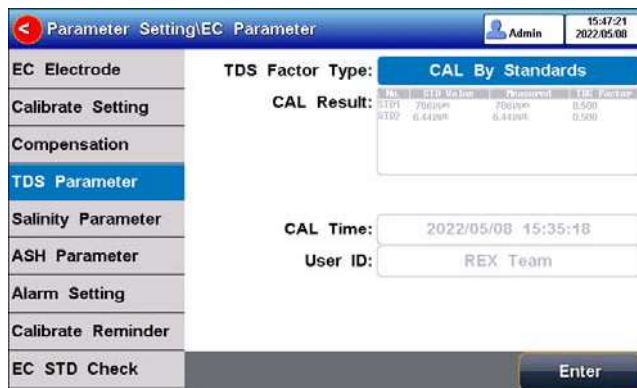
Figure 5-21 TDS measurement settings

### 5.4.8.7 Salinity parameter

The meter supports both default salinity and seawater salinity measurement types.

Sea salinity mode: it indicates the nominal salinity value used to measure seawater salinity correction. The standard seawater salinity concentration adopted here is 3.500%, also written as 35psu.

Default Salinity Mode: the salinity of the common sample, which can be used to approximate the salinity of the measured solution.

Figure 5-22 Salinity measurement settings

### 5.4.8.8 Conductivity Ash parameter

The meter supports the measurement of conductivity ash generally calculating the ash content using an ash conversion coefficient.

Ash Conversion Factor: Represents the conversion coefficient K of the conductivity of the solution to ash, which has a small value of  $10^{-4}$ .

Blank Conductivity: It allows for input manually or measures in the blank.

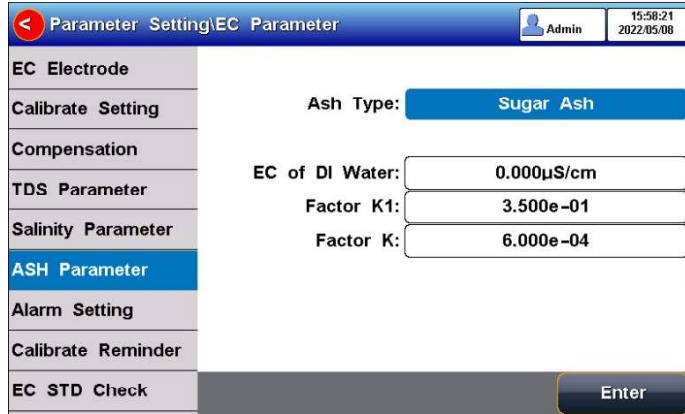
Figure 5-23 Conductivity Ash measurement parameter settings

### 5.4.8.9 Alarm setting

The meter support measurement and calibration result monitoring alarm. The parameter includes the conductivity, TDS, salinity, resistivity, sugar ash, and HFCS Ash.

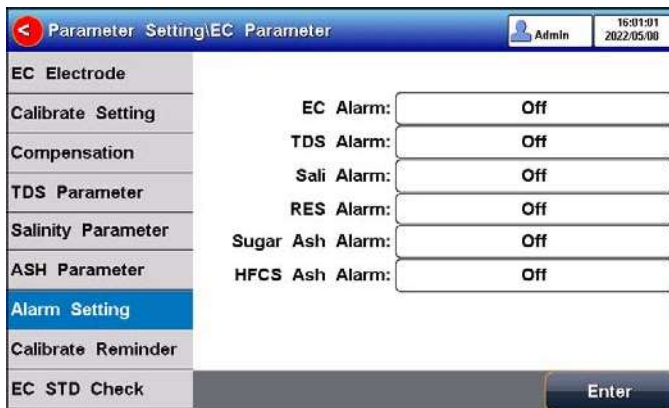
Figure 5-24 Alarm setting information

### 5.4.8.10 Calibration reminder settings

The meter provides a calibration prompts function. From the setting, the user can select calibration reminder (recommended) and calibration reminder (mandatory) for future calibration.

Calibration reminder (recommended): When enabling the calibration reminder, the instrument verifies whether the calibration data of the conductivity electrode is within the validity period. If calibration information expires, a striking calibration reminder will appear at "Calibrate" on the main interface, but it will not affect the measurement and data saved.

Calibration reminder (mandatory): When enabling the mandatory calibration, the instrument verifies whether the calibration data of the conductivity electrode is within the validity period. If calibration information expires, a striking calibration reminder will appear at "Calibrate" on the main interface, the measurement data cannot be saved and output.

Figure 5-25 Calibration reminder setting information

### 5.4.8.11 Verification Reminder settings

The meter provides a verification prompts function. The cell constant of conductivity electrodes drifts slightly over time, and this change will have a direct impact on the measurement results. So, the verification of standards is important for accurate results.

Verification reminder (mandatory): When enabling the verification reminder, the instrument verifies whether the verification data of the EC electrode is within the validity period. If verification information expires, a striking verification reminder will appear at "Calibrate" on the main interface, the measurement data cannot be saved and output.

It allows the user to input the conductivity standard value and the allowable deviation. If the deviation amount is set to 0, it means that no verification is required.

Figure 5-26 Conductivity verification information

## 5.4.9 Dissolved Oxygen Parameter Settings

### 5.4.9.1 DO Electrode management

The electrode information includes the name, serial No., cal result, cal time, and user ID.

No.	Name	Measured	Temp
1	DO Electrode	6.898	25.0°C

Figure 5-27 DO electrode information

### 5.4.9.2 DO salinity compensation settings

Salinity is the amount of sodium chloride dissolved in 1 L of water, in g/L. The dissolved oxygen concentration of water is highly affected by salinity. Typically, for every 1 g/L increase in salinity, the saturated dissolved oxygen of water decreases by 0.0559 ppm.

Users can set salinity compensation on the setting page. The salinity compensation range of measurement is (0.0~50.0) g/L.

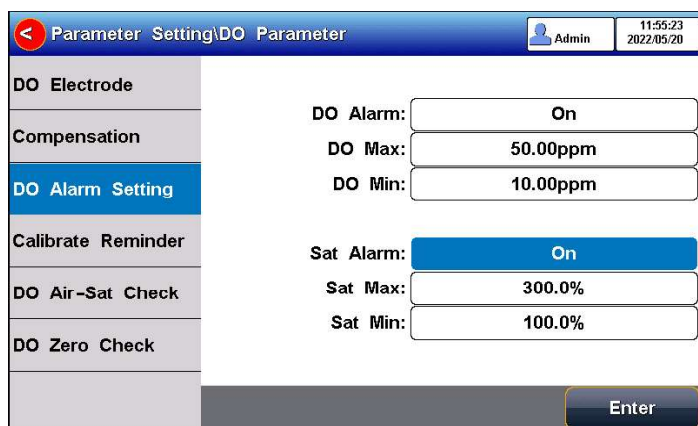
### 5.4.9.3 DO barometric compensation

Atmospheric pressure is an important factor for dissolved oxygen concentration and dissolved oxygen saturation measurement. The meter supports automatic barometric compensation and manual barometric compensation modes in the range of (600~1100) mbar. The default atmospheric pressure is 1013 mbar in manual compensation mode.

Users can select the necessary compensation mode and pressure unit for measurement in the setting.

### 5.4.9.4 Alarm setting

The meter support DO measurement and calibration result monitoring alarm. The setting includes the saturation value and monitoring options. When the measurement results out of the selected range, the meter alarms.

The screenshot shows a mobile application interface for 'Parameter Setting\DO Parameter'. The top bar includes a back arrow, the title 'Parameter Setting\DO Parameter', a user icon labeled 'Admin', and the time '11:55:23' and date '2022/05/20'. A vertical menu on the left lists: 'DO Electrode', 'Compensation', 'DO Alarm Setting' (highlighted in blue), 'Calibrate Reminder', 'DO Air-Sat Check', and 'DO Zero Check'. The main area contains the following settings:

DO Alarm:	On
DO Max:	50.00ppm
DO Min:	10.00ppm
Sat Alarm:	On
Sat Max:	300.0%
Sat Min:	100.0%

An 'Enter' button is located at the bottom right of the screen.

Figure 5-28 Alarm setting information

### 5.4.9.5 Calibration reminder settings

The meter provides a calibration prompts function. From the setting, the user can select calibration reminder (recommended) and calibration reminder (mandatory) for future calibration.

Calibration reminder (recommended): When enabling the calibration reminder, the instrument verifies whether the calibration data of the DO electrode is within the validity period. If calibration information expires, a striking calibration reminder will appear at "CAL&Check" on the main interface, but it will not affect the measurement and data saved.

Calibration reminder (mandatory): When enabling the mandatory calibration, the instrument verifies whether the calibration data of the DO electrode is within the validity period. If calibration information expires, a striking calibration reminder will appear at "CAL&Check" on the main interface, the measurement data cannot be saved and output.

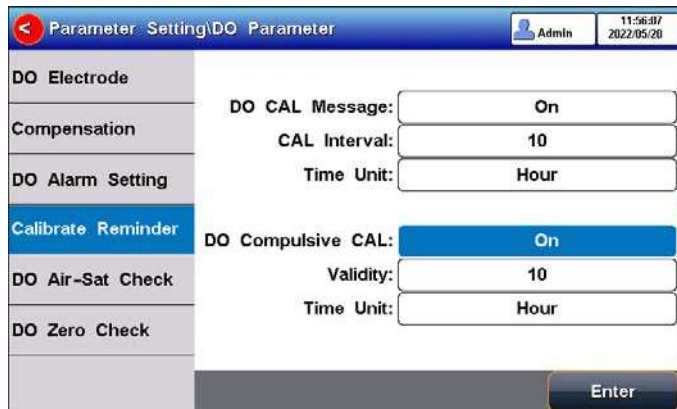
Figure 5-29 Calibration reminder setting information

#### 5.4.9.6 Verification Reminder settings

Verification reminder (mandatory): When enabling the verification reminder, the instrument verifies whether the verification data of the DO electrode is within the validity period. If verification information expires, a striking verification reminder will appear at "CAL&Check" on the main interface, the measurement data cannot be saved and output.

It allows the user to input the DO saturation/Zero standards and the allowable deviation. If the deviation amount is set to 0, it means that no verification is required.

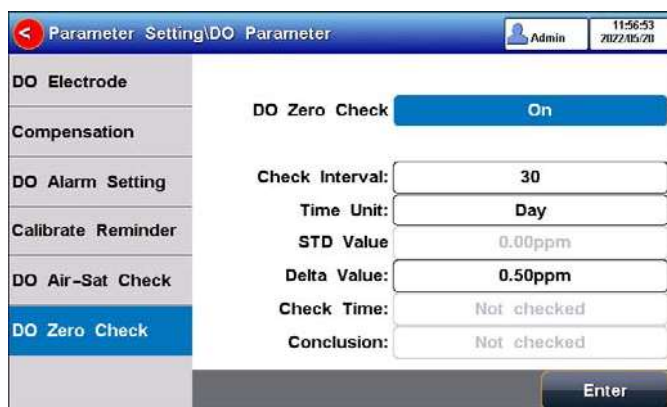
Figure 5-30 DO verification information

#### 5.4.10 Temperature Parameter Settings

The temperature unit of the meter is selectable in °C and °F. Temperature compensation mode: ATC and MTC.

ATC means automatic compensation. MTC means manual compensation. It allows the user to input the temperature.

#### 5.4.11 Data Management Settings

##### 5.4.11.1 Sample ID type

The instrument supports three setting methods of Sample ID: number order, time order, and manual.

- Number order: The sample ID No. is increasing with series number.
- Time order: The sample ID No. is increasing with sample measuring time. Format: Year/Y, Month/M, Day/D, Hour/H, Minutes/M, Second/S
- Manual: Manually set the sample ID No. It allows samples to manually enter the sample ID when saving or printing data.

##### 5.4.11.2 Autosave setting

When this function is enabled, the meter saves the results when the reading is stable in the auto-reading and interval timed reading mode.

### 5.4.11.3 Data Overwrite

The meter provides 1000 sets of measurement results storage space. When this function is enabled, the results data that exceeds capacity will overwrite the old results data.

### 5.4.12 Output option

Output to: Printer or U flash driver.

Output content and format: The number of characters per line and the title name can be set. The data formats are GLP, STD, and Custom.

### 5.4.13 User management

Figure 5-31 User management

The instrument supports user management and password security management, allowing to set the user ID and select the user type. Admin users can create sub-user for various operations.

Users can be divided into three types: Admin User, method manager user, and operator user. Create up to 8 users and support password security management. The first user is the default Admin user, and there is no default password. System administrators have permission to create users, and method administrators and operators are prohibited from creating users.

Password security management: The user password is unique. When GMP is applied, a password is required.

### 5.4.14 System Parameter Settings

#### 5.4.14.1 System Date & Time

Settings of system date and time.

#### 5.4.14.2 Buzzer setting

Users can set the key sound by this setting.

#### 5.4.14.3 Brightness setting

Users can adjust the screen brightness by this setting.

#### 5.4.14.4 Auto Power off

The meter provides an auto shutdown function. When the meter is not used, the meter switches off automatically.

#### 5.4.14.5 Restore Default

The meter supports "Restore Default" and "Restore Parameters". "Restoring Default" will restore all meter parameters to the factory state. "Restoring parameters" will restore the measurement parameters to the factory state.

### 5.4.14.6 Software version

Users can find the software version information on the general setting page.

### 5.4.15 GMP Mode

The meter supports the GMP mode, to achieve strict control of the measurement process and data traceability. In the mode, the meter enables login password protection, three-level user rights, data security protection, system log and run log, and other security control.

Press the "Setting"- "GMP Mode" to set the mode.

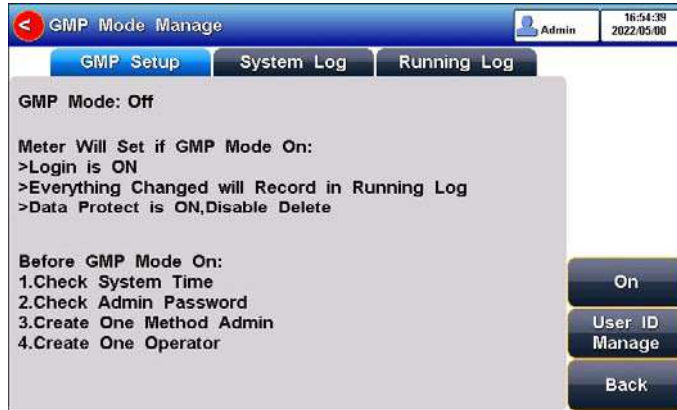
Figure 5-32 GMP mode setting

#### 5.4.15.1 System Log

Access the System Log to view the log, such as GMP mode open log, GMP mode close log, system time modification log, and creation log.

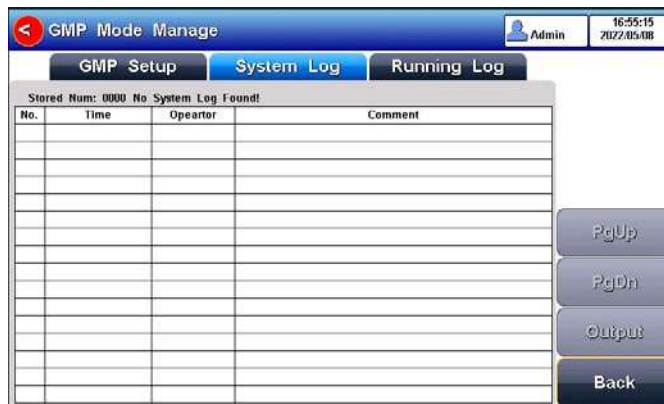
Figure 5-33 System log in GMP mode

#### 5.4.15.2 Running Log

Access the Operation Log to view the log, such as the data log and calibration log, print the data log, print the calibration log, and log in or log out of an existing account.

Figure 5-34 Operation log in GMP mode

## 5.5 pH Measurement

### 5.5.1 Calibration Preparation

The electrode slope and zero potential of pH electrodes drift slightly over time. To accurately measure pH, it is recommended to calibrate the pH electrode before use, the instrument supports 1-6 points calibration.

One-point calibration is a calibration process with a single standard solution, commonly applied in a quick test. The calibration slope is 100% here.

Two-point calibration is to use two pH standard buffer solutions to calibrate the electrode and construct a linear calibration curve through two points. Two-point calibration is the most commonly used calibration method, and it is usually recommended that the pH value of the solution be measured lies between the two standard buffer solutions. Two-point calibration can improve pH measurement accuracy.

Multi-point calibration is a calibration process with more than one standard solution. It is recommended to calibrate between two standard buffer solutions at the pH of the solution to be tested. Multi-point calibration covers a wider measurement range for accurate pH measurement. Before starting calibration, please prepare one or more pH standard buffer solutions.

### 5.5.2 Standards group selection

Before starting calibration, please prepare one or more pH standard buffer solutions. The meter has a standards recognition function. Please set the Standard Group before the measurement.

You can also set the identification mode to "Manual Mode" and manually enter the nominal value during the calibration process.

### 5.5.3 pH Calibration

Figure 5-35 pH electrode calibration information

The calibration process is as follows:

1. Select a method. The method includes the parameter (e.g., pH), NIST standard solution group, pH 4.01, pH 7.00, and pH 10.01, auto Mode recognition.
2. Connect the ATC probe or enter the temperature manually.
3. Press the "CAL & Check" -"pH Calibration".
4. Put the cleaned electrode into a pH 4.01 standard solution.
5. Wait for the instrument to display "Auto Mode Matched".
6. When the pH and temperature reading is stable, press the "Start".
7. If only 1-point calibration is required, after 1-point calibration is completed, press the "End" key to complete the calibration.

- If multi-point calibration is required, please replace the pH7.00 and pH10.01 standard buffer solutions. After cleaning the electrode, put the electrode into the standard solution. After the instrument recognizes it successfully, the instrument reads stably, press the "Next Point" to complete the calibration.
- After completing the calibration, press the "Enter" key to complete the calibration, save the calibration results and end the calibration, and directly enter the start interface. If the checked standard solution group is 8, automatically end the calibration after eight points of calibration.

### 5.5.4 pH Verification

The verification process is as follows:

- Select a method. The method includes the parameter (e.g., pH).
- Set the verification pH standard value (e.g., pH 4.01) and auto Mode recognition.
- Connect the ATC probe or enter the temperature manually.
- Press the "CAL & Check" -"pH Check".
- Put the cleaned electrode into a pH 4.01 standard solution.
- Wait for the instrument to display "Auto Mode Matched".
- When the pH and temperature reading is stable, press the "Start".
- If the pH standard value is in the allowable range, the verification result shows a pass. If the pH standard value is not in the allowable range, the verification result shows no pass.
- Press the "Save" key to save the verification results. Press the "End" key complete the verification.

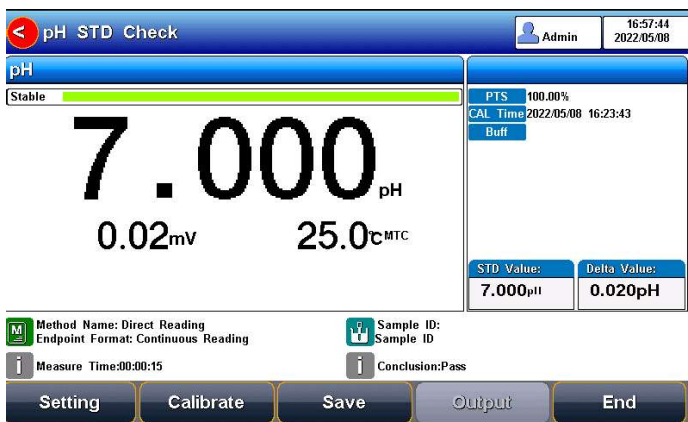
Figure 5-36 pH verification information

### 5.5.5 pH Measurement

The measurement process is as follows:

- Select a method. The method includes the parameter (e.g., pH), and reading mode (e.g., continuous reading, auto-reading, or timed format).
- Connect the ATC probe or enter the temperature manually.
- Rinse the pH electrode with DI water, and dry out.
- Put the electrode into the test solution under test.
- Put the measurement end of the electrode into the sample solution.
- Press "Measure" to enter into measurement status.
- When the reading is stable, read the results.
- Press the "Save" to save the measurement results and press the "Output" to print the result.

9. Between measurements, stored pH electrode in distilled or deionized water.
10. After measurement, rinse the pH electrode with deionized water thoroughly.

Figure 5-37 pH measurement information

## 5.6 Ion measurement

### 5.6.1 Preparation

The slope and zero potential value of the ion electrode will change with time, so the ion electrode needs to be calibrated before use, and the instrument supports up to 6-point calibration. For the specific use of the electrode, please refer to the electrode manual.

### 5.6.2 Ion-selective electrode

The ion-selective electrode is based on the ion-selective membrane, which can be divided into a single crystal membrane, salt membrane, glass membrane, and PVC ion-selective membrane. Ion-selective electrodes usually have single electrodes and composite electrodes. A single electrode can be used with different reference electrodes and has better measurement performance for some low-concentration ions. Composite electrodes are more convenient and simpler in the operation. You could be flexible to choose according to the requirements.

#### 5.6.2.1 Ionic strength adjustment buffer

The use of ion electrodes to measure ion concentration requires the addition of an ionic strength adjustment buffer.

The ionic strength of a solution has an important influence on the measurement of ion concentration. On the one hand, the ion-selective electrode directly measures the activity of the ion,  $\alpha = \gamma c$ . Wherein,  $\alpha$  is the activity of the ion,  $\gamma$  is the activity coefficient of the ion, and  $c$  is the ion concentration. Usually, the activity coefficient  $\gamma$  is affected by the ionic strength in the solution. By adding an ionic strength adjustment buffer to the standard solution and the test solution, the measured solution has a similar ionic strength to the standard solution, thereby having a similar activity coefficient  $\gamma$ . On the other hand, if a solution with low ionic strength, the potential of the reference electrode will show instability. The addition of the ionic strength adjustment buffer can help stabilize the reference electrode.

Various ion measurement needs various ionic strength adjustment buffer. Common ionic strength adjustment buffers are recommended in the following table

Table 5-4 Recommended ionic strength adjustment buffer

Ion category	ionic strength adjustment buffer
Na <sup>+</sup>	0.2 mol/L disopropylamine
F <sup>-</sup>	0.1 mol/L NaCl or TISAB
Cl <sup>-</sup>	0.1 mol/L KNO <sub>3</sub>
Br <sup>-</sup>	0.1 mol/L KNO <sub>3</sub>
I <sup>-</sup>	0.1 mol/L KNO <sub>3</sub>
Ag <sup>+</sup>	0.1 mol/L NaNO <sub>3</sub>
Cu <sup>2+</sup>	0.1 mol/L NaNO <sub>3</sub>
Pb <sup>2+</sup>	0.1 mol/L KNO <sub>3</sub>
S <sup>2-</sup>	0.1 mol/L KNO <sub>3</sub>
K <sup>+</sup>	0.05 mol/L MgAc <sub>2</sub>
Ca <sup>2+</sup>	0.1 mol/L KCl
NO <sub>3</sub> <sup>-</sup>	0.1 mol/L NaH <sub>2</sub> PO <sub>4</sub>
BF <sub>4</sub> <sup>-</sup>	0.1 mol/L Na <sub>2</sub> SO <sub>4</sub>
ClO <sub>4</sub> <sup>-</sup>	0.1 mol/L NaCl

\* The final concentration of the ionic strength modifier in the standard or sample.

### 5.6.2.2 Standard solutions preparation

The best way to prepare standards is to use serial dilutions. Sequential dilution refers to diluting an initially prepared standard using a volumetric flask to obtain a second standard. Dilute the second standard to prepare the third standard. And so on until the required standard solution is obtained.

In general, the concentration between two adjacent levels is a 10-fold relationship.

### 5.6.2.3 Activation of the ISE electrodes

When the electrode is used for the first time or has not been used for a long time, an activation is recommended. The electrode has better measurement performance after activation.

Table 5-5 Ion electrode activation solution and activation time recommendation

Ion category	Activation solution	Activation time
Na <sup>+</sup>	10 <sup>-3</sup> mol/L NaCl	2h
F <sup>-</sup>	10 <sup>-3</sup> mol/L NaF	2h
Cl <sup>-</sup>	10 <sup>-3</sup> mol/L KCl	2h
Br <sup>-</sup>	10 <sup>-3</sup> mol/L NaBr	2h
I <sup>-</sup>	10 <sup>-3</sup> mol/L NaI	2h
Ag <sup>+</sup>	10 <sup>-3</sup> mol/L AgNO <sub>3</sub>	2h
Cu <sup>2+</sup>	10 <sup>-3</sup> mol/L Cu(NO <sub>3</sub> ) <sub>2</sub>	2h
Pb <sup>2+</sup>	10 <sup>-3</sup> mol/L Pb(NO <sub>3</sub> ) <sub>2</sub>	2h
S <sup>2-</sup>	10 <sup>-3</sup> mol/L AgNO <sub>3</sub>	2h
K <sup>+</sup>	10 <sup>-3</sup> mol/L KCl	2h

Ca <sup>2+</sup>	10 <sup>-3</sup> mol/L CaCl <sub>2</sub>	2h
NO <sub>3</sub> <sup>-</sup>	10 <sup>-3</sup> mol/L NaNO <sub>3</sub>	2h
BF <sub>4</sub> <sup>-</sup>	10 <sup>-3</sup> mol/L NaBF <sub>4</sub>	2h
ClO <sub>4</sub> <sup>-</sup>	10 <sup>-3</sup> mol/L NaClO <sub>4</sub>	2h

### 5.6.2.4 Stirrer setting

The flow state of the solution influences the electrode potential of the ion-selective electrode. To improve the stability and repeatability of the measurement, it is recommended to use a stirrer to keep the flow rate of the solution stable during calibration and measurement.

### 5.6.3 Direct reading measurement

Direct Reading Measurement is a commonly used ion concentration measuring method. The Direct Reading mode uses the following Nernst formula to calculate the concentration:

$$E_x = E_0 + S \times \log(C_x + C_b)$$

Wherein,

$E_x$  ~ Equilibrium potential of the sample solution, in mV.  $E_0$  ~ zero potential value, in mV.

$S$  ~ Electrode slope (%).

$C_x$  ~ Concentration value of the sample solution, in mol/L.  $C_b$  ~ Blank concentration value in mol/L.

The electrode slope and zero potential value can be known from the calibration. When measuring the sample solution, the sample concentration result can be calculated from the formula. The direct-reading method is fast and suitable for quick test measurement of common samples.

#### 5.6.3.1 pX/ISE Calibration

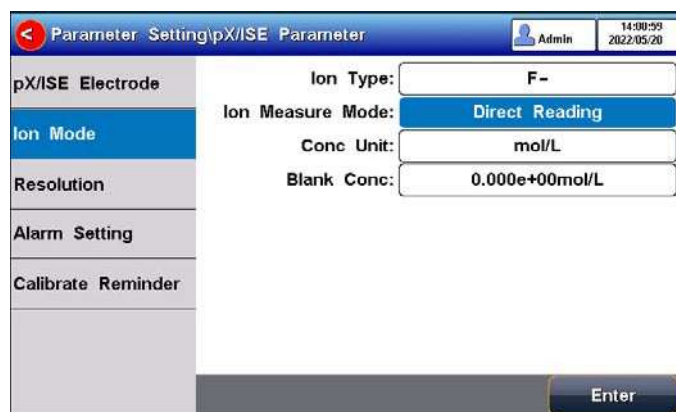
Figure 5-38 Ion measurement mode

The pX/ISE calibration process is as follows:

1. Select the "Direct Measure Method". The method includes:
  - 1) Set the parameters (e.g., pX).
  - 2) Set the ion mode (e.g., F-).
  - 3) Set the direct Reading as ion measure mode.
  - 4) Set the concentration unit (e.g., mol/L).
  - 5) Set the blank concentration (e.g., 0.000 mol/L).
2. Add an appropriate amount of standard solution (usually 100 ml) to the beaker, then add an ionic strength adjustment buffer. Adjust the stirring speed of the solution for measurement.

3. Press the F2 "CAL & Check" -"pX Calibration".
4. Put the cleaned electrode into a standard solution.
5. Press the F2 "STD value" to input the standard value of the standard solution.
6. Wait for the reading to be stable, and press the F4 "Start".
7. If only 1-point calibration is required, after 1-point calibration is completed, press the  " to complete the calibration.
8. If choosing multi-point calibration (up to 6), press "Next Point" to repeat the operation.
9. If the checked standard solution group is 6, automatically end the calibration after six points of calibration.

Figure 5-39 pX/ISE calibration information

### 5.6.3.2 Direct Reading Measurement

The measurement process is as follows:

10. Select the "Direct Measure Method". The method includes the parameter (e.g., pX), the ion type (e.g., F-), and reading mode (e.g., continuous reading, auto-reading, or timed format).
11. Or setting.
  - 1) Set the parameters (e.g., pX).
  - 2) Set the ion type (e.g., F-).
  - 3) Set the direct Reading as ion measure mode.
  - 4) Set the concentration unit (e.g., mol/L).
  - 5) Set the reading mode (e.g., continuous reading, auto-reading, or timed format).
12. Add an appropriate amount of standard solution (usually 100 ml) to the beaker, then add an ionic strength adjustment buffer. Adjust the stirring speed of the solution for measurement.
13. In the idle status, press F4 "Measure" to enter into measurement status.
14. When the reading is stable, read the results.
15. Press the "Save" to save the measurement results.
16. Press the "Output" to print the measurement result when connect to the printer.
17. Between measurements, stored ISE electrode in distilled or deionized water.
18. After measurement, rinse the ISE electrode with deionized water thoroughly.

Figure 5-40 PX/ISE measurement information

### 5.6.4 Standard addition method

Also known as the known additive method, it is mainly used to measure samples with tiny content. First, the equilibrium potential value of the system is determined, then a standard solution of known concentration is added to the system and the equilibrium potential value of the system is determined. From the potential change, the sample concentration result can be calculated by the formula as follows:

$$C_x = \frac{\rho \times C_s}{(1 + \rho) \times 10^{(E_2 - E_1)/S} - 1} + \frac{\rho \times C_b}{(1 + \rho) \times 10^{(E_{b2} - E_{b1})/S} - 1}$$

Wherein,

$C_x$  ~ the concentration value of the sample solution, in unit mol/L.  $C_s$  ~ concentration value of the standard (addition solution), in mol/L.  $S$  ~ electrode slope (%).

$C_b$  ~ blank standard concentration value in mol/L.

$E_1$  ~ The potential value before the standard addition, in mV.  $E_2$  ~ The potential value after the standard addition, in mV.

$\rho$  ~ Standard add volume ( $V_s$ ) / volume of sample to be tested ( $V_x$ ).

$E_{b1}$  ~ The potential value before the standard of calibration, in mV.  $E_{b2}$  ~ The potential value after standard of calibration, in mV.

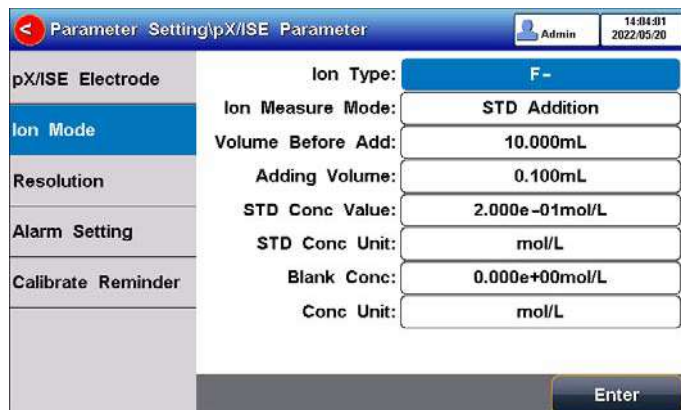
Figure 5-41 Standard addition method ion measurement mode The standard addition measurement process is as follows:

1. Select the "STD Addition Method". The method includes:
  - 1) Set the parameters (e.g., Ion Conc).
  - 2) Set the ion type (e.g.,  $F^-$ ).
  - 3) Set the Standard Addition as ion measure mode.

- 4) Set "Volume Before Add", "Adding Volume", "STD Conc Value", "STD Conc Unit", "Blank Conc" and "Conc Unit".
  - 5) Set the concentration unit (e.g., ppm).
  - 6) Set the reading mode (e.g., continuous reading, auto-reading, or timed format).
2. Add an appropriate amount of sample solution (usually 100 ml) to the beaker, then add an ionic strength adjustment buffer. Adjust the stirring speed of the solution for measurement.
  3. Put the cleaned electrode into a standard solution.
  4. Press the F3 "Measure".
  5. Press "Next" after the reading is stable.
  6. Follow the prompts to add the preset standard.
  7. Press "Next" to measure the potential value of the solution.
  8. After the reading is stable, press "Next" to end the measurement.
  9. The results are shown by auto calculation.
  10. Press the "Save" to save the measurement results.
  11. Press the "Output" to print the measurement result when connect to the printer.

### 5.6.5 Sample addition method

The method is similar to the standard addition method and is mainly used to measure samples with high content sample solutions. The standard is replaced by samples to add to the sample solution. The calculation formula is as follows:

$$C_x = C_s \times [(1 + \rho) \times 10^{(E_2 - E_1)/S} - \rho]$$

Wherein,

$C_x$  ~ the concentration value of the sample solution (addition solution), in mol/L.

$C_s$  ~ concentration value of the standard, in mol/L.

$\rho$  ~ Volume of the standard (Vs)/ Volume of the sample to be tested (Vx).  $E_1$  ~ The potential value before the sample addition, in mV.

$E_2$  ~ The potential value after the sample addition, in mV.

$S$  ~ Electrode slope (%).

Parameter Setting/pX/ISE Parameter		Admin	14:04:49 2022/05/20
pX/ISE Electrode	Ion Type:	F-	
Ion Mode	Ion Measure Mode:	Sample Addition	
Resolution	Volume Before Add:	10.000mL	
Alarm Setting	Adding Volume:	0.100mL	
Calibrate Reminder	STD Conc Value:	2.000e-01mol/L	
	STD Conc Unit:	mol/L	
	Blank Conc:	0.000e+00mol/L	
	Conc Unit:	mol/L	
		Enter	

Figure 5-42 Sample addition method ion measurement mode The sample addition measurement process is as follows:

1. Select the "STD Addition Method". The method includes:
  - 1) Set the parameters (e.g., Ion Conc).
  - 2) Set the ion type (e.g., F-).
  - 3) Set the sample addition as ion measure mode.
  - 4) Set "Volume before add", "Adding Volume", "STD Conc Value", "STD Conc Unit", "Blank Conc", and "Conc Unit".

- 5) Set the concentration unit (e.g., mol/L).
- 6) Set the reading mode (e.g., continuous reading, auto-reading, or timed format).
2. Add an appropriate amount of standard solution (usually 100 ml) to the beaker, then add an ionic strength adjustment buffer. Adjust the stirring speed of the solution for measurement.
3. Put the cleaned electrode into a standard solution.
4. Press the F3 "Measure".
5. Press "Next" after the reading is stable.
6. Follow the prompts to add the preset sample.
7. Press "Next" to measure the potential value of the solution.
8. After the reading is stable, press "Next" to end the measurement.
9. The results are shown by auto calculation.
10. Press the "Save" to save the measurement results.
11. Press the "Output" to print the measurement result when connect to the printer.

### 5.6.6 GRAN method

The measurement process is like the standard addition method, mainly for measuring samples with low content. According to the mathematical principle of GRAN mode, the ion concentration results of the sample can be calculated using the following equation:

$$\frac{(V_s + V_x) \times 10^{E/S}}{s} = 10^{E_0/S} \times \frac{(C_x V_x)}{x} + 10^{E_0/S} \times \frac{(C_s V_s)}{s}$$

Wherein,

$C_x$  ~ the concentration value of the sample solution (addition solution), in mol/L.

$C_s$  ~ concentration value of the standard, in mL.

$V_s$  ~ volume of standard, in mL.

$V_x$  ~ Volume of the sample to be tested in mol/L.

$E$  ~ The equilibrium potential of the sample solution (sample), in mV.

$E_0$  ~ zero potential value, in mV.  $S$  ~ electrode slope (%).

The screenshot shows a software interface for setting parameters for a pX/ISE measurement. The title bar reads "Parameter Setting pX/ISE Parameter" and includes a user icon for "Admin" and a timestamp of "14:05:22 2022/05/20". On the left, there is a vertical menu with options: "pX/ISE Electrode", "Ion Mode", "Resolution", "Alarm Setting", and "Calibrate Reminder". The "Ion Mode" option is currently selected. The main area on the right contains several input fields for configuration:

- Ion Type:** F-
- Ion Measure Mode:** GRAN Method (highlighted in blue)
- Volume Before Add:** 10.000mL
- Adding Volume:** 0.100mL
- STD Conc Value:** 2.000e-01mol/L
- STD Conc Unit:** mol/L
- Blank Conc:** 0.000e+00mol/L
- Conc Unit:** mol/L

An "Enter" button is located at the bottom right of the screen.

Figure 5-43 Ion measurement mode of GRAN measurement method The GRAN measurement process is as follows:

1. Select the "STD Addition Method". The method includes:
  - 1) Set the parameters (e.g., Ion Conc).
  - 2) Set the ion type (e.g., F-).
  - 3) Set the GRAN as ion measure mode.
  - 4) Set "Volume Before Add", "Adding Volume", "STD Conc Value", "STD Conc Unit", "Blank Conc", and "Conc Unit".

- 5) Set the concentration unit (e.g., mol/L).
  - 6) Set the reading mode (e.g., continuous reading, auto-reading, or timed format).
2. Add an appropriate amount of sample solution (usually 100 ml) to the beaker, then add an ionic strength adjustment buffer. Adjust the stirring speed of the solution for measurement.
  3. Put the cleaned electrode into a standard solution.
  4. Press the F3 "Measure".
  5. Press "Next" after the reading is stable.
  6. Follow the prompts to add the preset standard.
  7. Press "Next" to measure the potential value of the solution.
  8. After the reading is stable, repeat step 6 and step 7 for 3 or 8 times.
  9. The results are shown by auto calculation.
  10. Press the "Save" to save the measurement results.
  11. Press the "Output" to print the measurement result when connect to the printer.

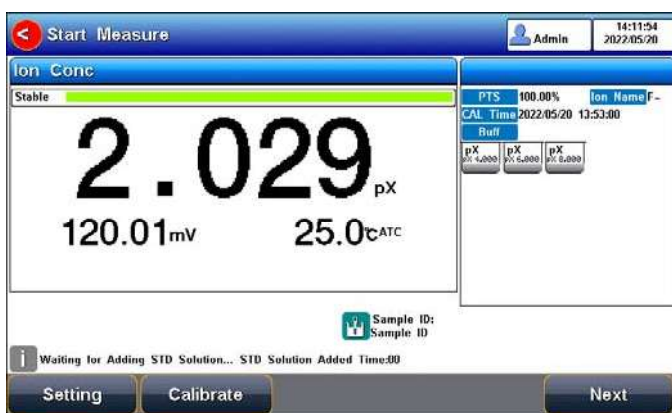
Figure 5-44 GRAN measurement information

## 5.7 ORP Measurements

### 5.7.1 Calibration Preparation

In general, ORP electrodes need few calibrations. When the electrode is used for the first time or has not been used for a long time, a calibration is needed.

The instrument supports ORP 1-point calibration, the electrode can be calibrated using the ORP calibration solution.

The instrument automatically calculates the offset and compensates for the measurement.

Before the calibration, prepare the ORP calibration standard solution.

### 5.7.2 ORP Calibration

The calibration process is as follows:

1. Select the "ORP Measure Method". The method includes:
  - 1) Set the parameters (e.g., ORP).
2. Connect the ATC probe or enter the temperature manually.
3. Press the "CAL & Check" -"ORP Calibration".
4. Put the cleaned electrode into ORP standard solution (e.g., 462mV ORP standard).
5. Press the "STD value" to input the ORP standard value.
6. When the reading is stable, press the "Start" to complete the first point calibration, and the instrument displays and stores the calibration results.

7. Press the "End" key to complete the calibration.

Figure 5-45 ORP electrode calibration information

### 5.7.3 ORP Measurement

The measurement process is as follows:

1. Select the "ORP Measure Method". The method includes:
  - 1) Set the parameters (e.g., ORP).
  - 2) Set the reading mode (e.g., continuous reading, auto-reading, or timed format).
2. Connect the ATC probe or enter the temperature manually.
3. Rinse the pH electrode with DI water, and dry out.
4. Put the electrode into the test solution under test.
5. Put the measurement end of the electrode into the sample solution.
6. Press "Measure" to enter into measurement status.
7. When the reading is stable, read the results.
8. Press the "Save" to save the measurement results and press the "Output" to print the result.
9. Between measurements, stored ORP electrode in distilled or deionized water.
10. After measurement, rinse the ORP electrode with deionized water thoroughly.

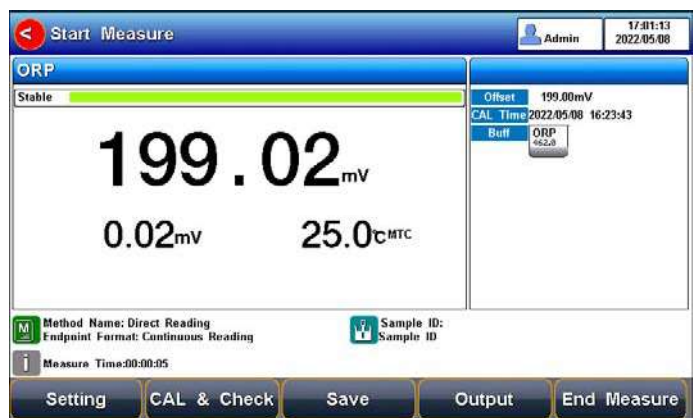
Figure 5-46 ORP measurement information

## 5.8 Conductivity Measurement

### 5.8.1 Cell Constant Input

Conductivity electrodes are precisely calibrated at the time of manufacture and marked with the exact cell constant. Before the measurement, by "Setting"- "EC Parameter" to enter the electrode cell constant.

## 5.8.2 Calibration Preparation

In general, conductivity electrodes need few calibrations. When the user gets an unexpected result, an electrode calibration is considerable.

Usually, a single standard solution is required for calibration. For accurate measurement of sample conductivity above 50mS/cm, a two-point calibration is required. Two standards are required, a low conductivity standard and a conductivity standard close to the sample.

The meter provides various Standards Group including Universal Standard Group and GB group. And allows the user to prepare the customized Standard groups.

## 5.8.3 Conductivity Calibration

Figure 5-47 Electrode cell constant calibration information

For conductivity electrodes with different cell constants, it is recommended to use the following conductivity standard solutions for calibration.

Table 5-6 KCl standards to electrode cell constants

Cell constant (cm <sup>-1</sup> ).	0.1	1	10
KCl solution Concentration (mol/L).	0.001	0.01 or 0.1	0.1 or 1

The calibration process is as follows:

1. In the idle status, press "Setting" -"ECParameter".
2. Press the "Constant type" to select the "1".
3. Press "Cal. Type" to select the "Cal by Standards".
4. Prepare one or more standard conductivity solutions (e.g., 1413μS/cm conductivity solution).
5. Prepare a thermostatic bath, and set the temperature to (25.0±0.1) °C.
6. Place a standard conductivity solution in a thermostatic bath, and set the temperature to (25.0±0.1) °C.
7. Place the conductivity electrode into a standard solution.
8. When the conductivity and temperature reading (e.g.,1413μS/cm, 25.0°C) are stable, press the "Start".
9. If choosing one-point calibration, press  to end the calibration.
10. If choosing multi-point calibration (up to 5), press "Next Point" to repeat the operation.
11. The meter saves calibration data automatically and turns to idle status.

## 5.8.4 Conductivity Measurement

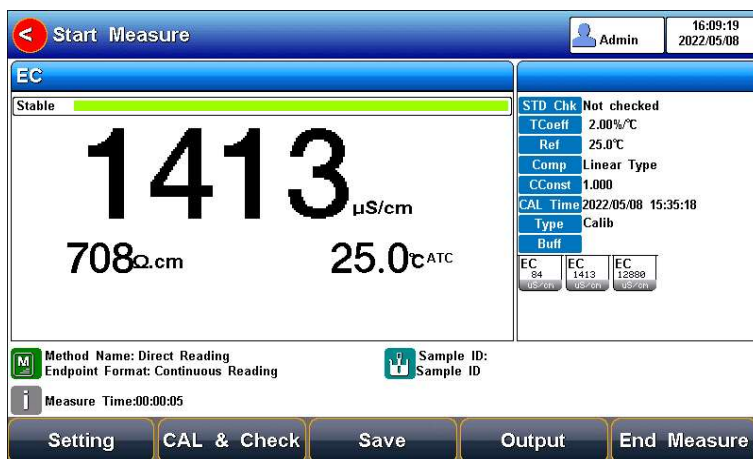
Figure 5-48 Conductivity measurement information The measurement process is as follows:

1. Setting.
  - 1) Set the parameters (e.g., conductivity).
  - 2) Set the reading mode (e.g., continuous reading, auto-reading, or timed format).
  - 3) Set the temperature compensation (e.g., Linear compensation, temperature compensation coefficient 2.00%/°C).
  - 4) Set the reference temperature (e.g., 25°C).
2. Rinse the conductivity electrode with DI water, and dry out.
3. Put the measurement end of the electrode into the sample solution.
4. In the idle status, press "Measure" to enter into measurement status.
5. When the reading is stable, read the results.
6. Press the "Save" to save the measurement results.
7. Press the "Output" to print the measurement result when connect to the printer.
8. Between measurements, stored EC electrode in distilled or deionized water.
9. After measurement, rinse the EC electrode with deionized water thoroughly and put it on the electrode protection cap.

## 5.8.5 Conductivity Verification

In the process of storage and use, the value of the conductivity standard solution may change, which will have a direct impact on the measurement results. Therefore, the verification of the standard substance is one of the important components of laboratory method development, validation, and quality control.

This meter supports the verification of conductivity electrodes. The measurement process is as follows:

1. Press "Calibrate"- "Calibration check" - "Conductivity verification"
2. Set the STD value and Delta value as the user requirement.
3. Rinse the electrode with DI water, and dry out.
4. Put the measurement end of the electrode into the sample solution.
5. When the reading is stable, read the results.
6. Press the "Save" to save the verification results.
7. Press the "Output" to print the measurement result when connect to the printer.
8. Between measurements, stored EC electrode in distilled or deionized water.
9. After measurement, rinse the EC electrode with deionized water thoroughly and put it on the electrode protection cap.

Figure 5-49 The conductivity verification information

## 5.9 TDS Measurements

### 5.9.1 TDS conversion factor

#### 5.9.1.1 Low Concentration TDS Sample

For samples with relatively simple composition and low concentration, the TDS of the solution can be estimated by conductivity. Compared with the weighing method, TDS estimation by conductivity is relatively simple and convenient with quite a good accuracy. For potassium chloride and sodium chloride solutions below 5000 µS/cm, the TDS coefficient is approximately 0.5. Therefore, 0.5 can be used as the TDS coefficient for approximate estimation in most situations.

The conversion factor adjust process is as follows:

1. Press "Setting" -"TDS Parameter".
2. Select the TDSF CAL Type as the set TDS Factor.
3. Input the TDS factor as the desired TDS coefficient. Table 5-7 Conductivity to TDS Standard Solution

Conductivity µS/cm	TDS standards		
	KCl(mg/L)	NaCl(mg/L)	442(mg/L)
23	11.6	10.7	14.74
84	40.38	38.04	50.5
447	225.6	215.5	300
1413	744.7	702.1	1000
1500	757.1	737.1	1050
2070	1045	1041	1500
2764	1382	1414.8	2062.7
8974	5101	4487	7608
12880	7447	7230	11367
15000	8759	8532	13455
80000	52168	48384	79688

1. 442 indicated the solution contains 40% Na<sub>2</sub>SO<sub>4</sub>, 40% NaHCO<sub>3</sub>, 20% NaCl.
2. The values listed in the table are values at 25°C.

For KCl and NaCl solutions below 5000 µS/cm, the TDS coefficient is about 0.5, so 0.5 can be used as an approximation in most cases.

### 5.9.1.2 High Concentrations TDS Sample Measurement

For samples with simple components and higher concentrations, such as high concentrations of NaCl solution, TDS factor re-calibration is needed.

For TDS measurements, the user may need to correct the TDS conversion factor by the TDS standard.

The conversion factor calibration process is as follows:

1. Setting.
  - 1) Set the parameters (e.g., TDS).
  - 2) Press "Setting" -"TDS Parameter".
  - 3) Select the TDSF CAL Type as the set Cal by STD.
  - 4) Set the reference temperature (e.g., 25°C).
2. Prepare TDS Standard.
3. Place a standard conductivity solution in a thermostatic bath, and set the temperature to (25.0±0.1) °C.
4. Rinse the conductivity electrode with DI water, and dry out.
5. Put the measurement end of the electrode into the sample solution.
6. Press the F2 "Calibrate"-"TDS Calibration".
7. Set the STD value as the sample STD value.
8. When the TDS and temperature reading (708ppm, 25.0°C) are stable, press the "Start".
9. If choosing one-point calibration, press  " to end the calibration.
10. If choosing multi-point calibration (up to 3), press "Next Point" to repeat the operation.
11. The meter saves calibration data automatically.

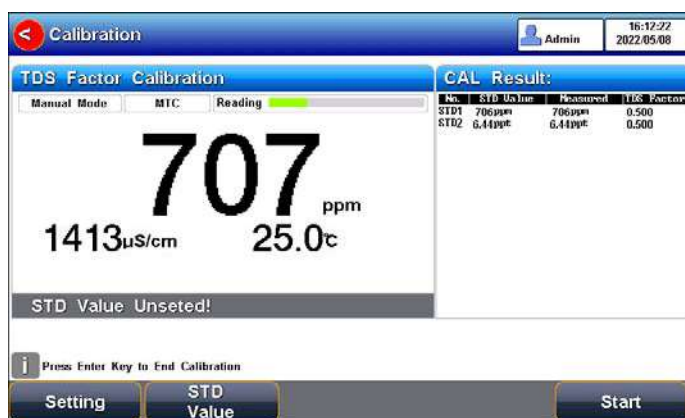
Figure 5-50 TDS coefficients calibration

### 5.9.1.3 Complex TDS Sample Measurement

For samples with complex compositions, the accuracy of TDS measurements can be improved by re-determination by laboratory methods and manual input of TDS coefficients. When the composition or concentration of the sample to be measured changes significantly, it is recommended to recalibrate the TDS coefficient.

The conversion factor calibration process is as follows:

1. Rinse the electrode with DI water. Put the measurement end of the electrode into the sample solution and set the temperature at (25.0±0.1) °C.
2. Using a weighing method to determine the TDS.
3. Calculate the TDS coefficient.
4. Press "Setting" -"TDS Parameter".

5. Select the TDSF CAL Type as the set TDS Factor.
6. Input the TDS factor as the desired TDS coefficient.

### 5.9.2 TDS Measurements

Users can switch the measurement parameter to TDS measurement by pressing the Conductivity/TDS measurement box on the screen.

## 5.10 Salinity Measurement

The instrument can be used to determine the salinity of sodium chloride. The salinity of sodium chloride can be used to approximate the salinity of the solution being measured. By measuring the conductivity of the sample, the mass percentage of the corresponding sodium chloride solution can be calculated to convert the sodium chloride salinity.

Users can switch the measurement parameter to salinity measurement by pressing the Conductivity/TDS measurement box on the screen. The detail refers to the measurement method of conductivity salinity measurement.

## 5.11 Resistivity Measurement

Resistivity and conductivity are reciprocal to each other, and conductivity can be measured at the same time when measuring resistivity.

Users can switch the measurement parameter to resistivity measurement by pressing the Conductivity/TDS measurement box on the screen. The detail refers to the measurement method of conductivity measurement.

## 5.12 Ash Measurement

### 5.12.1 White Sugar Ash

The ash content of sugar refers to the percentage of the weight of the sample that remains after the organic matter is completely removed after burning in the sugar. Ash is an important indicator of white sugar. If the ash content exceeds the standard, the grade of white sugar will drop.

The conductivity ash measurement method is to dissolve a certain quality of white sugar in a certain volume of water to measure the conductivity value of the solution. Conductivity ash is calculated at 20°C as follows:

$$C = 6 \times 10^{-4} (C_1 - 0.35C_2)$$

Wherein,

C: Conductivity ash (%)

C<sub>1</sub>: 31.3 g/100 ml of the conductivity of the sugar solution at 20 °C, μS/cm.

C<sub>2</sub>: The conductivity of the distilled water at 20°C for sugar dissolving, μS/cm.

When the temperature is lower than 20 °C, the following formula can be used to compensate, generally applied at 20±5 °C.

$$C_{20^\circ\text{C}} = C_t / [1 + 0.026(t - 20)]$$

Wherein,

C<sub>20°C</sub>: Conductivity value at 20°C, μS/cm.

C<sub>t</sub>: The actual measured conductivity value, μS/cm.

### 5.12.2 Fructose Glucose Syrup Ash

Fructose syrup Ash is measured similarly to the white sugar ash measurement method. Fructose syrup Ash calculation formula is as follows:

$$C=K*(C_1-K_1*C_2)$$

Wherein:

C: Fructose glucose syrup ash, %(g/100g).

K: The conversion coefficient of the solution conductivity converted to ash; the default is  $7.93 \times 10^{-4}$ .

K1: The correction coefficient of the solution with deionized water, default 0.39.

C1: 25% conductivity of the sugar solution at 25°C,  $\mu S/cm$ .

C2: Conductivity of dilution with deionized water at 25 °C,  $\mu S/cm$ .

### 5.12.3 Blank Measurement

There are 2 ways to measure blank values. It allows to input the blank value manually and measures the blank value in the calibration menu.

Manually enter the blank value for DI water in Parameter Settings - Conductivity Settings-Setting Ash Measurement Parameters.

The blank measurement process is as follows:

1. Select the Ash parameter.
2. Press "Calibrate"- "Calibrate ash blank conductivity".
3. After the reading is stable, press "Start" to complete the measurement of the blank conductivity.

### 5.12.4 White Sugar Ash Measurement

The white sugar ash measurement process is as follows:

1. Calibrate the EC electrode.
2. Blank Measurement.
3. Select the "White Sugar Ash Method" in "Method Management".
4. Dissolve a certain amount of white sugar in a certain volume of water (according to the ratio of 31.3g/100ml of sugar liquid).
5. Press "Measure" to start a measurement.

### 5.12.5 Fructose Glucose Syrup Ash Measurement

The white sugar ash measurement process is as follows:

1. Calibrate the EC electrode.
2. Blank Measurement.
3. Select the "Fructose Slurry Ash Method" in "Method Management".
4. Dissolve a certain amount of white sugar in a certain volume of water (according to the ratio of 31.3g/100ml of sugar liquid).
5. Press "Measure" to start a measurement.

## 5.13 Dissolved Oxygen Measurement

### 5.13.1 Calibration Preparation

Zero oxygen is calibrated using a freshly formulated 5% sodium sulphite solution as oxygen-free water.

#### 5.13.1.1 Preparation

Zero oxygen is calibrated using a freshly formulated 5% sodium sulphite solution as oxygen-free water.

### 5.13.1.2 Electrodes preparation

When the DO electrode (polarographic) is used for the first time or has not been used for a long time, the filling solution needs to be replaced.

Follow the below instruction to replace the filling solution:

1. Take the cap off the electrodes, rinse the cap with DI water and dry out.
2. Rinse the inner electrode with DI water and dry the electrode.
3. Add the filling solution (electrolyte) into the membrane cap up to 3/4.
4. Install the cap onto the electrode.

### 5.13.1.3 Electrode Polarization

Polarographic DO electrodes need to be polarized before every use.

Connect the DO electrode to the meter and turn on the meter, wait for 1 hour and the electrode are auto polarized. When the electrodes are unplugged from the meter for no more than 1 hour, measurements are allowed after 25 minutes of polarization.

### 5.13.2 DO Calibration

After the measurement parameters are selected, press "CAL&Check"- "DO Calibration".

The calibration process is as follows:

1. Rinse the electrode with DI water, and place it into an oxygen-free solution.
2. When the reading is stable, press the "CAL Zero" to complete the zero calibration. Users can change the calibration parameters by pressing the parameter setting during the calibration.
3. Rinse the electrode with DI water again, and place the probe in the upper part of a bottle filled with air-saturated (well shaken) water.
4. When the reading is stable, press the "Air-Sat" to complete the air calibration.

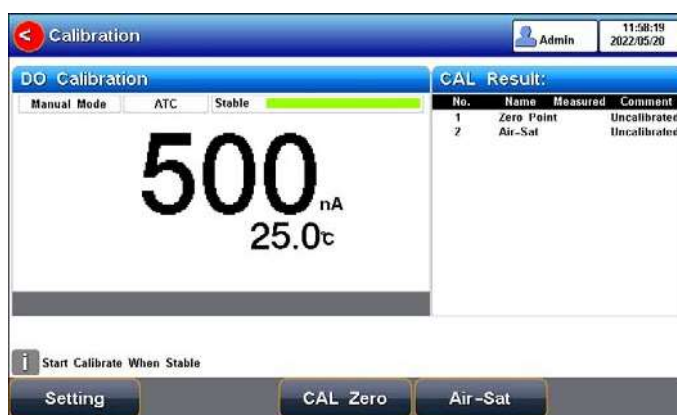
Figure 5-51 Dissolved oxygen calibration information

### 5.13.3 DO Measurement

After the calibration, press "Measure" to start a measurement. The measurement process is as follows:

- Put the electrode into the test solution under test and shake the electrode gently in a circle, in a circular motion, or use a stirrer to avoid air bubbles during the process.
- When the reading is stable, end the measurement and save the results.

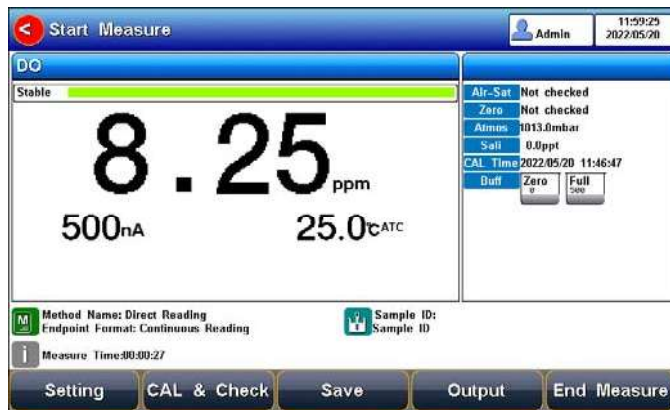
Figure 5-52 Dissolved oxygen measurement information.

### 5.13.4 DO standards verification

DO standard solutions are changed during storage and usage, and this change will have a direct impact on the measurement results. So, the verification of standards is important for accurate results.

#### 5.13.4.1 Air-Saturation standards preparation

This meter provides air-saturation standard verification. Before the measuring, the user can do a standard verification by "CAL&Check" - "Air-Satu standard verification", the process is as follows:

- Rinse the electrode with DI water. Put the measurement end of the electrode into the verification solution.
- Set the preset standard value and the allowable deviation.
- When the reading is stable, end the verification and save the results

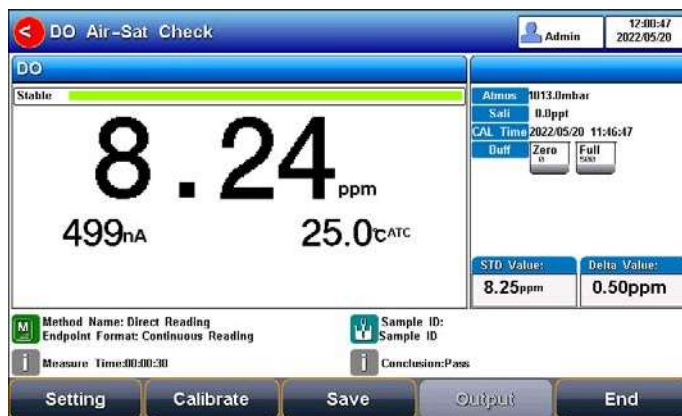
Figure 5-53 Air-Saturation standard verification

#### 5.13.4.2 Zero oxygen standards preparation

This meter provides zero-oxygen standard verification. Before the measuring, the user can do a standard verification by "CAL&Check" - "DO Zero standard verification", the process is as follows:

- Rinse the electrode with DI water. Put the measurement end of the electrode into the verification solution.
- Set the preset standard value and the allowable deviation.
- When the reading is stable, end the verification and save the results.

Figure 5-54 Zero-oxygen standard verification

## 5.14 Saturation Measurement

Dissolved oxygen saturation refers to the ratio of the dissolved oxygen concentration to the saturated dissolved oxygen concentration under the same conditions.

Users can switch the measurement parameter to Saturation measurement by pressing the DO measurement box on the screen.

## 5.15 Batch Sample Measurement

Batch sample measurement is suitable for multiple sample measurements to improve work efficiency. This function needs to match the automatic sampler to achieve multi-sample measurement.

Enter the batch sample measurement function by "Sample List".

## 5.16 Data Management

Press "Data" to view the detail of the results.

The meter stores the measurement results independently according to the measured parameters. The meter provides data Storage of 1000 sets for each parameter (pH/pX/ISE/ORP), 100 sets of calibration data, and 1000 sets of verification data.

### 5.16.1 Filter

The data type includes the measurement data and verification data, and calibration data. The parameter can be selected by pH, pX, ISE, and ORP. By the lookup filter, the user can view the result data by storage number, time, operator, method name, sample ID, and electrode Name. The filter data shows in a graph or curve. By the filter setting, press "Start Search" to look up the data.

Figure 5-55 Data setting view

### 5.16.2 Result

In the result page, press "Page Up", "Page Down", "<<" and ">>" to choose data. Users can press "Delete" to delete the current result.

#### 5.16.2.1 Measurement Data

In the measurement data, the result is displayed as a graph or a sheet, and the coordinates correspond to the relevant information.

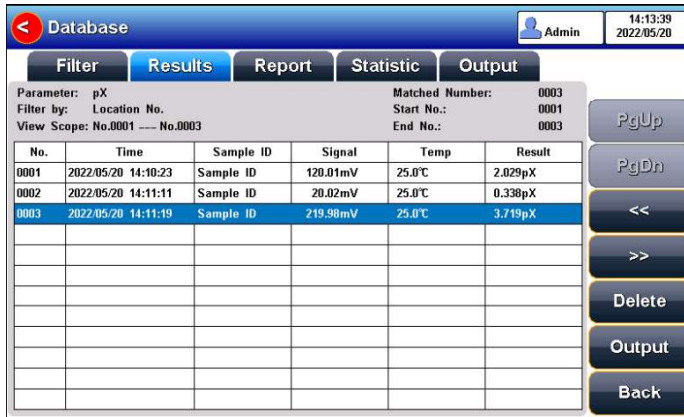
Figure 5-56 Matched measurement data in a sheet

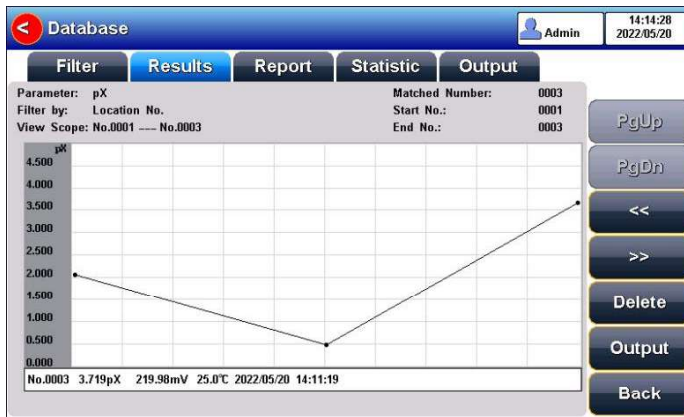
Figure 5-57 Matched result data results in a graph

#### 5.16.2.2 Verification Data

In the verification data, the pH verification result is displayed as a sheet, the coordinates correspond to the relevant information.

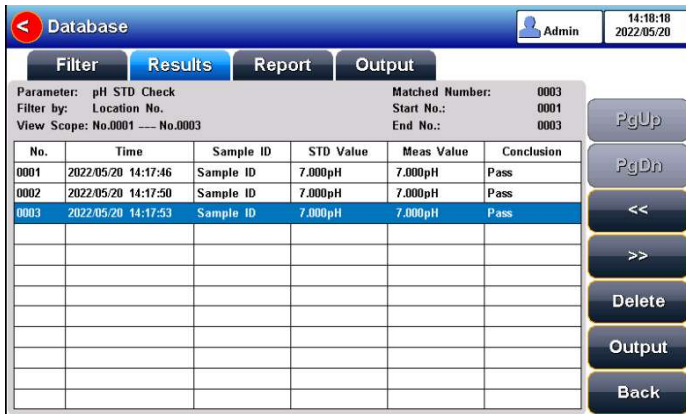
Figure 5-58 Matched verification data results in a sheet

### 5.16.2.3 Calibration Data

In the calibration data, the pH calibration result, pX calibration result, or ORP calibration result is displayed as a sheet, and the coordinates correspond to the relevant information.

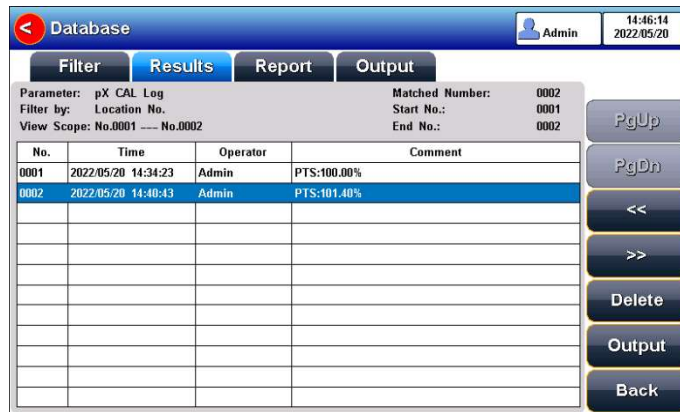
Figure 5-59 Matched calibration data results in a sheet

### 5.16.3 Report

The report shows relevant information of the data in detail. It includes the result, calibration result, and basic information.

Figure 5-60 Measurement result report

### 5.16.4 Statistic

The meter supports the statistical analysis function. By pressing "Statistic", the meter calculates basic statistical results, including maximum, minimum, average, standard deviation, relative standard deviation, and related statistical information of the results.

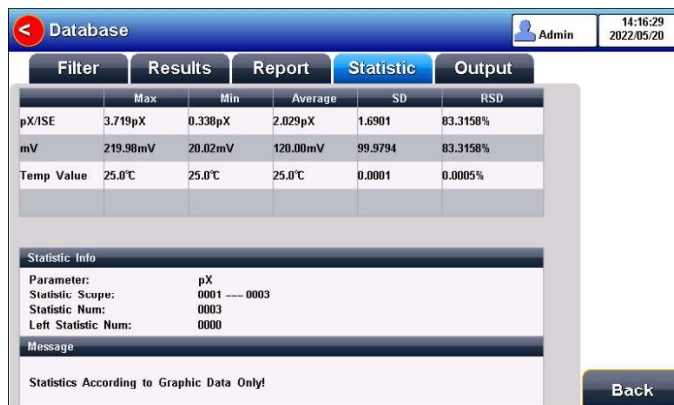
Figure 5-61 Statistical analysis

## 5.16.5 Output Settings

It is the same as the output setting.

The output device can be a printer or U disk.

Output the measurement results by selecting the data format. Format: Standard, GLP, and custom.

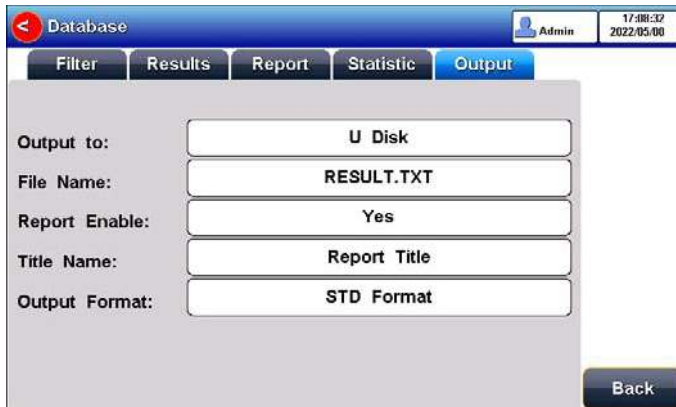
Figure 5-62 Output settings

The output format is as follows:

```
*****
Report Title
-----
Measure Time:2021/01/19 12:27:28
Operator: Operator 1
Model: M500T Multi-parameter Analyzer
Serial Number:
SW Version: Ver 1.00
-----
..... MATCHED INFO
Stored Num:                28
Matched Num:               1
Stored No. :               15
-----
..... CALIB INFO
Calib Operator:           REX Team
Calib Time:               2020/05/13 08:30:00
Calib Num:                3
-----
..... CALIB RESULT
STD 1:  4.01pH    176.8mV    25.0c
STD 2:  7.00pH     0.0mV    25.0c
STD 3: 10.01pH   -178.1mV    25.0c
pH Slope 1:                100.00%
pH E0 1:                   0.0mV
pH Slope 2:                100.00%
pH E0 2:                   0.0mV
-----
..... BRIEF INFO
Reading Mode:              Timed Reading
Stable Type:              Medium
Temp Comp Type:           ATC
-----
..... SAMPLE INFO
Sample ID:                 Sample 1
-----
..... RESULT
Result:                   7.000pH
Signal Value:             -0.0mV
Temp Value:               25.0c
-----
Signature:
```

## 6 Maintenance/Troubleshooting

### 6.1 Meter Maintenance

The correct use and maintenance of the instrument can ensure the accurate and reliable performance of the instrument. Additionally, exposure to chemicals or harsh use environments can affect performance.

- If the meter is not used for a long time, please disconnect the power supply.
- The electrode socket of the instrument must be kept clean and dry, and should not be in contact with acid, alkali, and salt solutions.
- Keep the meter and accessories clean and away from acids, alkalis, and any corrosive solutions/gases.
- Users can clean the meter surface with clean water and detergent.
- When the meter is transported, please follow the instructions:
  - please remove all connected cables.
  - Please remove the electrode holder.
  - Please use the original packaging in the long-distance transport to avoid damage.

### 6.2 Electrodes Maintenance

For more detailed information, please refer to the electrode instruction manual.

### 6.3 Troubleshooting

Table 6-1 Troubleshooting

Phenomenon	Probable reasons	Solutions
1. No Display	Not powered on. Damage to the meter.	Connect the adapter and press the power key to turn it on. Replace or repair as required.
2. Incorrect mV measurement is	1. The electrode is out of service life 2. The electrode plug is in poor contact	1. Replace the electrodes 2. Connect the protection plug, if the potential is not 0mV, please contact the after-sales service.
3. Incorrect pH measurement	1. Refer to as 2.2 2. Refer to as 2.2 3. The electrodes are not calibrated or are calibrated incorrectly	1. Refer to as 2.2 2. Refer to as 2.2 3. Recalibrate the electrode or replace the standard solution
4. Incorrect pX/ISE measurement	1. Refer to as 2.2 2. Refer to as 2.2 3. The electrodes are not calibrated or are calibrated incorrectly 4. Incorrect ISE probe	1. Refer to as 2.2 2. Refer to as 2.2 3. Recalibrate the electrode or replace the standard solution 4. Buy the correct ISE probe. Add ionic strength adjustment buffer.
5. Incorrect conductivity measurement	1. The electrode is out of service life 2. The electrodes are not calibrated or are calibrated incorrectly	1. Replace the electrodes 2. Recalibrate the electrode or replace the standard solution
6. Incorrect DO measurement	1. The electrode is out of service life 2. The electrodes are calibrated incorrectly 3. Stirring speed is not good	1. Replace the electrodes 2. Recalibrate the electrode 3. Keep the bubbles in the stirring state consistent

If the meter still does not work, please contact your local dealer for further assistance.

## 7 Technical Supports

### Accessories

Please refer to the accessories table for purchasing recommendations.

Table 7-1 Meter accessories

Name	Description
E-301-QC 3 in 1 pH composite electrode	pH Measurement Probe
501 ORP composite electrode	ORP Measurement Probe
PF-2-01 Fluoride ion electrode	Measure the fluoride ion content
232-01 Reference electrode	Use with ion electrode
T-818-Q Temperature electrode	Measure temperature value
DJS-1VTC conductivity electrode	Conductivity, TDS Measurement Probe
REX-6 Electrode stand	Place electrodes during the measurement
pH 4.01/7.00/10.01 standard sachets	To prepare the standard
462mV/296mV/419mV ORP calibration reagent	Calibrate ORP Electrodes
Conductivity solution 1413 $\mu\text{s}/\text{cm}$	Standard solution

## 8 Appendixes

### Appendix 1

pH-Temperature Relationship Table of pH Standard Solutions

Temperature(°C)	1.68	4.01	7.00	10.01
5	1.67	4.00	7.09	10.25
10	1.67	4.00	7.06	10.18
15	1.67	4.00	7.04	10.12
20	1.68	4.00	7.02	10.06
25	1.68	4.01	7.00	10.01
30	1.68	4.01	6.99	9.97
35	1.69	4.02	6.98	9.93
40	1.69	4.03	6.97	9.89
45	1.7	4.04	6.97	9.86
50	1.71	4.06	6.97	9.83

## Appendix 2

Approximate concentrations of KCl solutions and their conductivity values relationship

T(°C)	84μS/cm	1413μS/cm	12.88mS/cm
5	53.02	896	8.22
10	60.34	1020	9.33
15	67.61	1147	10.48
20	75.80	1278	11.67
25	84.00	1413	12.88
30	92,19	1552	14.12
35	100.92	1696	15.39

### Notice to customers:

The product is made up of various components and various materials that must be recycled or, failing that, deposited in the corresponding debris removal sites when the product's life has been completed or when otherwise it is necessary to dispose of it. To do this, the end user who acquires the product must know the current regulations of each municipality and / or locality based on the waste electrical and electronic equipment. The user who acquires this product must be aware of and responsible for the potential effects of the components on the environment and human health as a result of the presence of hazardous substances. Never place the product in a conventional container of citizen scope if a previous dismantling and knowledge of the components that incorporates. If you do not know the procedure to follow, consult with the city council for more information.