CENTRÍFUGA CENTRIFUGE

**7001485
7001486
7001476
7001477**

**MACROTRONIC BLT
MACROFRIGER BLT
MEDITRONIC BLT
MEDIFRIGER BLT**

Indice

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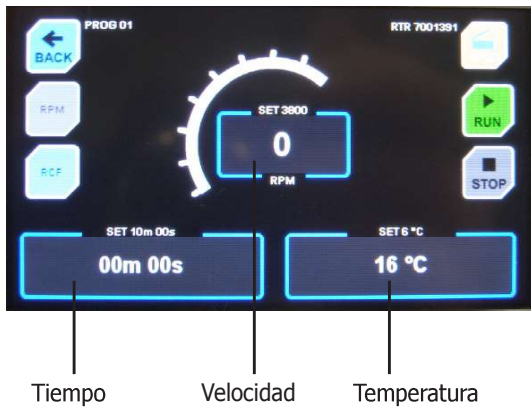
ENGLISH VERSION

1 Safety

The equipment includes appropriate security measures.
Risky situations that must be respected are indicated along this manual.

1.1 Safety icons

The safety icons identify the risky situations and the safety measures that must be taken.

Risk of danger
Electrical risk

Respect the indicated instructions to perform the operation.

Electrical risks

Risk of electrical accident when accessing the areas indicated by this sign or when performing the operations indicated by this icon in this manual.

Respect the indicated instructions to perform the operation.

Important information

- Important information for good results or for an optimum equipment operation.
- Important information to extend life of the equipment or to prevent degradation of any of the components.

1.2 Personnel qualification

This equipment can only be used by personnel who have been properly qualified.

This equipment can only be used by personnel who have read and understood these instructions or who has been properly qualified in the operation of this equipment.

2 General information

Handle the package with care. Unpack and check that the contents coincides with the "Packing List". If any part is damaged or missing, please advise the distributor immediately.

Do not install or use the equipment without previously reading this instructions manual. They are an inseparable part of the equipment and should be available to all users of the machine. Any questions can be clarified by contacting J.P. SELECTA's technical service.

Any modification, deletion or lack of maintenance of the device could cause damages. The manufacturer is not liable for the damages that may arise.

Do not use the equipment with fluids which can give off vapors or form explosive or flammable mixtures.

Maintenance

Manipulation of electronic circuits by unauthorized persons may cause damages difficult to repair. Check with J.P. SELECTA's technical service.

To clean the accessible parts, use cleaner or a similar product. Do not use aggressive solvents.

Guarantee

This product is guaranteed for one year. The warranty does not cover damages caused by an improper use or causes beyond J. P. SELECTA, s.a.u.

Any manipulation of the equipment by an unauthorized personnel cancels the warranty benefits.

3 Technical specifications

Voltage supply 115/230V 50/60 Hz as indicated in the machine's features plate.

Code:	7001476	7001477	7001485	7001486
Dimensions (mm): (Height/Width/Depth)	380x460x520	380x680x520	450x600x660	970x600x660
Power (W):	520	1230	720	1230
Max speed (rpm):	11000	15000	10000	12000
Max RCF:	25470	(Depending on the rotor)	11066	15935
Weight (kg):	44	70	76	156

4 Packing List

4.1 Standard package content

The standard equipment consists of the following components:

- | | |
|--|-----------------|
| - Centrifuge MEDIFRIGER/MACROFRIGER BLT | 7001477/7001486 |
| - Centrifuge MEDITRONIC/MACROTRONIC BLT | 7001476/7001485 |
| - Network connection cable | 7140 |
| - Rotor nut wrench | 5377 |
| - Drainage nipple in refrigerated models | |
| - Instructions manual | 80316 |

5 Introduction

The centrifuge has an induction motor maintenance free and a maximum noise level of 60 dBA. It has an internal security camera and a bucket, stainless steel cover.

The electronic circuit is controlled by a microprocessor that controls the operating parameters: speed, acceleration, braking, RCF, operating time and temperature on models 7001486/7001477. It can store up to 10 operating programs.

6 Installation

6.1 Location

- Place the equipment near a power socket suitable for consumption and over a flat level surface and strong enough to support the equipment weight and possible vibrations.
- The surface must allow the adherence of the centrifuge legs.
- Leave a free space of 30 cm. around it (see danger indication).
- This equipment can cause vibrations that could affect other nearby equipments.
- Do not install the centrifuge near heat sources or in places where it is exposed to direct sunlight, as this may affect temperature regulation.
- Place the nipple provided in outlet (16) and connect to a hose to drain condensates.
- Equipments with wheels, ensure that they are braked prior to start-up.

FREE SPACE AROUND CENTRIFUGE

Safety zone:

Leave a free space of 30 cm. around the centrifuge.

With the centrifuge running:

- Do not place hazardous substances in the security area.
- The staff should be outside the safety area.

DO NOT INSTALL THE EQUIPMENT IN EXPLOSIVE ATMOSPHERES OR PLACES WHERE SUBSTANCES LIABLE TO FORM EXPLOSIVE OR FLAMMABLE MIXTURES ARE USED OR STORED.

6.2 First lid opening and emergency opening

Upon receiving the centrifuge, the network connection cable is inside the tank and therefore the only way to open the equipment is following these instructions:

- Remove the equipment side plastic caps.
- Gently push from both caps, and the lid will open.

USE THIS OPENING SYSTEM FOR EMERGENCY OPENING ONLY.

NEVER use the device emergency opening when the engine is rotating.

6.3 Connection to the mains

Make sure that the equipment is connected to the mains and that the voltage coincides with the indicated on the features plate.

Do not use an equipment without being earthed.

Only use the supplied cable for network connection.

6.4 Rotor positioning

Before starting the centrifuge, ensure that you have placed the rotor in a proper way, otherwise it could cause damages on the equipment.

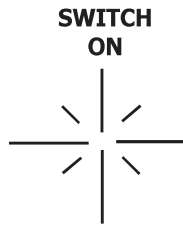
Make sure to fit the pin of the motor shaft in its housing.

Firmly tighten the nut of the rotor with the provided key.

Place the adapters and tubes evenly.

OPERATING TIME:	00004h 10m
ROTOR TIME:	00004h 10m

Example of accumulated working times of components with limited service life

7 Operating menus

Press the ON/Off switch.

The hour counter of the centrifugal elements appears with limited service life.

Operating time

Rotor time*

** It is not possible to use the centrifuge if there is no mounted head. The device has a head reader that identifies it to select the correct parameters. Without head is locked.*

After a few seconds, the main menu screen appears.

You can select from the main menu, by pressing on the touch screen:

SETTINGS: To update date, time, language, data download, software updates, life counters initialization.

PROGRAMMING: To define the programs for samples treatment and to select the active program:

- Times
- Speed
- Temperature
- Rotor

WORKING: To operate the selected program and see its evolution.

Main menu screen

7.1 Settings

Set the following settings in this screen::

LANGUAGE: Spanish, English & French.

DATE & TIME.

TEMPERATURE UNITS in °F or °C.

ENTER / CHANGE PASSWORD: Password access for software updates or restart life counters.

ROTORS: View the list of rotors available for this equipment. Rotor in use will be marked with an (*) next to the code. The rotor type is selected in the work program.

Return to main menu

Access authorization for firmware update.

7.2 Programming

In this menu the program parameters are set. Select the program to edit.

- Rotor key: Enter the rotor identification. This limits the programming to avoid dangerous situations.
- Temperature key: To set the working temperature (only refrigerated models)
- Save key: To save the data.
- Home key: To return to the main menu.
- Edit key: To edit "Phases".

Press "Edit" to edit the "Phases". There must be at least 2 phases:

1- Acceleration (slope)

Adjust the working speed and acceleration time to reach that speed (slope).

Set the time that will be maintained at the programmed speed (rpm).

Up to 4 additional steps can be added, and up to the braking phase (total 6 steps).

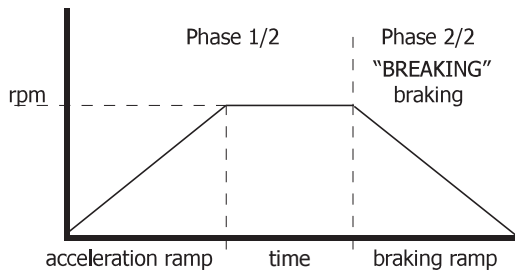
Adjust as in the acceleration phase, speed, time and acceleration time. To add steps, press the + key.

2- Brake phase

The last phase of the program is always the braking phase. In this phase it is only allowed to adjust the braking time (since the final speed will be 0 rpm).

Press the "SEL" icon to select the current program as the active one.

Note: All programs are modifiable by the user.

7.3 Run a program (WORKING)

From the main menu, press on «WORKING».

Program selected in the former paragraph will be running.

Press on «RUN» to start the program.

The program starts and displays the following information:

- Program and rotor number.
- Current and target speed.
- Elapsed (minutes, seconds) and target time.
- Temperature in the bucket and target temperature.

Press on «RPM» so that the display shows «RCF».

RPM: Revolutions per minute.

RCF: Number of «g» (acceleration of gravity) to which the sample is subjected to.

The program notifies the end with a long beep.

Press "STOP" to end the program execution.

Press "OPEN" to open the lid.

Press "BACK" to return to the main screen.