**DRYGLASS**

**2000381**

**ESTUFA PARA SECAR VIDRIO  
GLASS DRYING OVEN**

**INDICE**

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## GENERAL INFORMATION

- 1) Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.
- 2) Do not install or use the equipment without reading this handbook before.
- 3) This handbook must always be attached to the equipment and it must be available for all users.
- 4) If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.
- 5) IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 2009/104/UE, the manufacturer is not responsible for the damage that can occur.
- 7) Do not use the equipment with liquids which can give off vapours capable of making explosive mixtures.

## PACKING LIST

The standard equipment consists of the following components:

| Piece               | Reference | Quantity |
|---------------------|-----------|----------|
| Dryglass            | 2000381   | 1        |
| Glass doors         |           | 2        |
| Shelf               | 2000091   | 2        |
| Shelf runner        | 2000081   | 4        |
| Instructions Manual | 80025     | 1        |

## ACCESSORIES

(The following components, at user's request, must be set in factory)

### Denomination

### Code

TIMER: 0-120 minutes.

2000002

Installed in the control panel, it disconnects the oven once the preset time has finished.

PROGRAMMABLE TIMER OF 24 HOURS 2000009

For continuous cycle with "ON-OFF" switch levers that can be operated every 15 minutes.

## TECHNICAL FEATURES

Voltage supply 115-230 V 50/60 Hz as indicated on the characteristics plate.

| Code    | Capacity (L) | Internal measures (cm) |       |       | External measures (cm) |       |       | Self runner device | Consumption (W) | Weight Kg. |
|---------|--------------|------------------------|-------|-------|------------------------|-------|-------|--------------------|-----------------|------------|
|         |              | Alto                   | Ancho | Fondo | Alto                   | Ancho | Fondo |                    |                 |            |
| 2000381 | 126          | 45                     | 70    | 40    | 66                     | 94    | 54    | 8                  | 3000            | 65         |

## EQUIPMENT DESCRIPTION

Glass drying ovens DRYGLASS are designed with fan circulation for adjustable temperatures from 40°C up to 170°C.

Hydraulic thermostat for temperature control and air circulation by turbo fan.

Removable sliding doors of tempered glass.

## INSTALLATION

Put the oven on a flat, horizontal, level surface leaving a space of about 10cm at the back side of the equipment.

For a better oven function, do not situate it in places where it can be directly exposed to the sun radiation or heat sources.

For an optimum function, the ambient temperature must not be over 30°C.

**CAUTION!!! IMPORTANT FOR YOUR SAFETY**

**Do not use the oven to dry or for thermal treatment of products which can give off vapours capable of making an explosive mixture.**

**Be sure that the voltage supply is the same as the one indicated on the characteristics plate.**

**Do not use the equipment if it is not earthed.**

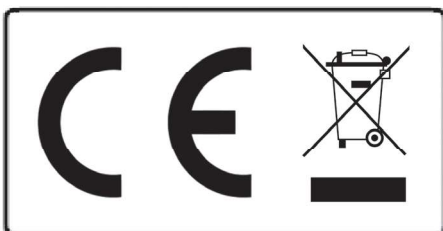
**If you change the plug, bear in mind the following**

|                            |                 |
|----------------------------|-----------------|
| <b>Blue cable:</b>         | <b>Neutral.</b> |
| <b>Brown cable:</b>        | <b>Phase.</b>   |
| <b>Yellow/green cable:</b> | <b>Earth.</b>   |

**CAUTION!!! DUE TO  
HOT WORKING TEMPERATURES  
OF THE OVEN, BE CAREFUL TO AVOID BURNS  
WHEN USING THE EQUIPMENT.****GUARANTEE**

This product is guaranteed for one year. The guarantee does not cover damages caused by an incorrect use or causes beyond the control of J.P. SELECTA,s.a.u.

Any manipulation of the equipment by personnel not authorized by J.P. SELECTA, s.a.u. automatically cancels the guarantee.

**Notice to customers:**

The product is made up of various components and various materials that must be recycled or, failing that, deposited in the corresponding debris removal sites when the product's life has been completed or when otherwise it is necessary to dispose of it. To do this, the end user who acquires the product must know the current regulations of each municipality and / or locality based on the waste electrical and electronic equipment. The user who acquires this product must be aware of and responsible for the potential effects of the components on the environment and human health as a result of the presence of hazardous substances. Never place the product in a conventional container of citizen scope if a previous dismantling and knowledge of the components that incorporates. If you do not know the procedure to follow, consult with the city council for more information.

## OPERATION

### CONTROL PANEL:

1. Main switch
2. Temperature regulator thermostat
3. Mains indicator lamp (green).
- 4a. Heater operation indicator lamp (amber) 50%.
- 4b. Heater operation indicator lamp (amber) 100%.
5. [OPTIONAL] 24 hour timer or timer control.
6. Safety thermostat adjustment knob.
7. Safety thermostat indicator lamp.
8. Analog thermometer temperature reader.
9. Locking gear. Thermostat control immobilization device.
10. [OPTIONAL] End of cycle signaling lamp.
11. Power selector 50% / 100%.
12. Safety thermostat reset push-button.
13. Cyclical or one-time programming selector.

**NOTE:** The panel shown includes an optional timer supplied as an accessory.

### STARTING UP:

1. Press the mains switch (1). The green indicator lamp will light up (3).
2. Turn the temperature control knob (2) clockwise until the working temperature coincides with the pointer. The amber indicator lamp (4) will light up.
3. Block the temperature control knob (2) with the blockage device (9) to avoid accidentally moving the knob.
4. After a certain time, the heater operation indicator lamp (4) will start operating intermittently. When these intermittences become uniform, the temperature of the oven is stabilized.
5. The thermostat scale is intended as a guide, the real temperature level inside the oven is the one that the thermometer (8) shows.

\* This oven is equipped with a power selector (11) which permits the use of the oven at 50% or at 100% of its power.

### ADJUSTING THE SAFETY THERMOSTAT:

To adjust the safety temperature thermostat, do as follows:

1. With the oven connected to the mains, turn the safety thermostat control knob (6) clockwise to the end of the scale.
2. Ensure that the safety thermostat is reset by pressing the reset button (12). The red indicator lamp (7) must be off.
3. Select the desired safety temperature with the temperature control knob (2). This temperature has to be minimum at 5°C above the working temperature.
4. With the door closed, turn on the oven with the switch (1).

5. When the temperature of the oven is stabilized, slowly turn the safety thermostat knob (6) to the left until it clicks, the red indicator lamp (7) lights up and the heating turns off permanently.
6. Block the temperature control knob (2) with the blockage device (9) to avoid accidentally moving the knob.
7. Stop the oven with the switch (1). Open the door and leave it cool with the door open until the temperature goes down 50% of the working temperature (approx).
8. Start the oven once again by switching it on (1).
9. Reset the safety thermostat by pressing the reset button (12) and the red pilot (6) must turn off.
10. Select the working temperature with the thermostat control knob (2) (minimum 5°C below the desired safety temperature).
11. Now the oven is ready to work.

**NOTE: The adjusting process of the safety thermostat is long. For that reason, the safety thermostat is useful when the users normally work at the same temperature.**

OVENS WITH TIMER (option):

1. Put the switch selector (1) into the  position.
2. Choose the time you want the oven to stop automatically with knob (5). At the end of the cycle, the indicator lamp "END" (10) will light up.
3. If you want to work to an indefinable time, put the timer selector (1) in the "ON" position.

OVENS WITH PROGRAMMABLE TIMER (option):

1. Put the switch selector (1) into the  position.
2. Set the starting up and stop time with the levers of the programmable timer.
3. Use the selector (13) to select whether you want it to only do the first cycle marked by the levers (a single cycle) by pressing to the right or to be cyclical (which will repeat the sequence until the user stops the oven) by pressing to the left.
4. If you want to work to an indefinable time, put the timer selector (1) in the "ON" position.

COMMENTS:

\* Do not use the oven at a lower temperature than ambient because the thermostat will not operate correctly.

\* All the temperature regulations must always be done in ascent. Before choosing any value, the knob must be turned left to the end of the scale. Then, a more exact regulation is obtained, as it compensates for the thermostat spindle inevitable movement.

## MAINTENANCE

**Before removing the casing, disconnect the equipment from the mains.**

**The manipulation of internal electronic circuits of the oven by unauthorized personnel can cause irreparable damages. Take it to one of the J.P.SELECTA, s.a.u. authorized technical services.**

CLEANING:

For the cleaning of the different parts, use the following products:

Cleaning of stainless steel: alcohol.

Cleaning of plastic: alcohol with cotton duster.

## SPARE PARTS

**To guarantee the safety of the equipment, all spare parts must be purchased from J.P.SELECTA, s.a.u.**

| Description            | Código / Code |
|------------------------|---------------|
| Model                  | 2000381       |
| Control knob           | 6010          |
| Thermostat setting     | 43030         |
| Termometer             | 42053         |
| Indicator lamp (green) | 22017         |
| Indicator lamp (amber) | 22020         |
| Switch 3 positions     | 20126         |
| Switch 2 positions     | 20061         |
| Heating element        | 39047         |
| Turbine                | 24301         |
| Safety thermostat      | 43019         |
| Glass door             | 47164         |