VACUO-TEMP

4000474

DESECADOR TERMOSTÁTICO AL VACÍO
THERMOSTATIC VACUUM DRYER

Indice

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GENERAL INFORMATION

- 1) Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.
- 2) Do not install or use the equipment without reading this handbook before.
- 3) This handbook must always be attached to the equipment and it must be available for all users.
- 4) If you have any doubts or enquiries, please contact your supplier or J.P. Selecta's technical service.
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.
- 7) Do not use the equipment with liquids which can give off vapours capable of making explosive mixtures.

PACKING LIST

The standard equipment consists of the following components:

Description	Code	Qty.
"Vacuo-temp" vacuum dryer	4000474	1
Glass bell	4000475	1
Silicone gasket	4000476	1
Instruction manual	80047	1

ACCESSORIES

Denomination	Code
VACUUM PUMP 2XZ-1G 10 Pa Rotary blade pump with built-in oil return. Motor with thermic protection device and starting up switch. Maximum working temperature: 60°C.	5900621

EQUIPMENT DESCRIPTION

VACUO-TEMP thermostatic vacuum dryer is designed for drying at adjustable temperatures from +5°C up to 170°C with a temperature fluctuation of $\pm 2^\circ\text{C}$. Electronic temperature regulation and digital reading (3 digits) with 1 digit resolution.

It includes Pt100 probe, shielded heating element, vacuum pump rear connection, vacuum valve and tempered glass bell with silicone gasket.

The upper plate is made of special aluminium alloy with surface grinding and sealing gasket. Outer casing is made in AISI 304 stainless steel.

WORKING OF THE SAFETY THERMOSTAT

If a bad function of regulation produces an abnormal increase of the temperature, the safety thermostat starts working blocking the heating of the machine, and the overtemperature indicator lamp on the frontal panel lights up.

To reset the thermostat, stop the machine, leave it cool and start up again. In case that the failure continues, call the authorized technical service.

TECHNICAL FEATURES

Voltage supply 115-230V 50/60 Hz according to the features plate indications.

Code		4000474
Maximum vacuum		10 ⁻² mm Hg.
Capacity litres		--
Useful volumen litres		3
Internal dimensions cm.	Ø	23,5
	Depth	--
External dimensions cm.	Height	17
	Length	28
	Depth	34
Consumption W		540
Time range		1 to 999 min
Weight Kg		9

INSTALLATION

Put the equipment on a flat, horizontal and level surface leaving a space of 10 cm at the back and on each side.

For maximum efficiency, do not place the equipment in direct sunlight or other sources of heat.

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Do not use the equipment for drying or for treatment of products which can give off gases or make explosive or inflammable mixtures.

Be sure that the voltage supply is the same as the one indicated on the equipment features plate.


Do not use the equipment if it is NOT earthed.


If you have to change the plug, bear in mind the following:

Blue cable:	Neutral
Brown cable:	Phase
Yellow/green cable:	Earth

OPERATION

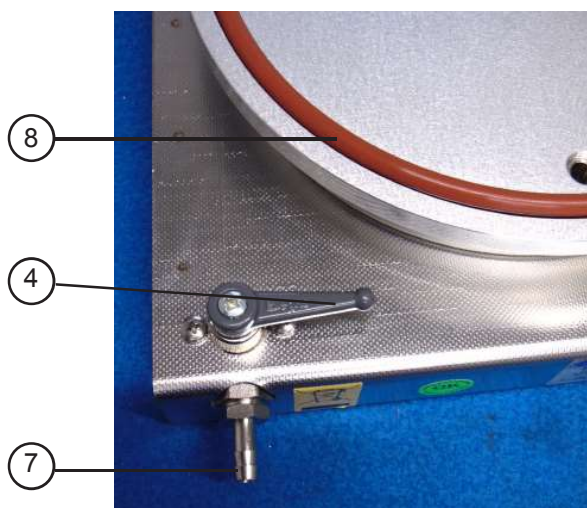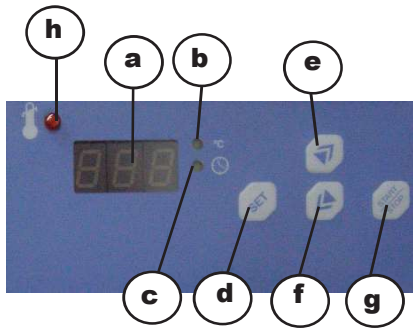1. Main switch.
2. Temperature control thermostat.
3. Vacuum gauge.
4. Ventilation valve.
5. Tempered glass bell.
6. Hotplate.
7. Nipple.
8. Silicone gasket

CONTROL PANEL

- a. Digital display indicating speed and time in minutes.
- b. LED indicator for speed selection.
- c. LED indicator for time selection.
- d. Button for operation speed and time parameters selection.
- e. Button for increasing parameter value.
- f. Button for reducing parameter value.
- g. Start/Stop button.
- h. Overtemperatura pilot lamp indicator.

STARTING UP

1. Press the main switch (1).
2. Select the temperature led (b) to be lit with key (e). Then press key SET (d) and keep it pressed to select temperature with buttons (e) and (f). This will be shown in the display (a).
3. Select the time led (c) to be lit with key (f). Then press key SET (d) and keep it pressed to select time with buttons (e) and (f). This will be shown in the display (a).
4. Check that the gasket is adjusted and place the glass bell.
5. Open the valve (4) and connect the vacuum pump in to the nipple (7).
6. Do the vacuum, and when the vacuum gauge shows the desired vacuum, close the valve (4).
7. Disconnect the vacuum pump from the nipple (7).
8. Press the Start key (g) in order that the equipment begins to operate with the temperature and time parameters already changed. Temperature (b) or time (c) led will begin blinking, which means that the equipment is warming up. The temperature will increase till reaching the programmed temperature. Once reached, the time programmed begins to discount. When the time arrives to zero, the led will remain fixed on and the machine will stop warming up.
9. When it is necessary to remove the glass bell, open the valve (4).

Nota: If the material inside the glass bell can produce gases that can damage the membrane of the vacuum pump, it is recommended to place a water jet pump between the vacuum pump and the Vacuo-temp.

CAUTION! *Serious burns can be produced, with the hotplate surface, even after switching off the equipment.*

MAINTENANCE

Before removing the cover, disconnect the equipment from the mains.

The manipulation of the dryer internal electronic circuits by unauthorized personnel can cause irreparable damage. Take it to one of the J.P.SELECTA's authorized technical services.

CLEANING:

For the cleaning of the different parts use the following products:

Cleaning of stainless steel: Alcohol

Cleaning of plastic: Alcohol with cotton duster.

SPARE PARTS

To guarantee the safety of the equipment, the spare parts must be purchased from J.P.SELECTA, s.a.u.

Code	4000474
Vacuum gauge	16006
Spherical valve	34086
Heating element	39021
Electronic circuit	29329
Temperature limiter circuit	29159
Pt100 probe	43073
Tempered glass bell	4000475
Silicone gasket	4000476

GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damage caused by an incorrect use or causes beyond the control of J.P. SELECTA, s.a.u.

Any manipulation of the equipment by unauthorized personnel by J.P. SELECTA, s.a.u. cancels the guarantee automatically.

Notice to customers:

The product is made up of various components and various materials that must be recycled or, failing that, deposited in the corresponding debris removal sites when the product's life has been completed or when otherwise it is necessary to dispose of it. To do this, the end user who acquires the product must know the current regulations of each municipality and / or locality based on the waste electrical and electronic equipment. The user who acquires this product must be aware of and responsible for the potential effects of the components on the environment and human health as a result of the presence of hazardous substances. Never place the product in a conventional container of citizen scope if a previous dismantling and knowledge of the components that incorporates. If you do not know the procedure to follow, consult with the city council for more information.