VB-1423

3001423

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ENGLISH VERSION

GENERAL INFORMATION

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.*
- 5) ***IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.***
- 6) *If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the equipment with liquids which can give off vapours capable of making explosive mixtures.*

PACKING LIST

The standard equipment consist of the following components:

Descripción / Description	Codigo / Code	Cant. / Qty.
VB-1423	3001423	1
Aro tapa / Lid	2037+27028	3
Manual de instrucciones / <i>Instruction manual</i>	80154	1

ACCESSORIES

(They must be installed in the factory)

Denomination	Code
<i>Universal support</i>	1001453
<i>Digital Chronometer "Diamond"</i>	5900501
<u><i>Thermometers:</i></u>	
<i>At 38,6°C to 41,4°C Div. scale 0,05°C</i>	1001454
<i>At 98,6°C to 101,4°C Div. scale 0,05°C</i>	1001455
<i>At 20,0°C to 50,0°C Div. scale 0,1°C</i>	1001456
<i>At 40,0°C to 70,0°C Div. scale 0,1°C</i>	1001457
<i>At 60,0°C to 90,0°C Div. scale 0,1°C</i>	1001458
<i>At 80,0°C to 110,0°C Div. scale 0,1°C</i>	1001459

EQUIPMENT FEATURES

Temperature sensor; Pt100 thermo-resistor.

Stainless steel AISI 304 lid with three viscometer locations ports, three independent lids and an additional location port for the control thermometer.

The main body of the bath is made of a 20 litre borosilicate glass tank. A white plate is located at the back to help optimise and read the viscometers.

Control temperature with an electronic controller at 25°C up to 100°C. With an adjustable hydraulic safety thermostat supplied.

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TECHNICAL FEATURES

Voltage supply 115V 50/60Hz or 230V 50/60Hz as indicated on the characteristics plate.

Rango de regulación / <i>Regulation range</i>		25°C ... 100°C
Temperatura máxima / <i>Maximum temperature</i>		100°C
Estabilidad / <i>Stability</i>		0.05°C
Resolución / <i>Resolution</i>		0.1°C
Error de consigna / <i>Set error value</i>		0.09°C
Capacidad litros / <i>Capacity litres</i>		20
Medidas Recipiente <i>Tank dimensions (mm)</i>	Alto / <i>Hight</i>	320
	∅	300
Medidas totales / <i>Total dimensions (mm)</i>	Alto / <i>Hight</i>	470
	∅	300
Potencia calefacción / <i>Heating power</i>		1000W
Potencia agitación / <i>Stirring power</i>		60W
Peso / <i>Weight</i>		8 Kg.

INSTALLATION

Place the "VB-1423" in the tank in his placement.

Fill the tank until the liquid level desired, always is necessary to cover 2 or 3 cm over the heating element.

Connect "VB-1423" plug to the mains, press the maing switch, the controller display will lights on and it will be prepared to configure the set working temperature..

Do not use the equipment in explosive atmospheres.

Be sure that the voltage supply is the same as the one indicated on the characteristic plate of the equipment.

Do not use the apparatus if it is NOT earthed.

If you have to change the plug bear in mind the following:

Blue cable: Neutral.


Brown cable: Phase.

Yellow/green cable: Earth.

OPERATION


CONTROL PANEL:

1. Main switch.
2. Temperature regulator.
3. Display.
4. Increase value push button.
5. Decrease value push button.
6. Configure operation push button.
7. Safety thermostat indicator lamp.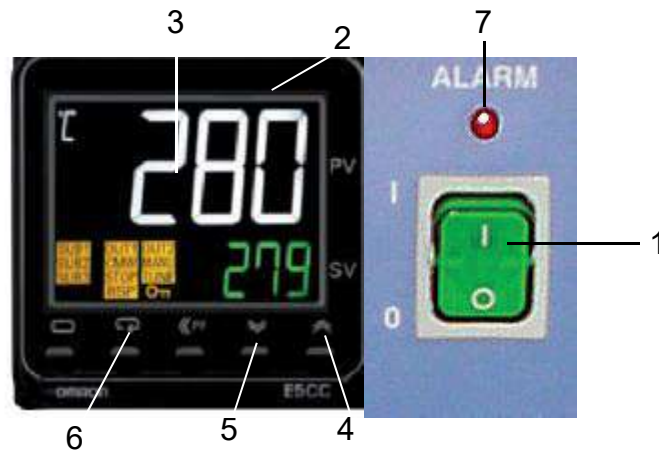

REAR PART:

8. Radiator.
9. Fuse.
10. Safety thermostat.
11. Main supply.

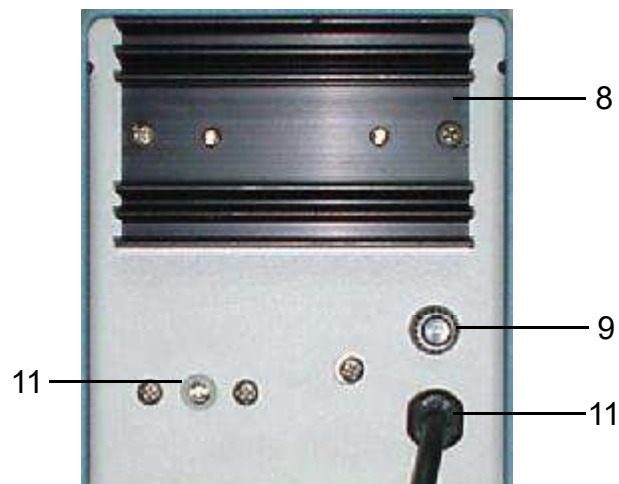SAFETY THERMOSTAT:

To adjust the safety thermostat (10) temperature do as follows:

1. Turn the safety thermostat control knob (10) clockwise to the end of the scale.
2. Select the desired safety temperature with the electronic regulator.
3. When the temperature is stabilized, turn the safety thermostat control knob (10) slowly to the left until it clicks, the red indicator lamp (7) lights up.
4. Stop the incubator with the main switch 1 and leave it to cool.
5. Start the apparatus and with the regulator select the working temperature (minimum at 5°C below the previous selected temperature).

The safety thermostat is useful when the user works habitually at the same temperature.

MAINTENANCE

Before removing the cover disconnect the equipment from the mains.

The manipulation of the internal electronic circuits of the equipment by unauthorized personnel can cause irreparable damage. Take it to one of the J.P.SELECTA authorized technical services.

CLEANING:

For the cleaning of the different parts use the following products:

Cleaning of stainless steel: Alcohol


Cleaning of plastic: Alcohol with cotton duster.

SPARE PARTS

To guarantee the safety of the equipment, all spare parts must be purchased from J.P.SELECTA, s.a.u.

Descripción / Description	Código / Code
Regulador de temperatura / Temperature controller	16123
Relé estático / Static relay	36027
Motor / Motor	24167
Termostato de seguridad / Safety Thermostat	43031
Sonda Pt100 / Pt100 probe	43059
Resistencia 3000W / Heating element 3000W	39018

STARTING UP

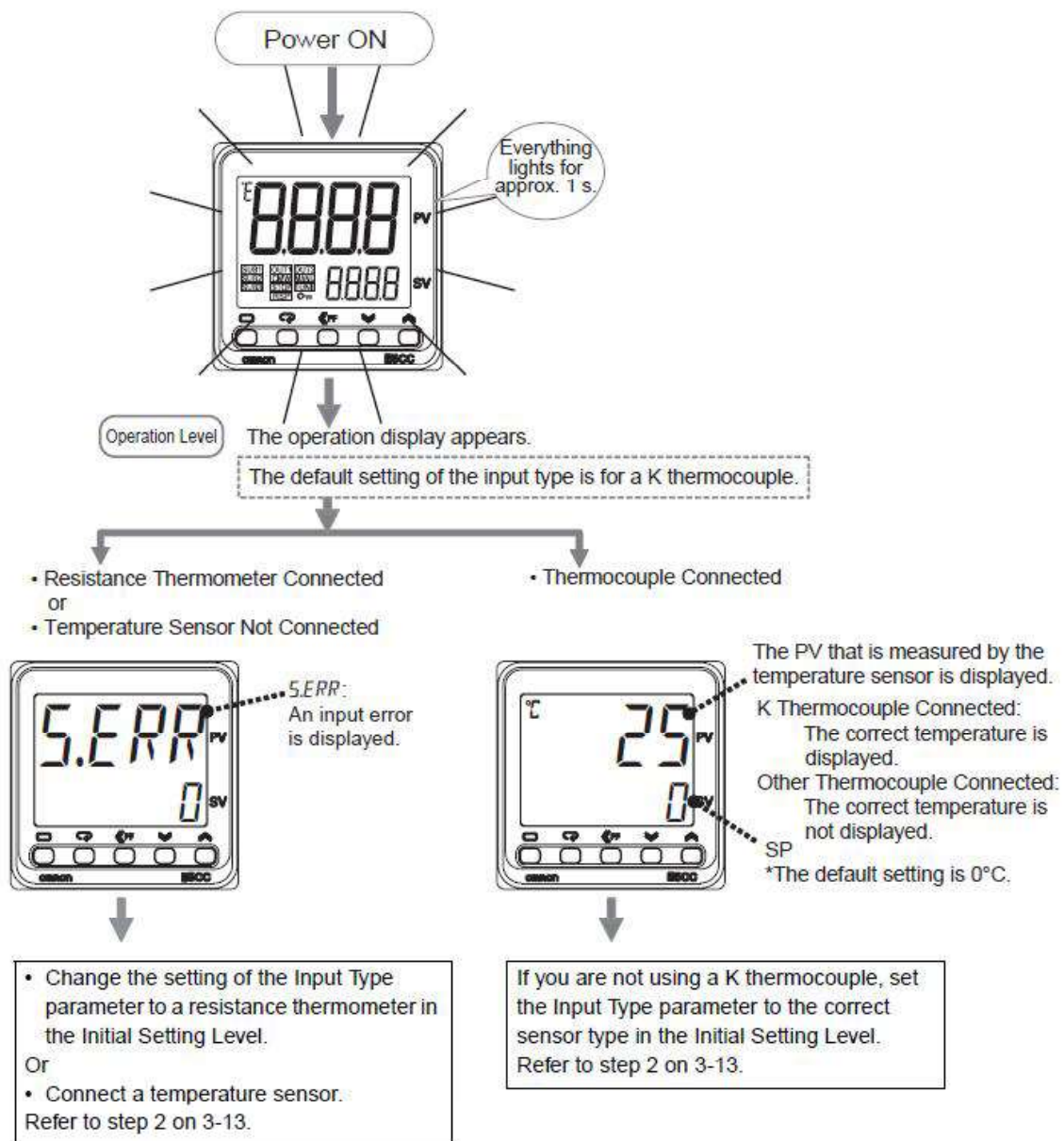GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damage caused by incorrect use or causes beyond the control of J.P. SELECTA, s.a.u.

Any manipulation of the equipment by personnel not authorized by J.P. SELECTA, s.a.u. automatically cancels the guarantee.