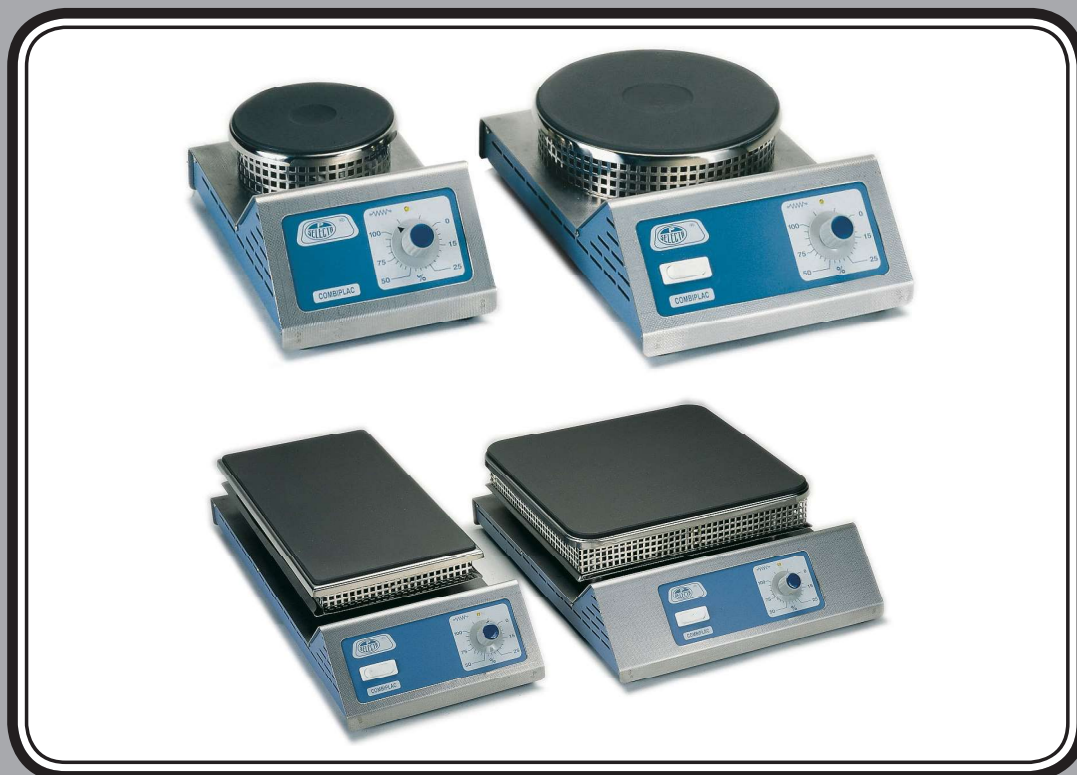
COMBIPLAC	3000156
COMBIPLAC	3000718
COMBIPLAC	3002301
COMBIPLAC	3002302
COMBIPLAC-SAND	6000708
COMBIPLAC-SAND	6000709

PLACAS CALEFACTORAS RECTANGULARES Y CIRCULARES
RECTANGULAR AND CIRCULAR HOTPLATES

ÍNDICE

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ENGLISH VERSION

GENERAL INFORMATION

- 1) Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please inform the distributor immediately.
- 2) Do not install or use the equipment without reading this handbook before.
- 3) This handbook must always be attached to the equipment and it must be available for all users.
- 4) If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.
- 7) Do not use the equipment with liquids which can give off vapours capable of making explosive mixtures.

PACKING LIST

The standard equipment consists of the following components:

Piece	Code	Quantity
Combiplac	3000156	1
Combiplac	3000718	1
Combiplac	3002301	1
Combiplac	3002302	1
Combiplac-Sand	6000708	1
Combiplac-Sand	6000709	1
Instruction manual	80030	1

ACCESSORIES

Denomination

Code

* Only for hotplate of 12 cm Ø

3002301

- Tank in AISI 304 stainless steel, 140 mm Ø and 75 mm high, with reducing rings. Suitable for water, oil and sand baths.

7002510

- Adaptor in aluminium for 500-800ml Kjeldahl flasks and 250 to 750 ml round-bottomed flasks.

1002280

- Adaptor in aluminium for 30-250ml Kjeldahl flasks and 50-250 ml round-bottomed flasks.

1002290

- Stand rod 12mm Ø and 500mm high adjustable to casing

1000725

* Electric contact thermometers

Adjustable temperatures by magnetic rotor with fixture.

Inert gas-filled capillary.

Code	Total length mm	Immersion stem length mm	Regulation range °C	Scale divi- sion °C
3521159	465	150	0 / 100	1 / 1
3522159			0 / 200	2 / 1
3523159			0 / 300	2 / 1

Supplied with connection cable.

* Electronic contact thermometer "SENSOTERM TFT". With digital reading and separated probe, permits regulation away from the area of heat or steam, eliminating the danger of glass breakage.

1001009

* Electronic temperature regulator "ELECTEMP TFT".

Digital temperature regulator P.I.D. operation controlled by a microprocessor, adaptable to every user's requirements.

3000897

Code	Overall dimensions cm			Maximum load W (at 230 V)	Weight Kg
	Height	Length	Width		
3000897	10,5	13,5	21	2000	1250

Pt100 probe

1000893

K-type probe

1000895

* Sand bath attachments

For plate code 3000156: 5.5cmx20cmx40cm (HxLxW)

6000719

For plate code 3000718: 5.5cmx31cmx31cm (HxLxW)

6000720

TECHNICAL FEATURES

Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

Code	Ø Plate cm	Plate dimensions cm.			Overall dimensions cm.			Con- sump. W.	Weight Kg.
		Height	Length	Width	Height	Length	Width		
3000156	---	---	20	40	12	22	55	2300	9
3000718	---	---	30	30	13	33	46	2800	13
3002301	12	---	---	---	10	16	27	550	2
3002302	19	---	---	---	13	21	30	950	3
6000708	---	5	18	38	18	22	41	2300	13
6000709	---	5	30	30	21	32	35	2800	15

EQUIPMENT DESCRIPTION

COMBIPLAC heating plates are equipped with electronic regulation of heaters electrical power, distributed around the surface of the plate, made of non-deformable steel, which allows temperature to be distributed over the plate (at maximum power it can reach a minimum of 400°C).

INSTALLATION

Place the equipment on a flat, horizontal and level surface, trying to leave a space of about 10 cm at the back and on each side.

For a longer life of the equipment provided with hotplate, it is recommended that its surface will be covered as much as possible by the beaker to be heated. This avoids overheating and deformation of the plate by absorbing the heat more.

Keep the plate turned off when not in use, to prolong the life of the equipment.

CAUTION!!! IMPORTANT FOR YOUR SAFETY

Do NOT use the plate for thermic treatment of products which made form explosive gases, not even in special protection area.

Although the plate is strong, it should be treated as fragile and any hard knocks should be avoided.

Be sure that the voltage supply is the same as the one indicated on the characteristics plate of the equipment.

Do not use the equipment, if it is NOT EARTHED.

If you have to change the plug bear in mind the following:

Blue cable: Neutral.

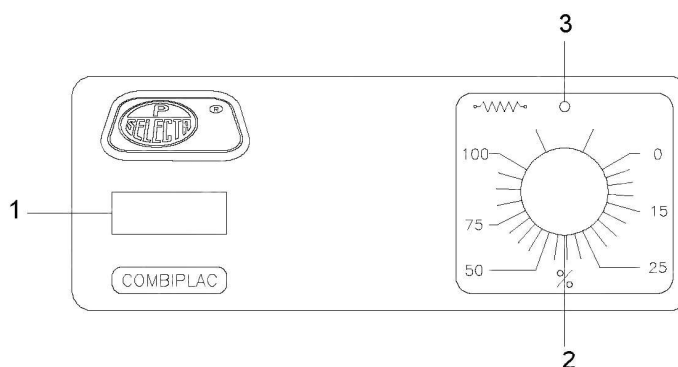
Brown cable: Phase.

Yellow/gree cable: Earth.

OPERATION

CONTROL PANEL

1. Main switch (except **3002301**).
2. Control knob of the heating power.
3. Heating operation indicator lamp.

STARTING UP:

1. Turn on the main switch (1) to start up. Plate code **3002301** has this switch at the rear side.
2. Using the control knob of the heating power, adjust the operation cycle of the hotplate to the desired working power.
3. Amber indicator lamp shows that the heating element is working.
4. If temperature precision is required, it can be connected to a contact thermometer or "ELECTEMP TFT" regulator described in the accessories section, at the base of connection at the back of the equipment.

WARNING!! Take all necessary precautions to avoid burns, the plate may be at a high temperature even when the equipment is off.

**MAINTENANCE**

Before removing the cover, disconnect the equipment from the mains.

The manipulation of the plate's internal electronic circuits by unauthorized personnel can cause irreparable damage. Take it to one of the J.P.SELECTA's authorized technical services.

CLEANING:

For the cleaning of the different parts, use the following products:

Cleaning of stainless steel: alcohol.

Cleaning of plastics and covers: alcohol with cotton or a soft duster.

GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damages caused by an incorrect use or causes beyond the control of J.P. SELECTA, s.a.u.

Any manipulation of the equipment by unauthorized personnel by J.P. SELECTA, s.a.u. cancels the guarantee automatically.

SPARE PARTS

To guarantee the safety of the equipment, all spare parts must be purchased from J.P. SELECTA, s.a.u.

Description	Code					
	3000156	3000718	3002301	3002302	6000708	6000709
Schuko cable	7001		7002		7001	
Electronic P.C.B.	11043	11044	11045	11046	11043	11044
Switch	20100					
Cable retainer	28004					
Plate	32013	32018	32004	32005	32013	32018

Notice to customers:

The product is made up of various components and various materials that must be recycled or, failing that, deposited in the corresponding debris removal sites when the product's life has been completed or when otherwise it is necessary to dispose of it. To do this, the end user who acquires the product must know the current regulations of each municipality and / or locality based on the waste electrical and electronic equipment. The user who acquires this product must be aware of and responsible for the potential effects of the components on the environment and human health as a result of the presence of hazardous substances. Never place the product in a conventional container of citizen scope if a previous dismantling and knowledge of the components that incorporates. If you do not know the procedure to follow, consult with the city council for more information.