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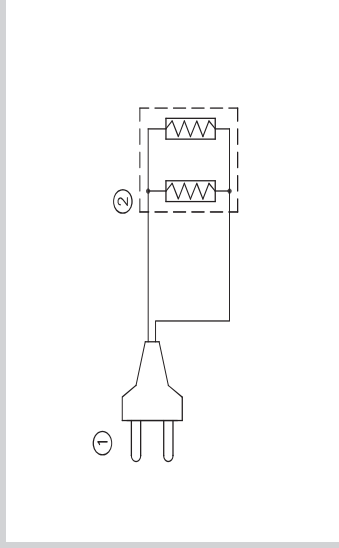
**CAUTION:**

Read these instructions carefully before installing and starting up the equipment.

Do not touch the heating element. Burns can be caused due to its high temperature.

Do not use the equipment if it is not earthed.

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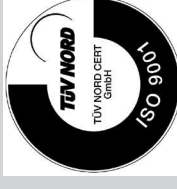
*This product is guaranteed for one year. It does not cover damages caused by an incorrect use or causes beyond the control of J.P. SELECTA, s.a.u.*

*Any manipulation of the equipment by personnel not authorized by J.P. SELECTA, s.a.u. automatically cancels the guarantee.*

J.P. SELECTA, s.a.u. Autovía A2 Km 585,1 ABRERA 08630 (Barcelona) SPAIN  
Tel.: (34) 93 770 08 77 Fax.: (34) 93 770 23 62

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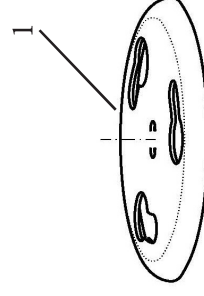
### FEATURES:

- Temperature up to 400°C on the mantle surface.
- For spheric flasks from 250 to 2.000ml.
- Woven mineral glass fiber, fully free of asbestos.
- Heating element evenly distributed throughout the body of the mantle.
- With attachment for AGIMAN magnetic stirrer.

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
### INSTRUCTIONS:

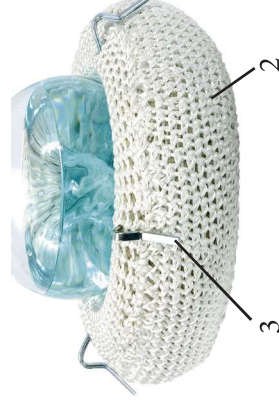- During the first starting up, you will observe that the heating mantle gives off some fumes and a characteristic smell; it is due to the dressing of the heating material. These circumstances are completely normal and will disappear in a short period of time.
- It is recommended to use flasks suited to the mantle size.
- Avoid spilling liquids on the heating material surface, since it means the immediate damage of the mantle.
- Make sure that the mains power corresponds to the one indicated on the description plate.
- For your safety, do not switch on the equipment if it is not earthed.
- Mount the mantle on top of the AGIMAN magnetic stirrer. The stirrer adapter should be located in the center of the mantle.
- Fix the mantle with the fixing plate, turning the plate a few degrees.
- Start the stirrer. Adjust speed and electrical power by using the agitator controls.

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