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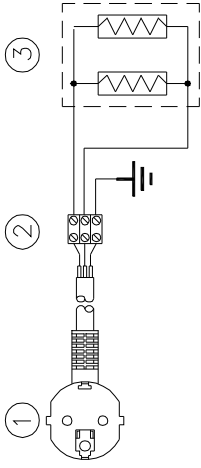
CAUTION:

Read these instructions carefully before installing and starting up the equipment.

Do not touch the heating element. Burns can be caused due to its high temperature.

Do not use the equipment if it is not earthed.

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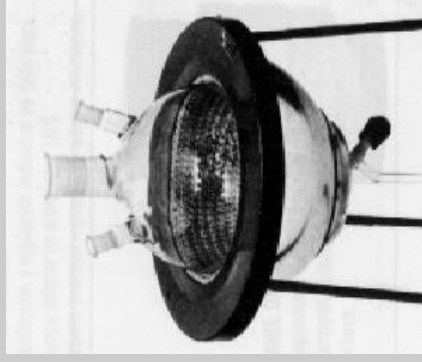
This product is guaranteed for one year. The guarantees does not cover damaged caused by misuse of the equipment or outside causes to J.P. SELECTA, s.a.

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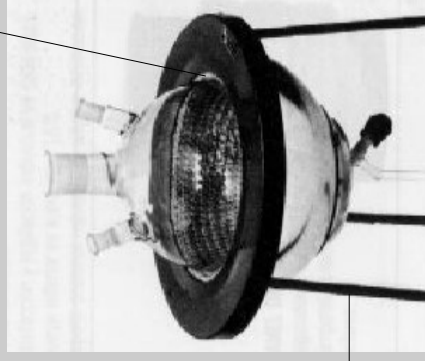
FEATURES:

- *Temperature up to 450°C on the mantle surface.*
- *For spheric flasks from 100ml. to 2.000ml.*
- *Woven mineral glass fiber, fully free of asbestos.*
- *Heating element evenly distributed throughout the body of the mantle.*
- *Three power selector.*
- *Vent for reactors with lower outlet.*

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INSTRUCTIONS:

- *During the first starting up, you will observe that the heating mantle gives off some fumes and a characteristic smell; it is due to the dressing of the heating material. These circumstances are completely normal and will disappear in a period of time.*
- *It is recommended to use flask suited to mantle size.*
- *Avoid spilling liquids on the heating material surface, since it means the immediate damage of the mantle.*
- *Make sure that the mains power corresponds to the one indicated on the description plate.*
- *For your safety, do not switch on the apparatus if it is not earthed.*
- *Turn on the main switch «POWER» (1) on selecting the desired power.*