PLACA CIRCULAR	1000718
PLACA CIRCULAR	1000442
PLACA CIRCULAR	1000443

PLACFIN	7000497
----------------	----------------

UNIPLAC	6002314
UNIPLAC	6002316

**PLACAS CALEFACTORAS
HOTPLATES**

INDICE

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GENERAL INFORMATION

- 1) Handle the parcel with care. Unpack and check that the contents coincide with the packing list. If any part is damaged or missing, please advise the distributor immediately.
- 2) Do not install or use the equipment without reading this handbook before.
- 3) This handbook must always be attached to the equipment and it must be available for all users.
- 4) If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) Any modification, elimination or lacking in maintenance of any part of the equipment transgress the directive 2009/104/CE, the manufacturer is not responsible for the damage that can occur.
- 7) Do not use the equipment with liquids which can give off vapours capable of making explosive mixtures.

PACKING LIST

The standard equipment consist of the following components:

Description	Code	Quantity
Circular hotplate	1000718	1
Circular hotplate	1000442	1
Circular hotplate	1000443	1
Rectangular hotplate "Placfin"	7000497	1
Digital thermometer	42056	1
Circular hotplate unit "Uniplac"	6002314	1
Vertical rods Ø12x60	12103	2
Horizontal rods Ø12x77	12101	1
Bosshead for rods	7008133	2
Head key	9100005	1
Circular hotplate unit "Uniplac"	6002316	1
Vertical rods Ø12x60	12103	2
Horizontal rods Ø12x105	12107	1
Bosshead for rods	7008133	2
Head key	9100005	1
Instruction manual	80061	1

ACCESSORIES

	Circular hotplate unit UNIPLAC		Rectangular hot- plate PLACFIN
	6002314	6002316	7000497
Plate code	6002314	6002316	7000497
Tank suitable for water, oil and sand baths	7002510	7002510	No
Bosshead for clamps	1000069	1000069	No
Clamp	7000403	7000403	No
Stand bar	6023141	6023161	No
Glass fume tubes	1000104	1000106	No
Clamp for steam collector tubes	6000008	6000008	No
Adaptor for slides	No	No	7000498
Adaptor for flasks from 500 up to 800 ml	1002280	1002280	No
Adaptor for flasks from 30 up to 250 ml	1002290	1002290	No

EQUIPMENT DESCRIPTION

The hotplates are designed to work with a big temperature range, from 130°C (Placfin) up to 400°C (Circular hotplates and Uniplac).

Temperature regulation is done by a hydraulic thermostat (Placfin) or impulse regulators (Uniplac) independent for each plate.

TECHNICAL FEATURES

Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

		Circular hotplates			Rectangular hotplate PLACFIN	Circular hotplate units UNIPLAC	
Code		1000718	1000442	1000443	7000497	6002314	6002316
N° of plates		1	1	1	1	4	6
Plate dimensions cm	Length	--	--	--	23	--	--
	Width	--	--	--	22	--	--
Ø Plate cm		12	15	19	--	12	
Overall dimensions cm	Height	8			8	23	
	Length	20	22	26	22	77	105
	Width	20	22	26	30	27	
Consumption W		800	1500	2000	550	2200	3300
Maximum temperature		400°C			130°C	400°C	
Weight Kg		2.1	3	4.5	3.7	15	22

INSTALLATION

Place the hotplate on a flat horizontal, level surface, leaving a space of 10 cm at the back and both sides of the equipment.

For a longer life of the equipment provided with a hotplate, it is recommended to cover its surface as much as possible by the container to be heated. This avoids overheating and deformation of the plate by absorbing more heat.

Because of that, when using a round-bottomed, Soxhlet or Kjeldahl types, it is necessary to use adapters in aluminium described in the "HOTPLATES" section of our catalog.

CAUTION!!! IMPORTANT FOR YOUR SAFETY 

Be sure that the voltage supply is the same as the one indicated on the characteristics plate.

Do not use the equipment if it is not earthed.


If you change the plug bear in mind the following:

- Blue cable: Neutral**
- Brown cable: Phase**
- Yellow/green cable: Earth**


OPERATION

CIRCULAR HOTPLATE CONTROL PANEL:

These heaters include a selector of 6 heating powers settings and a stop. In the table below the power in watts is given according to the selector position.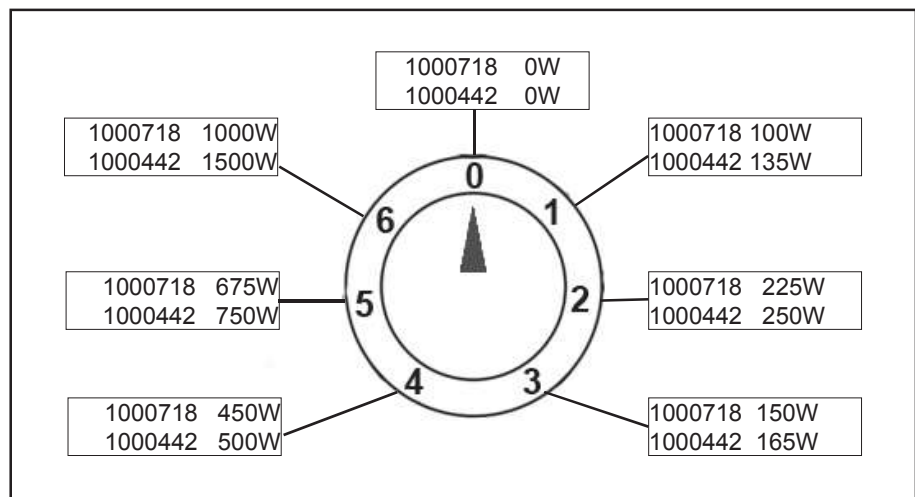

CIRCULAR HOTPLATE UNIPLAC UNITS CONTROL PANEL: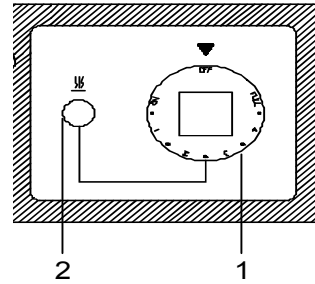

CIRCULAR HOTPLATE UNIPLAC UNITS STARTING UP:


1. Move the control knob of temperature (1) from the "OFF" position (disconnected) to the "FULL" position (maximum power).
2. It will be noted that according to the control position, the ON-OFF cycle switch of the indicator lamp for the working of the heater (2) will be longer or shorter, which will permit adjusting to the desired working temperature.
3. To stop, move the control button to "OFF" position.

RECTANGULAR HOTPLATE PLACFIN:

Rectangular hotplate "PLACFIN" is provided with a temperature regulator thermostat whose sensitive part is in the center of the plate in order to register the temperature of the container to be heated.

Move the thermostat button numbered from 1 to 12 to the right and fix it at one point, bearing in mind that the higher the point is fixed, the higher the temperature is. If after a reasonable time the desired temperature has not been reached, move the button of the thermostat to more or less, as desired.

To disconnect turn the button to 0.

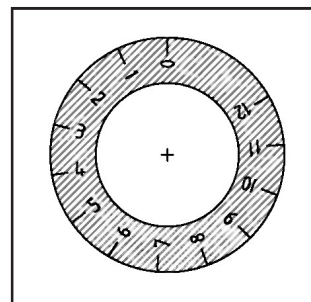WARNING!!! Caution with the hotplate surface, serious burns can be produced, even after switching off the equipment.

MAINTENANCE

Before removing the cover, disconnect the equipment from the mains.

The manipulation of the internal electronic circuits of the equipment by unauthorized personnel can cause irreparable damage. Take it to one of the J.P.SELECTA's authorized technical services.

CLEANING:

For the cleaning of the different parts, use the following products:

Cleaning of stainless steel: Alcohol

Cleaning of plastic: Alcohol with a cotton duster.

SPARE PARTS

To guarantee the safety of the equipment, the spare parts must be purchased from J.P.SELECTA, s.a.u.


Model	6002314	6002316	7000497
Reading thermometer	No	No	42056
Control knob	06001		No
Simerstat regulator	23054		43020
Indicator lamp	22020		No

GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damages caused by an incorrect use or causes beyond the control of J.P. SELECTA, s.a.u.

Any manipulation of the equipment by unauthorized personnel by J.P. SELECTA, s.a.u. cancels the guarantee automatically.

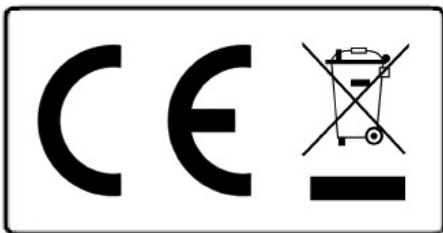Notice to customers:

The product is made up of various components and various materials that must be recycled or, failing that, deposited in the corresponding debris removal sites when the product's life has been completed or when otherwise it is necessary to dispose of it. To do this, the end user who acquires the product must know the current regulations of each municipality and / or locality based on the waste electrical and electronic equipment. The user who acquires this product must be aware of and responsible for the potential effects of the components on the environment and human health as a result of the presence of hazardous substances. Never place the product in a conventional container of citizen scope if a previous dismantling and knowledge of the components that incorporates. If you do not know the procedure to follow, consult with the city council for more information.