**PLX-IN**

**3000958**

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**GENERAL INFORMATION**

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.*
- 5) ***IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.***
- 6) *If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the equipment with liquids which can give off vapours capable of making explosive mixtures.*

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**CONTENTS**

GENERAL INFORMATION  
 CONTENTS  
 PACKING LISTS  
 TECHNICAL FEATURES  
 EQUIPMENT DESCRIPTION  
 INSTALLATION  
 OPERATION  
 MAINTENANCE

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**PACKING LIST**

The standard equipment consists of the following components:

<b>Pieza</b>	<b>Piece</b>	<b>Código / Code</b>	<b>Cant. / Qty.</b>
Cámara	<i>Chamber (with closed bottom base)</i>	3000958	1
Manual de instrucciones:	<i>Instruction manual</i>	80283	1

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**TECHNICAL FEATURES**

**Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.**

Temperature range:	10°C to 60°C
Temperature variation:	±0.5%
Temperature fluctuation:	±0.5%
Consumption:	200W
Weight:	11Kg
Overall dimensions (H x L x W)	61x51x51

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**EQUIPMENT DESCRIPTION**

The chamber is made from methacrylate that allows to view inside the chamber during incubation. Its fan circulation system ensures the homogeneity of temperature. The temperature regulation is made by an electronic circuit board.

The chamber can be fitted onto the platelet shaker APL-54 catalogue number 3000980.

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## INSTALLATION

To install the incubation chamber PLX-IN on the platelet shaker APL-54 follow the steps below:

- Remove the lateral screws from the mobil upper part of the shaker
- Place the chamber PLX-IN on the shaker APL-54
- Lock the set with the same screws

Place the incubator on a level surface keeping a separation of about 25 cm from the wall to allow ventilation and equipment movement.

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Make sure that the voltage of the power supply is the same as the one indicated on the features label attached on the equipment.

Do not use the equipment if it is NOT earthed.

If you have to change the plug, bear in mind the following:

**Blue cable: Neutral.**  
**Brown cable: Phase.**  
**Yellow/green cable: Earth.**

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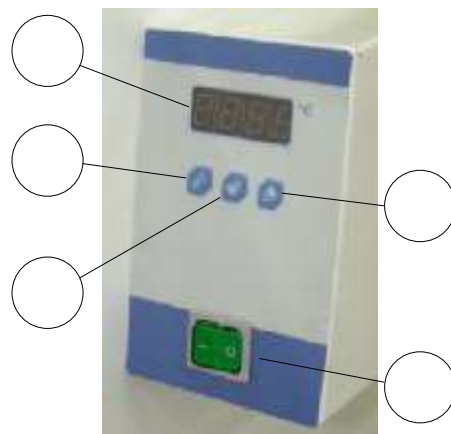
CONTROL PANEL:

1. *Main switch (POWER).*
2. *SET push-button.*
3. *Push-button for temperature increase.*
4. *Push-button for temperature decrease.*
5. *Temperature display.*

## OPERATION

To start up the chamber, press the main switch. The chamber starts at the preset temperature in the temperature regulator.

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**STARTING UP:**

1. Turn on the main switch (1).
  2. Press simultaneously SET key (2) and then the key (3) to increase or (4) to decrease the working temperature. The selected temperature will be shown on the display (5).
- \* During the working, the chamber temperature will be shown on the display (5).
  - \* By pressing the SET key (2), the set temperature will be shown on the display (5).

**WORKING WITH THE PLATELET SHAKER «APL-54»**

When the incubator works on the platelet shaker must be connected by the wire on the rear part shown on the pictures below. In these conditions, on opening the incubator door the shaker stops in order to make easier the working.

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**MAINTENANCE****CLEANING:**

To clean the different parts it is advised to use:

Cleaning of stainless steel: alcohol

Cleaning of plastic: alcohol with cotton or a soft piece of cloth.

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