AGIMATIC-S
AGIMATIC-N
AGIMATIC-E

7000270
7000271
7000273

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GENERAL INFORMATION

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing list. If any part is damaged or missing, please advise the distributor immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.*
- 5) ***IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.***
- 6) *If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the equipment with liquids which can give off vapours capable of making explosive mixtures.*

PACKING LIST

The standard equipment consist of the following components:

DESCRIPTION	CODE	QUANTITY
AGIMATIC S	7000270	1
AGIMATIC N	7000271	1
AGIMATIC E	7000273	1
Power supply wire	7140	1
Safety ring	2049	1
Teflon magnet 8Øx39mm	1000039	1
Instruction manual	80002	1

ACCESSORIES

Denomination	Code
<i>Magnetic stirring-bar 7Øx9mm</i>	1000009
<i>Magnetic stirring-bar 4Øx12mm</i>	1000012
<i>Magnetic stirring-bar 7Øx19mm</i>	1000019
<i>Magnetic stirring-bar 8Øx32mm</i>	1000032
<i>Magnetic stirring-bar 8Øx39mm</i>	1000039
<i>Magnetic stirring-bar 10Øx45mm</i>	1000045
<i>Magnetic stirring-bar 12Øx75mm</i>	1001061
<i>Magnetic stirring-bar 16Øx102mm</i>	1001062
<i>Magnetic stirring-bar 16Øx127mm</i>	1001063
<i>Magnetic stirring-bar 19Øx150mm</i>	1001064
<i>Triangular magnetic stirring-bar TRIMAG 10 (sidex35 length)</i>	1000536
<i>Triangular magnetic stirring-bar TRIMAG 14 (sidex39 length)</i>	1000537
<i>Triangular magnetic stirring-bar TRIMAG 14 (sidex45 length)</i>	1000538
<i>Triangular magnetic stirring-bar TRIMAG 14 (sidex55 length)</i>	1000539
<i>Magnetic stirring-bar ROTOR-PAT 10Øx32 mm</i>	1000001
<i>OVAL magnetic stirring-bar 16 Øx30 mm length</i>	1000788
<i>OVAL magnetic stirring-bar 20 Øx17 mm length</i>	1000789
<i>Crosshead magnetic stirring bars 25Øx15 mm height</i>	1000790

<i>Floating magnet IMANFLOTTER. Structure 50 mm. Magnet 39 mm</i>	<i>1000897</i>
<i>Floating magnet IMANFLOTTER. Structure 60 mm. Magnet 45 mm</i>	<i>1000898</i>
<i>Concentric magnet CENTRIMAN (length 32 mm)</i>	<i>1000900</i>
<i>Concentric magnet CENTRIMAN (length 50 mm)</i>	<i>1000901</i>
<i>Concentric magnet CENTRIMAN (length 50 mm)</i>	<i>1000901</i>
<i>Concentric magnet CENTRIMAN (length 55 mm)</i>	<i>1000902</i>
<i>Stirring bar retriever</i>	<i>1000020</i>
<i>Box with assorted magnetic stirring bars</i>	<i>1000487</i>
<i>Pt100 probe with handle which allows read and control liquid temperature (model E)</i>	<i>1001610</i>
<i>Digital contact thermometer SENSOTERM -40-270°C (models N & E)</i>	<i>1001009</i>
<i>Electric contact thermometer 0-100°C (models N & E)</i>	<i>3521159</i>
<i>Electric contact thermometer 0-200°C (models N & E)</i>	<i>3522159</i>
<i>Electric contact thermometer 0-300°C (models N & E)</i>	<i>3523159</i>
<i>Stay rod 12Øx500mm length</i>	<i>1000725</i>
<i>Clamp to hold thermometer</i>	<i>7000025</i>
<i>Stainless steel round tray 2 litres</i>	<i>1001059</i>
<i>Reducing discs set with cover</i>	<i>3000917</i>
<i>Bosshead</i>	<i>1000069</i>

EQUIPMENT DESCRIPTION

The magnetic electronic stirrers described in this manual comprise three models with stainless steel plate.

*The stirrers **AGIMATIC S, N and E** is a group of stirrers with analogic technology for the control of speed and temperature. This technology allows setting parameters in a precise and repetible way.*

S model allows digital reading and control of speed.

N model allows setting a heating level as a rate of total heating power. To regulate the temperature of the stirred liquid in a better way, use an electronic controller Sensoterm II as an accessory.

E model allows setting the heater plate temperature. Including 1001610 probe as an accessory, it allows controlling liquid temperature at same time that it is stirring.

In some application it is important bearing in mind that heater plate temperature and the liquid one are very different.

POWER	AGIMATIC S	25W
	AGIMATIC N	550W
	AGIMATIC E	
Maximum temperature (N & E models)		350°C
Temperatura resolution (E model)		± 2°C
Heating control (N & E models)		N:0...100% E:50...350°C
External temperature control connection (N & E models)		Yes
Speed		60...1600rpm
Stirring power		10.000ml
Dimensions		14,5x19x26cm
Weight (S model)		3,0Kg
Weight (N & E models)		3,2Kg
Temperatura limiter (N & E models)		370...390°C

INSTALLATION

Place the stirrer on a flat, horizontal, level surface near the mains.

Connect the connection cable at the rear part of the stirrer base and then plug it into the mains.

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Be sure that the voltage supply is the same as the one indicated on the characteristics plate.


Do not use the equipment if it is not earthed.

If you change the plug, bear in mind the following:

Blue cable:	Neutral
Brown cable:	Phase
Yellow/green cable:	Earth


OPERATION

CONTROL PANEL: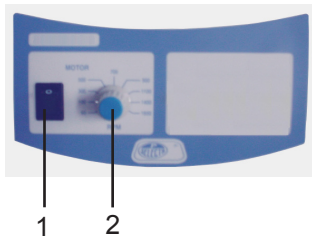


S MODEL

1. Agitation switch
2. Speed set button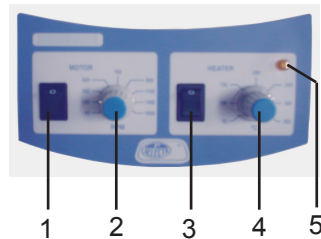




N MODEL

1. Agitation switch
2. Speed set button
3. Heating switch
4. Set % heating power
5. Heating indicator lamp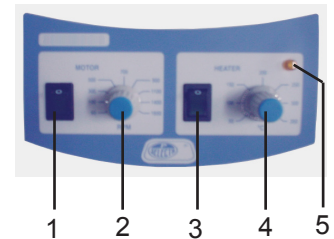

E MODEL

1. Agitation switch
2. Speed set button
3. Heating switch
4. Set temperature
5. Heating indicator lamp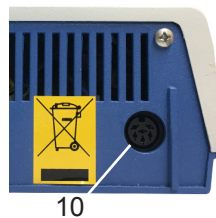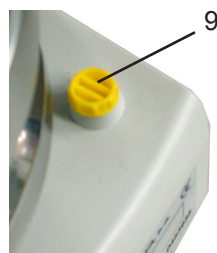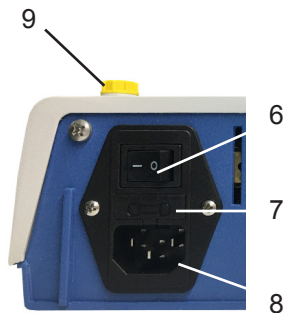

REAR PART

6. Main switch
7. Fuse carrier
8. Mains plug
9. Thermometer support rod housing
10. External control connection or liquid probe (N & E models)

STARTING UP :**AGIMATIC-S:**

Once the indications in the **Installation** section have been made, switch on the main switch on the rear part, place the recipient to be stirred on the top plate, let the teflon-coated magnetic stirring bar drop into the recipient, turn on the "MOTOR" switch and adjust speed with "RPM" knob.

AGIMATIC-N:

For its starting up operation and to select the speed, do as the above model, AGIMATIC-S. There is no heating until motor is working.

Then switch on the HEATER switch, and with the regulator for heating power, select the required "%". The yellow lamp will show the working of the heater.

AGIMATIC-E:

To start it working and to adjust speed, follow the above indications. With the regulator "HEATER" the suitable working temperature may be selected with the "°C" knob.

**LIQUID TEMPERATURE CONTROL**

(N & E models)

1. Connection to work without external connection of a temperature control device.

2. Connection for contact thermometer.

- Put the thermometer with the bulb submerged into the liquid. Connect the thermometer in the stirrer rear part plug.
- Set the temperature turning the superior part of the thermometer. The stirrer must be configured to work at maximum of power rate or temperature.
- The connection is as shown on figure 1.

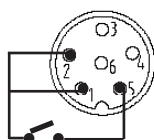
Figure 1

3. Connection for regulator "SENSOTERM II" (1001009)

- Put the probe submerged into the liquid. Connect the probe in the stirrer rear part plug.
- Set the temperature on the «SENSOTERM II».
- Set the power rate or temperature on the stirrer at maximum.
- The connection is as shown on figure 2.

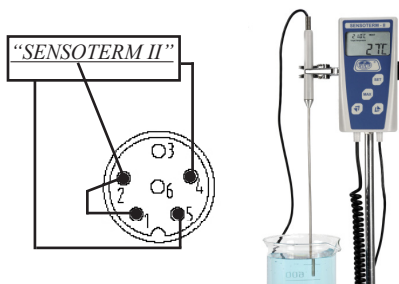
Figure 2

MAINTENANCE

Before removing the stirrer casing, disconnect the equipment from the mains.

The guarantee will not have any effect if the equipment is manipulated by somebody who has not been expressly authorized by J.P.SELECTA, s.a.u.

CLEANING:

For the cleaning of the different parts, use the following products:

Cleaning of stainless steel: Alcohol

Cleaning of plastic: Alcohol with a cotton duster.

WARRANTY

This product is guaranteed for one year. The guarantee does not cover damages caused by an incorrect use or causes beyond the control of J.P. SELECTA,s.a.u.

Any manipulation of the equipment by not authorized personnel by J.P. SELECTA,s.a.u. automatically cancels the guarantee.

SPARE PARTS

Heating plate	0032026
Control circuit:	
for model S	0038515
for model N	0038514
for model E	0038513
Pattern magnet	0015449
Motor	0024283
Pt100 probe	0043028 (model E only)