
**AGIMATIC ED
AGIMATIC ED-C**

**7000280
7000281**

**AGITADORES MAGNÉTICOS DE CONTROL DIGITAL
CON PLATO CERÁMICO O DE ACERO INOXIDABLE
DIGITAL CONTROL MAGNETIC STIRRERS
WITH VITROCERAMIC OR STAINLESS STEELTOP PLATES**

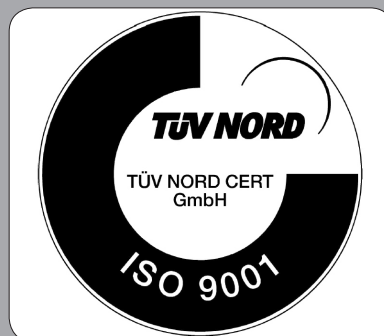INDICE

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GENERAL INFORMATION

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.*
- 2) *Do not install or use the equipment without reading this handbook first.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.*
- 5) ***IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.***
- 6) *If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the equipment with liquids which can give off vapours capable of making explosive mixtures.*

PACKING LIST

The standard equipment consists of the following components:

DESCRIPTION	CODE
AGIMATIC ED	7000280
AGIMATIC ED-C	7000281
Power Supply Wire	0007140
Safety Ring	0002049
Teflon-coated magnet 8Øx39mm	1000039
Instructions manual	M.80148

ACCESSORIES

Denomination	Code
<i>Magnetic stirring-bar 7Øx9mm</i>	1000009
<i>Magnetic stirring-bar 4Øx12mm</i>	1000012
<i>Magnetic stirring-bar 7Øx19mm</i>	1000019
<i>Magnetic stirring-bar 8Øx32mm</i>	1000032
<i>Magnetic stirring-bar 8Øx39mm</i>	1000039
<i>Magnetic stirring-bar 10Øx45mm</i>	1000045
<i>Magnetic stirring-bar 12Øx75mm</i>	1001061
<i>Magnetic stirring-bar 16Øx102mm</i>	1001062
<i>Magnetic stirring-bar 16Øx127mm</i>	1001063
<i>Magnetic stirring-bar 19Øx150mm</i>	1001064
 <i>Triangular magnet 10 x35</i>	 1000536
<i>Triangular magnet 14 x39</i>	1000537
<i>Triangular magnet 14 x45</i>	1000538
<i>Triangular magnet 14 X55</i>	1000539
 <i>Magnet ROTOR-PAT 10 Øx32 mm</i>	 1000001
 <i>OVAL magnet 16 Øx30 mm length</i>	 1000788
<i>OVAL magnet 20 Øx17 mm length</i>	1000789
 <i>Crosshead magnet 25Øx15 mm height</i>	 1000790
<i>Crosshead magnet 40Øx17 mm height</i>	1000791
 <i>IMANFLOTTER 50 mm. Magnet 39 mm</i>	 1000897
<i>IMANFLOTTER 60 mm. Magnet 45 mm</i>	1000898
 <i>Magnet CENTRIMAN (length 32 mm)</i>	 1000900
<i>Magnet CENTRIMAN (length 50 mm)</i>	1000901
<i>Magnet CENTRIMAN (length 55 mm)</i>	1000902
 <i>Stirring bar retriever</i>	 1000020
 <i>Box with 17 assortment magnets</i>	 1000487
 <i>Pt100 probe with handle which allows read and control liquid temperature</i>	 1001610
 <i>Stay rod 12Øx500mm length</i>	 1000725
<i>Clamp to hold thermometer</i>	7000025
<i>Stainless steel round tray 2 litres*</i>	1001059
<i>Reducing discs set with cover *</i>	3000917
<i>Bosshead</i>	1000069

* Except ED-C model

EQUIPMENT DESCRIPTION

The magnetic digital stirrers described in this manual comprise one model with stainless steel plate (ED) and another one with ceramic plate (ED-C).

Stirrers **AGIMATIC ED** and **ED-C** are a group of stirrers with digital technology for the control of the speed and temperature. This technology allows setting parameters in a precise and repeatable way.

ED and ED-C models allow setting the heater plate temperature. Including 1001610 probe as an accessory, it allows controlling liquid temperature at same time that it is stirring.

ED-C model includes a ceramic heater plate that allows a better liquid temperature control.

In some application it is important being in mind that heater plate temperature and the liquid one are very different.

TECHNICAL FEATURES

Voltage supply 115-230 V 50/60 Hz as indicated on the characteristics plate.

Power	550W
Maximum Temperature	350°C
Temperature Resolution	1°C
Heating Control	50 ... 350°C
Electronic Control Plug	Yes
Speed	60 ... 1600 rpm
Stirring Power	10.000ml
Dimensions	14.5 x 19 x 26
Weight	3.2Kg
Temperature Limiter	370 ... 390°C

* 50 ... 200°C WITH THE EXTERNAL PROBE PT100

INSTALLATION

Place AGIMATIC on a flat, horizontal, level surface near mains.

Be sure that the voltage supplied is the same as the one indicated on the characteristics plate.

Do not use the equipment if it is not earthed.

If you change the plug, bear in mind the following:
Yellow/green cable: Earth

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
IMPORTANT:

Before any operation inside stirrer disconnect it from mains

OPERATION

CONTROL PANEL: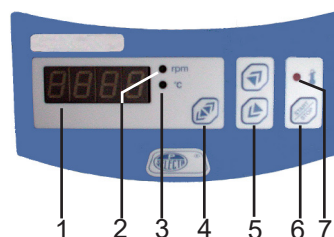

1. Speed display (rpm) / temperature display °C)
2. Speed indicator led (rpm)
3. Temperature indicator led (°C)
4. Temperature / speed selection
5. Up/Down parameters.
6. Start / Stop
7. Over temperature alarm

STARTING UP :

- Switch on rear part main switch.
- Place the flask centred on the stirrer's plate.
- Leave the magnet into the flask.
- To adjust a temperature value, press key (4) until the temperature indicator (3) is on.
- Press keys (5) to select the desired temperature value.
- To adjust a speed value, press key (4) until the speed indicator (2) is on.
- Press keys (5) to select the desired speed value.
- To start up stirring, press "START/STOP" (6) key.
- Press «START/STOP» (6) again to stop stirring.

(*) NOTE1: During operation it is possible to change the set values in the same way as in the waiting mode.

(*) NOTE 2: Processor memory retains the last operating values used before pressing «START / STOP» (6).

Over temperature protective device:

ED and ED-C models have an over temperature device which cuts the power to the heating element when the temperature reaches over 390°C (approx.). When this occurs, pilot lamp "7" lights up. If this failure usually repeats, please contact your technical service.

Special caution with ED-C model:

ED-C model includes a ceramic heater plate which allows a better temperature control but has, as disadvantage, its greater fragility.

To avoid electric risks do not use the stirrer if ceramic plate surface shows any crack.

LIQUID'S TEMPERATURE CONTROL


1. Connection to work without external connection of a temperature control device.

This configuration is the default one.

2. Connection for Pt100 probe (1001610)

ATTENTION: for safety reasons, the temperature regulation by the external probe is limited from 50 to 200°C.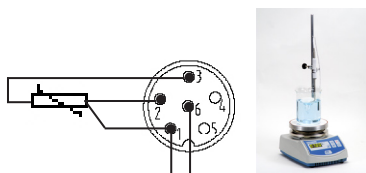

- Put the probe submerged in the liquid. Connect the probe in the stirrer rear part plug.
- Set the temperature on the stirrer control panel.

SPARE PARTS

To guarantee the safety of the equipment, all spare parts must be purchased from J.P.SELECTA, s.a.u.

DESCRIPTION	CODE
Main Board	0038511
Display Board	0038512
Motor	0024283
Heating Element (ED-C)	0039248
Ceramic Disc (ED-C)	0032048
Heater Plate (ED)	0039221
Probe	0043028

MAINTENANCE

The guarantee will not have any effect if the electronic circuit boards are manipulated by personnel not expressly authorized by J.P. SELECTA, s.a.u.

CLEANING:

For the cleaning of the different parts use the following products:

Cleaning of stainless steel: Alcohol

Cleaning of control panel and ceramic plate: Alcohol with cotton duster.

WARRANTY

This product is guaranteed for one year. The guarantee does not cover damages caused by an incorrect use or causes beyond the control of J.P. SELECTA, s.a.u.

Any manipulation of the equipment by personnel not authorized by J.P. SELECTA, s.a.u. automatically cancels the guarantee.