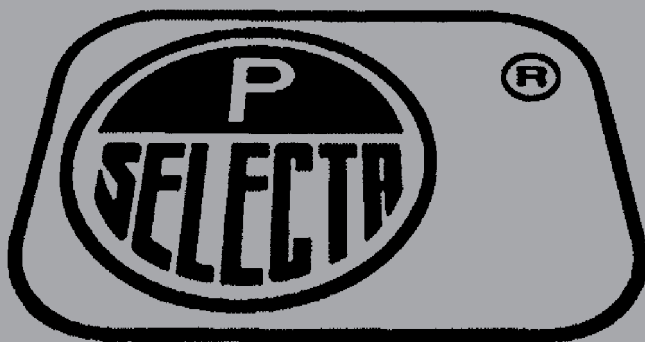
Agimatic-Rev-TFT 7000837
Agimatic-Rev-TFT-C 7000838

Agitadores magnéticos de giro reversible
Reversible magnetic stirrers

Indice

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Index

1. General Information	8
2. Packing List	8
3. Technical Specifications	8
4. Information about safety in the use of the equipment	9
4.1 Risks the operator is subjected to	9
4.2 Staff qualification	9
4.3 Use with hazardous substances	9
4.4 Maintenance	9
4.5 Warranty	9
Menu Structure	10
5. Equipment Description	11
6. Installation	11
6.1 Location	11
6.2 Connection to the mains	11
7. Operation	11
7.1 Information screen and Settings	11
7.2 Quick start for stirring	12

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1. General Information

- Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributors immediately.
- Do not install or use the equipment without reading this handbook first. This instruction manual must be attached to the equipment and must be available for all users.
- If you have any doubts or enquiries, contact with J.P. SELECTA's technical service.
- Do not use the equipment in explosive, flammable or corrosive atmospheres.

2. Packing List

Model:	7000837	7000838
Content:	Stirrer, power cord, centering ring, instructions manual, pen for touch screen.	

3. Technical Specifications

Model:	7000837	7000838
Rotational speed: (rpm)	de 100 a 1600	
Maximum volume: (L)	1	
External: high/width/depth (mm)	145/190/260	
Power supply:	230 VAC 50Hz	
Electrical power:	550	
Weight: (Kg)	3.2	
External fuse:	Replaceable (Type: 10x5mm)	
Temperature range:	From... 30°C ... to ... 350°C	
Display resolution:	1°C	
Timing operation:	Minutes, seconds	
Temperature profiles:	Up to 10 profiles of 6 segments, from 1m to 99 hours. (5940 minutes)	

«External probe» accessory

The external probe allows a better regulation of the sample temperature.

pT 100 temperature probe	1000893
Measures	Ø4mm x 136mm

4. Information about safety in the use of the equipment

Hazard risk

Electric risk

Risk of high temperature contact

4.1 Risks the operator is subjected to

- Risk of burns
- Electrical risk

4.2 Staff qualification

This equipment may only be used by properly trained personnel who have read and understood these instructions or have been properly trained in this equipment operation.

4.3 Use with hazardous substances

Do not treat flammable or explosive materials or those which chemically react with violence.

ATTENTION!

Burning risk

Take proper precautions
(use of thermal gloves, clamps,...)
to extract samples from inside the stirrer.

It can be at a high temperature.

4.4 Maintenance

It does not need maintenance, only cleaning the heating plate.

Please disconnect the equipment before handling inside.

The manipulation of the equipment internal electronic circuits by unauthorized personnel can cause irreparable damage.

Take it to one of J.P.SELECTA's authorized technical services.

It is recommended to clean the accessible parts of the oven with a cloth and a cleaner, or a similar solution. Do not use aggressive solvents.

4.5 Warranty

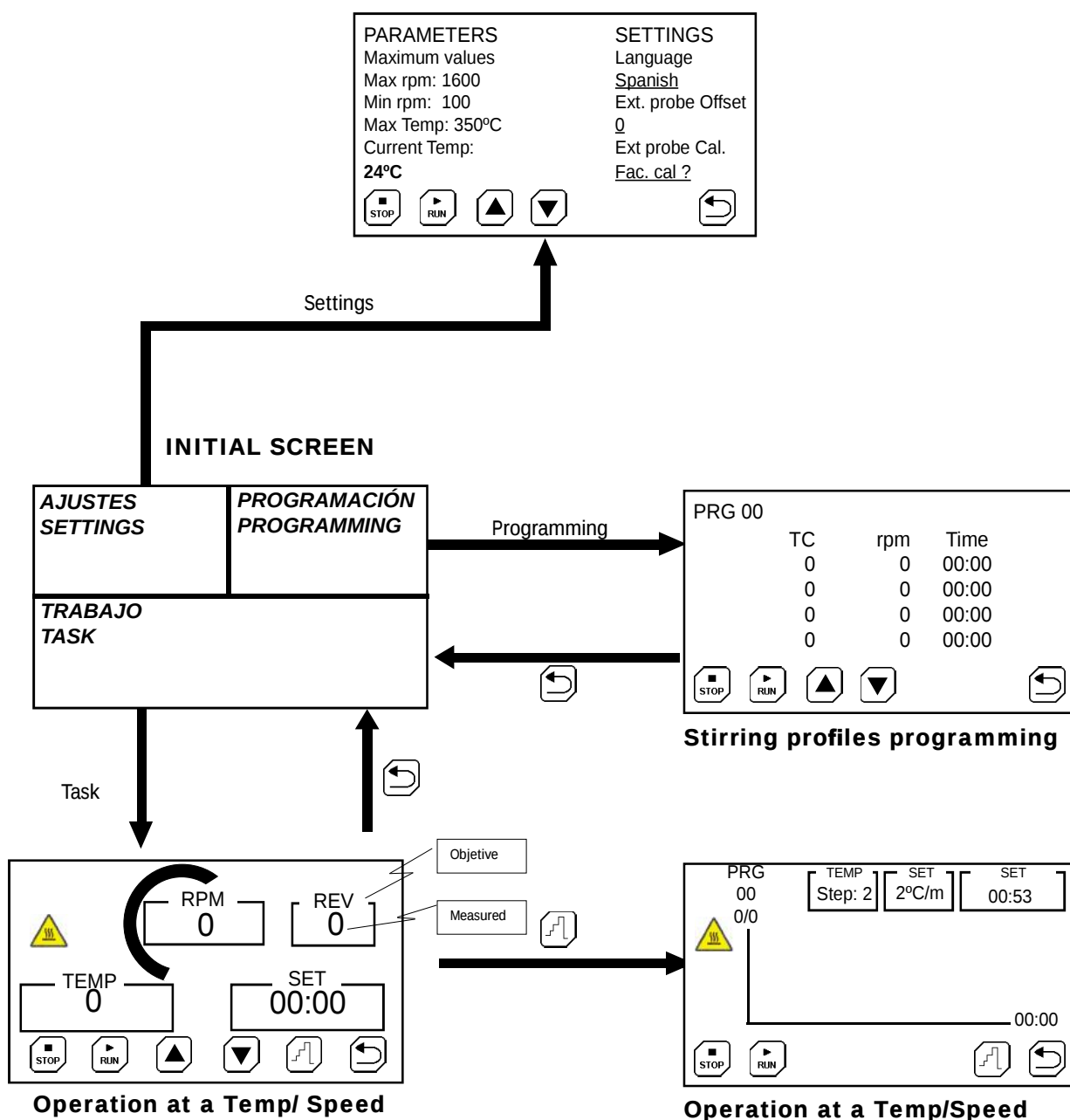
This product is guaranteed for one year. It does not cover damages caused by an incorrect use or causes beyond the control of J. P. SELECTA, s.a.u.

Any manipulation of the equipment by personnel not authorized by J.P. SELECTA, s.a.u. automatically cancels the guarantee.

Menu Structure

All functions and parameters of the stirrer are selected by clicking on the touch screen icons.

It is recommended to use the pointer delivered with the equipment.

5. Equipment Description

The stirrer provides the preparation of solutions through a combination of temperature and stirring speed.

A temperature profile is a succession of several steps of heating and stirring speed.

6. Installation

Unpack the equipment and visually check that the contents are in good conditions. Keep the package for some days.

6.1 Location

Place the equipment on a flat level stable surface, suitable for the equipment weight.

To work in a comfortable way, please leave a free space of 20cm around the equipment.

6.2 Connection to the mains

Plug the equipment to a near socket suitable for the equipment power. Use the connection cable supplied or a similar one. For your safety, the equipment must be earthed. Before plugging the machine, check that the voltage supply and the frequency corresponds to the one indicated in the equipment characteristics plate.

7. Operation

All the equipment functions are set using the touch screen. Gently press with your finger in the center of each icon.

The computer has two operation modes:

- Simple stirring
- Several steps stirring program

When starting the equipment, starting screen in Fig 7.1 will be shown.


- SETTINGS:**
 - Language selection and calibration of external probe.
- PROGRAMMING**
 - Select the temperature profile of the external probe.
- TASK**
 - Starts the stirrer by selectiong the speed and the heating or starts an stirring profile.

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7.1 Information screen and Settings

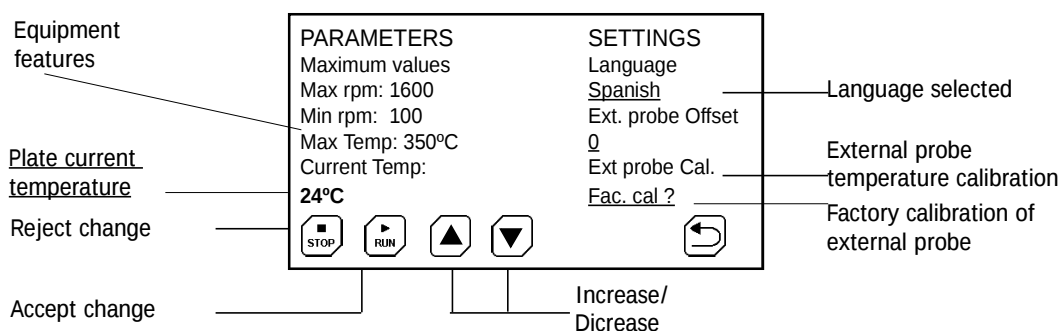Icons meaning:

7.1.1 Calibrating the temperature and the alarms

With a certain periodicity (annually, for instance), it is recommended to check the temperature calibration, for example, by using a thermometer or a temperature pattern.

If as a result of calibration a reading correction is needed, a correction must be entered.

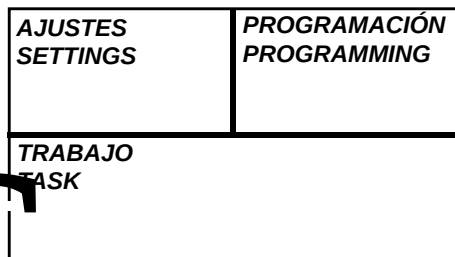
In case of doubt about the calibration done, factory calibration can be recovered.

7.1.2 Set Language


Click on the Language icon. Successively click on the box to change from Spanish, French or English.

Set by pressing on the Increase and Decrease icons.


7.2 Quick start for stirring


1 Press on «TASK».

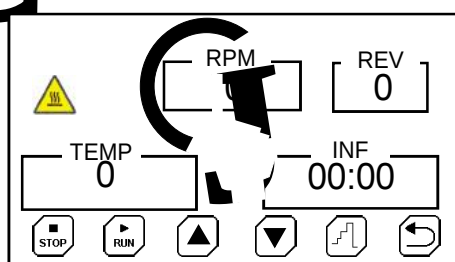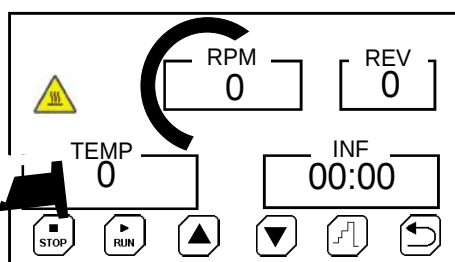2 Press on the speed selector.
Press to select speed.

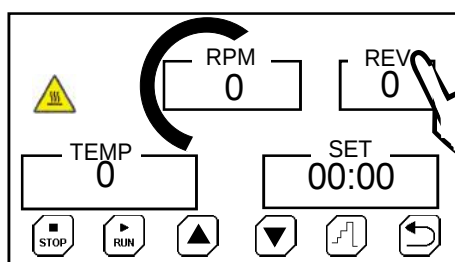
3 Press on the temperature selector.
Press and select the temperature.

4 Press on «REV» to change the rotational direction every 10, 20 or 30sec.

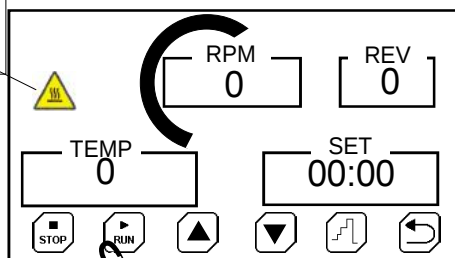Warning sign:
The plate is at a higher temperature than 50°C.

5 Press on «SET» to enter an operating time.
If left 00:00 means a continuous operation.

6 Press on «RUN» to start stirring.

7 Alternatively press on «profile» to run an stirring profile previously programmed.