**AGITADOR ROTATIVO «MOVIL ROD II»**

**TUBE SHAKER «MOVIL ROD II»**

**7001726**

**Aviso a los clientes:**

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**Notice to customers:**

The product is made up of various components and various materials that must be recycled or, failing that, deposited in the corresponding debris removal sites when the product's life has been completed or when otherwise it is necessary to dispose of it. To do this, the end user who acquires the product must know the current regulations of each municipality and / or locality based on the waste electrical and electronic equipment. The user who acquires this product must be aware of and responsible for the potential effects of the components on the environment and human health as a result of the presence of hazardous substances. Never place the product in a conventional container of citizen scope if a previous dismantling and knowledge of the components that incorporates. If you do not know the procedure to follow, consult with the city council for more information.

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**INDEX**

1 Packing list.....4  
 2 Description and installation .....4  
 3 Technical specification.....4  
 4 Operation .....4  
 5 Warranty .....4

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## 2 Description and installation

### Unit description:

Shaker with 5 rollers with double motion action: rotation and swing, this provide a suitable homogenize effect. Suitable for use in liquid suspension, biology and haematology.

### Application

Suitable for mixing blood in a sample tube, for mixing dry powder and lyophilized agent, and also for mixing liquid in a tube. This equipment has a higher tilt angle. Its mixing method is simple, safe and reliable.

### Location:

Place the equipment on a flat and levelled surface, suitable for the shaker's weight, and leave about 20cm free around it.

### Connection to power supply:

- Check that power socket have suitable voltage, according unit plate.
- Power socket have earth connector. Unit must be grounded.
- Unit could be dangerous is not properly grounded.

## 3 Technical specification

Rolling and vibrating speed:	36 rpm
Tilting angle:	±11°
Power consumption:	6W
Fuse:	F1A1250V
Voltage:	220V AC ±10% (50Hz)
Weight:	3Kg
Dimensions:	260 x 155 x 140 mm
Working shift:	Continuous

## 4 Operation

- Place the tubes to stirr.
- Press the main switch
- Adjust the speed.

About once a year, lightly lubricate both ends of the rollers, to prevent the bearings from rubbing against each other.

## 5 Warranty

The warranty covers the defective components or faulty assembling of the unit for 1 year since the purchasing date. Keep purchasing bill as the only valid document for the purchasing date.

The warranty does not covers the damages caused for a unit malfunction.

Return goods are not accepted without previous J.P. SELECTA, s.a.u. approval.