FLOCUMATIC 4 PLAZAS/SAMPLES LED 3000820
FLOCUMATIC 6 PLAZAS/SAMPLES LED 3000822

FLOCULADORES DE LABORATORIO
LABORATORY FLOCCULATOR

INDICE

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ENGLISH VERSION

GENERAL INFORMATION

1) Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.

2) Do not install or use the equipment without reading this handbook before.

3) This handbook must always be attached to the equipment and it must be available for all users.

4) If you have any doubts or enquiries, please contact your supplier or J.P. Selecta's technical service.

5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.**

6) If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.

7) Do not use the equipment with liquids which can give off vapours capable of making explosive mixtures.

PACKING LIST

The standard equipment consists of the following components:

DESCRIPTION	CODE
FLOCUMATIC 4 SAMPLES LED	3000820
Illuminating element Flocumatic 4 samples	3000821
FLOCUMATIC 6 SAMPLES LED	3000822
Illuminating element Flocumatic 6 samples	3000823
Schuko cable	7140
Instruction manual	80385

ACCESSORIES

Denomination	Code
Additional illuminating element, for simultaneous illumination.	
- For model code 3000820 4 samples	3000821
- For model code 3000822 6 samples	3000823

TECHNICAL FEATURES

Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

Code	No samples	Speed regulation r.p.m.	Overall dimensions (cm)			Total consumption (W)	Illumination power (W)	Weight (Kg)
			Height	Length	Width			
3000820	4	15-200	52	70	23	160	40	24
3000822	6			99		170	50	36

EQUIPMENT DESCRIPTION

FLOCUMATIC LED are stirring equipments which fit beakers up to 1000 ml (high form) or 2000 ml (low form).

Stirring speed is adjustable from 15 to 200 r.p.m. Noiseless operation and a synchronized movement of the agitation elements.

Fixed blade impellers, easily changeable and adjustable height.

INSTALLATION

Place **FLOCUMATIC LED** on a flat, horizontal steady surface.

Level **FLOCUMATIC LED** by adjusting the screwed legs supplied for this purpose.

To install the stirring bars, raise the anti-slip base and insert them into the holes.

To obtain the required illumination, proceed as follows:

- a. To adjust the height, loosen the lateral screws, slide the element to the desired height, and tighten the screws again.
- b. To change the lighting angle, loosen the lateral screws completely until you can see that the pivot of the washer is free, move the light to the desired angle, and tighten the screws again, checking that the pivots fit back into their housing.

Place **FLOCUMATIC LED** mains suitable with the equipment consumption.

Do not install the equipment in areas where inflammable liquids are stored or in zones with special protection.

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Do not use the equipment if it is not earthed.

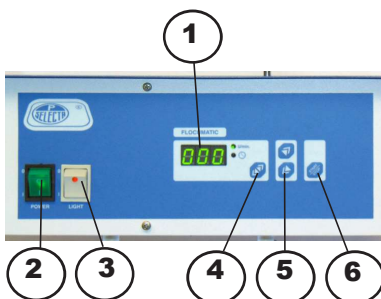
Be sure that the voltage supply is the same than the one indicated on the characteristics plate.

If you change the plug, bear in mind the following:

Blue cable:	Neutral
Brown cable:	Phase
Yellow/green cable:	Earth

OPERATION

CONTROL PANEL:

1. Digital display indicating speed and time in minutes.
2. **POWER** button
3. **LIGHTing** button
4. Selection key. Time / rpm
5. Key for reducing and increasing parameter value.
6. **Start/Stop** key

STARTING UP:

1. Loosen the bar supports and slide the bars to the upper part.
2. Place the beakers.
3. Adjust the bars to the desired height.
4. Connect the equipment to the mains.
5. Turn on the main switch **POWER** (2).
6. To set speed parameters, press key (4) and select the speed with key (5). The speed selected will be shown on the digital display (1).
7. To set time parameters, press key (4) pressed and select the time with key (5). The time selected will be shown on the digital display (1).
8. To start the machine, press the **Start/Stop** key (6). Stirring will finish when pressing the **Start/Stop** key (6) again or when finishing the set time.
9. Connect the power cable of the lighting element to the back of the flocculator.
10. If necessary, please switch on the **LIGHT** (3).

MAINTENANCE

Before removing the cover, disconnect the equipment from the mains.

The manipulation of the internal electronic circuits by unauthorized staff can cause irreparable damage. Take it to one of the J.P.SELECTA authorized technical services.

CLEANING:

For the cleaning of the different parts use the following products:

Cleaning of stainless steel: Alcohol

Cleaning of plastic: Alcohol with cotton duster.

GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damage caused by incorrect use or causes beyond the control of J.P. SELECTA, s.a.u.

Any manipulation of the equipment by unauthorized people by J.P. SELECTA, s.a.u. cancels the guarantee automatically.

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Notice to customers:

The product is made up of various components and various materials that must be recycled or, failing that, deposited in the corresponding debris removal sites when the product's life has been completed or when otherwise it is necessary to dispose of it. To do this, the end user who acquires the product must know the current regulations of each municipality and / or locality based on the waste electrical and electronic equipment. The user who acquires this product must be aware of and responsible for the potential effects of the components on the environment and human health as a result of the presence of hazardous substances. Never place the product in a conventional container of citizen scope if a previous dismantling and knowledge of the components that incorporates. If you do not know the procedure to follow, consult with the city council for more information.