**DESTILADOR KJELDAHL AUTOMÁTICO**  
**AUTOMATIC KJELDAHL DISTILLER**

**PRO - NITRO A 4002430**

# Indice

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# Table of contents

<b>1 Packing list &amp; accessories .....</b>	<b>36</b>
1.1 Standard package content.....	36
1.2 Accessories .....	36
<b>2 Safety .....</b>	<b>37</b>
2.1 Safety symbols.....	37
2.2 Risks which the operator is put under .....	38
2.3 Personnel training.....	38
2.4 Proper uses of the equipment.....	38
2.5 Modifications.....	38
<b>3 Overview .....</b>	<b>39</b>
<b>4 Installation .....</b>	<b>41</b>
4.1 Suitable location .....	41
4.2 Connection to power supply .....	41
4.3 Connection to tap water and waste.....	41
4.4 Water for steam generator .....	41
4.5 Connection to reagents tanks .....	41
<b>5 Operation.....</b>	<b>42</b>
5.1 Menu for function selection .....	42
5.2 Alarms .....	45
5.3 Insert/ Remove the sample tube.....	45
<b>6 Sample analysis.....</b>	<b>46</b>
6.1 Check levels of tanks .....	46
6.2 Unit warm-up and cleaning.....	46
6.3 Distillation. (Sample analysis) .....	48
6.4 Analysis interruption .....	48
6.5 Protein contents (%) calculation. ....	48
6.6 Printing results. ....	48
6.7 Storing results on a computer.....	48
<b>7 Reagents .....</b>	<b>50</b>
7.1 Prepared reagents .....	50
7.2 Ammonia fixative solution preparation .....	50
7.3 Fixative solution checking and adjusting. ....	51
<b>8 Checking the distiller (OQ).....</b>	<b>52</b>
8.1 Functional checking .....	52
8.2 Checking Nitrogen Recovery. ....	54
8.3 Blank test.....	54
8.4 Checking Nitrogen recovery with Acetanilide. ....	55
<b>9 Maintenance .....</b>	<b>56</b>
9.1 Cleaning the unit: .....	56
9.2 Hose inspection (monthly).....	56
9.3 Preparing the unit for shipping or storing. ....	57
9.4 Level sensor cleaning.....	58
<b>10 Technical specification .....</b>	<b>59</b>
<b>11 Technical service .....</b>	<b>60</b>
11.1 Spare parts .....	60
11.2 Hydraulic tubing diagram.....	61
11.3 Wiring diagram. (1 of 2).....	62
11.4 Diagnostic menu.....	64
<b>12 Kjeldahl analysis step by step.....</b>	<b>65</b>
12.1 Sample preparation.....	65
12.2 Digestion.....	65
12.3 Dilution .....	65
12.4 Distillation .....	65
12.5 Protein % contend calculation.....	65
<b>13 Guidelines for IQ y OQ.....</b>	<b>66</b>
<b>14 Warranty .....</b>	<b>67</b>
<b>15 Troubleshooting table .....</b>	<b>68</b>
15.1 General operation .....	68
15.2 Alarm anomalies.....	68

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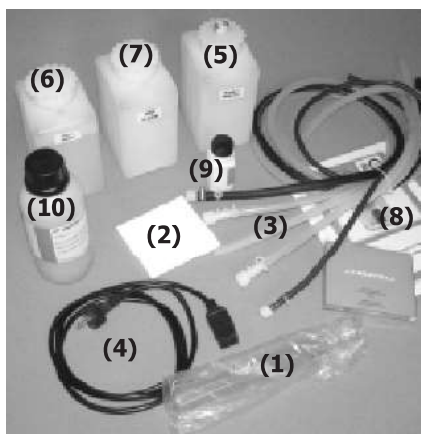
# 1 Packing list & accessories

## 1.1 Standard package content

- Distilling unit (1)
- MACRO size sample tube Ø 42 (2)
- Tray (2)
- Tube 1.5m incoming water. High pressure, coupling 3/8" (3)
- Tube 1.5m water outlet. Silicone 8x14, coupling 3/8" (3)
- Tube 1.5m remaining steam. Silicone 8x14, coupling 3/8" (3)
- Tube 1.5m generator emptying. Silicone 8x14 (3)
- Tube 1.5m waste outlet. ISOVERSINIC 7x12 (3)
- Power supply wire (4)
- 2 litre tank for NaOH (5)
- 2 litre tank for Fixative solution (6)
- 2 litre tank for HCl (7)
- Instructions Manual (8)
- Flask 100g of Ammonia sulphate (9)
- 250ml bottle of mixed indicator 4.8 (10)

## 1.2 Accessories

- Ink printer 4120113  
56mm paper wide roll  
Serial connection wire included

### Tubes, racks and tongs:

- MACRO sample tube Ø 42 4042300
- MICRO sample tube Ø 26 4001045

### Racks:

- Rack MACRO 6 tubes. 4005071
- Rack MACRO 12 tubes. 4005081
- Rack MACRO 20 tubes. 4005091
- Rack MICRO 12 tubes. 4001053
- Rack MICRO 24 tubes. 4001054
- Rack MICRO 40 tubes. 4001055

### Safety goods:

- Tongs for tubes 7000532
- Protection gloves 1001255
- Anti-acid gloves 5000044
- Safety glasses 5028192

### Complete digestion sets

- Digestion unit MACRO 6 tubes. 4000629
- Digestion unit MACRO 12 tubes. 4000630
- Digestion unit MACRO 20 tubes. 4000631
- Digestion unit MICRO 12 tubes. 4001047
- Digestion unit MICRO 24 tubes. 4001048
- Digestion unit MICRO 40 tubes. 4001049

### Smokes removal aids:

- Scrubber. 4001611
- Water circulation vacuum pump. 4001612

## 2 Safety

PRO-NITRO "A" unit has the suitable measures for its use in laboratory. However, the nitrogen analysis by the Kjeldahl method can be hazardous and the user must be informed of some of those risks.

In order to be respected, across this manual, the risk of danger situations are remarked.

### 2.1 Safety symbols

It identifies risky situations and the safety actions to be applied.

The symbols make reference to the paragraph marked with a grey line.

#### **Warning**

Warning of a dangerous operation.

Indicated instruction must be followed.

#### **Hazard of contact with corrosive substances**

Hazard of contact with sodium hydroxide and strong acid, substances that attack skin, clothes and other materials easily.

Follow the instructions for the describe operation.

#### **Hazard of electrical shock**

Electric hazard for entering in areas marked with this signal, or for following the manual instruction signalled by this icon.

Follow the instructions for the described operation.

#### **Hazard of burns to be in contact with hot surfaces**

Temperature in this area can exceed 60°C. Use heat protected gloves for the described action.

Follow the instructions for the described operation.

#### **Wear gloves mandatory**

The use of gloves is obligatory to run the operation.

In the section «Accessories» an advised model is indicated.

#### **Mandatory use of safety glasses**

The use of safety glasses or complete face protection is obligatory to conduct the described operation.

In the section «Accessories» an advised model is indicated.

### Important information

- Information important to obtain good results or for an optimal operation of the equipment.
- Information important to extend the life of the equipment or to avoid the degradation of some of its components.

### 2.2 Risks which the operator is subjected to

The process of determination of Kjeldahl Nitrogen with the described equipment exposes to the operator to the following situations of risk:

- Manipulation of concentrated sulphuric acids and Hydrochloric.
- Sodium manipulation Hydroxide 40%.
- Manipulation of glass parts.
- Possibility of touching parts to a temperature above to 60°C
- Possible exhibition to irritating acid steam.
- Electrical risk.

### 2.3 Personnel training

This equipment must only be operated by personnel trained and suitable qualified to work in a chemical laboratory and its common hazards.

This equipment must only operated by personnel who have read and understood this manual or have been qualified on its manipulation and operation.

### 2.4 Proper uses of the equipment

This equipment is for laboratory use, for sample distillation by steam pulling. Mainly in Kjeldahl nitrogen analysis.

The improper use of this unit (not following manual instructions) can endanger the protection prepared by the equipment.

Do not use to distillate explosive or flammable substances.

This equipment is not designed and not manufactured to work in an explosive ambient area.

### 2.5 Modifications

The operation modification or the manipulation of the safety systems unauthorized by the manufacturer could expose the operator to hazards not indicated in this manual.

### 3 Overview

The PRO-NITRO "A" is a steam distillation unit that includes a colorimetric titration unit, designed, mainly for organic nitrogen determination according to direct Kjeldahl method.

This highest automation degree assures more reliable and robust results.

It is very easy to use but its operation must be perfectly known in order to get good results and avoid losing valuable samples.

Before starting to analyse valuable samples, do some training sessions running the operations related on the chapter 8, until the operator will be familiar with the unit.

Also the situations when the collector could be contaminated with not neutralized rests must be understood to avoid wrong results.

Please, start identifying the name of the analyser components:

#### The steam generator

Generates steam from the water reservoir placed at the top of the unit.

The steam generator includes two safety devices: **Safety thermostat**: which cuts the electrical power to the heater when overtemperature is detected on the steam generator (probably for poor level of water), and the: **Pressure switch**: which cuts the electrical power to the heater when there is an overpressure on the steam generator.

Also includes a capacitive probe to keep a constant level of water in the steam generator.

#### The distillation circuit

It is made, mainly, by two glassware parts: The **splash protector**: its special design avoids that NaOH splash reaches the cooler glassware, and the **cooler**: it condensates the steams that are leaking from the sample. This glassware is cooled by a water circuit that comes from an external tap. An internal valve opens and closes this circuit automatically.

#### NaOH and ammonia fixative solution addition system

Composed by a properly specified pumps to dispense NaOH directly to sample tube when distillation is running and fixative solution (Boric acid with mix of indicators) directly to collector before the distillation.

#### Emptying systems

It consists in some pumps and valves to empty the sample tube and the collector once the analysis is finished.

#### Colorimetric automatic titration unit

It performs the titration at the same time that the distilled is obtained.

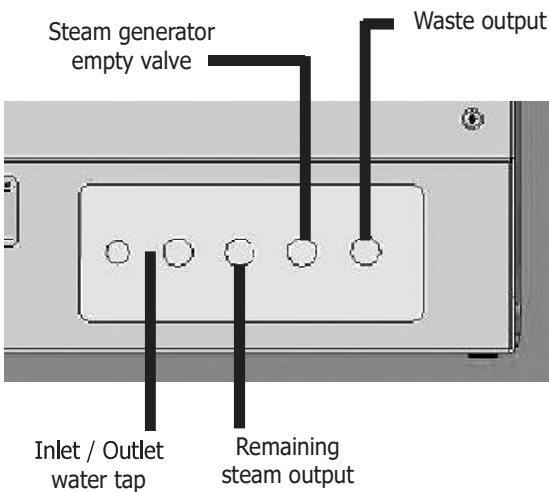
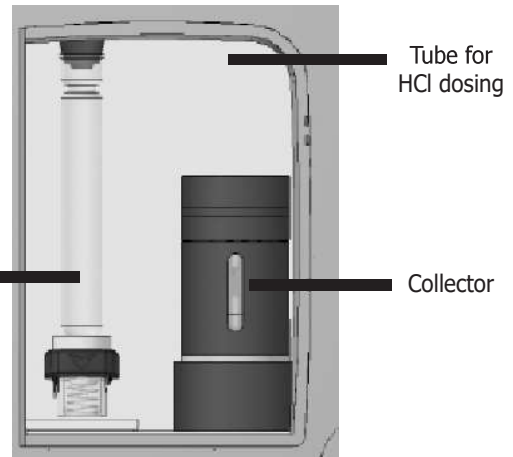
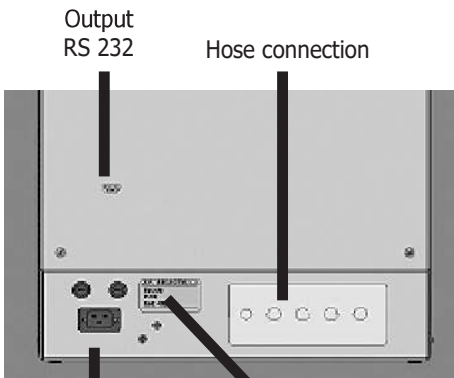
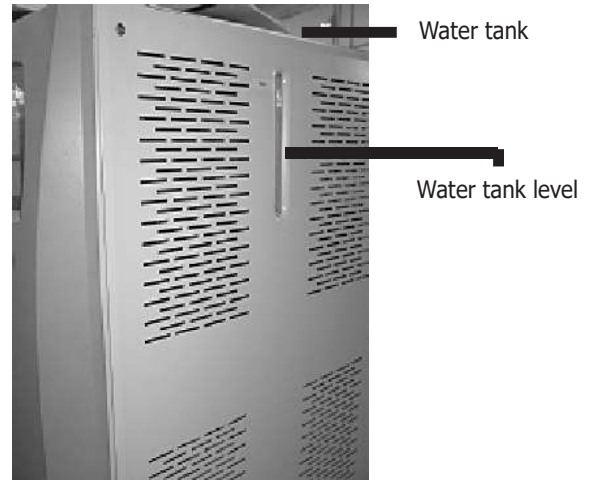
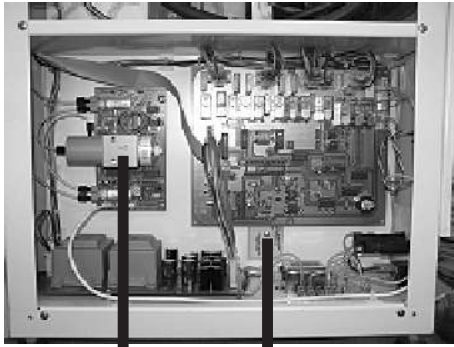
It has two basic components: The **digital dispenser**: which dispenses the titration reagent (usually HCl 0.25N) in volumes of 10 and 50 microlitre, and the **colorimeter** which consists in a LED light source and a photodiode.

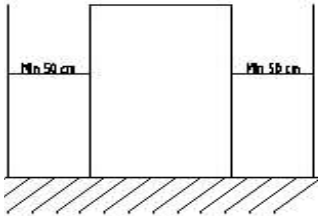
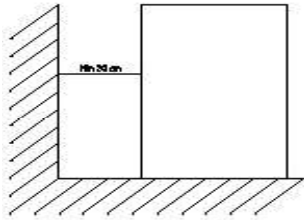
#### Software for recording and reporting preparation

PRO-NITRO «A» includes a RS-232 serial output to send data to a printer or to a computer. This last option allows storing results on the computer and preparing a report from the batch analysed, according GLP practices.

Computer must have WINDOWS 95 or higher version installed on the PC.

The digital dosifier builds components from **Lee Company U.S.A.** ([www.TheLeeCo.com](http://www.TheLeeCo.com)) who pioneered the miniature fluid control components.

## 4 Installation

Once the unit is unpacked, visually check, the packing list (See 1.1) and that there are no damaged parts during transportation. Keep the pack for some days after receiving it.

### 4.1 Suitable location

Place the unit over a flat level and stable surface, suitable for the unit weight.

No longer that 1.5m from a water tap, waste output and power supply plug.

To work comfortably, let a free space of 50cm on each side and not least than 20cm on the rear part.

Prepare to refill the water reservoir from the top of the unit.

### 4.2 Connection to power supply

Locate a power supply socket (rated to 16A) near the unit.

Use the power cord supplied or another of similar specifications.

For safety reason the socket must have ground terminal, since the unit must work grounded.

Before to plug on the socket, check that its voltage match the voltage marked on the unit plate.

### 4.3 Connection to tap water and waste.

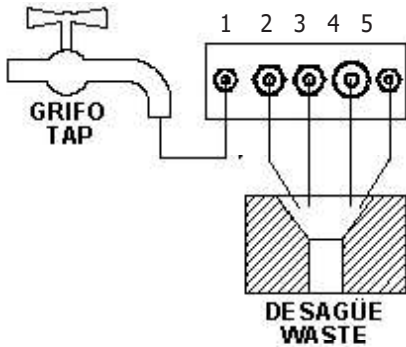
Connect the water entry hose to a water tap (1). Connect the water output hose from (2) to waste.

The hoses «remaining steam» (3) and «steam generator empty» (4) and "waste" (5) must go to waste.

Do not forget the white rubber white nuts on the hose's connections.

Press plastic connections by hand. Do not use any tool.

The tap water must have, at least, 1 bar pressure.

**Note:** It is critical in order that the unit works properly, the hose «Remaining steam» (3) MUST NOT BE IMMERSED under water, or blended or blocked by any reason.

### 4.4 Water for steam generator

The unit includes a 6 litres water reservoir, at the top, to store the water for the steam generator.

As the system used to control the level of water on the steam generator works by a capacity probe, water of at least 20 microSiemens must be used. Water, also, must have a low content of calcium to increase the life of the steam generator.

Water with this conductivity could be obtained from a demineralised unit. If the conductivity is lower, add about 50mg of salt or sodium carbonate. (per 10L water).

Reagents tanks

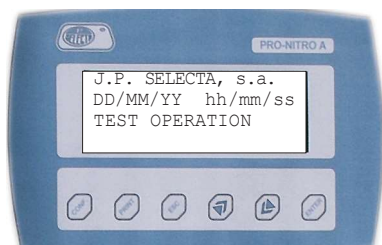
### 4.5 Connection to reagents tanks

At the left hand side, there is a door to access reagent tanks, NaOH, fixative solution and HCl. Connect each one to the hose indicated with his name.

Software version

«MAIN MENU»

## 5 Operation

The control panel allows function selection and shows the distiller alarms status.

Once the unit is plugged in, put it on, by pressing the main switch. The unit goes to stand by status (ON).

**Note:** If main switch is pushed consecutively with an interval of less than 3 seconds, the unit could start in abnormal mode. So wait few seconds to restart the unit.

First the display shows the software version installed on the unit.

Wait until the display shows the «MAIN MENU» before to push any key.

On starting the unit for the first time, the water pump starts to work, automatically, to fill the steam generator.

This operation usually takes near 40 seconds and it will stop, also automatically when the steam generator is filled. At 20s the «STEAM» alarm sounds, but it will also stop when the steam generator is filled.

If this operation is not finished after approx. 40 seconds, please check that the water tank has some water (visual check on the right hand side level), also be sure that the water has the suitable conductivity. (See 4.4)

### 5.1 Menu for function selection

All functions will be selected from the control panel. As a general rule, select a function and point the function by the arrows and then press «ENTER».

#### 5.1.1 Configuration menu:


From the «MAIN MENU» pressing the «CONF» key this menu is get. Use this menu to configure some parameters:

- **Date and time:** Set date and time with keys «▲» «▼» and «ENTER».
- **NaOH:** Pump factor: Adjust the precision on the NaOH dispensing. Range 1 to 5. Factory set at 3.
- **Boric:** Pump factor: Adjust the precision on the Boric dispensing. Range 1 to 5. Factory set at 3.
- **Recovery:** Sets the titration sensibility: Range 1 to 10. Factory set at 5.


Value 1: Low sensibility. Low recovery, fast titration.  
Value 10: High sensibility. High recovery, Slow titration.

**Note:** To modify this value could produce some errors on the titration.

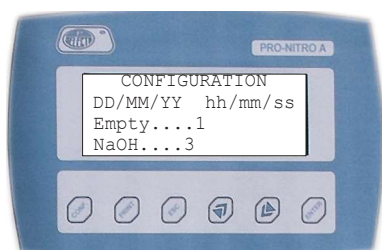- **Power:** Steam power factor: Range: 10 to 5. Factory set at 8.
- **Emp.col:** This number sets the working time of the emptying collector pump. This allows adjust increasing or decreasing it the exact time for emptying the collector completely. Range: 1 to 5. Factory set at 1.
- **LS:** Tritation parameter. Range: 0 to 10. Factory set at 5.
- **Light:** Tritation parameter. Range: 0 to 10. Factory set at 0.
- **Prn:** Configures the mode for data sending to printer or computer.  
**Not Continuous:** (Default) Send data after press «PRINT».

«CONFIGURATION MENU 1»


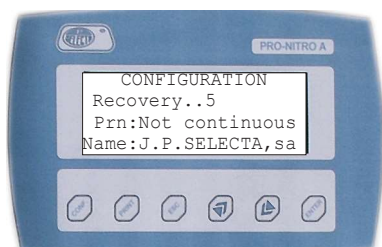
«CONFIGURATION MENU 2»

- **Continuous:** Send data, automatically, on finishing the analysis.
- **Name:** Laboratory name that will appear on the printed reports. Using keys «▲» «▼» and «ENTER» entry 14 Characters.

- **Emp sample:** Configures the collector empty function.

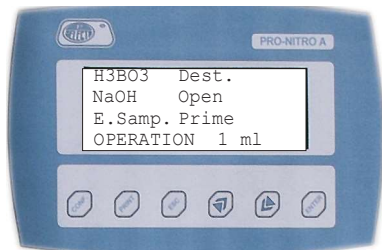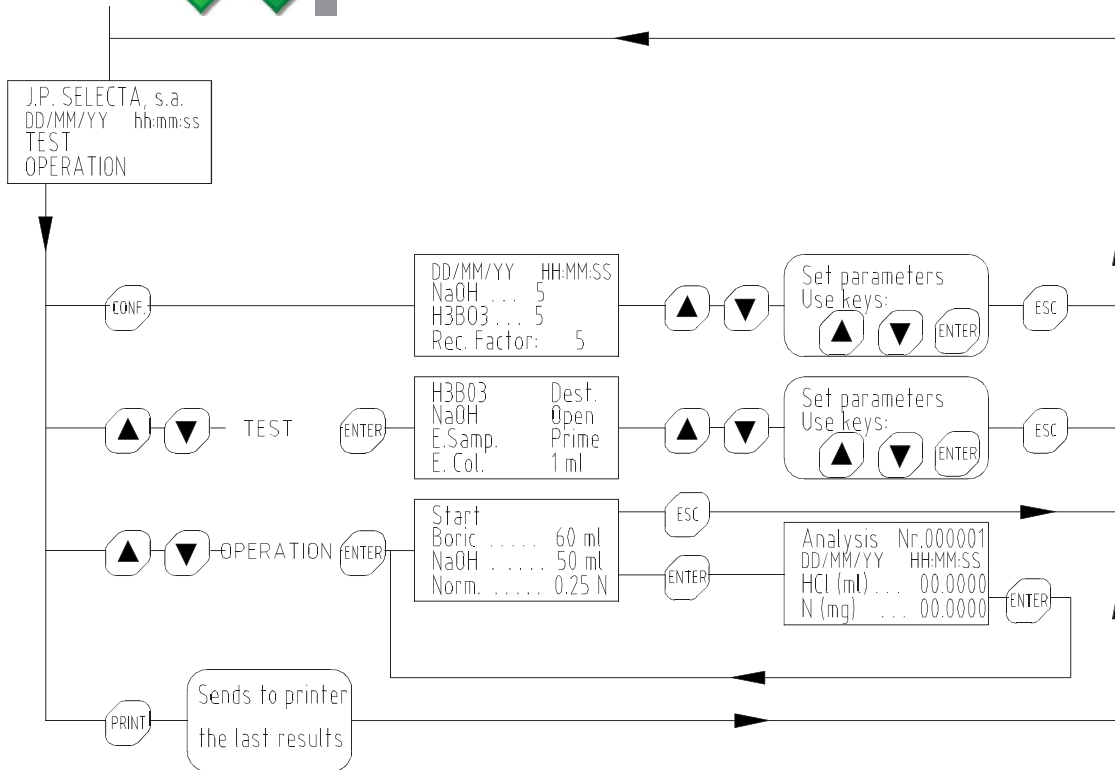To set the default settings, press "CONF" key for 5 seconds in the display.

«MENU TEST»

### 5.1.2 «TEST» Menu

This menu allows execute essential functions for checking, maintenance and setting up operation.

Please achieve a good knowledge of these functions and when it is necessary use each one in order to get optimum performance of the analyser.

As a general rule, select a function by pointing with the arrows and then press «ENTER», and «ESC» to stop it.

The essential functions are the following:

**H3BO3:** Starts the pump to dispense ammonia fixative solution (Boric acid + indicator mixed 5) to the collector. This function is used to prime the dispensing tubing or for its cleaning.

**NaOH:** Starts the pump to dispense NaOH to the sample tube. This function is used to prime the dispensing tubing or for its cleaning.

**E. Sample:** Starts the pump to empty the sample tube.

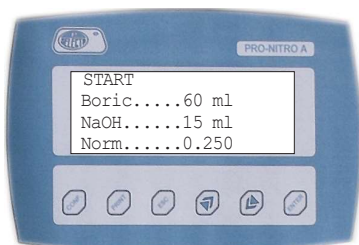
**E. Col.:** Starts the pump to empty the collector. This function is used to clean the distillation circuit.

**Dest:** Initiates the distillation (without automatic titration). This function is used to clean the distillation circuit.

**Open:** Opens the coolant water entry valve (Tap water). This function is used to empty the coolant circuit.

**1 ml:** It gives order to the digital dispenser to dispense 1ml. This function is used for checking or validation purposes.

**Prime:** It gives order to the digital dispenser to dispense continuously until the key «ESC» be pressed. This function is used to prime the dispensing tubing or for its cleaning.

«OPERATION MENU»

### 5.1.3 «OPERATION» Menu

This menu allows programming the analysis parameters and initiating the analysis.

By selecting the «START», the analysis is initiated with the parameters that are shown on the display.

The parameters can be modified pointing it by the arrow keys and pressing «ENTER».

Analysis parameters:

«**Boric**»: Sets the fixative solution volume to be dispensed on the collector. Value factory set is 55 ml and, usually, it will never be modified.

«**NaOH**»: Sets the NaOH to be dispensed on the sample tube. Usually a volume from 50 to 75 ml should be used for sample analysis.

**Note:** This volume depends on the sulphuric acid volume used on the digestion step.

«**Norm.**»: Set the titration reagent concentration (Normality) (usually HCl 0.25N). Values from 0.01N to 1N could be set.

The parameters programmed remain stored on the unit for later use.

**Note:** The precision of the results are affected directly by the precision of the acid normality used for titration. **This is a common cause of error**, because, after some time, opened bottles tend to increase its normality.

Also, often there are errors in concentration calculation to prepare the acid for titration.

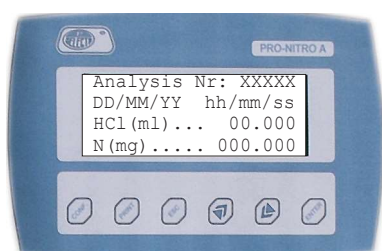
Check often the real value of the normality and update its value on the analyser.

On selecting «START» the distillation & titration are initiated.

On this distillation & titration process, four steps could be identified:

1. Fixative solution (Boric) and NaOH dispensing. The display shows the reagent that is being dispensed.
2. Distillation: for 1.5 minutes the sample is distilled without titration.
3. Distillation & titration: The digital dispenser starts the distilled titration until the colour change is detected.
4. Titration: Just in case that distillation was very long (sample with high nitrogen contents o titration reagent with low Normality) the distillation is stopped but titration continues.

During the steps 2, 3 and 4 the display shows the «RESULTS SCREEN» where the consumed HCl volume (ml) and detected nitrogen (mg) is indicated.

«RESULTS SCREEN»

## 5.2 Alarms

In order to avoid dangerous situations or to protect the analyser life, some alarm systems are provided on the unit:

### «Steam» alarm

It indicates that a low level of water was detected on the steam generator for a long time. The microprocessor stops the distillation.

Probably the water reservoir is empty or the water does not have the suitable conductivity.

### «Tube/Door» alarm

Safety door not closed or sample tube is not inserted.

This alarm blocks the distillation and the NaOH adding.

### «Overtemp» alarm

An excess temperature on the steam generator has switched the safety thermostat.

Block the distillation function.

Once switched, it must be left to get cool for some minutes. Solve the root problem that causes the switching and reset the safety thermostat by pushing the central button. The safety thermostat is located on the electrical box, on the right hand side. A cover must be removed.

The safety thermostat switching is due to low level of water on the steam generator, probably because the water reservoir is empty or water pump malfunctions.

### «Overfill» alarm

Overfill of distilled on the collector.

Blocks the distillation function.

From the menu «TEST» select «E. col» to empty the collector.

### «Low / High signal level» alarm

The initial colour of the ammonia fixative solution present on the collector is wrong. Probably there were some residuals from previous sample not neutralized. Abort the analysis (press «ESC») and proceed with the cleaning operation.

Safety thermostat

**CAUTION**  
Wear Anti-acid,  
Anti-thermal gloves.

The white PTFE tube could have rest of  
Sodium hydroxide or Acid.

The sample tube's temperature is  
above 60°C after the distillation.

## 5.3 Insert/ Remove the sample tube

- Open the safety door.
- Insert the tube starting to introduce the PTFE tube.
- Push down the fixing lever.
- Close the safety door.

Proceed similar to remove the sample tube.

## 6 Sample analysis

To get good analysis' results and a good performance of the distiller it is suggested to follow a working **routine**.

Check water level

Before starting to analyse valuable samples do some training sessions running the operations related on the chapter 8, until the operator will be familiar with the unit.

### 6.1 Check levels of tanks

Before starting a batch analysis, check the following levels:

#### Steam generator water tank level

Refill it with demineralised water.

Assure its conductivity adding about 50mg of sodium carbonate or NaCl.

The water reservoir is 6 litres capacity, enough for analysing more than 20 samples.

#### Level of NaOH tank

Refill with NaOH 40% from the 2 litres flask, placed inside the unit.

Get the NaOH flask from the door placed on the left hand side cover.

The NaOH solution must be free of ammonia salts, to avoid its interference on the result.

#### Level of Ammonia fixative solution tank

Refill with Fixative solution (Boric acid 1% + mixed indicator 4.8) 2 litres flask. It is placed inside the unit.

**Note:** See chapter 7.2 for ammonia fixative solution preparation.

#### Level of Titration reagent

Refill the 2 litres flask placed inside the unit with titration reagent, usually HCl or  $H_2SO_4$  with a concentration of 0.25 N. (or 0.1N)

**Note:** See chapter 7 about reagents.

**Note:** The precision of the results are affected directly by the precision of the acid normality used for titration. **This is a common cause of error**, because bottles opened for a long time tend to increase its normality.

#### Priming the dispensers

The tubing for NaOH, fixative solution and titration reagent must be properly primed, air free, to provide a good and exact dispensing.

Carry on the refill operations to avoid air entry on the tubing. See 5.1.2 and 8.1 to know how to carry on the prime operations.

### 6.2 Unit warm-up and cleaning.

In order to get the best results with your Kjeldahl distiller, you have to run a warm-up step before starting samples analysis.

To do so, just insert a sample tube with 25ml of clean water.

From the «TEST» select «DEST.» and push «ENTER» distillate for 5 minutes and push «ESC» to stop the distillation.

Check the reagents tanks levels before starting a batch analysis.

Then select «E.COL.» and push «ENTER» and observe how the collector is emptied and push «ESC» to stop the emptying function.

Run this warm-up operation, also, if the unit is 2 or 3 hours without distillation.

This procedure is also used for cleaning the distillation circuit.

### 6.3 Distillation (Sample analysis)

Please follow these steps: (on the same order as indicated)

- Open the safety door.
- Check visually that on the collector do not remain rests of the previous analysis. If necessary, empty from the «TEST» menu. (See 5.1.1 to adjust emptying time)
- Insert the sample tube. See 5.2
- From the «MAIN MENU» select «OPERATION». If necessary modify the NaOH volume and the HCl normality.
- Select «START» and push «ENTER»
- The PRO-NITRO "A" run the following steps automatically:
  - Dispense fixative solution on the collector.
  - Initiates sample distillation.
  - NaOH addition to sample.
  - After 1.5 minutes the automatic titration is initiated.
  - The digital dispenser adds HCl on the collector until the colour change is produced (from green to red). The detected nitrogen is shown on the display.
    - Once the red colour is permanent on the collector, the analyser stops the analysis.
      - It empties the collector.
      - It empties part of the sample tube.
- Open the safety door.
- Remove the sample tube. See 5.3
- Check, visually, that the collector remains completely emptied after the analysis. If some rests remains, push the arrow key «▼» for empty operation. (See 5.1.1 to adjust emptying time)
- The unit is completely prepared to analyse the following sample. Any cleaning operation is needed, unless any malfunction has occurred during the analysis.

At the end of the analysis check, visually, through the collector window, that it is completely empty.

#### Visual Checking during the analysis:

Some visual checking can be applied from the colour that takes the solution on the collector or in the sample tube.

This checking is useful at the beginning to set the optimal parameters or to solve unit malfunctions.

#### Visual Checking on the sample tube:

When the NaOH volume is sufficient, the sample becomes blue. (If not, program more NaOH on the following analysis)

#### Visual Checking on the collector:

The initial colour must be the fixative solution colour (Red). If it is grey or green, it is because there are not neutralized rests of the previous sample remaining on the collector. In this case, stop the analysis and run a cleaning operation. See 6.2

## 6.4 Analysis interruption

Press «ESC» to abort the analysis. The process is stopped. Press «ESC» to confirm or «ENTER» to follow with the analysis.

**Note:** The analysis interruption due to not neutralized rests left on the collector will corrupt the following analysis. Run a cleaning operation (See 6.2) before running new analysis.

### Common protein factors

Almonds	5,18
Nuts	5,30
Nut- peanuts	5,41
Jello	5,55
Soya	5,71
Barley, trenches, rye	5,83
wheat, whole flour	5,83
Flours (not whole one)	5,70
Rice	5,95
Corn	6,25
<b>General</b>	<b>6,25</b>
Saved	6,31
Milk and milky products	6,38

## 6.5 Protein contents (%) calculation.

From the HCl volume (ml) consumption on the titration the Kjeldahl Nitrogen (mg) is calculated:

$$\text{Nitrogen (mg)} = 14 \times \text{Volume}_{\text{HCl}} \text{ (ml)} \times \text{Normality}_{\text{HCl}}$$

And, from the nitrogen we calculate the sample protein content:

$$\% \text{ Protein} = \frac{P_2}{P_0} \times 100 \times F$$

P2: Nitrogen (mg).

P0: Sample weight (mg).

F: Protein factor.

## 6.6 Printing results.

The PRO-NITRO «A» includes a serial output for printer connection as printer code 4120113 (See accessories).

Printer must be configured as: 2400, No parity, 8 bits, 1 bit stop,

Once the printer is connected, press "PRINT" for the analysis result to be printed.

See 5.1.1 to set the printer operation configuration.

## 6.7 Storing results on a computer

The PRO-NITRO «A» includes a software to record results on the computer and allows batch reporting.

After each analysis the PRO-NITRO «A» sends the results to the computer: Analysis number, date, time, HCl consumption (ml) and the nitrogen detected (mg).

The user can configure the data sending mode:

- Automatic mode: Once the analyser finishes the analysis, it sends the results to the computer. According to 5.1.1, configure the printer as "Continuous mode".
- Manual mode: Once the analyser finishes the analysis, the user must press the «PRINT» button to send the results to the computer. According to 5.1.1, configure the printer as "Not continuous mode".

## Installation

Connect a serial wiring (See figure) on the RS-232 output from PRO-NITRO «A» to a computer.

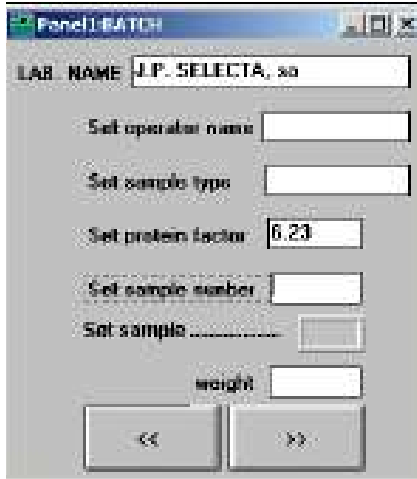
Install the software running the SETUP application from the provided CD.

## COM Port

Each time that the software is opened, choose the COM port where the PRO-NITRO «A» is connected.

Wiring diagram for serial connection from PRO-NITRO «A» and computer serial port

PRO-NITRO «A»	PC
2	3
3	2
5	5
(9 pin female)	(9 pin female)

**The software panels:**

**The batch panel:**

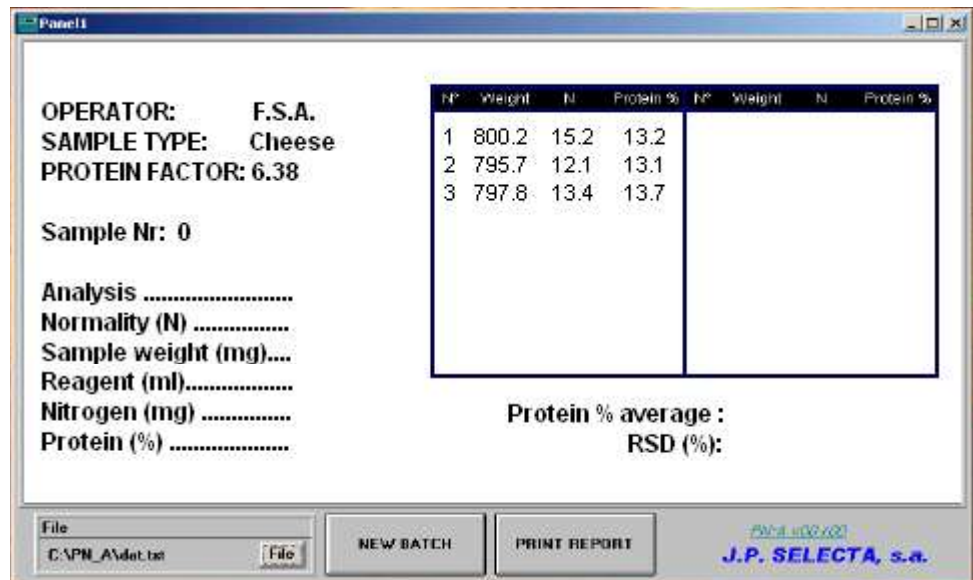
Enter the batch general data on this panel:

- **Lab. Name:** Enter the laboratory name that will appear on the report.
- **Set operator name:** Enter the batch operator name.
- **Protein factor:** Enter the protein factor according to sample type.
- **Sample size:** Enter the quantity of samples that the batch has.
- **Sample:** Choose, using the << and >> keys, the sample and enter the weight (It can be modified).
- **Weight:** Enter sample's weight.

Once all samples' weights are entered, press "OK" to return to the main panel.

**The main panel: (batch analysis)**

- **General data:** (upper left corner) General batch data.
- **Sample results:** (left side) Results of last sample analysed:
  - **Sample Nr:** Number of sample in the batch.
  - **Analysis:** Analysis number.
  - **Normality:** Number of sample in the batch.
  - **Sample weight (mg):** Sample weight.
  - **Reagent (ml):** Reagent volume consumption on the titration.
  - **Nitrogen (mg):** Detected nitrogen.
  - **Protein (%):** Sample protein contents in %.
- **Batch results:** (upper right table) this table summarizes the batch samples results.
- **Global batch results:** (down right corner) results of the batch:
  - **Protein % average:** Batch average protein contents.
  - **RSD (%):** Relative standard deviation (over average)

## 7 Reagents

The reagents to be used must be free of ammonia.

### 7.1 Prepared reagents

The best thing is to use reagents factory prepared, mainly the HCl for titration. Any error on its preparation has a direct influence on the result.

As an example, the following reagents could be used, or the equivalent one in other brands:

- |  |                |
|--|----------------|
| • Boric acid in powder 99.5%   | PANREAC 141015 |
| • Mixed indicator 4.8 (or 5) RV<br>(Red of Metile + Green Bromocresol) | PANREAC 283303 |
| • Mixed indicator 4.4 RV<br>(Red of Metile + Blue of Metilene)         | PANREAC 282430 |
| • HCl 0.1N SV  | PANREAC 171023 |
| • HCl 0.25N SV   | PANREAC 182318 |
| • H <sub>2</sub> SO <sub>4</sub> 0.1N SV                               | PANREAC 181061 |
| • H <sub>2</sub> SO <sub>4</sub> 0.2N SV                               | PANREAC 182011 |
| • Sodium Hydroxide 40% RE<br>(For N determination)                     | PANREAC 171220 |
| • Acetanilide 99% (Validation standard)                                | PANREAC 151005 |
| • Ammonia sulphate (Validation standard)                               | PANREAC 131140 |
| • Kjeldahl Catalyst 6.25% Cu tabl. 8g                                  | PANREAC 174428 |

### 7.2 Ammonia fixative solution preparation

The preparation of the ammonia fixative solution **is critical** to get a good operation with PRO-NITRO «A».

This solution must be prepared, working accurately, and using material clean of strong acids or alkalis.

#### Note:

- The use of different fixative solutions to the specified ones do not warranty the equipment good operation.

- Leave the ammonia fixative solution between 24h and 48h before checking, adjusting and using it.

#### 7.2.1 Preparation for Mixed indicator 4.8 (or 5) (Easy)

Prepare according to this formula for 1 litre of solution:

- |  |                |
|--|----------------|
| - Weight 10g of Boric acid (in powder) | PANREAC 141015 |
| - Solve in 1 litre of distilled water. |                |
| - Add 15ml of mixed indicator 5        | PANREAC 283303 |

#### 7.2.2 Preparation for Mixed indicator 4.4

Prepare according to this formula for 1 litre of solution:

- |  |                |
|--|----------------|
| - Weight 10g of Boric acid (in powder) | PANREAC 141015 |
| - Solve in 1 litre of distilled water. |                |
| - Add 10ml of mixed indicator 4.4      | PANREAC 282430 |

#### 7.2.3 Preparation for indicator according to ISO 5663

Prepare according to standard ISO 5663 / EN 25663. For water analysis.

Modify on the «CONFIGURATION» menu the following parameters:

- LS: Change to value 10
- Light: Change to value 10

**Note:** If this solution is not prepared with sufficient accuracy, it may be necessary to adjust the assessment parameters with several tests.

### 7.3 Fixative solution checking and adjusting

Due to the differences existing between the water used for ammonia fixative solution preparation (deionised, distilled from different assay,...) it is necessary to check and adjust this solution to achieve a correct operation of the equipment.

#### 7.3.1. Material

- Spectrophotometer suitable to work at 520 nm.
- 10 mm glass cuvette.
- Flask with 0.01 ml resolution.

#### 7.3.2 Using Boric acid solution 1% + 15 ml indicator: (4.8 mixed, Metihyl red - Bromocresol green)

##### 7.3.2.1 Checking:

- Select 520 nm in the spectrophotometer.
- Adjust blank with distilled water.
- Measure the fixative solution absorbance value.
- The value must be approximately around 0.650 (between 0.630 and 0.670).

##### 7.3.2.2 Adjusting:

If the value is lower than 0.630:

Add HCl 0.1M till you get a value between 0.630 and 0.670 according to the following formula:  $V = (0.650 - L) \times 1515$

where:

V = HCl 0.1M volume to be added per dissolution litre (ul).

L = Solution absorbance at 520 nm.

If the value is higher than 0.670:

Add NaOH 0.1M till you get a value between 0.630 and 0.670 according to the following formula:  $V = (L - 0.650) \times 1515$

where:

V = NaOH 0.1M volume to be added per dissolution litre (ul).

L = Solution absorbance at 520 nm.

##### Note:

- These formulas are orientative. Add a lower quantity than calculated one, and check again the result.
- The expression "ul" should be read as "microlitre".

##### 7.3.2.3 Example:

We have a solution whose absorbance (at 520nm) is: **L = 0.559**

- According to 7.3.2.2, we have to add:  
 $V = (0.650 - 0.559) \times 1515 \text{ ul} = 137.86 \text{ ul}$   
(of HCl 0.1 M per solution litre)
- Instead of adding **137.86 ul**, we add **100 ul** of HCl 0.1 M.
- We repeat the reading.
- The new reading is: L = **0.611**
- We add 50 ul more and repeat the reading L = **0.640**
- Now we have the solution prepared to work.

## 8 Checking the distiller (OQ)

Once the unit is properly installed and the operator has acquired the knowledge to work with it, the unit must be checked.

This process is known as Operation Qualification.

There are four basics checking to do:

### 8.1 Functional checking

Run this checking after the installation and also for some maintenance.

The «TEST» menu is used for components checking.

#### 8.1.1 Checking the steam generator filling

When starting the unit for the first time, the water pump starts to refill automatically the steam generator. This operation usually takes around 40 seconds and then the unit stops also automatically. After 20 seconds, the «STEAM» alarm is activated, which will also stop by itself.

If, approx. after these 40 seconds this operation is not finished, switch the equipment off. Check that maybe the water tank has no water or that the water has not the suitable conductivity.

If so, check, empty the steam generator by opening the drain and repeat the operation.

#### 8.1.2 Checking the ammonia fixative solution

Insert a MACRO tube.

From «TEST» menu, select «H3BO3», press «ENTER», and this starts the dispensing pump, it takes liquid from the 2 litres tank to collector flask.

Check, visually that the liquid gets the collector and press «ESC» to stop the pump.

Run this operation when the dispensing tube must be primed. If tubing is not primed the dispensed volume will be wrong.

#### 8.1.3 Checking the NaOH pump and tubing

Insert a MACRO tube.

From «TEST» menu, select «NaOH» and press «ENTER», this starts the NaOH dispensing pump, it takes liquid from the 2 litres tank to sample tube.

Check visually that the liquid gets the collector and press «ESC» to stop the pump.

Run this operation when the dispensing tube must be primed. If tubing is not primed, the dispensed volume will be wrong.

#### 8.1.4 Checking and priming the digital dispenser

Run this operation to check the digital dispenser.

From «TEST» menu, select «PRIME», and press «ENTER», the digital dispenser starts to pump HCl, in a continuous way from the 2 litres tank to collector flask.

Press «ESC» to stop the operation.

Run this operation when the dispensing tube must be primed. If tubing is not primed, the dispensed volume will be wrong.

### 8.1.5 Check the precision of the digital dispenser

This operation can be used to measure the digital dispenser accuracy.

To get an accurate dispensing the tubing must be primed. (8.1.5)

Remove the dispensing tube (see figure) from the collector to dispense into a calibrated burette.

From «TEST» menu, select «1ml» and press «ENTER».

The dispensed volume must be 0.97ml and 1.03ml.

**Note:** This operation can be done in a more accurate way by using a scale.

After operation, place the dispensing tube into the collector. Check visually, across the collector window, that the tube is properly inserted.

### 8.1.6 Checking distillation «Cleaning»

This operation is often used when working with the PRO-NITRO «A» because it performs a cleaning of the distillation glassware. It is also used for unit warm-up.

Place a MACRO sample tube with near 25ml of distilled water.

From the «TEST» menu, select «Dest», press «ENTER», this starts the steam production and the coolant water circulation. This is the distillation.

After the warming-up, the distillation rate will be between 30 - 40 ml / minute.

Left 5 minutes distilling and then press «ESC» to stop distillation.

After this operation, the collector must be emptied. See 8.1.9

Run this operation when the distillation circuit need to be cleaned.

Run this operation for warming-up the unit before a batch analysis or after more than 2 hours of inactivity.

### 8.1.7 Checking sample tube empty function.

Place a MACRO sample tube with around 100ml distilled water.

From «TEST» menu select «E.Samp», press «ENTER». This starts the pump to empty the sample tube to the waste output.

### 8.1.8 Checking collector empty function.

This operation is used together with 8.1.7 for cleaning purposes of the equipment.

After 5 minutes of water distilling (See, 8.1.7) the collector will collect around 100ml of water.

From «TEST» menu, select «E.col», press «ENTER». This start the pump to empty the collector to waste output.

### 8.1.9 Collector contaminated

The titration will be wrong if at the beginning of the analysis the collector has any drop of HCl.

The titration will be also wrong if at the beginning of the analysis the collector has any drop of distilled from previous sample not neutralized by titration.

The equipment detects this situation by indicating: "Low/High Signal Level", which is normally solved by cleaning the collector.

So you must be sure the collector is clean before any analysis, see 8.1.7.

Normally, a cleaning operation must be run after running operations 8.1.3 and 8.1.5.

### 8.2 Checking Nitrogen Recovery.

This checking is widely used to certify that the unit works properly.

Run some of these tests for training purposes, before starting to analyse valuable samples, until the operator will be familiar with the unit.

It consists on preparing some samples that have a known Nitrogen content. It is distilled, titrated and the nitrogen detected is calculated.

The rate (%) of nitrogen detected versus the sample Nitrogen content gives us the **Nitrogen Recovery**.

To allow this checking, a 100gr ammonia sulphate flask is delivered together with the distiller.

- Weight near 100mg (50mg reagent 0.1N) of ammonia sulphate. The exact weight will be  $P_0$

- The exact quantity of Nitrogen is:  $P_1 \text{ (mg)} = P_0 \times 0,212$

- Add around 25ml of  $H_2O$

- Distillate adding 15ml of NaOH

- Use HCl 0.25N (0.1N) as a titration reagent.

- Once distillation and titration are finished, the PRO-NITRO «A» gives the detected nitrogen (mg):  $P_2$

- The recovery is calculated:  $R(\%) = P_2 / P_1 * 100$

- Acceptable values are from 99.0 to 101%.

**Note:** Some literature sets that the acceptable values for recovery must be above 99.5%. To get this result you must work with high precision and use reagents with high purity degree.

### 8.3 Blank test

This checking must be run on each batch in order to quantify the absence of ammonia on the reagents.

It consists on distilling and titrating a sample of 25ml distilled water.

Add 15ml of NaOH on the distillation.

Run some blank tests in order to get a constant value of blank. If this value is repeated, it could be subtracted from the detected Nitrogen.

#### How many ammonia sulphate must I use on the test sample?

The HCl volume ( $Vol_{HCl}$ ) for a quick analysis must be between 6 and 12 ml.

From the expresion:

$$P_1 \text{ (mg)} = 14 \times Vol_{HCl} \text{ (ml)} \times N_{HCl}$$

We arrange it as:

$$P_0 \text{ (mg)} = \frac{14 \times Vol_{HCl} \text{ (ml)} \times N_{HCl}}{0.212}$$

And get the table:

Normality	Sample amount suphate amm.
0,05	20 ... 40 mg
0,1	40 ... 90 mg
0,25	100 ... 200 mg
0,5	200 ... 400 mg

The volume consumption of HCl ( $Vol_{HCl}$ ) on the titration can be forecasted from the expression:

$$P_1 \text{ (mg)} = 14 \times Vol_{HCl} \text{ (ml)} \times N_{HCl}$$

thus we obtain that HCl must be:

$$Vol_{HCl} \text{ (ml)} = 14 / [P_1 \text{ (mg)} \times N_{HCl}]$$

### 8.4 Checking Nitrogen recovery with Acetanilide

This test is widely used to certify the whole Kjeldahl analysis process, including the digestion, distillation and titration steps.

It consists on preparing some samples that have a known Nitrogen content. The samples are digested, distilled, titrated and the nitrogen detected is calculated.

Run this test some times to get familiar with the equipment and the Kjeldahl analysis before analysing valuable samples.

The rate (%) of nitrogen detected versus the sample Nitrogen content give us the **Nitrogen Recovery**.

Acetanilide is used to run this test.

#### Preparing the sample:

Nitrogen content of an acetanilide sample:

Formula:  $C_8H_9NO$

Molecular weight: 135.17

Factor:  $14 / 135.17 = 0.1035$ mg of Nitrogen =  $0.1035 * \text{mg Acetanilide}$ .

- Weight around 250mg (100mg) of acetanilide. The exact weight  $P_0$
- The Nitrogen content is:  $P_1 \text{ (mg)} = P_0 \times 0,1035$

#### Sample digestion:

• Place the acetanilide on a MACRO tube, add 10ml of Sulphuric acid 98% (for example PANREAC 173163) and one tablet (or 8gr) of Kjeldahl catalyser.

• Digest at 400°C for 1h. (The result must be a transparent blue liquid)

• Let cool at ambient temperature. Add 25ml of distilled water on each tube. Operate with caution to avoid splash. Water reaction over the sulphuric acid could be violent.

#### Distillation:

• Distillate by adding 75ml of NaOH. Use HCl 0.25N (0.1N) for tritration.

• Once sample is distilled and titrated following chapter 6, the detected Nitrogen is calculated as:  $P_2$

• Recovery is calculated as:  $R(\%) = P_2 / P_1 * 100$

• Acceptable recovery is from 99.0 to 101 %.

#### Note:

If MICRO size tube is used, weight 0.3g of acetanilide, add 5ml of sulphuric acid, and one tablet of Kjeldahl catalyser. On the dilution add 10ml water.

Distillate for 6 minutes adding 35ml of NaOH.

#### Note:

Some literature sets that the acceptable values for recovery must be above 99.5%. To get this result, you must work with high precision and use reagents with high purity degree.

## 9 Maintenance

The distiller PRO-NITRO «A» uses NaOH and concentrated acids for the analysis process. These substances are very aggressive and will degrade some parts of the unit.

To avoid derating and increase distiller's life a suitable cleaning routine should be done.

### IMPORTANT NOTE

Read the preventive maintenance instructions at the last pages of this instruction manual.

### 9.1 Cleaning the unit:

#### 9.1.1 Clean the distillation circuit:

After any analysis session, distillate for 5 minutes in order to remove any rest of NaOH on the distillation glassware. See 6.2

#### 9.1.2 Cleaning the chamber:

The chamber is protected with a chemical resistant coating. But it should be daily cleaned to increase its life.

Use a well damp cloth to clean the chamber and its components to remove any rests of NaOH.

Clean the tray with abundant water.

#### 9.1.3 Cleaning the NaOH and Boric acid hose connection

Approximately every 3 months, clean the NaOH and the fixative solution from the tank hose connection with a wet cloth.

#### 9.1.4 Cleaning the steam generator (weekly)

Switch off the equipment. Open the «Steam generator drainage valve» and wait till it will be completely empty. Close the valve and start the equipment. See how the steam generator is filled in chapter 5 of this manual.

### 9.2 Hose inspection (monthly)

Visually check the hose condition and its connections to components.

To avoid premature derating of the unit due to liquid leakage, replace as soon as possible any hose that has any leakage or damage.

#### 9.2.1 Hose replacing:

Take these cautions to hose replacing:

- Unplug the unit from the electrical power supply.
- Check, as you can, (See 9.3) that the hose to remove is empty.
- Use only tubes material as indicated on 11.2.
- Use clamps as is factory assembled.
- Once the hose is replaced, it must be checked for leakages.

Clean periodically the NaOH tank and fixative solution tank hose connector.

### 9.3 Preparing the unit for shipping or storing.

Once the unit has been used with NaOH, proper actions should be taken for its transportation or storing.

Initially clean the unit according to chapter 9.1

#### 9.3.1 Remove completely the NaOH from the unit:

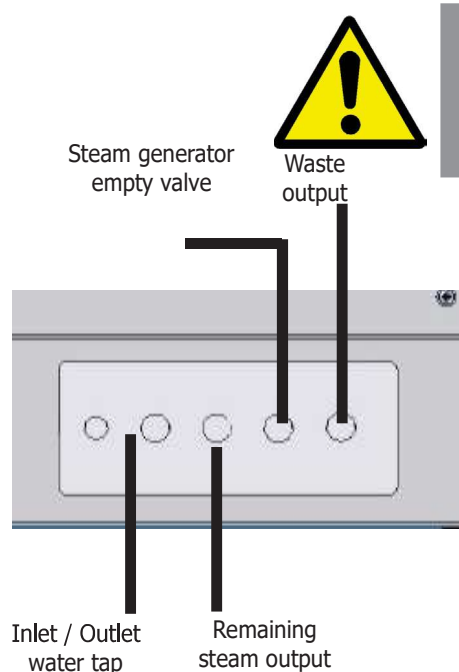
- **Clean the NaOH tank:**

Remove the NaOH 2 litres tank from the unit.  
Clean it with abundant water for some minutes.

- **Cleaning the NaOH dosing circuit:**

Fill the NaOH 2 litres tank with water.  
Connect it to NaOH circuit.  
Insert a MACRO tube on the unit. From «TEST» menu, select «NaOH» and press «NaOH» several times till the sample tube is full.

Repeat the operation with the Fixative solution tank. (On the menu «TEST» select «H3BO3» and later «E.Col»)

Cooling circuit valve

**CAUTION**  
The NaOH is very corrosive for the skin  
Use chemical resistant gloves.  
Take proper cautions procedures for NaOH handling.

- **Emptying the HCl tubing:**

1.- Remove the HCl dosing tube from the collector. (See 8.1.6) Place a flask at the HCl tube output to collect the HCl.

2.- Remove the HCl tank from the unit.

3.- From «TEST» menu, select «PRIME». The digital dispenser starts to dispense all the HCl that is on its tubing. Press «ESC» when no more HCl outputs from the tubing.

#### 9.3.2 Remove completely the water from the unit:

- **Emptying the water reservoir and the steam generator:**

1. Open the tap "Steam generator empty valve".

2. Start the unit, and wait that the pump empties all the water from the water reservoir. The «STEAM» alarm will be activated but you do not have to do anything.

- **Emptying the water from the cooling circuit.**

1. Open the "cooling circuit valve".

**CAUTION**  
Do not ship off the unit if the chemicals (NaOH, H3BO3 and HCl) and water have not been completely removed.

During transportation, chemicals or water leakages could fall over important components and cause severe damages.

During transportation, chemicals or water leakages could fall outside of the package.

### 9.4 Level sensor cleaning

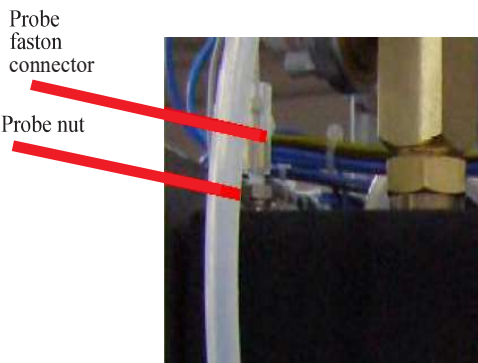
Inside the steam generator a constant water level is necessary.

To provide this water level there is a conductivity based level probe (a steel rod) that when it loses contact with the water, it shows a high resistance to the electronics and then the microprocessor orders the water pump to feed the steam generator with water from the reservoir up to the probe level.

After some time working with this unit (some months), some residues are settled on the probe causing a malfunction.

Effects of this problem:

The residues short-circuit the path between the probe and the steam generator walls (connected to earth) and so the sensor shows a LOW resistance to the electronics. It is the same signal as if the steam generator has the correct water level. This makes that the microprocessor does not supply water and this will probably cause the safety thermostat action.

This means that the probe must be removed and clean. You will need a cross screwdriver and a 12mm spanner and a 19mm one.

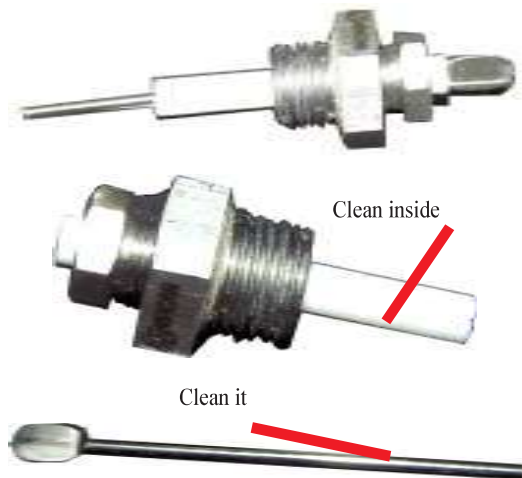
The cleaning frequency depends on two factors:

- The hardness of the water used.
- The usage of the unit.

As a general rule, please do it every **6 months**.

- Remove the left hand side cover (6 screws).
- Locate the steam generator.
- Unplug the probe connector (white cable with FASTON connector).
- Unscrew the sensor with the 19mm spanner.
- Unscrew the probe with a 12mm spanner.
- Remove the steel rod from the PTFE (white part).
- Clean the steel rod and the white PTFE part from the calcium deposits. Use diluted HCl or any home cleaner product for calcium deposits.
- Assemble it again.

## 10 Technical specifications

Mains voltage:	230V 50/60Hz (See plate)
	Other power supply voltages could be available (ask to your dealer)
Electrical power:	1900W (See plate)
Maximum consumption:	8.5 A
Fuses:	10 A fast. dimensions: 20x5 mm
Power supply cord:	Phase, Neutral, Earth (section 1mm <sup>2</sup> )
Weight (Net):	35 Kg
Dimensions: W x L x H	500 x 490 x 740 mm
Water reservoir volume:	6 litres
NaOH, H3BO3 and HCl tank volume:	2 litres
Nitrogen recovery:	>99.5%
Detection range:	0.2 to 200 mg
Reproducibility: relative standard deviation	< 1%
Normality range	0.010 to 1.000 N
Detected nitrogen resolution:	0.0001
Distillation rate:	approx. 30 / minute
Analysis time for 50mg of N:	6 minutes approx.
Analysis time for 100mg of N:	9 minutes approx.
Cooling water entry connector:	3/4"
Cooling water output connector:	3/4"
Remaining steam output connector:	3/4"
Dosification for NaOH pump:	0 to 250 ml $\pm$ 5%
Dosification for H3BO3 pump:	50 to 65 ml $\pm$ 5%
Sound level:	Lower than 60 dBA
Ambient conditions:	Inner use. Altitude up to 2000m Temperature from 5°C to 40°C. Relative maximum damp 80% for above 31°C temperatures, decreasing lineally to 50% for 40°C.
Overvoltage category:	Category II
Pollution degree:	2
Glass material:	Borosilicate 3.3
Door and front material:	ABS injected
Carriage material:	Iron with overlap
Chamber material:	Aluminium with RILSAN coating chemical resistant.
Steam generator material:	Stainless steel.
Hosepipe material:	Silicone, synthetic fluorated elastomer.
RS-232 output:	Connector 9 pins male. Pin 2 and 3 interchanged, pin 5 common. Baud rate 2400, no parity, 8 bit data, 1 bit stop.

# 11 Technical service

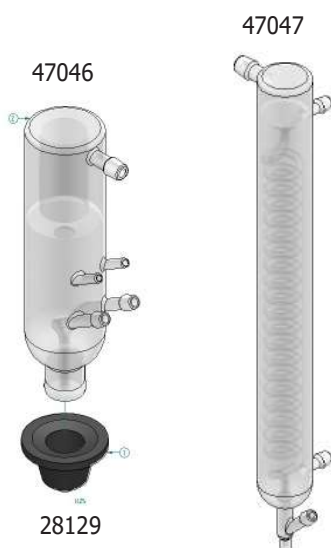
For repair and substitution of deteriorated parts please contact with J.P. SELECTA s.a.u. or its distributor. They will inform you about the technical services authorized in your area.

Find a spare parts list for easy buying and identifying.

## 11.1 Spare parts

32052

47046

47047

28129

24218

- 16053** Electrovalve N. C. 24V PRO-NITRO A Collector
- 16087** Electrovalve N. A.
- 16067** Water inlet electrovalve
- 24218** Water, collector emptied and sample emptied
- 24278** KNF pump (NaOH and fixative solution)
- 24120** Stirrer motor
- 28129** Lock black rubber
- 28404** Shelf
- 32052** Steam generator 1800W
- 47021** Colorimeter flask
- 47046** Splash protector
- 47047** Refrigerant
- 47259** Anti feedback valve
- 29503** PRO-NITRO A Colorimeter PCB
- 29506** PRO-NITRO A LED's PCB
- 29508** Lee digital dispenser PCB
- 29515** PRO-NITRO A Keyboard PCB
- 29516** PRO-NITRO A Power supply PCB
- 29517** CPU board and PRO-NITRO A relay
- 29530** Level

Hoses:

- |              |                                 |                              |
|--------------|---------------------------------|------------------------------|
| <b>46030</b> | <b>Teflon tube 8x10</b>         | Tube immersed on the sample. |
| <b>46041</b> | <b>Silicone tube 8x14</b>       | For water and steam.         |
| <b>46042</b> | <b>Silicone tube 6x10</b>       | For water and steam.         |
| <b>46143</b> | <b>Pressure hose</b>            | Tap water entry.             |
| <b>46084</b> | <b>ISO VERSINIC Black 10x14</b> | Distilling glassware joint.  |
| <b>46139</b> | <b>ISO VERSINIC Black 6x10</b>  | NaOH                         |


- Water (24218)
- Collector emptied (24218)
- Sample emptied (24218)
- Boric (24278)
- NaOH (24278)


**WARNING**  
 Before any work inside the equipment  
**UNPLUG** it from the MAINS.

**11.2 Hydraulic tubing diagram**


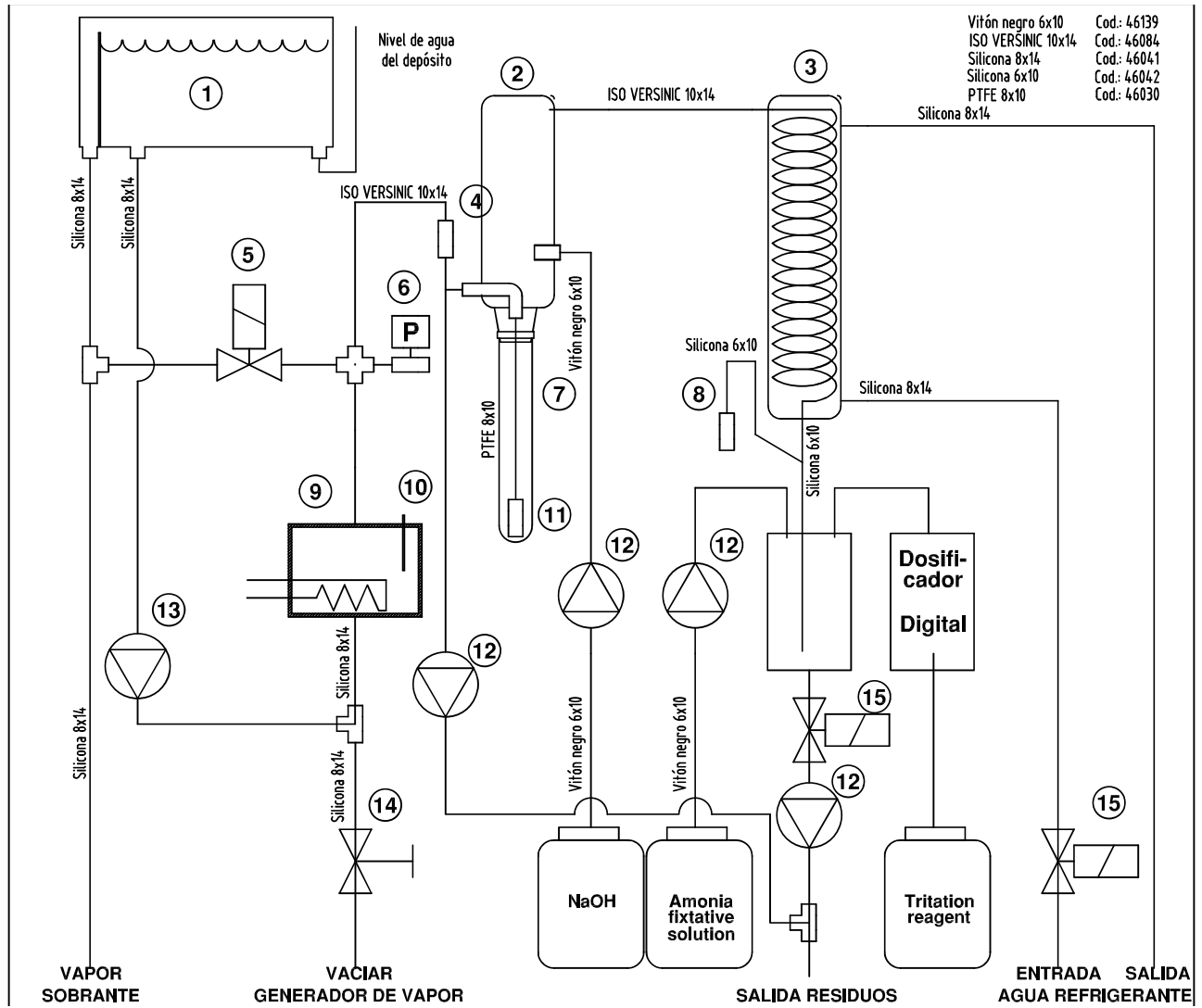
### 11.3 Wiring diagram (1 of 2)




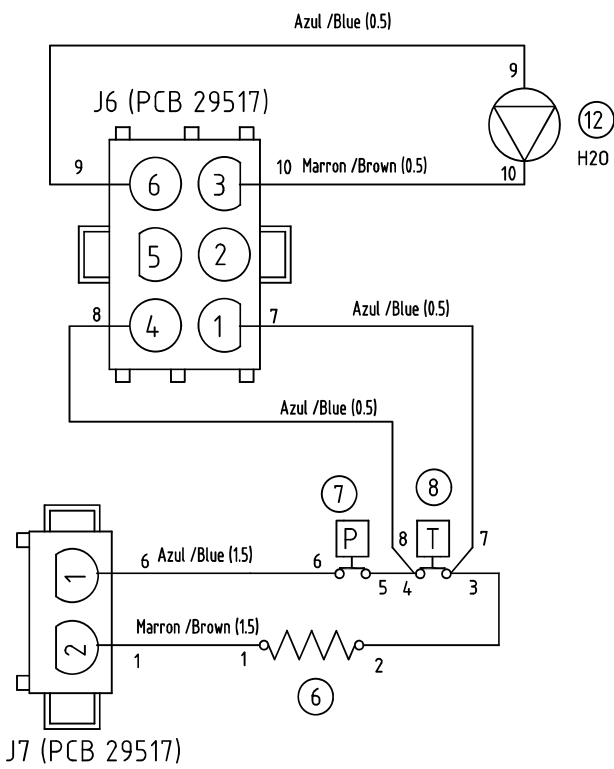
J12 (PCB 29517)

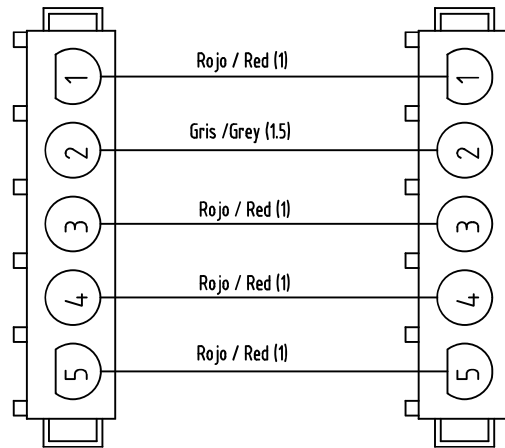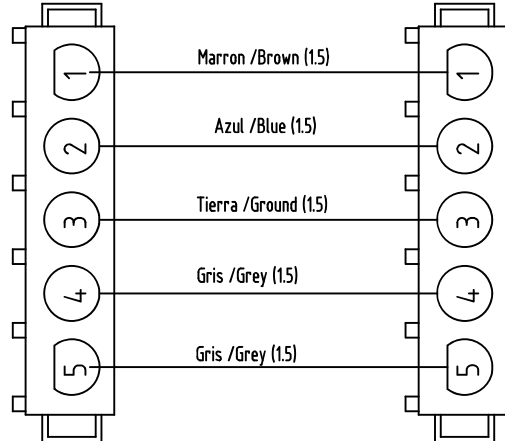J2 (PCB 29516)





J13 (PCB 29517)

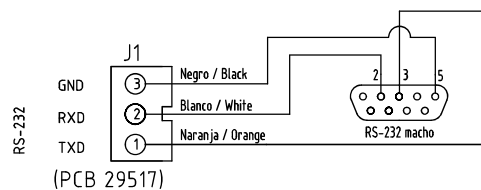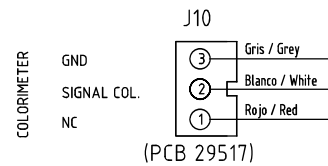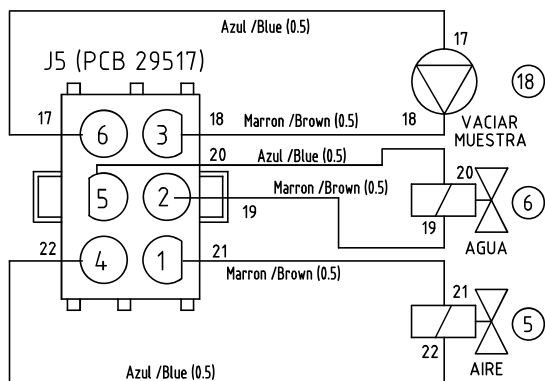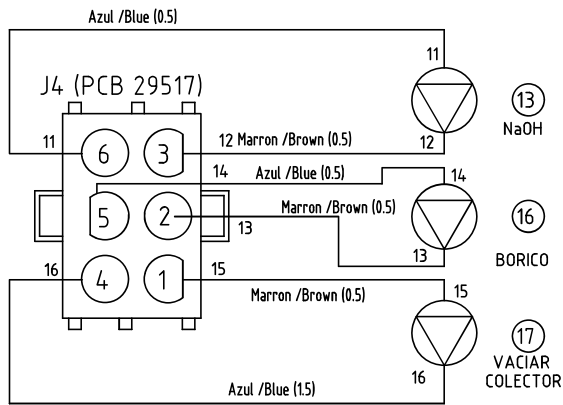J3 (PCB 29516)

### 11.3 Wiring diagram (1 of 2)
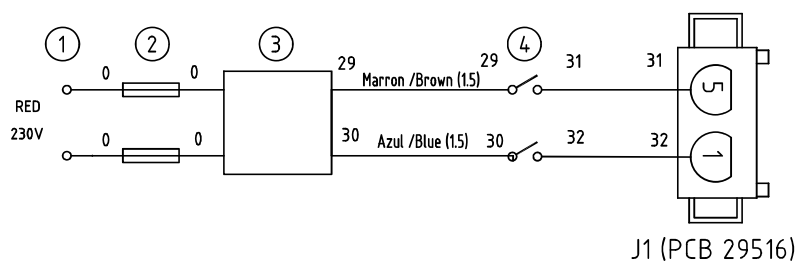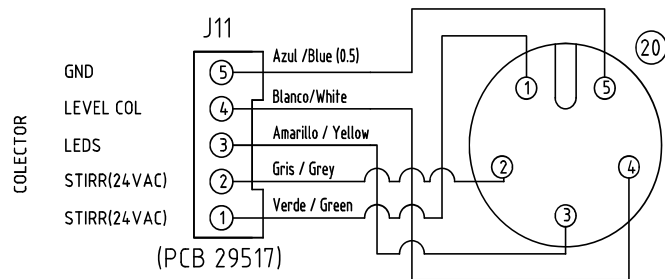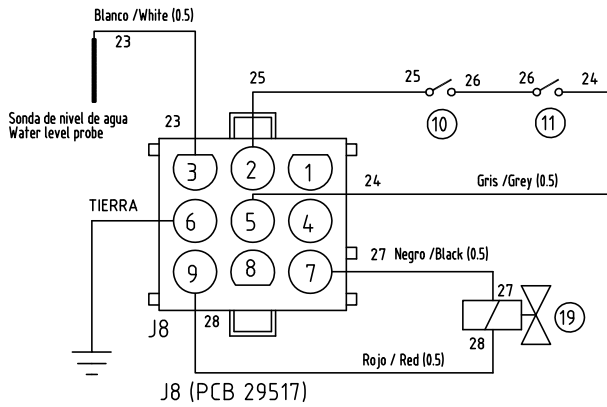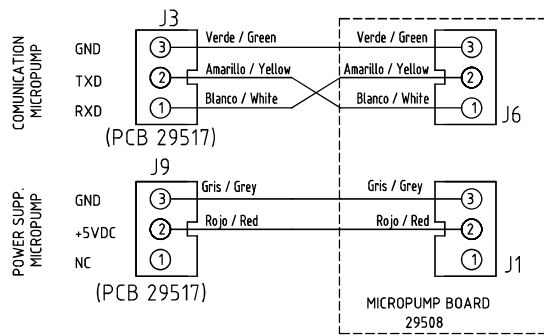
### 11.4 Diagnostic menu

The analyser PRO-NITRO «A» includes an utility for checking and adjusting some important parameters for properly working. It is mainly used for:

- Checking light level.
- Checking functions 8.1 without alarms.
- Checking colorimeter output signal level during titration and detecting problems.

#### Accessing to diagnostic utility:

1. The PRO-NITRO «A» must be OFF (Mains switch OFF).
2. Press «ENTER» key. Keep pressed and press mains switch ON.
3. After a few seconds «STIRRING» menu appear on the display. See attached diagram.

#### Diagnostics utility operation:

For function selection, use the keys:

1. The keys ▲«» and ▼«» to point any function.
2. Press «ENTER» to start or stop the pointed function.
3. Press «ESC» to output from the diagnostics utility.

#### Diagnostic functions

Some of the functions on this menu do the same as others from the «TEST» menu, except that here the alarms are disabled.

#### CAUTION

When the diagnosis functions are running the alarm systems are **DISABLED**

#### 1. Check the light level

The display shows the light level: C0 = Level from 200 ±75.

#### 2. Diagnosis of basic functions

<b>STIRRING:</b>	Stop / run of collector stirrer.
<b>EMPTY COOLING:</b>	Open/Closes the coolant (Tap) water.
<b>TITRATION:</b>	See. 3 (below)
<b>ADD 1ml:</b>	Orders to digital dispenser a 1ml volume.
<b>PRIME:</b>	Stop / run digital dispenser.
<b>EMPTY COLLECTOR:</b>	Stop / run collector empty system (pump + valve).
<b>EMPTY SAMPLE:</b>	Stop / run sample tube empty system (pump).
<b>STEAM COOLANT:</b>	Stop / run distillation (without titration)
<b>BORIC:</b>	Dispenses 50ml of fixative solution (Boric).
<b>NaOH:</b>	Dispenses 50ml of NaOH.

#### 3. Colorimetric titration diagnostics

Using keys «▲» y «▼» point «TITRATION».

To run this diagnostic, a sample of ammonia sulphate must be prepared (See 8.2). During the analysis the following figures appears on the display:

«Ls»	Upper limit.	«Li»	Lower limit.
«Rd»	Colorimeter reading.	«D»	Dispensed volume of HCl.

Diagnostic: (normal values): Ls: [100 ... 250] , Li: [75 ... 150]

- The reading Rd, with green colour must be **above** «Ls»
- The reading Rd, with red colour must be **under** «Li»

## 12 Kjeldahl analysis step by step

### 12.1 Sample preparation

- Triture, homogenize and mix the sample.
- Weight between 1 and 2 grams of the sample.
- In low nitrogen content samples, (waste waters, etc...) take enough sample to get 5mg nitrogen.

### 12.2 Digestion

- Add between 10-15ml (MACRO tube) of  $H_2SO_4$  96-98% and one catalyser tablet (8gr) (for MICRO tube maximum  $H_2SO_4$  is 5ml).
- Rig up a system for smoke extraction or scrubber with  $Na_2CO_3$ .
- Make this digestion in 3 steps:

1. According to the level of water in the sample, begin digestion at 150°C for 20 or 60 minutes.
2. Make a second step by 280°C for 30 minutes to reduce white smoke production.
3. Continue the digestion with 400°C for 60-90 minutes.

**Visual checking:** the result is a neat transparent liquid with a clear blue coloration or green or yellow depending on the catalyser used. No black rest must be stuck on the tube's wall.

**Note:** During digestion you must control foam production in the samples. If there is too much foam, length step 1.

### 12.3 Dilution

- Take out the sample tube off the digester block and let it cool to ambient temperature.
- Add about 25ml of water in each tube (10ml for MICRO tube).
- Slow add water, moving the tube to not let solidify the sample. Heat the tube if necessary. (e.g. introducing the tube into the digester block)
- Let it cool again.
- To avoid nitrogen loss and violent reactions do not introduce the hot tube inside the distiller.

### 12.4 Distillation

- Check the tank level of NaOH and Boric acid. Check the dispensing pumps are primed and give the correct volume.
- Put a 250ml Erlenmeyer in the outlet cooler with 50ml Boric acid and some indicator drops.
- Program a 50 to 75 ml NaOH dosage (35ml for MICRO tube).
- Program the HCl normality.
- Insert the tube with the sample in the distiller.
- Initiate distillation / titration.
- The PRO-NITRO «A» shows the nitrogen detected at the end of the analysis

**Visual checking:** Once the NaOH has been added, the sample must get a blue coloration, if not add more NaOH. It can be added more NaOH by pressing «▲» before titration display appears.

### 12.5 Protein % content calculation

- Apply the protein factor, according the sample type. Use:

$$\% \text{ Protein} = \frac{P_2}{P_0} \times 100 \times F$$

P2: Nitrogen (mg).  
 P0: Sample weight (mg).  
 F: Protein factor.  
 (6.25 default)

## 13 Guidelines for IQ y OQ

Some laboratories need the installer make an IQ document to certificate the good installation of this equipment. As well as in OQ document to certificate the good working of this equipment.

The installer can elaborate those documents following the examples:

### Checking list for IQ (Installation Qualification):

- Installer name.
- Manufacturer name.
- Unit identification: Code, model and serial number.
- Unit location: Laboratory, person in charge.
- Checking packing list
- Unit location matches the requirements?

Water and electricity requirements are suitable?

- Electrical supply  
Socket:
  - Ground terminal?
  - Suitable to power consumption (16A)
  - Voltage according plate  $\pm 10\%$
- Water connection.
  - Water pressure bigger than 1bar.
  - Hose to waste are properly fixed and not blocked.
- Water for steam generator: Low level of calcium, conductivity above 20 Micro-Siemens.
- Personnel training: Operator personnel understand the instruction manual?
- Date and name of who has made the installation (Signed).

### Checking list for OQ (Operation Qualification):

- Installer name.
- Manufacturer name.
- Unit identification: code, model and serial number.
- Unit location: laboratory, person in charge.
- Check the equipment distillate correctly. See: 8.1
- Check nitrogen recuperation with ammonia sulphate sample. See 8.2.  
Note : ammonia sulphate weight, HCl normality, HCl consumption and obtained recovery.
- Check the white test See: 8.3  
Note: Ammonia sulphate weight, HCl normality, HCl consumption and obtained recovery.
- You can check the nitrogen recovery with acetanilide sample. See 8.4.  
Note: Acetanilide weight, HCl normality, HCl consumption and obtained recovery.
- Date and name of who has made the installation (Signed).

## 14 Warranty

The warranty covers the defective components or faulty assembling of the unit for 1 year since the purchasing date.

Keep purchasing bill as the only valid document for the purchasing date.

The warranty does not cover the damages caused for a unit malfunction. The unit must be suitably qualified before to be used for valuable samples processing.

Return goods are not accepted without previous J.P. SELECTA, s.a.u. approval.

## 15 Troubleshooting table

### 15.1 General operation:

- **The green indicator «ON» does not light.**

Check fuses and wiring connections cables.

Check the main switch.

- **The equipment still adding water till it breaks out off the sample.**

The equipment does not detect the water level for low conductivity. Add about 250ml of tap water in the water tank.

Check the cable (white) of the level probe of the steam generator.

- **At the end of distillation the sample tube content is absorbed to the steam generator.**

Check that the steam hose is not obstruct.

Locate the aerated electrovalve:

Check it actives, check that the voltage with a tester (230V).

Check (detaching it) that it is not obstructed.

### 15.2 Alarm anomalies:

- **The alarm (2) for hose presence and lock door is always active and block the unit running.**

Check the door is correctly closed.

Check the microswitches are o.k. (you will hear a click) when you close the door or lower the lever.

Check the cable (grey) continuity.

- **The refrigerating water does not enter.**

Check if the tap is open.

Locate the refrigerating water input valve, check the voltage with a tester(230V). If there is voltage and the valve does not open, knock it. That is due to a small lime rest blocking the water pass. If knocking is not enough, detach the valve to clean the water pass.

- **Security thermostat alarm activates and the equipment is blocked.**

This alarm activates when there is an excessive temperature in the steam generator. In this case, the heating element stays disconnected by the thermostat and the control electronics block all the unit functions.

To rearm the thermostat, wait some minutes to let the temperature of the generator fall down to a more normal level, and press the rearm switch.

The cause of the thermostat disconnection is due to the steam generator's lack of water. Check the hydraulic wiring to generate water, tank, pump, hoses, level probe.