## DESTILADOR CABINA CABINET DISTILLER

**L-4 CABINET**

**4903002**

## 1.1. Tabla de contenidos

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## 1.2. General information

- 1) Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please, notify to your supplier immediately.
- 2) Do not install or use the equipment without reading this handbook before.
- 3) This handbook must always be attached to the equipment and it must be available for all users.
- 4) **IMPORTANT! WE WILL NOT ACCEPT ANY APPARATUS TO BE REPAIRED IF IT IS NOT DULY CLEANED**
- 5) If any modification, elimination or lack of maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.
- 6) Do not use the apparatus with liquids which can give off vapours capable of making explosive mixtures.

## 1.3. Packing list

The standard equipment consists of the following components:

Description	Qty.
Heater	2
Filter	2
Level sensor	1
Feed tube	1
Support for water still	1
Set of silicone tubes	1

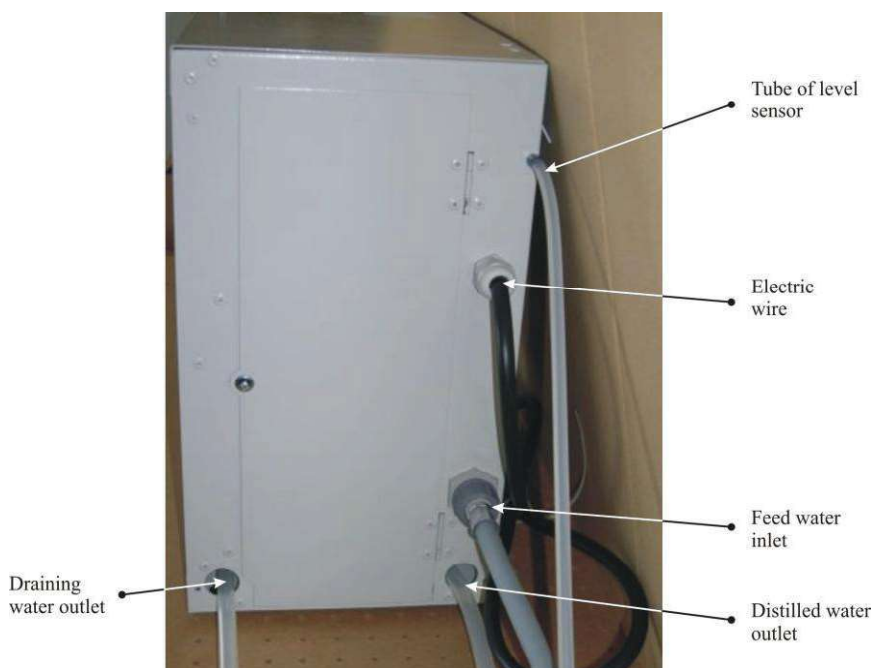
## 1.4. Recommendations

- Electric:
  - Electrical wiring must be, at least, of 4 mm<sup>2</sup> section.
  - It must have ground.
  - Power point must let intensity of 13A pass.
- Outlet:
  - The tap water should have screw mouth of 3/4".

### 1.5. Technical features

Code	Length (cm)	Width (cm)	Height (cm)	Weight (Kg)
L-4 CABINET	55	22	40	12

Code	Output (lts/h)	Conductivity at 20 °C	Water consumption (l/min)	Consumption (W)	Voltage (V)	Intensity (A)
L-4 CABINET	4	2 $\mu$ S / cm	1,4	2 x 1500	230	13

#### ATTENTION

Since distilled water is in contact with the atmosphere before the conductivity measurement can be carried out, its value is actually much less (up to 10 times less) than what it is actually measured. The measurement of any ultrapure water in contact with atmosphere has a conductivity of 1-2  $\mu$ S/cm and a pH of 5, due to a small amount of CO<sub>2</sub> (0.05 ppm) absorbed from the atmosphere.

## 1.6. Assembly

1. Be sure that heater elements (washer of silicone rubber and PBTP nut) are in the order and position that the photo shows.
2. Insert the heaters through the screw mouth.
3. Connect the end wiring of the heaters to the corresponding sockets.
4. Put the water inlet tube on the inlet still.
5. Be sure of putting the filter, of tightening and connecting the other end to the tap.
6. Be sure that the silicone tubes hold suitable position to let water pass and to avoid water pressures.
7. Be sure that the level sensor is connected.

## 1.7. Operating

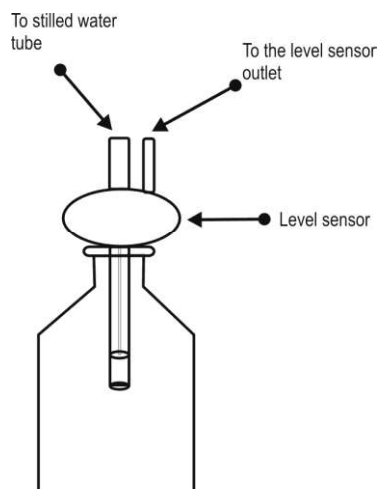
1. Connect the feed water inlet to the tap water.
2. Connect the water still to the mains.
3. Switch on the switch.
4. Turn the tap on slowly. When the water supply is enough, the pilot of “NOT ENOUGH WATER” will turn off and the distillation will start.

## 1.8. Manteinance

- You should clean the still regularly. So, in this way, the results will be better and you can get a still more durable.

## 1.9. Storage of distilled water

- If you wish to fill in a container without worrying when it finished, insert the level sensor in the chosen container.

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**Notice to customers:**

The product is made up of various components and various materials that must be recycled or, failing that, deposited in the corresponding debris removal sites when the product's life has been completed or when otherwise it is necessary to dispose of it. To do this, the end user who acquires the product must know the current regulations of each municipality and / or locality based on the waste electrical and electronic equipment. The user who acquires this product must be aware of and responsible for the potential effects of the components on the environment and human health as a result of the presence of hazardous substances. Never place the product in a conventional container of citizen scope if a previous dismantling and knowledge of the components that incorporates. If you do not know the procedure to follow, consult with the city council for more information.