Electemp Basic

3000896 Regulador de temperatura
Temperature controller

3000896 Termómetro
Thermometer

Indice

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1. General Information

- Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributors immediately.
- Do not install or use the equipment without reading this handbook first. This instruction manual must be attached to the equipment and must be available for all users.
- If you have any doubts or enquiries, contact with J.P. SELECTA's technical service.
- Do not use the equipment in explosive, flammable or corrosive atmospheres.

2. Packing list

Process Controller	3000896
pT100 probe with wire and handle	1000893

3. Technical specification

Model:	3000896 (temp. controller)	3000896 (thermometer)
Power supply:	110VAC 230 VAC 50/60Hz	110VAC 230 VAC 50/60Hz
Electrical power (W):	1700W (110VAC) 2500W (230VAC)	5W
Weight (Kg):	1,1	1,1
External Fuse:	Replaceable (Type: 10x5mm) 15A	Replaceable (Type: 10x5mm) 15A
Timed operation:	Hour, minute	--
Temperature range:	0 to 310°C	0 to 310°C

4. Information about safety in the use of the equipment

Hazard risk

Electric risk

Risk of high temperature contact

4.1 Risks the user is subjected to

- Risk of burns
- Electrical risk

4.2 Staff qualification

This equipment may only be used by properly trained personnel who have read and understand these instructions or have been properly trained in this equipment operation.

4.3 Use with hazardous substances

Do not treat flammable or explosive materials or those which chemically react with violence.

ATTENTION!

Burning risk

Take proper precautions
(use of thermal gloves, clamps,...)
to extract samples from inside the equipment.

It can be at a high temperature.

5. Installation

Unpack the equipment and visually check that the contents are in good conditions. Keep the package for some days.

5.1 Location

Place the equipment on a flat level stable surface, suitable for the equipment weight.


To work in a comfortable way, please leave a free space of 50cm around the equipment and a minimum of 20cm at the rear part.

5.2 Connection to the mains

Plug the equipment to a near socket suitable for the equipment power. Use the connection cable supplied or a similar one.

For your safety, the equipment must be earthed.

Before plugging the machine, check that the voltage supply and the frequency corresponds to the one indicated in the equipment characteristics plate.

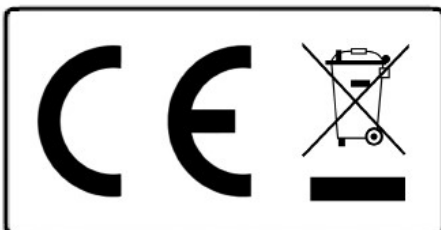Notice to customers:

The product is made up of various components and various materials that must be recycled or, failing that, deposited in the corresponding debris removal sites when the product's life has been completed or when otherwise it is necessary to dispose of it. To do this, the end user who acquires the product must know the current regulations of each municipality and / or locality based on the waste electrical and electronic equipment. The user who acquires this product must be aware of and responsible for the potential effects of the components on the environment and human health as a result of the presence of hazardous substances. Never place the product in a conventional container of citizen scope if a previous dismantling and knowledge of the components that incorporates. If you do not know the procedure to follow, consult with the city council for more information.

6. Operation (temp. controller model)

CONTROL PANEL:

1. Adjustment (change between temperature / time)
2. Increase Value key
3. Reduce Value key
4. Stop button / Start cycle

STARTING UP:

1. Press the main switch (rear) and the screen will light up at the same time.
2. Press keys (2)/(3) until you select the desired temperature value in the green display (the upper one).
3. Press key (1) if you want to select an operating time.
4. Press key (4) to start the cycle.
5. The heating symbol  will light and blink.
6. Press the key (4) again to stop the process. Note: during operation, the temperature or time adjustment value can be changed.

COMMENTS:

- * Do not use the equipment at a lower temperature than ambient +5°C, because the thermostat might have difficulties in regulating correctly.
- * When using the time function, the green display (the upper one) will indicate the remaining time. If you want to see the target temperature value, press key (1).
- * Before pressing the START / STOP key (4), make sure that the temperature probe is connected to the equipment and mounted on the point to be regulated.

6. Operation (thermometer model)

1. Press the main switch (rear) and the screen will light up at the same time.
Make sure that the temperature probe is connected to the equipment and mounted on the point to be measured. The device will show the temperature by means the lower digits at the screen.

7. Error messages (temp. controller)

P.FAIL Power failure. The controller continues with the process.

Err will appear in green on the top line of the display.

SAFE/

TSEG Safety thermostat is open.

SENS Open temperature probe. Control does not detect the temperature probe.

HEAT The heating element does not work. The equipment detects that the temperature does not increase.

8. Settings

The offset (A) and the °C/°F unit (B) can be adjusted.

A. Adjust the probe reading correction (offset)

Press keys (2) and (3) continuously and at the same time (about 5 seconds). The **OFFs** message will appear.

Use keys (2) / (3) to set the actual temperature reading value. Press key (4) to confirm.

B. Presentation unit: °C / °F

Press key (1). The message **Unit** will appear. Using keys (2) / (3), select the unit: °C / °F. Confirm with key (4).

9. Maintenance

Before removing the casing disconnect the equipment from the mains.

The manipulation of the internal electronic circuits of the equipment by unauthorized personnel can cause irreparable damage. Take it to one of the J.P. SELECTA, s.a.u. authorized technical services.

CLEANING:

For the cleaning of the different parts, use the following products:

Cleaning of stainless steel: Alcohol.

Cleaning of plastic: Alcohol with cotton duster.

10. Guarantee

This product is guaranteed for one year. The guarantee does not cover damages caused by incorrect use or causes beyond the control of J.P. SELECTA, s.a.u.

Any manipulation of the equipment by personnel not authorized by J.P. SELECTA, s.a.u. automatically cancels the guarantee.