3000897

Electemp-TFT

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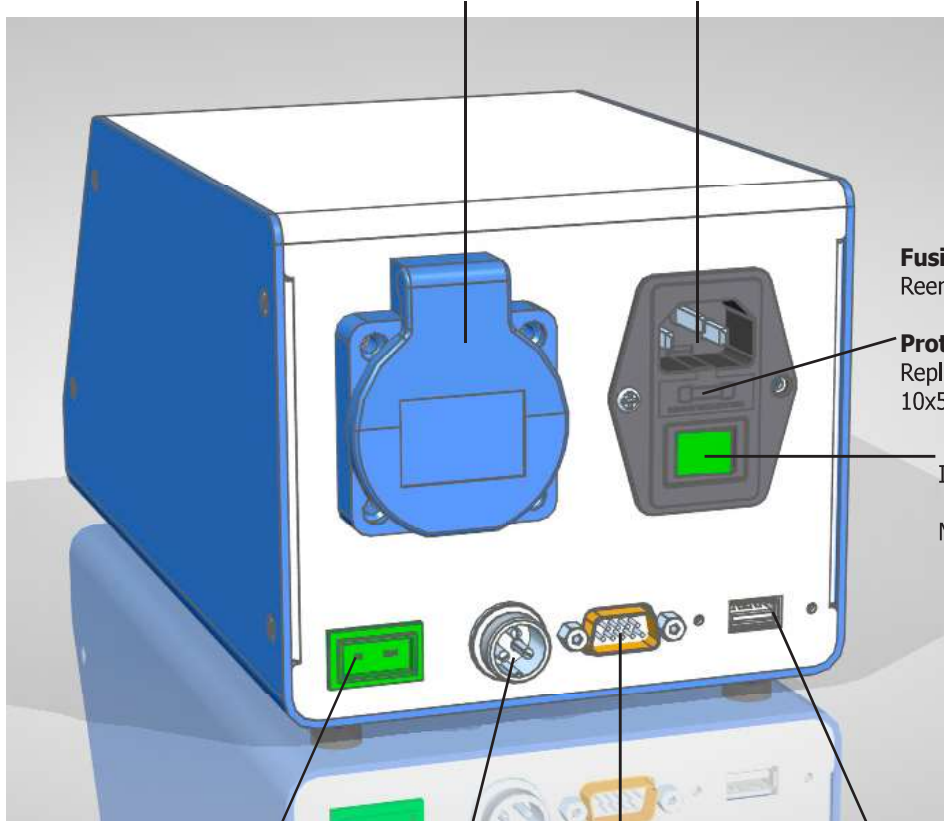
Index

1. General Information	14
2. Packing list	14
3. Accesories	14
4. Technical specification	14
5. Information about safety in the use of the equipment	15
5.1 Risks the operator is subjected to	15
5.2 Staff qualification.....	15
5.3 Use with hazardous substances	15
5.4. Maintenance.....	15
5.5. Warranty.....	15
6. Equipment Description	16
7. Installation	16
7.1 Location	16
7.2 Connection to the mains.....	16
8. Operation	16
8.1 Quick start to work at a temperature.....	18
8.2 SETTING screen	19
8.3 TASK screen	21
8.4 PROGRAMMING screen	22

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Mains plug for the device to be controlled (Heating mantle, bath, ...)
Max 3100W (230V)
Max 1650W (110V)

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Protective fuse
 Replaceable
 10x5 15A

On/Off
 Interruptor principal.
 Main switch.

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K type Thermocouple
 Temperature probe

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Serial connector for data logging during process.

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1. General Information

- Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributors immediately.
- Do not install or use the equipment without reading this handbook first. This instruction manual must be attached to the equipment and must be available for all users.
- If you have any doubts or enquiries, contact with J.P. SELECTA's technical service.
- Do not use the equipment in explosive, flammable or corrosive atmospheres.

2. Packing list

Process Controller **3000897**

3. Accesories

Probe pT100 **1000893**

Probe termocouple K type **1000895**

4. Technical specification

(voltage as indicated on the equipment label)

Model:	3000897	
Power supply:	230 VAC 50/60Hz	110 VAC 50/60Hz
Electrical power:	3100W	1650W
Weight (Kg):	1.25	
External Fuse:	Replaceable (Type: 10x5mm) 15A	
Timed operation:	Weekly, hour, minute	
Temperature Profiles:	Up to 10 profiles of 6 segments. Segment from 1minute to 99hours.	
Data recording:	By USB memory device and Serial RS-232	
External dimensions H x W x D (cm):	10.5 x 13.5 x 21	
Temperature reading range	-150°C to + 1300°C * (depends on the type of probe connected)	
Stability 0-100°C / other ranges:	+ -0,1°C / + -1,0°C	
Resolution:	+ -0,1°C PT100 probe/ + - 1°C K probe	
Working mode:	Cold / heat	

5. Information about safety in the use of the equipment

Hazard risk

Electric risk

Risk of high temperature contact

5.1 Risks the operator is subjected to

- Risk of burns
- Electrical risk

5.2 Staff qualification

This equipment may only be used by properly trained personnel who have read and understand these instructions or have been properly trained in this equipment operation.

5.3 Use with hazardous substances

Do not treat flammable or explosive materials or those which chemically react with violence.

ATTENTION!

Burning risk

Take proper precautions
(use of thermal gloves, clamps,...)
to extract samples from inside the equipment.

It can be at a high temperature.

5.4. Maintenance

Unplug the equipment from the mains before removing the equipment's cover to handle inside.

The manipulation of the equipment internal electronic circuits by unauthorized personnel can cause irreparable damage.

Take it to one of J.P.SELECTA's authorized technical services.

It is recommended to clean the accessible parts of the equipment with a cloth and a cleaner, or a similar solution. Do not use aggressive solvents.

5.5. Warranty

This product is guaranteed for one year. It does not cover damages caused by an incorrect use or causes beyond the control of J. P. SELECTA, s.a.u.

Any manipulation of the equipment by personnel not authorized by J.P. SELECTA, s.a.u. automatically cancels the guarantee.

6. Equipment Description

This equipment is designed to control equipments with heating or compressor and to adjust the heating or cooling power value by means of a temperature probe. To do this, connect the power supply of the equipment to be controlled in the rear base of the ELECTEMP - maximum admissible power 3100W (230V) / 1650W (110V). It will give some electrical pulses to reach the desired temperature value. It is recommended that if the equipment to be controlled has power adjustment, it should be placed in its maximum position. On the other hand, place the temperature probe in contact with the control point. Make sure the probe has good mechanical contact so that the reading is correct.

7. Installation

Unpack the equipment and visually check that the contents are in good conditions. Keep the package for some days.

7.1 Location

Place the equipment on a flat level stable surface, suitable for the equipment weight.

To work in a comfortable way, please leave a free space of 30cm around the equipment and a minimum of 40cm at the rear part. Note that all connections are made on the back of the computer.

7.2 Connection to the mains

Plug the equipment to a near socket suitable for the equipment power. Use the connection cable supplied or a similar one.

For your safety, the equipment must be earthed.

Before plugging the machine, check that the voltage supply and the frequency corresponds to the one indicated in the equipment characteristics plate.

8. Operation

All the equipment functions are set with the touch screen. Gently press with your finger in the center of each icon.

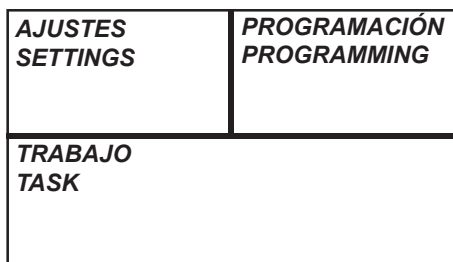
Fig 8.1. Home screen

When starting the computer, the home screen will be shown (Fig 8.1)

- **SETTINGS:**
 - Time and date: menu to set the date and time.
 - PID and Autotuning: menu to set the parameters of temperature control.
 - Downloading recorded data.
 - Temperature correction.
 - Temperature alarm limits.
- **PROGRAMMING:**
 - Allows you to program the temperature profile, by defining temperatures, times and ramps up or down, if you work in cold mode.
- **TASK:**
 - Allows you to select and launch the operating temperature or a temperature profile.

The possibility of giving a heat treatment following a "temperature profile" allows to avoid cracks, breaks, changes of color, etc.

A temperature profile is a succession of "ramps" and "dells".

A ramp is specified by a slope (always positive in heat) of °C/minute.

A dell is specified by a temperature and a time.



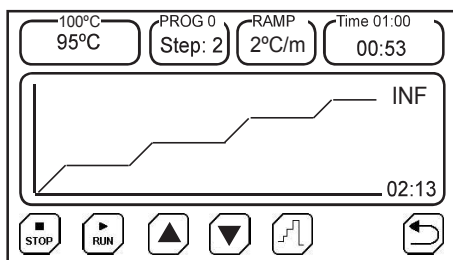
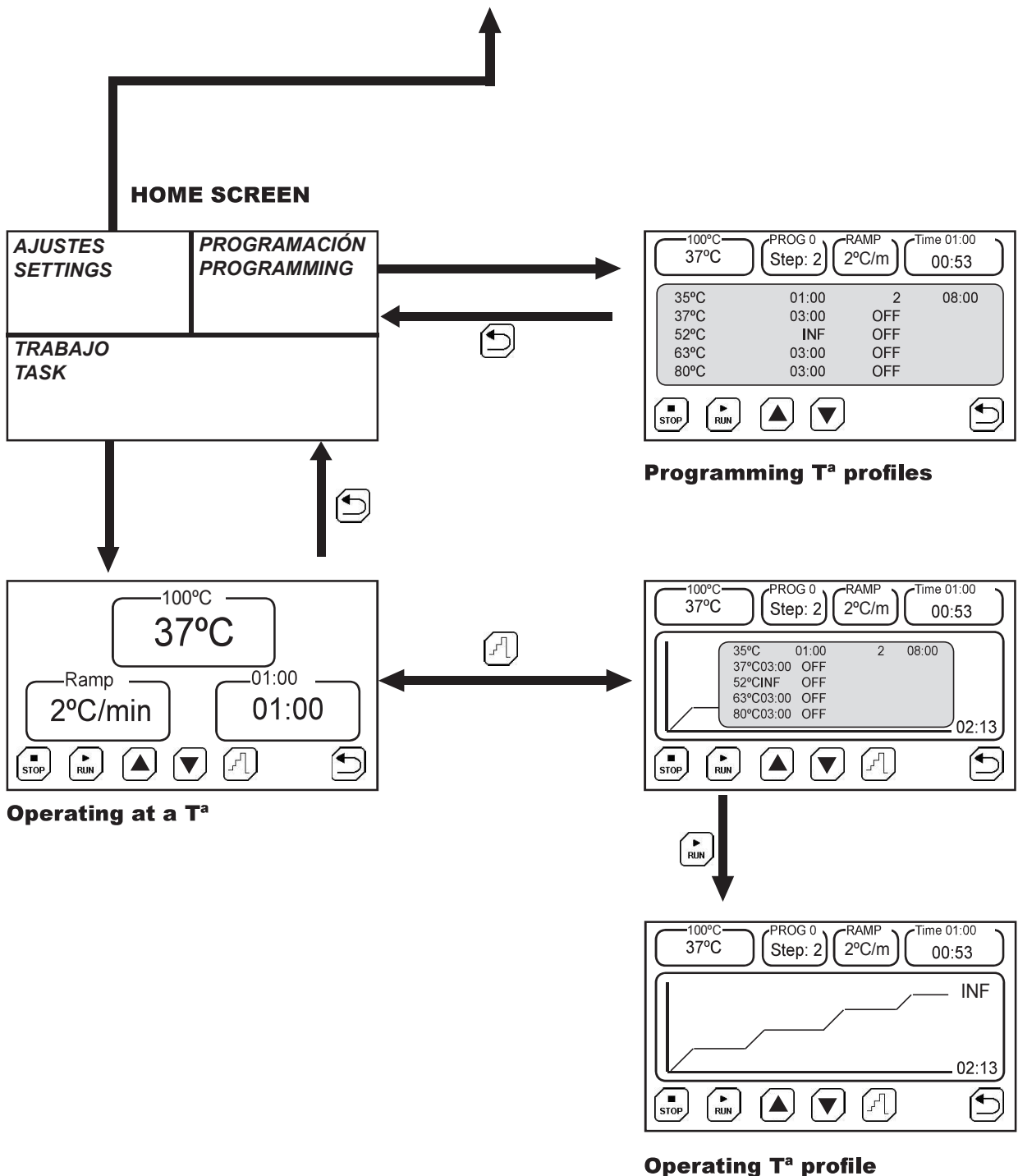
Fig 6. Temperature profile

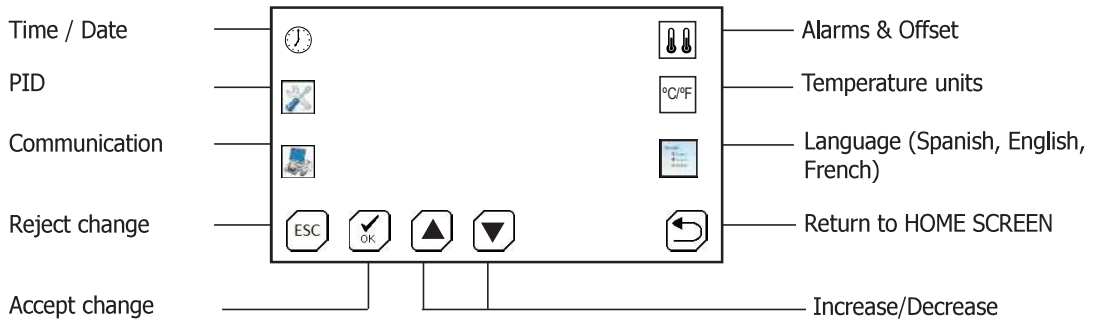
Menus diagram

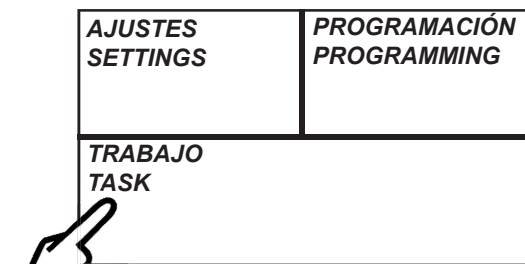
8.1 Quick start to work at a temperature


1 Press on «TASK».

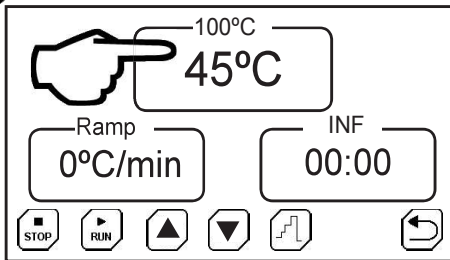
2 Press over the temperature panel.

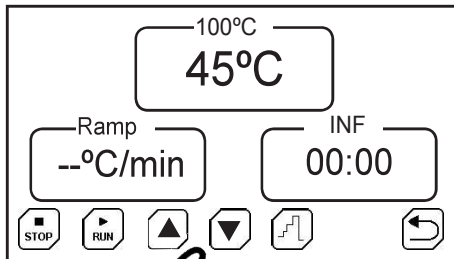3 Press   to set the temperature.

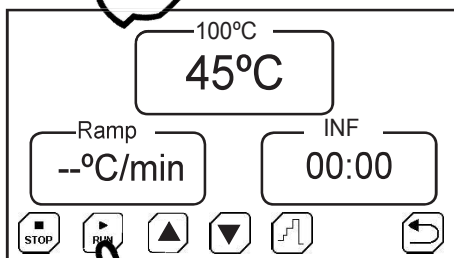
4 Press on  to start.

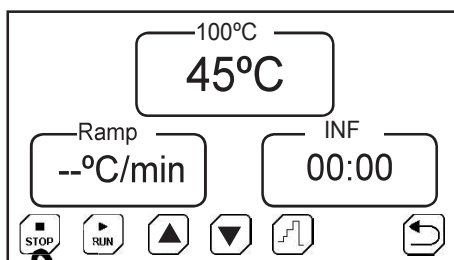5 Press on STOP to stop the cycle.

8.2 SETTING screen

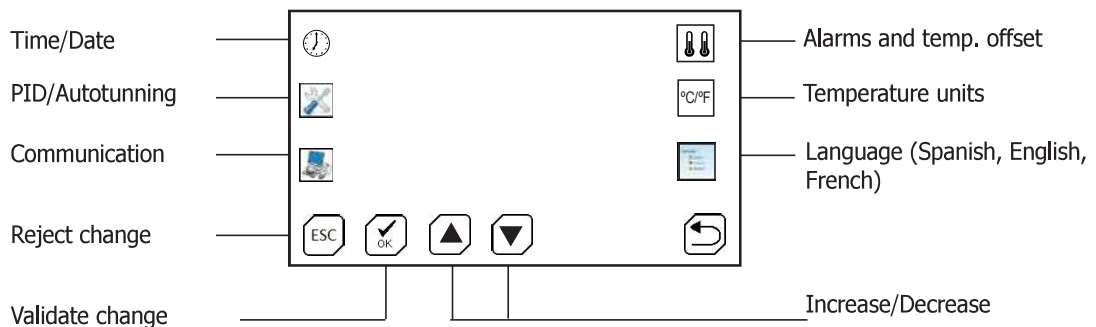Icons meaning:

8.2.1 Set time and date

Press on the time/date icon. Press successively on the box to select the day, month, year, minute and second adjustment.

Adjust by pressing on the icons Increase and Decrease.

8.2.2 Set parameters of temperature control, «Autotuning», Probe type & Ramp units

For an optimal use of the equipment, P, I, D parameters of temperature control should be the appropriate values.

With a value of the appropriate P, I, D parameters, the equipment will reach the temperature in the shortest possible time and without exceeding the set temperature. However, those values can be quite different at high and low temperatures.

The equipment comes with some generic values which are appropriate for most jobs.

These values can be modified, but it's advisable to perform an «Autotuning» when you need a better accuracy.

The «Autotuning» automatically adjusts the optimal parameters for a given temperature and load.

Press on the PID/Autotuning icon. Press successively on the box to select the P, I, D values adjustment and modify the value by pressing on the Increase and Decrease icons.

Attention: It is not recommended to change these parameters. Manually use the «Autotuning» option to automatically adjust these values.

Press on the PID/Autotuning icon.

Select the «Autotuning» temperature with:  
(for example, the temperature most used)

Start «Autotuning» by pressing on the icon: 

If for some reason you want to stop, press on: 

Note: keep in mind that the «Autotuning» process is long because the equipment needs to perform several startup/stop processes to obtain the new PID values.

The «Autotuning» function has two variants: 40 and 100. The 40 variant is faster but less accurate and the 100 variant is slower but more accurate.

- Temperature probe type:
pT100: from -150.0°C to 250°C
K: from 10°C to 1300°C

Note: Connect only one type of probe. It can give erroneous values. When changing the probe, turn off/on the device to ensure a correct reading.

Working mode: heat/cold

HEAT: indicates to the equipment that the system regulates equipments which generate heat.

COLD: indicates to the equipment that the system regulates equipments which generate cold.

8.2.3 Communication and logs type

The equipment can store readings from time to time and send them to a computer or store them in the internal memory and then download them to a USB memory stick.

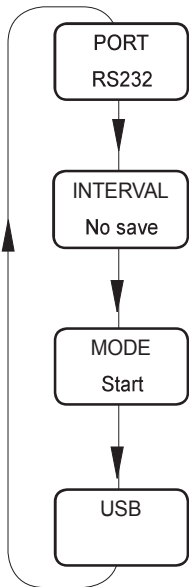
Fig 8.2.3 Communication

«PORT» Set the readings storing mode:

- RS232: Send the temperature readings to a computer by a serial cable (DB9 type).
- USB: Store the readings.

«INTERVAL» Select with   the register time interval and the data sending.

«MODE» Select the data to be stored:

- Start: Save all temperature readings.
- Set point: Save temperature readings, since the set point temperature has been reached.
- PC: Send data, it is not stored.

«USB» Update the equipment internal firmware through a file that can be requested from JP Selecta, s.a.u.

Note: the equipment only saves the last process. When you start a new process, the stored data will be deleted from memory.

If you have the USB mode activated, simply connecting the USB memory at the end of a process will be enough to download the data file. Follow the instructions that appear on the screen.

The generated file is a text file that can be opened with any text editor or spreadsheet.

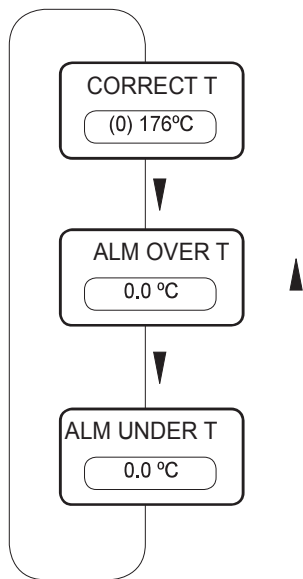
Fig 8.2.4 Temperature alarms

8.2.4 Set the temperature and the temperature alarms

It is recommended to verify calibration of temperature with some regularity (for example, annually), by using for instance a thermometer or a temperature pattern.

If it is necessary to correct the reading, as a result of calibration, one must enter a correction.

To do this, press on this icon:

A box indicating the current correction value appears.

Adjust with keys until you get the indicated value on the temperature pattern.

Applied correction will appear in brackets.

8.2.4.1 Temperature alarms

By pressing successively on the temperature correction box, one may enter an over-temperature and low-temperature limit. Both alarms activate a message on screen, but it does not control the process, it's just an alarm.

8.2.5 Temperature units

Press on the Temperature units icon. Press successively to change between Celsius or Fahrenheit degrees.

8.2.6 Set language

Press on the Language icon. Press successively on the box to select among Spanish, French or English.

Adjust by pressing on the Increase and Decrease icons.

8.3 TASK screen

There are two operation modes:

- Working at a unique temperature.
- Working according to a pre-programmed temperature profile.

Press on to switch between modes.

8.3.1 Operation at a temperature

- Select the set point temperature by pressing on the temperature indicator.
- Select a temperature increase ramp if you want to increase it in a slowly way.
The value "--" makes that the ramp increases as fast as possible (without ramp control)
- Select an operating time. This time starts to count down when the temperature reaches the set point.
- The 0°C value sets a continuous operation.
- Start by pressing on this icon:
The process will end if, the STOP icon is pressed or if it has a programmed time which comes to an end. When finished, the message "Process completed" will appear in a text box. Click on the text box to accept the completion.

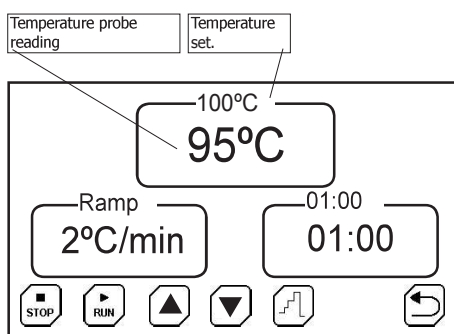
Fig 8.3.1. Run screen operating at one temperature.

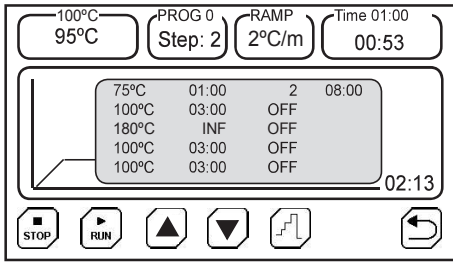
Fig 8.3.2 Run screen operating with temperature profile.


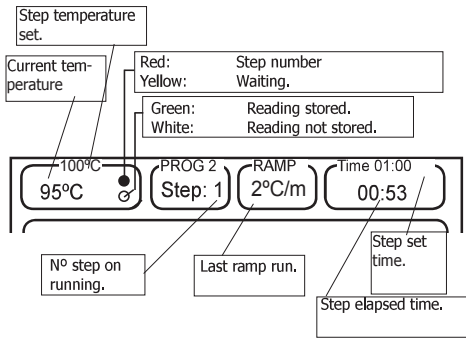
Fig 8.3.2b Working screen.

8.3.2 Operation with a temperature profile

- Temperature profiles must be previously programmed from the «PROGRAMMATION» display.
- Select a profile (from 0 to 9) by pressing on keys
- Start by pressing on the icon:
- In the temperature box (top-left corner) the current temperature and the set point are indicated.

- When RUN is pressed, the profile table disappears and the temperature evolution graph (in red) and the programmed temperature profile (in green) appear.

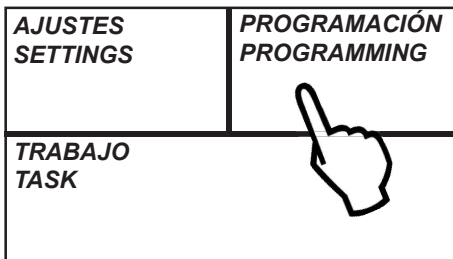The process will end if the STOP icon is pressed or at the end of the program. The message "Process completed" will appear in a text box. Click on the text box to accept the completion.

8.4 PROGRAMMING screen

Go to PROGRAMMING screen to:

- Set the temperature profiles.
- Set the timed operation for a profile.
(For instance, you can set a time an a week day to start a profile. As : start, profile n^{er} 1 on Sunday at 19h00')

Profile no. 1-10	Segment temperatur. °C	Step duration hh:mm	Ramp °C/minute	Profile start
1	75°C	01:00	OFF	--:--
	100°C	03:00	25	Mon
	180°C	02:00	OFF	Tue
	100°C	03:00	OFF	Wed
	100°C	03:00	OFF	Thu
	100°C	03:00	OFF	Fri
	100°C	03:00	OFF	Sat
	600°C	00:30	OFF	Sun

8.4.1 Programming a temperature profile

A temperature profile is a sequence of steps, each step could be a ramp or a dell.

To go from the end of a «dell» to the following «dell» could be done in two ways:

- With a «temperature ramp»:
(go from 100°C to 200°C following a ramp at 2°C per hour).
Value 2°C/h
- As fast as possible: value «--».

Profile parameters input:

Profile no. 1-10	Segment temperatur. °C	Step duration hh:mm	Ramp °C/minute	Profile start
1	75°C	01:00	OFF	--:--
	100°C	03:00	25	Mon
	180°C	02:00	OFF	Tue
	100°C	03:00	OFF	Wed
	100°C	03:00	OFF	Thu
	100°C	03:00	OFF	Fri
	100°C	03:00	OFF	Sat
	600°C	00:30	OFF	Sun

Special values:

- Select the profile number between 0 and 10 with  .
- Press on the parameter to be modified. (Temp, ramp, length ...)
- Press on these icons:   to enter the value.
- Modify another parameter.
- If you enter **OFF** value in the duration of the step, the equipment will end when it reaches this point. The following steps will be deactivated and will not be executed.
- If you enter **INF** in **Step length**, the equipment remains permanently at this temperature.
- If you enter «--» in the **ramp**, the equipment makes the temperature rise at the fastest speed possible.

Profile no. 1-10	Segment temperatur. °C	Step duration hh:mm	Ramp °C/minute	Profile start
1	75°C	01:00	OFF	--:--
	100°C	03:00	25	Mon
	OFF	02:00	OFF	Wed
	100°C	03:00	OFF	Thu
	100°C	03:00	OFF	Fri
	100°C	03:00	OFF	Sat
	600°C	00:30	OFF	Sun

Note: The INF or OFF values are reached by selecting the value 00:00 --> INF --> OFF

Starting of temperature profile

Profile no. 1-10	Segment temperatur. °C	Step duration hh:mm	Ramp °C/minute	Profile start
1	75°C	01:00	OFF	--:--
	100°C	03:00	25	Mon
	180°C	02:00	OFF	Tue
	100°C	03:00	OFF	Wed
	100°C	03:00	OFF	Thu
	100°C	03:00	OFF	Fri
	100°C	03:00	OFF	Sat
	600°C	00:30	OFF	Sun

- To set the profile starting time, press on «--:--»  .
- Enter time by pressing   in **hh:mm** format.
- Press on the days of the week when the profile should be run.
- Accept the programming by pressing on .

Note: if you program a day and a start time, you must press RUN, and the machine will remain in standby mode until the start of the process. This process will be cyclical, that is, each week will be repeated.