MEDILOW XS

2101286

**ARMARIO REFRIGERADOR
COOLING CABINET**

INDICE

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ENGLISH VERSION

GENERAL INFORMATION

1) Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.

2) Do not install or use the equipment without reading this handbook before.

3) This handbook must always be attached to the equipment and it must be available for all users.

4) If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.

5) IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.

6) If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.

7) Do not use the equipment with liquids which can give off vapours capable of making explosive mixtures.

PACKING LIST

The standard equipment consists of the following components:

Piece	Part number	Quantity
Medilow-XS cabinet	2101286	1
Shelves		8
Instructions manual	80366	1

TECHNICAL FEATURES

Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

Code	Cap. litres	Shelves number	Internal dimensions (cm)			Overall dimensions (cm)			Cons. (W)	Weight (Kg)
			Height	Length	Width	Height	Length	Width		
2101286	105	8	146,5	28	27	182,5	36	38,5	110	65

EQUIPMENT DESCRIPTION

MEDILOW XS refrigeration cabinets are designed to work with adjustable temperatures from +2°C to +8°C. They are especially indicated for samples conservation, incubation, cosmetics, botany, food, etc ...

The cabinet is epoxy coating while the inside is molded in ABS, durable and easy to clean.

Inside the cabinet there is a turbine for air circulation, and a hermetic and vibration-free compressor group.

Front foot supports are adjustable in height to achieve a perfect leveling.

TRANSPORTATION & STORAGE

Keep the unit upright when moving, the angle to the horizontal line must not exceed 50°, otherwise, it will damage the compressor and affect the normal operation of the business.

The unit should be placed on a solid and firm floor, remove the lower part of the package to reduce the impact and noise. Allow at least 20cm distance between the back of the unit and the wall for a good ventilation and leave a space with an open angle of 160° for the door.

Keep the unit away from heat sources and avoid direct sunlight.

Never place the unit in a place where water can reach it or which be very humid. Splashed water and dirt should be cleaned quickly using a soft cloth. Do not place the unit in an environment cold enough to freeze.

Adjust height with the four adjustable feet.

CAUTION!!! IMPORTANT FOR YOUR SAFETY

Be sure that the voltage supply is the same as the one indicated on the characteristic plate of the equipment.

Do not use the equipment if it is NOT earthed.

If you have to change the plug bear in mind the following:

Blue cable: Neutral.

Brown cable: Phase.

Yellow/green cable: Earth.

PREPARATION

Remove the box and the packaging foam (including the bottom foam) before using the equipment. Clean the unit before use. Adjust the feet so that the unit stabilizes according to the ground conditions to reduce noise.

After moving the unit, let it stand for about 30 minutes before turning it on.

Running: close the door and turn on the power, turn on the light switch at the back. The temperature should be adequate for normal use after half an hour. Once the power is turned on, the unit must remain empty for one hour.

CONFIGURATION & TEMPERATURE SETTINGS

Adjust the storage temperature through the temperature controller located at the back of the unit. Adjust the temperature according to the ambient temperature and the product to be stored.

The equipment has an adjustable control for the temperature, which ranges from 1 to 7. Set the storage temperature by turning the temperature knob clockwise. 7 shows the lowest storage temperature. The knob can rotate freely but cannot be turned clockwise from 7 to 1. The unit is factory set to 3.

Check if the knob is in the proper position before using the equipment. Frequent setting is not appropriate when using the equipment. Normally, in summer, the ambient temperature is high, so we will turn the knob counterclockwise to 7; while in winter the ambient temperature is low, therefore we will turn the control to the right until 1.

0 position means stop the equipment.

PRODUCTS STORAGE

Do not put too many products at the same time in the equipment, you should leave some space around the products so that cold air can flow into the chamber. Keep bottled items away from the inside walls of the chamber so they do not freeze as they might break.

Shelves can be adjusted according to different needs.

Do not place hot products directly on the equipment until it cools down to normal temperature, to avoid reducing the power consumption and prolonging life of the compressor.

MAINTENANCE

Disconnect power supply before performing maintenance on the machine.

Keep the door seal clean, remove it from the door and wash it. Dry and seal as before. Put some talcum powder on the magnetic surface to prolong its life.

If there is a lot of ice in the evaporator after prolonged use, remove things from inside the chamber, turn the power off, and automatically the frost will melt. Clean the water and turn the power on again.

To clean the unit, use a soft cloth and wipe it with a little soap and water (or a neutral, non-caustic detergent). Dry it after cleaning it to avoid oxidation.

Do not use organic liquids, reservoir water, detergent powder or acid to clean the unit. Do not use a hard or steel brush to clean the equipment.

If you are not going to use the equipment for an long period of time, disconnect it from the power supply, clean it by following the steps above, and dry the water inside the camera with a soft cloth.

TROUBLESHOOTING

Problem	Possible cause	Solution
The equipment does not work	Plug's bad connection	Unplug it and reconnect it to the power supply
Abnormal compressor operation or buzz	Supply voltage is outside the nominal range	Disconnect the power source immediately and reconnect it when the voltage is normal. It is necessary to have a voltage stabilizer in case of poor power.
The compressor has been running for too long and there is no frost on the evaporator's surface.	The cooling system is damaged (leaking or blocked)	Send the unit to the technical service for repair.
There is frost or ice on the camera's internal walls and it has a too low internal temperature. The compressor always runs, it never stops.	The thermostat does not work	Take it to the technical service to repair.
The internal temperature is too high and the compressor runs, it never stops.	Bad heat dissipation and ventilation of the housing, the room temperature is too high.	Improve ventilation and lower ambient temperature.
	Door frequently open and for a long time.	Decrease time and number of times of door openings.
Too much noise	Unit is not level	Adjust the adjustable feet
	Pipes touch the compressor.	Carefully separate the pipes that touch each other.
The light is not on	The light has melted	Replace it

Warning:

- Keep the ventilation openings unobstructed, inside the appliance or in the built-in structure.
- Do not use mechanical devices or other means to accelerate the defrosting process.
- Do not use electrical appliances inside the storage compartments of the equipment.
- It is forbidden to disassemble the unit by non-professional personnel, it can be dangerous.

PRECAUTIONS

This equipment is not suitable for blood storage neither for pharmaceutical and biological products.

Do not place heavy items on top of the unit.

In case of problems, request assistance from a professional, do not disassemble the unit without prior authorization to avoid risks.

ORDINARY PHENOMENA

If you hear a slight sound of water flowing in the unit, it is the sound of the coolant flowing inside the pipe.

The compressor, the condenser and part of the unit near the door might be hot or too hot.

In a place with humidity, frost may be formed on the housing or on the door. Dry it with a dry cloth.

GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damages caused by an incorrect use or causes beyond the control of J.P. SELECTA, s.a.u.

Any manipulation of the equipment by unauthorized personnel by J.P. SELECTA, s.a.u. cancels the guarantee automatically.

Notice to customers:

The product is made up of various components and various materials that must be recycled or, failing that, deposited in the corresponding debris removal sites when the product's life has been completed or when otherwise it is necessary to dispose of it. To do this, the end user who acquires the product must know the current regulations of each municipality and / or locality based on the waste electrical and electronic equipment. The user who acquires this product must be aware of and responsible for the potential effects of the components on the environment and human health as a result of the presence of hazardous substances. Never place the product in a conventional container of citizen scope if a previous dismantling and knowledge of the components that incorporates. If you do not know the procedure to follow, consult with the city council for more information.