**ARMARIOS REFRIGERADORES
COOLING CABINETS**

**MEDILOW-L
MEDILOW-LG**

**2101301
2101302**

INDICE

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INDEX

GENERAL INFORMATION	10
PACKING LIST	10
ACCESSORIES	11
TECHNICAL FEATURES	11
EQUIPMENT DESCRIPTION	11
INSTALLATION	12
OPERATION	12
MAINTENANCE	15
SPARE PARTS	16
GUARANTEE	16

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Notice to customers:

The product is made up of various components and various materials that must be recycled or, failing that, deposited in the corresponding debris removal sites when the product's life has been completed or when otherwise it is necessary to dispose of it. To do this, the end user who acquires the product must know the current regulations of each municipality and / or locality based on the waste electrical and electronic equipment. The user who acquires this product must be aware of and responsible for the potential effects of the components on the environment and human health as a result of the presence of hazardous substances. Never place the product in a conventional container of citizen scope if a previous dismantling and knowledge of the components that incorporates. If you do not know the procedure to follow, consult with the city council for more information.

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GENERAL INFORMATION

- 1) Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.
- 2) Do not install or use the equipment without reading this handbook before.
- 3) This handbook must always be attached to the equipment and it must be available for all users.
- 4) If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 2009/104/CE, the manufacturer is not responsible for the damage that can occur.
- 7) Do not use the equipment with liquids which can give off vapours capable of making explosive mixtures.

PACKING LIST

The standard equipment consists of the following components:

Piece	Code	Quantity
Cooling cabinet Medilow-L	2101301	1
Shelves		3
Cooling cabinet Medilow-LG	2101302	1
Shelves		4
Instruction manual	80432	1

ACCESSORIES

OXIGEN BIOLOGICAL DEMAND BOD_{5D} Code: 5906000
Manometrical method, without sample dissolution.

Code	Overall dimensions cm			Consump. W.	Weight Kg
	Height	Length	Width		
5906000	32	42	23	25	8

TECHNICAL FEATURES

Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

Code	Cap. litres	No. shelves	Internal dimensions cm			Overall dimensions cm			Con-sump. W.	Weight Kg.
			Height	Length	Width	Height	Length	Width		
2101301	610	3	145	60	56	191	68	80	330	73
2101302	544	4	145	60	56	164	75	95	330	82

* Maximum weight per shelf, 50Kg (evenly distributed)

EQUIPMENT DESCRIPTION

MEDILOW cooling cabinets are designed to work with adjustable temperatures from +1°C up to +40°C, with a stability of $\pm 1,5^\circ\text{C}$. They are specially appropriated for BOD_{5D} determination at 20°C, enzymatic measurement at 25°C, perservation of sample, incubator, cosmetics, botanics, food, etc.

Model 2101301 (MEDILOW-L) is made of easy-to-clean AISI -304 stainless steel, inside and outside.

Both models incorporate an internal turbine for air circulating, and also an hermetic compressor group without vibrations. Side hole for connections, probes, etc.

Height-adjustable front base supports. Lockable door, and it can be opened from left or right.

Note: model 2101302 (MEDILOW-LG) has the same features as model 2101301 but with glass door, external cabinet covered in epoxy and inside molded in ABS, resistant and easy to clean.

INSTALLATION

Place the equipment on a flat, horizontal and level surface, trying to leave a space of about 10 cms. at the back and on each side.

Avoid positioning the equipment in direct sunlight or exposed to similar sources of heat. Poorly ventilated rooms are also unsuitable.

The ventilation grille at the back of the equipment must NOT be blocked not even change its position.

CAUTION!!! IMPORTANT FOR YOUR SAFETY


Be sure that the voltage supply is the same as the one indicated on the characteristic plate of the equipment.

Do not use the equipment if it is NOT earthed.

If you have to change the plug bear in mind the following:

Blue cable: Neutral.
Brown cable: Phase.
Yellow/green cable: Earth.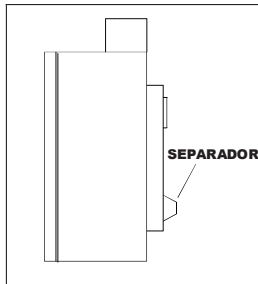


VERY IMPORTANT BEFORE THE FIRST START UP:


Place the buffers at the rear part of the equipment to avoid contact with the wall (see drawing).

OPERATION

CONTROL PANEL: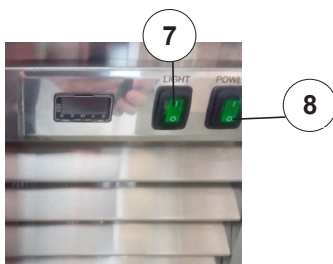

In case of int connection (plugs), activate the circuit breaker located in the rear part

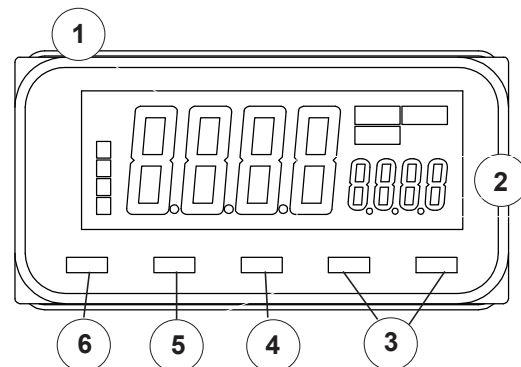1. Screen 1
2. Screen 2
3. Up/Down keys
4. Digit change key
5. Parameter change key
6. Parameter input key
7. Light switch
8. Main POWER switch

START UP:

- Connect to the mains.
- Press POWER (8)
- Select the temperature value with the UP & DOWN keys (3) and in display 2 (green) the value flashes. The value is automatically saved.

part number: 0013014

TEMPERATURE CORRECTION:

You can correct the temperature displayed on the screen if contrasted with a more accurate thermometer.

- Press the parameter input key to enter the setup menu.
- Press the parameter change key until the message iNS appears.
- Adjust the temperature value with the Up / Down keys.

Example 1:

If the standard thermometer indicates 10°C while the MEDILOW unit temperature control indicates 9 °C. The value to be entered is 1.

Example 2:

If the standard thermometer indicates 10°C while the MEDILOW unit temperature control indicates 11°C. The value to be entered is -1.

Note: For more details on regulator operation, use additional specific instructions for it.

MAINTENANCE

Before removing the covers disconnect the equipment from the mains.

The parameters of temperature regulator must not be manipulated by unauthorized personnel. It can cause damages of difficult solution.

When its moving is necessary, put the set vertically on a palet whose dimensions (at least) are the same as the ones of the base of the cabinet. Once the moving is finished, 12 hours must be passed before starting up the set.

CLEANING:

The cleaning of the cabinet freezer frequently is recommended. For it, disconnect the equipment from the power. Clean the chamber and casing with lukewarm water with a small quantity of detergent, rinse with clear water and dry well. **Do not use abrasive, alkaline, acid or base cleaning products.**


DEFROSTING:

The equipment is equipped with an automatic defrost system. However, when operating with the regulator in the coldest settings, frost may occur on the inner bottom wall. In this case, the equipment must be defrosted manually. • To defrost it, disconnect the equipment by removing the plug from the socket!
• Remove the products from inside the unit and store them in a cool place. • During the defrosting process, leave the door open.

Collect the defrost water with an absorbent wipe and clean the equipment. Make sure that the water can circulate freely through the drain located at the back of the cooling chamber. If necessary, disengage the drain with a pointed object. **To defrost, never use mechanical devices or other means, as it can deteriorate the cabinet!**

CHANGING THE SIDE (HINGES) OF THE DOOR: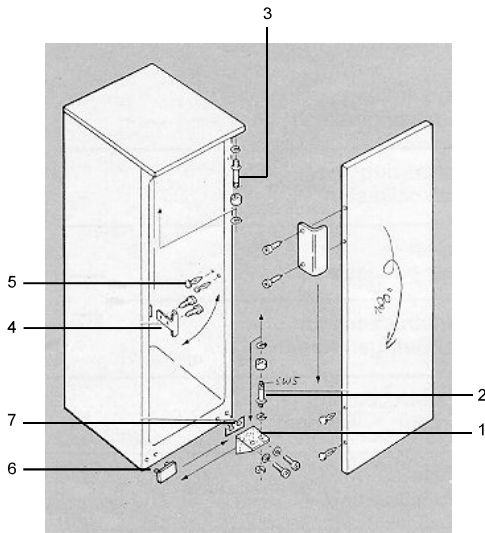

1. Unscrew the edge of the hinge (1) with the door closed. Remove the door downwards.
2. Remove the hinge-bolt (2) from the edge of the hinge and screw it into the next support.
3. Unscrew the upper bolt of the hinge (3) and screw it to the opposite side.
4. Separate the covering parts (4, 5 and 6) from the left side of the equipment and cover the support holes left open on the right.
5. Hang the door from the hinge-bolt (3) and close it.
6. Screw up the edge of the hinge (1 and 7) onto the left side of the equipment (don't forget the small plates under the hinges).

SPARE PARTS

To guarantee the safety of the equipment, the spare parts must be purchased from J.P.SELECTA, s.a.u.

Description	Code	
	2101301	2101302
Fan	24055	
Heating element 150W 220V	39125	
Temperature regulator	16131	
Pt100 probe	43071	

GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damages caused by an incorrect use or causes beyond the control of J.P. SELECTA, s.a.u.

Any manipulation of the equipment by unauthorized personnel by J.P. SELECTA, s.a.u. cancels the guarantee automatically.