**REFRICUB**

**4100519**

**CUBETA REFRIGERADORA  
COOLING CABINET**

**INFORMACIÓN GENERAL**

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**GENERAL INFORMATION**

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please notify to your J.P. Selecta supplier immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubt or enquiries, please contact with your supplier or J.P. Selecta's technical service.*
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY APPARATUS TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) *If any modification, elimination or lack of maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the apparatus with liquids which can give off vapours capable of making explosive mixtures.*

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**CONTENTS**

GENERAL INFORMATION
CONTENTS
PACKING LISTS
TECHNICAL FEATURES
EQUIPMENT DESCRIPTION
INSTALLATION
OPERATION
MAINTENANCE
PROGRAMMING
ELECTRIC DIAGRAM
COOLER CIRCUIT DIAGRAM
GUARANTEE
"EC" CONFORMITY DECLARATION

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*The standard equipment consist of the following components:*

DESCRIPTION <i>DESCRIPTION</i>	CODIGO <i>CODE</i>	CANTIDAD <i>QUANTITY</i>
Refricub	4100519	1
Manual de instrucciones/ <i>instruction manual</i>	80086	1

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*Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.*

Capacidad (litres)/ <i>Capacity (litres)</i>		18
Medidas útiles <i>Tank dimensions (cm)</i>	Alto/ <i>Height</i>	20
	Ancho/ <i>Length</i>	30
	Fondo/ <i>Width</i>	30
Medidas exteriores <i>Overall dimensions (cm)</i>	Alto/ <i>Height</i>	88
	Ancho/ <i>Length</i>	50
	Fondo/ <i>Width</i>	48
Potencia (CV)/ <i>Power (HP)</i>		3/8
Consumo (W)/ <i>Consumption (W)</i>		333
Peso (kg)/ <i>Weight (kg)</i>		64

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*The cooling tank REFRICUB is especially designed to keep liquids and objects at low temperature.*

*REFRICUB allows work and control temperatures until -30°C, by means of an electronic controller and digital reading.*

*Outer casing in epoxy and inner tank and lid in AISI 304 stainless steel. Hermetically sealed compressor is installed on an anti-vibration mounting with ventilated evaporator, regulating valve and condenser.*

*The control panel is equipped with main switch with indicator lamp, analogue electronic temperature regulator and digital reading with NTC probe and resolution 1 degree.*

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## INSTALLATION

Place the apparatus on a flat horizontal, level surface, leaving a space of about 10 cm at the back and on each sides of the apparatus.

Fill inner tank with approximately 8 litres of methanol. Methanol is not supplied with equipment so it must be acquired from a chemical supplier

To fill this tank you must do it in following way:

- Open the screw top placed on top part using an allen key nummer 6.
- Fill the tank through the hole.
- Put on the screwed top again using the same key.

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*Do not use the apparatus with inflammable liquids which can give off vapours capable of making explosive mixtures or in special protection zones.*

*Be sure that the voltage supply is the same as the one indicated on the characteristics plate.*

*Do not use the apparatus if it is not earthed.*

*If you change the plug be in mind the following:*

**Blue cable: Neutral.**  
**Brown cable: Phase.**  
**Yellow/green cable: Earth.**

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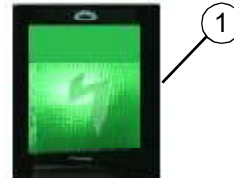
## OPERATION

### CONTROL PANEL:

1. Main switch with indicator lamp.
2. Temperature reading display.
3. Temperature selection buttons.

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### FUNCIONES DEL FRONTAL:

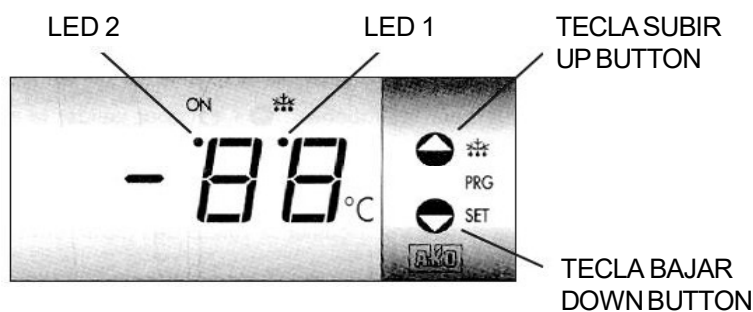
### STARTING UP:

1. Connect the apparatus into the mains.
2. Press the main switch (1).
3. Set the temperature.

Before introducing the first charge, it is advisable to leave the chest working for about two hours at the pre-set temperature.

NOTE: The apparatus do not start up until a minute later user switch on the main switch.

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*By pressing for 5 seconds, a manual defrosting will start. REFRICUB does not have this function enabled.*

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Pushing for a 5 seconds, present set point is shown.

Pressing  and  simultaneously users can fix the new value

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## MAINTENANCE

*Before removing the casing disconnect the apparatus from the mains.*

*The manipulation of the internal electronic circuits of the apparatus by unqualified personnel can cause irreparable damage. Be sure that apparatus is repaired by a J.P.SELECTA authorized technical service.*

### CLEANING:

*For the cleaning of the different parts use the following products:*

*Cleaning of stainless steel: Alcohol*

*Cleaning of plastic: Alcohol with cotton duster.*

*From time to time clean the condenser wings of the cooling circuit using a wet cloth.*

*CAPACITY of the internal tank: 6 litres of methilic alcohol. (Not supplied)*

COMPRESSOR: 3/8 CV.  
Gas 404A. 300 gr

## PROGRAMMING

**VERY IMPORTANT:** Parameters only should be programmed or modified by personnel who know REFRICUB limitations.

These parameters are programmed in factory. If you need change them you would better contact to J.P. SELECTA, s.a.

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