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
1. Introduction

- 1.01 Electricity is very dangerous, please pay attention to safety, so this equipment must be installed by a professional person.
- 1.02 In order to prevent possible electric shock, THIS EQUIPMENT MUST BE GROUNDED / EARTHED.
- 1.03 Must double check the voltage and frequency of electricity and make sure it is suitable for the chiller.
- 1.04 Must place the chiller in a well-ventilated area indoor, and ambient temperature shall be lower 40°C.
- 1.05 Keep all power cords and hoses away the heating and moving parts.
- 1.06 Feeding the chiller with purified water/glycol is necessary for the chiller, the deionized water and distilled water are the best choose, don't feed the seawater for normal chiller.
- 1.07 If the target temperature less than 5°C is required, please DON'T FEED THE CHILLER WITH WATER, feed it with GLYCOL ONLY! Otherwise, the water will be iced and bust the circulating loop and damage the chiller.
- 1.08 Must empty the water inside the chiller for storage, otherwise, the water will be iced and bust the circulating loop and damage the chiller if the ambient temperature is too low.
- 1.09 Regular maintenance is necessary.
- 1.10 J.P. Selecta s.a.u. can't be responsible for anything that is caused by poor installation and use.

2. Working schematic of chiller

According to this working diagram, please install the chiller to your system first.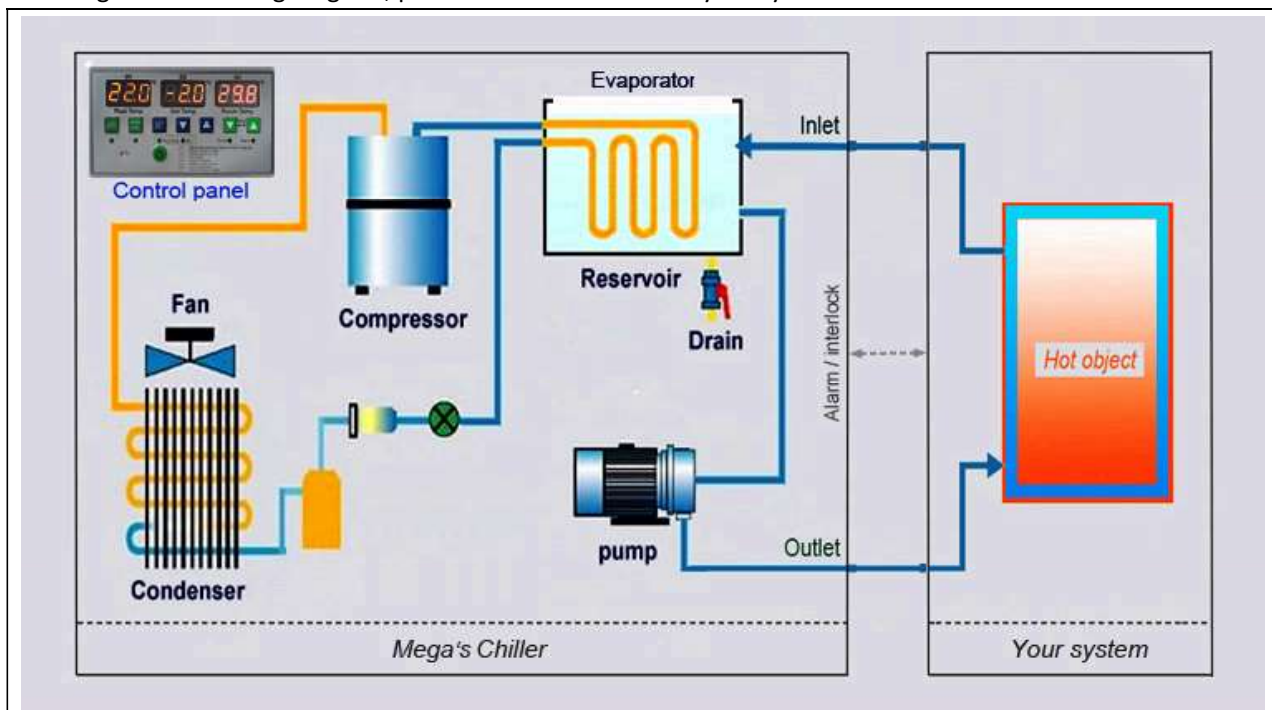

REMINING:

- * If the hot object is a closed loop, please use the open loop chillers like our standard items.
- * If the hot object is an open loop, please use the closed loop chillers or open loop chiller and extra heat exchanger.

3. Product specifications

MODEL	CH-1100	CH-1400
PART NUMBER	6001492	6001493
POWER SUPPLY	220-240V/50Hz	220-240V/50Hz
INPUT POWER	560W	610W
COMPRESSOR POWER	430W	480W
COOLING CAPACITY	1100W	1450W
REFRIGERANT	R134A	R134A
WATER TEMPERATURE RANGE	+5..+35°C	+5..+35°C
TEMPERATURE PRECISION	±1°C	±1°C
TANK VOLUME	6,7L.	6,7L.
PUMP FLOW	11L/min	19L/min
PUMP LIFT	8m	12m
COOLING MODE	Air Forced Cooling	Air Forced Cooling
INLET/OUTLET Ø	Ø ext: 11mm	Ø ext: 11mm
DIMENSIONS	28,6(W)x57(L)x42,2(H)cm	28,6(W)x57(L)x42,2(H)cm
CUSTOMIZATION	Available	Available
PROTECTIVE DEVICES	Compressor overheating, over current, high & low pressure, overtemperature, water flow & level, antifreezing	Compressor overheating, over current, high & low pressure, overtemperature, water flow & level, antifreezing

MODEL	CH-4200	CH-8200
PART NUMBER	6001494	6001496
POWER SUPPLY	220-240V/50Hz	3PH-380V, 50Hz
INPUT POWER	1235W	3570W
COMPRESSOR POWER	980W	2600W
COOLING CAPACITY	4200W	8200W
REFRIGERANT	R410a	R410a
WATER TEMPERATURE RANGE	5-35°C	5-35°C
TEMPERATURE PRECISION	±1°C	±1°C
TANK VOLUME	26L.	46L
PUMP FLOW	28L/min	50L/min
PUMP LIFT	40m	26m
COOLING MODE	Air Forced Cooling	Air Forced Cooling
INLET/OUTLET Ø	Ø ext: 19mm	Ø ext: 19mm
DIMENSIONS	48(W)x55(L)x89(H)cm	61(W)x75(L)x90(H)cm
CUSTOMIZATION	Available	Available
PROTECTIVE DEVICES	Compressor overheating, over current, high & low pressure, overtemperature, water flow & level, antifreezing	Compressor overheating, over current, high & low pressure, overtemperature, water flow & level, antifreezing

3.1 Packaging Contents:

- Chiller according reference
- Silicone hose Ø14x8mm (5mts). Cod. 46041 for models 6001492 & 6001493
- Reinforce silicone hose Ø15x24. Cod.46156 for models 6001494 & 6001496
- Accessories kit (only model 6001496): code 21281 (4pcs); code 34551 (2pcs); code 34560 (2pcs); code 34594 (2pcs)
- Cooling fluid code 100028. Models 6001492 & 6001493 (2x5L); 6001494 (5x5L); 6001496 (10x5L)
- Instruction Manual M.80433.00
- EU Declaration of Conformity

4. Control touch panel

4.1 D1 digital display (Real Temp):

- *It displays the real water temperature.
- *Or displays the setting item(s) C# under the setting mode.

4.2 D2 digital display (Set Temp):

- *It displays the setting value under setting mode.
- *Or displays the error code during alarm.

4.3 D3 digital display (Room temperature):

- *It displays the ambient/room temperature.

4.4 AUTO button (Smart mode):

- *Press once to enter the AUTO mode.
- *Press it again to display the room temperature, and automatically return to the display real water temperature after 3 seconds.

4.5 CONSTANT TEMP button (constant temperature mode):

- *Press once to enter the CONSTANT TEMP mode.
- *Press it again to display the room temperature, and automatically return to display the real water temperature after 3 seconds.

4.6 SET (Settings):

- *Press SET button to enter the setting mode,
- *Press it again or no operation 10 seconds later to save the setting and exit the setting mode.
- * **✘ RESRT**: Press and hold the SET button for 4 seconds to reset the chiller. When system setting is too mess or crash that reset the system is need, the parameter values of default settings: C0:15 C1:1.5 C2:60 C3:1 C4:20 C5:30 C6:3 C7:50.

4.7 ▼ & ▲ (UP / DOWN) button on display D3:

- * All parameter C0-C8 can be selected to be modified.

4.8 ▼ & ▲ (UP / DOWN) button on display D2:

- *Click the up and down button to modify setting values from C0 to C8 under setting mode.

5. Working mode selection and target temperature setting

5.1 Under AUTO mode:

This is the intelligent temperature control mode. It is just for keeping the water/glycol temperature close to the ambient temperature. The Set temperature D2 represents the temperature deviation between the target water temperature and the ambient temperature, so the target water/glycol temperature would be the ambient temp/Room temperature D3 plus the Set temp/temp deviation. That is, the Set temporary value in -2 degrees and ambient temperature 22 degrees, so the target water/glycol temperature is in 20 (22-2) degrees.

Ambient temp/Room Temp on D3 (Degrees Celsius)	T0/D3	Take 22°C example
Setting temperature/Set Temp on D2 (Degrees Celsius)	T1	Take -2°C example
Target water temperature/Real Temp on D1 (Degrees Celsius)	$T=T0+T1$	20°C

How to set the temperature under AUTO mode?

According to the Set temperature, D2 displays the temperature deviation between the target water temperature and the ambient temperature.

So calculate the temporary deviation first after switch on the chiller---click the AUTO button to enter AUTO mode---Click or hold the UP and DOWN button to set the value of temperature deviation you calculated—OK.


(Refer to 5.3 if the SET Temp cannot be increased or reduced)

5.2 Under CONSTANT mode:

D2 displays the real-time water temperature, directly press the UP / DOWN button on D2 display to set the target temperature as your needs under CONSTANT mode.

How to set the temperature under CONSTANT mode?

For example, how to set the target temperature in +22 degrees: Switch on the chiller- Click the CONSTANT button to enter the CONSTANT mode- Click or hold the UP or DOWN button to get the target water temp as your needs.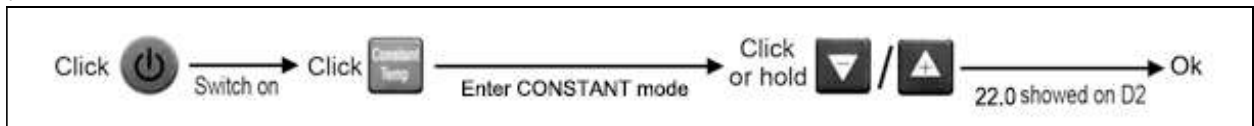


(Refer to 5.3 if the SET Temp cannot be increased or reduced)

5.3 How to unlock the temperature limitation?

Under any modes, if the Set temperature can't be reduced, please reduce the value of minimum of water temperature C4, or if the Set temperature can't be increased, please increase the value of maximum of water temperature C5 as below:

Click the SET button to enter the SET mode -click the UP or DOWN button to choose setting item C4/C5 at the **D3 display**. Click the UP/DOWN button to get the value of C4/C5 showed on **D2 display**, the value of C4 must be no higher than the target temp or the value of C5 what must be no lower than the target temp you expected- click the SET button to save the setting or wait for 10 sec to autosave the setting and return: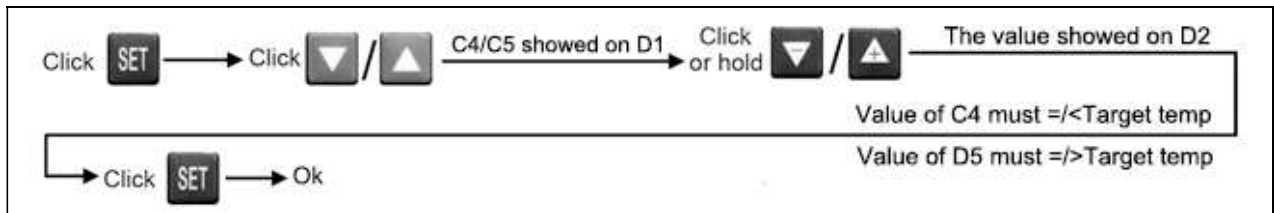

And then go back to 5.1 or 5.2 to set the temperature again.

6. Setting items and parameters


- C0: For setting the alarm of temp difference between the water temp and setting temp (5-15 degree), default (±)15 degree.
- C1: For setting the deviation of water temp (±0.1-2.0 degree), default ±1.5 degree.
- C2: For setting the alarm of customized temperature (1-100 degree) - Optional for extra temperature detection, default 60 degree.
- C3: For setting the delay time of alarm for lack of the water flow (0-100s), default 1 second, 0 for switch off, 1-100 is the delay time in second.
- C4: For setting and limiting the minimum of water temperature (1-25 degree), default 20 degree
- C5: For setting and limiting the maximum of water temperature (25-50 degree), default 30 degree
- C6: For setting the heating target temperature (1-20 degree)- Optional, default 3 degree
- C7: For setting the alarm of ambient temperature (42-50 degrees), default 50 degree.
- C8: For setting the boot mode, default 1 for powering on the chiller by the ON/OFF button, 0 for powering on the chiller by main switch.

REMINDING:

- *The temperature display range is in -20°C to 99.9°C, but it is invalid during setting temperature with buzzer three times changing the parameter value is required to unlock the limitation first if need, such as:
 - Reduce the value of C4 first if the temperature can't be reduced when setting the temperature.
 - Increase the value of C5 first if the temperature can't be increased when setting the temperature.
 - Modify any setting item(s) to solve the problem/error as below.

*How to modify the parameter value of setting items C#.

Click SET button- Click the UP/DOWN button **on D3 display** to get C#- Click the UP/DOWN button **on D2 display** to get the target value as your need- Click SET to save the setting.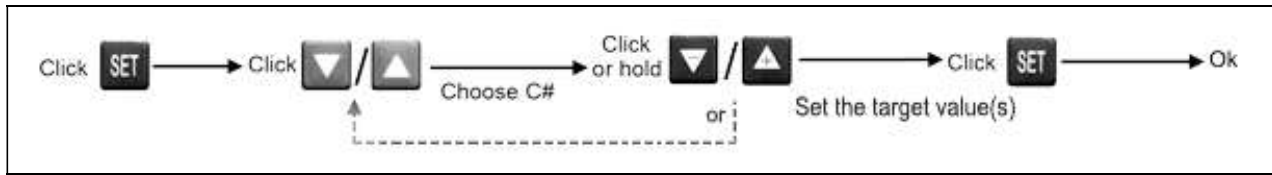

7. Error codes and simple solutions

- E0: Communication problem between the control panel and main board, check the cable or terminal is damaged or loose.
- E1: Over temperature protection, wait for the water temperature to drop, or appropriately increase the value of C0.
- E2: Low temperature protection, wait for the water temperature to rise, or appropriately reduce the value of C0.
- E3: Water flow problem, check the pump, hosing, water inlet & outlet, water volume or some things blocked inside.
- E4: Cooling system problem, check the fan, dedust the condenser and air inlet & outlet, and the space around the chiller.
- E5: High voltage protection.
- E6: Low voltage protection.
- E7: High room temperature alarm, low down the room temperature and replace the chiller in a well-ventilated area.

If problems persist, please seek professional help.

8. Equipment maintenance

- * Use the clean and soft fabric to clean the dust on the machine once a week.
- * Check the water capacity volume, tubes and connector, plug and cord once a week.
- * Thoroughly clean the anti-dust screen once a month.
- * Thoroughly clean the condenser and fan once a month.
- * Exchange the water in the tank of chiller once every 2 months at least.
- * Thoroughly clean the water-circulating loop if the water flow gets smaller and smaller.
- * Other regular maintenance is necessary.

To keep the machine working in good position, the maintenance period can be adjusted by user according to a specific working environment. For example, in a dusty place, maybe thoroughly clean the anti-dust screen, condenser and fan twice a month. However, in a dust-free plant, perhaps it would be cleaned one a year.

9. Warranty


Products manufactured and sold by JP SELECTA, s.a.u. are capital goods intended for industrial and nondomestic use. It is therefore that the guarantee applied is not regulated by the consumers and users law but by the commercial guarantee

laws. Thus, the guarantee of JP SELECTA, s.a.u. is **ONE YEAR** from the date of purchase and it covers, during the period stipulated in the commercial contract, any manufacturing defect or any hidden defect of the device.

The warranty does not cover damages caused by improper use or causes beyond J.P. SELECTA, s.a.u. Any manipulation of the device by personnel not authorized by J.P. SELECTA, s.a.u. automatically voids warranty benefits.

10. Warnings

1. To provide the clean cold recirculating water to the system and protect the circulating loop, it is best to install an extra water filter ahead of the inlet of chiller.
2. If the setting temperature can't be changed under any modes, or the default settings don't match your requirements, or the error codes occur more often, modifying the relevant parameters is needed.
3. To increase cooling efficiency and save energy, please make the hoses as short as you can, and sleeve/wrap the hoses with the thermal insulation tube/sheet, and place the chiller in the well-ventilated area.
4. When the first time to connect the chiller to your system, must double check the water/glycol volume to see if need to feed more water/glycol again after running the chiller for 3-5 minutes.
5. If you want some low temperature water or glycol to shock the hot object, please precool the water or glycol in the tank in advance.
6. When system setting is too mess or crash that reset the system is needed, press and hold SET Button 4 seconds, buzzer rings three times, reset the settings by default.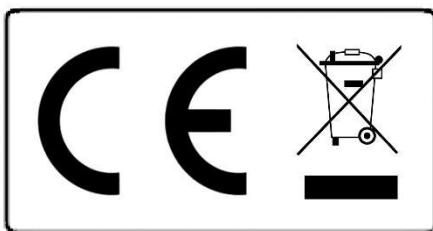

Notice to customers:

The product is made up of various components and various materials that must be recycled or, failing that, deposited in the corresponding debris removal sites when the product's life has been completed or when otherwise it is necessary to dispose of it. To do this, the end user who acquires the product must know the current regulations of each municipality and / or locality based on the waste electrical and electronic equipment. The user who acquires this product must be aware of and responsible for the potential effects of the components on the environment and human health as a result of the presence of hazardous substances. Never place the product in a conventional container of citizen scope if a previous dismantling and knowledge of the components that incorporates. If you do not know the procedure to follow, consult with the city council for more information.