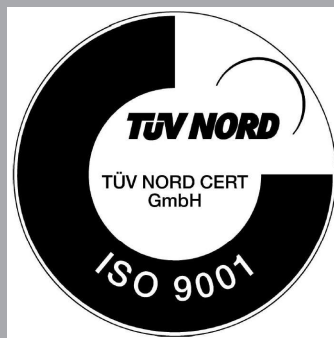
**STERILBIO****3000788**

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GENERAL INFORMATION

- 1) Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.
- 2) Do not install or use the equipment without reading this handbook before.
- 3) This handbook must always be attached to the equipment and it must be available for all users.
- 4) If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.
- 7) Do not use the equipment with liquids which can give off vapours capable of making explosive mixtures.

PACKING LIST

The standard equipment consists of the following components:

Piece	Code	Qty.
Sterilbio	3000788	1
Instruction manual	80106	1

TECHNICAL FEATURES

Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

Power consumption	Weight	Overall dimensions cm			Internal dimensions quartz tube cm	
		Height	Length	Width	Ø	Length
120 W	1.5 Kg	20	13	18	1.4	15

EQUIPMENT DESCRIPTION

Infra-red incinerators **STERILBIO** are made of ceramic parts with an internal tube of high quality quartz. No oxygen required and can be used in anaerobics environments. It eliminates problems of aerosols. It avoids bacterial contamination resulting from flame methods of sterilization. Side stand with four housing include for Kolle handles, needles, lances, dissecting forceps, etc...

Effective incineration of culture handles, needles, culture tubes and small instruments and of all organic material in 5/8 seconds at a temperature of 900°C.

INSTALLATION

It is not necessary any special installation, simply plug it to the mains.

Important:

During the first starting up, the equipment can gives off some fume because of the initial warming of the materials. That circumstance is completely normal and will disappear in a period of time.

CAUTION!!! IMPORTANT FOR YOUR SAFETY

Be sure that the voltage supply is the same as the one indicated on the characteristic plate of the equipment.

Do not use the equipment if it is NOT earthed.

If you have to change the plug, bear in mind the following:

Blue cable: Neutral.

Brown cable: Phase.

Yellow/green cable: Earth.

OPERATION

For starting up, you only have to press the main switch.

After 7 or 8 minutes, a temperature of 900°C will be reached inside the tube and the equipment is ready for use.

To sterilize an instrument, introduce it by the central tube, and after 5/8 seconds, the end of the instrument will be sterilized because of the high temperature.

The sterilizer can be on without stopping for about 3 or 4 hours, as temperature stops raising when it reaches 950°C. But if it is not necessary, we recommend to disconnect it after using and foresee the connection 10 minutes before the use. Doing that, the live of the sterilizer will be longer.

CAUTION
**The ceramic tube envelope can reach
temperatures of about 80°C**

GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damages caused by an incorrect use or causes beyond the control of J.P. SELECTA, s.a.u.

Any manipulation of the equipment by unauthorized personnel by J.P. SELECTA, s.a.u. cancels the guarantee automatically.

Notice to customers:

The product is made up of various components and various materials that must be recycled or, failing that, deposited in the corresponding debris removal sites when the product's life has been completed or when otherwise it is necessary to dispose of it. To do this, the end user who acquires the product must know the current regulations of each municipality and / or locality based on the waste electrical and electronic equipment. The user who acquires this product must be aware of and responsible for the potential effects of the components on the environment and human health as a result of the presence of hazardous substances. Never place the product in a conventional container of citizen scope if a previous dismantling and knowledge of the components that incorporates. If you do not know the procedure to follow, consult with the city council for more information.