CONTADOR DE COLONIAS DIGITAL-LED
DIGITAL CELL COUNTER-LED

DIGITAL-S 4905002

INFORMACIÓN GENERAL

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GENERAL INFORMATION

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.*
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) *If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the equipment with liquids which can give off vapours capable of making explosive mixtures.*

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CONTENTS

GENERAL INFORMATION
CONTENTS
PACKING LISTS
TECHNICAL FEATURES
EQUIPMENT DESCRIPTION
INSTALLATION
OPERATION
MAINTENANCE
GUARANTEE
SPARE PARTS

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PACKING LIST

The standard equipment consists of the following components:

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EQUIPMENT DESCRIPTION

Colony counter characteristics:

- Allows petri dishes up to 100mm.
- Working area with stable light.
- Counts by pressing with a pointed sharpe object as a pencil, pince, ...
- Selectable background black / white colour to get a better contrast
- Magnifier 120 mm diameter 2x. Adjustable height and rotation angle.
- Digital count up to 999 colonies.
- Bip sound indicator on any count.
- Reset count push button.
- Removable transparent dish.
- Selectable counting dish.
- External dimensions: 25 x 23 x 30 cms
- Weigth: 3,5 Kg
- Consumption: 10 W
- Power supply voltage from 110 to 230V.

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INSTALLATION

- Place the counter on a flat and leveled surface.
- A mains plug must be near.
- Screw the blade lent holder (3) into the screwed hole of the colony counter.
- Fix the holder bosshead (4) with lent (5) in the blade (3). The bosshead is adjustable, allowing the lent to be raised or lowered to your preference. This lent can rotate a full 360°.
- Place the transparent base into the holder capsule hole (6).
- Place the Wolffhuegel counting plate (7) on the holder capsule base (6).
- Plug the power supply cable into a 220V single-phase socket.

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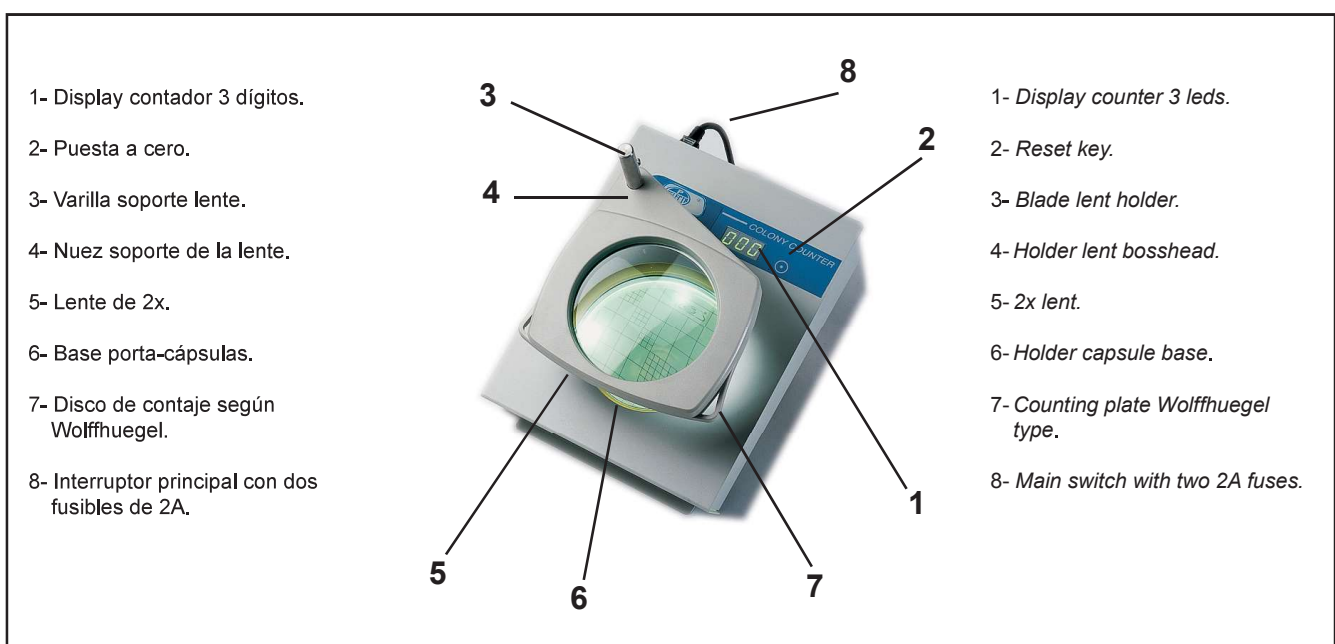
Be sure that the voltage supply match the one indicated on the characteristics plate.

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Do not use the equipment if it is NOT EARTHED.

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If you have to change the plug bear, in mind that the yellow/green cable is the protective earth and it is connected to the metallic envelope of the equipment

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OPERATION

- *Turn on the main switch (8) placed at rear part.*
- *The digital counter display (1) must show 000, if not, press the "reset" key (2).*
- *Colony counting is done by pressing the capsules (6) with any sharp-pointed tool (like a pen, puch, etc...).*
- *Once the counting is finished, press the "reset" key (2).*

MAINTENANCE

Before removing the covers, disconnect the equipment from mains.

The manipulation of the internal electronic circuits of the plate by unauthorized personnel can cause damage of difficult solution. Make sure to take it to one of J.P.Selecta, s.a.u. authorized technical services.

CLEANING:

Use alcohol or similar solvents applied with no abrasive wool to clean lens, plastic parts and envelope. Do not use inorganic solvers which can damage plastic parts

GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damages caused by an incorrect use or causes beyond the control of J.P. SELECTA, s.a.u.

Any manipulation of the equipment by unathorized personnel by J.P. SELECTA,s.a.u. cancels the guarantee automatically.

SPARE PARTS

To guarantee the safety of the equipment, all spare parts must be purchased from J.P.SELECTA, s.a.u.

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