**CC-1431**

**7121431**

**CONTADOR DE CÉLULAS  
CELL COUNTER**

## Indice

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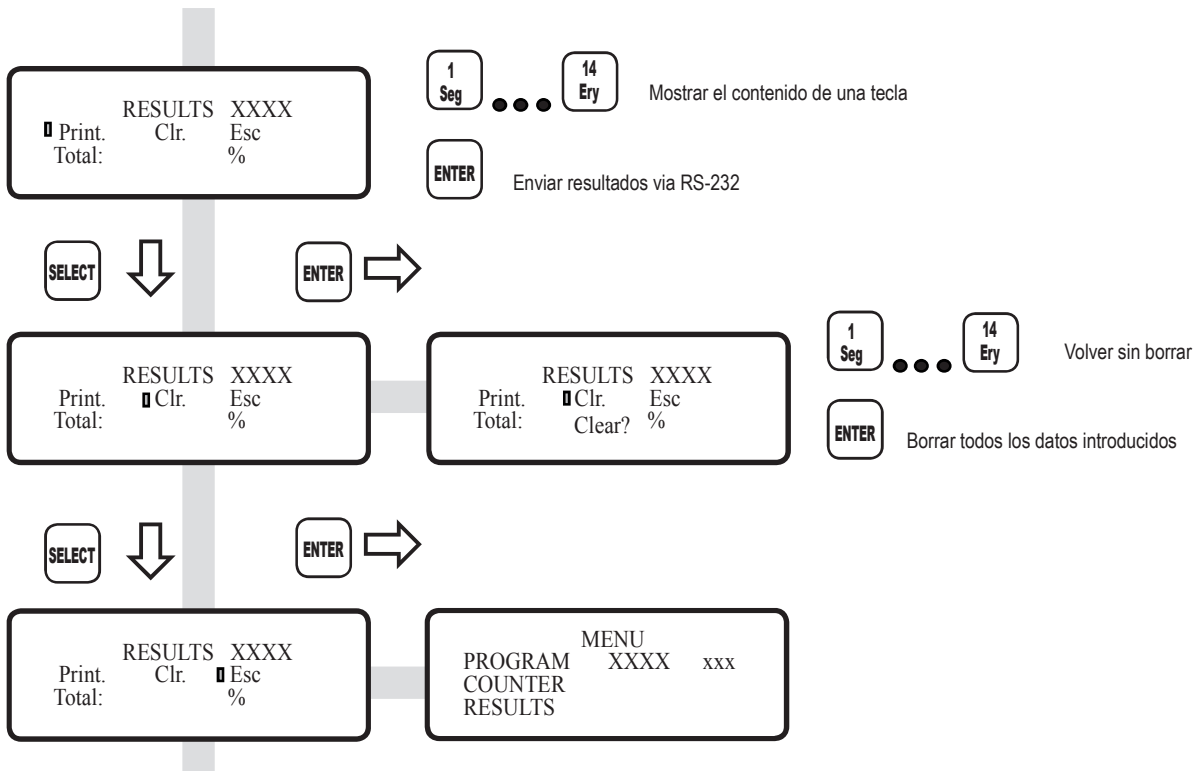
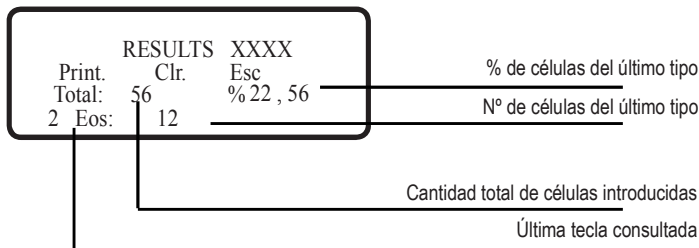
## 1 GENERAL INFORMATION

- 1) Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.
- 2) Do not install or use the equipment without reading this handbook before.
- 3) This handbook must always be attached to the equipment and it must be available for all users.
- 4) If you have any doubts or enquiries, please contact your supplier or J.P. Selecta's technical service.
- 5) IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.
- 6) If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.

## 2 PACKAGE LIST

Components supplied with package:

- Main unit.
- 12V power adapter
- USB serial wire connection
- Software (on demand)
- Instructions manual

## 3 TECHNICAL DATA

- USB output (ask for software to [service@jpselecta.es](mailto:service@jpselecta.es))
- 3 selectable languages: Spanish, English and French.
- 3 different acoustic beeps
  - «Key push»;
  - «End of count»;
  - «Disabled key»
- Four programs:
  - LEUCO: leucocytes counting
  - MYELO: MieloBlastes counting
  - RETI: Reticulocytes counting
  - FREE: Free counting
- Display LCD 16x4 characters.
- 16 keys (14 for cell counting).

## 4 DESCRIPTION

The Cell counter CC-1431 is designed to simplify the manual recount of sanguine cells.

Their ergonomic design allows a comfortable use during long spaces of time.

Basically the operation consists in that each pulsation of a key (that means a cell) adds to an accumulated total. When you reach the total (for example 100 cells) the counter emits a triple beep and it doesn't admit more pulsations. The result shows the subtotal of each cell and the percentage of each cell on the total.

In the program «RETI» «fields» are counted instead of cells, which contains a pre-set number of cells.

It includes four different programs:

**LEUCO:**

- Counting of Leucocytes.
- 14 types of different cells.
- 3 keys for cells whose counts are not accumulated to total («11. Normo» «12.Tox Gr» and «14. Ery»).
- Total of cells (end of the count) seleccionable from 100 to 1000 in steps of 100.

**MYELO:**

- Counting of Myelocytes
- 14 types of cells that accumulates to the total.
- Total of cells seleccionable from 100 to 1000 in steps of 100.

**RETI:**

- Counting of Reticulocytes.
- Number of cells for field configurable between 50 and 250.
- One key for reticulocytes fields.
- One key for eritrocytes that doesn't accumulate in the total.
- Total (end of count) fixed in 1000.

**FREE:**

- Configurable counting.
- Each key can be enabled or disabled for the counting.
- Total of cells (end of the count) seleccionable from 100 to 1000 in steps of 100.

In the LCD display of data entry, the following is displayed:

- The count of cells of the last entered type.
- The total of counted cells
- The «count end»

In any moment of the counting cells, the results screen can be selected without danger of losing data.

## 5 EXTERNAL WIRING

Plug the power adapter.

Pay attention with the plug from AC adapter.

To download data to PC computer:

- Install the software.
- Connect the serial wire.

To get the USB connection software, contact with JP Selecta, s.a.u. at [service@jpselecta.es](mailto:service@jpselecta.es)

If power adapter is replaced, be sure that it matches the following specifications:

Output voltage: 12 V  
Output current available: 400mA

## 6 WORKING WITH COUNTER

Power-up the cell counter by the rear panel switch.

Select some of the counter programs:

### LEUCO program:

The LEUCO program uses 14 keys. Eleven of them, when pressed, increase the total counting and partial counting corresponding key.

The other three ones «11. Normo» «12.Tox Gr» and «14. Ery» increase the value of the corresponding key but not the total one and therefore they don't have influence in the moment to calculate the percentage.

### MYELO program:

In this program all the keys, when being pressed, increase the value of the corresponding key and at the same time increase the total value.

### RETI program:

In this program there are only two active keys:  
«13. Reti» and «14. Ery».

Key «13. Reti» increases their own value in 1 when being pressed and it doesn't act on the total.

Key «14. Ery» increases their own value in 50,100,150,200 or 250 according to the elected configuration by the user in the SETTINGS MENU and, at the same time, it increases the total in this same quantity.

The «count end» is 1000.

### FREE program:

In the FREE program all the keys are active and the user can choose from the screen FREE PROGRAM which keys increase the total and which don't.

Abbreviations used in the keys:

Seg:	Segmentados o Granulocitos.	Segmented or Granulocytes
Eos:	Eosinofilos	Eosinophiles.
Lym:	Linfocitos	Lymphocytes.
Mono:	Monocitos	Monocytes.
Stab:	Bandas	Stab
Baso:	Basofilos	Basophiles
Met My:	MetaMielocitos	MetaMyelicocites
Myelo:	MieloBlastos	MieloBlastos
At Ly:	Linfocitos Atípicos.	Atypical Lymphocytes
My Bi:	Mielocitos Binucleados	Binucleated myelicocytes
Normo:	NormoBlastos	NormoBlastes
Tox. Gr.:	Granulocitos toxicos	Toxical granulocytes
Reti:	Reticulocitos	Reticulocytes
Ery:	Eritrocitos / EritoBlastos	Erythrocytes / Erythroblasts

## 7 MAIN MENU

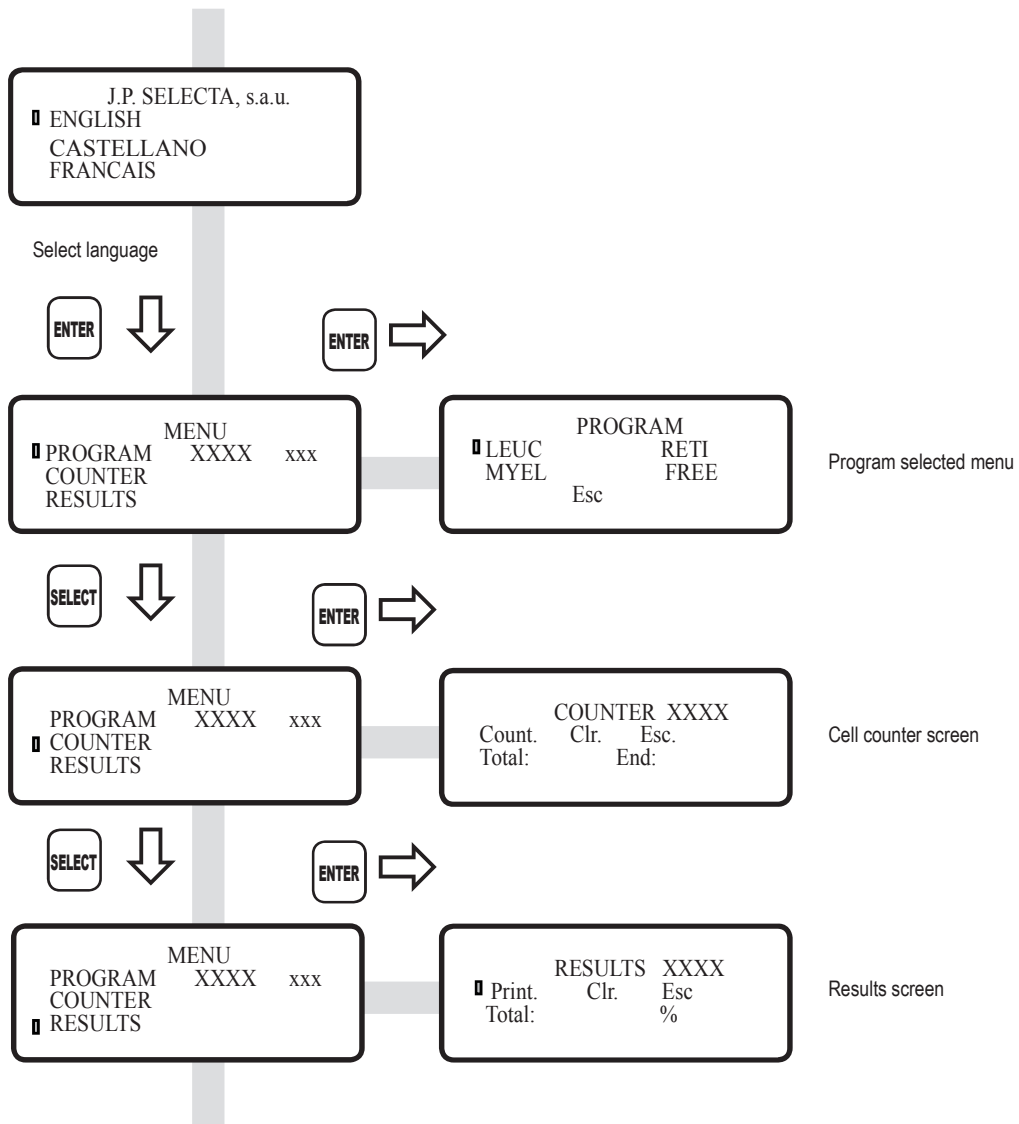
After switching-on the cell counter, the first choice is to select the language from Spanish, English or French.

Once language is selected, main menu appears. Select one of the following sub-menus:

- Program menu
- Counter screen.
- Results screen.

To choose one item from menu, select it first by pressing **SELECT**

and then press key **ENTER** to accept.

```

                J.P. SELECTA, s.a.
    ▣ ENGLISH
      CASTELLANO
      FRANCAIS
    
```

**Language selection screen:**

Select with **SELECT** and accept with **ENTER**

```

                MENU
    ▣ PROGRAM   XXXX   xxx
      COUNTER
      RESULTS
    
```

**Main menu screen:**

Select with **SELECT** and accept with **ENTER**

```

                PROGRAM
    ▣ LEUC      RETI
      MYEL     FREE
      Esc
    
```

**Program menu screen:**

Select with **SELECT** and accept with **ENTER**

```

                COUNTER XXXX
    ▣ Count.  Clr.  Esc.
      Total:   End:
    
```

**Counter screen:**

Select with **SELECT** and accept with **ENTER**

```

                RESULTS XXXX
    ▣ Print.  Clr.  Esc
      Total:   %
    
```

**Results screen:**

Select with **SELECT** and accept with **ENTER**

```

    SET   xxx
    
```

**«End counting» setting screen:**

On program RETI, select the number of cells for field Ery

```

                PROGRAM FREE
    1  2  3  4  5
    6  7  8  9  10
    11 12 13 14
    
```

**Free program setting screen:**

Select the key enabled to add in total counting

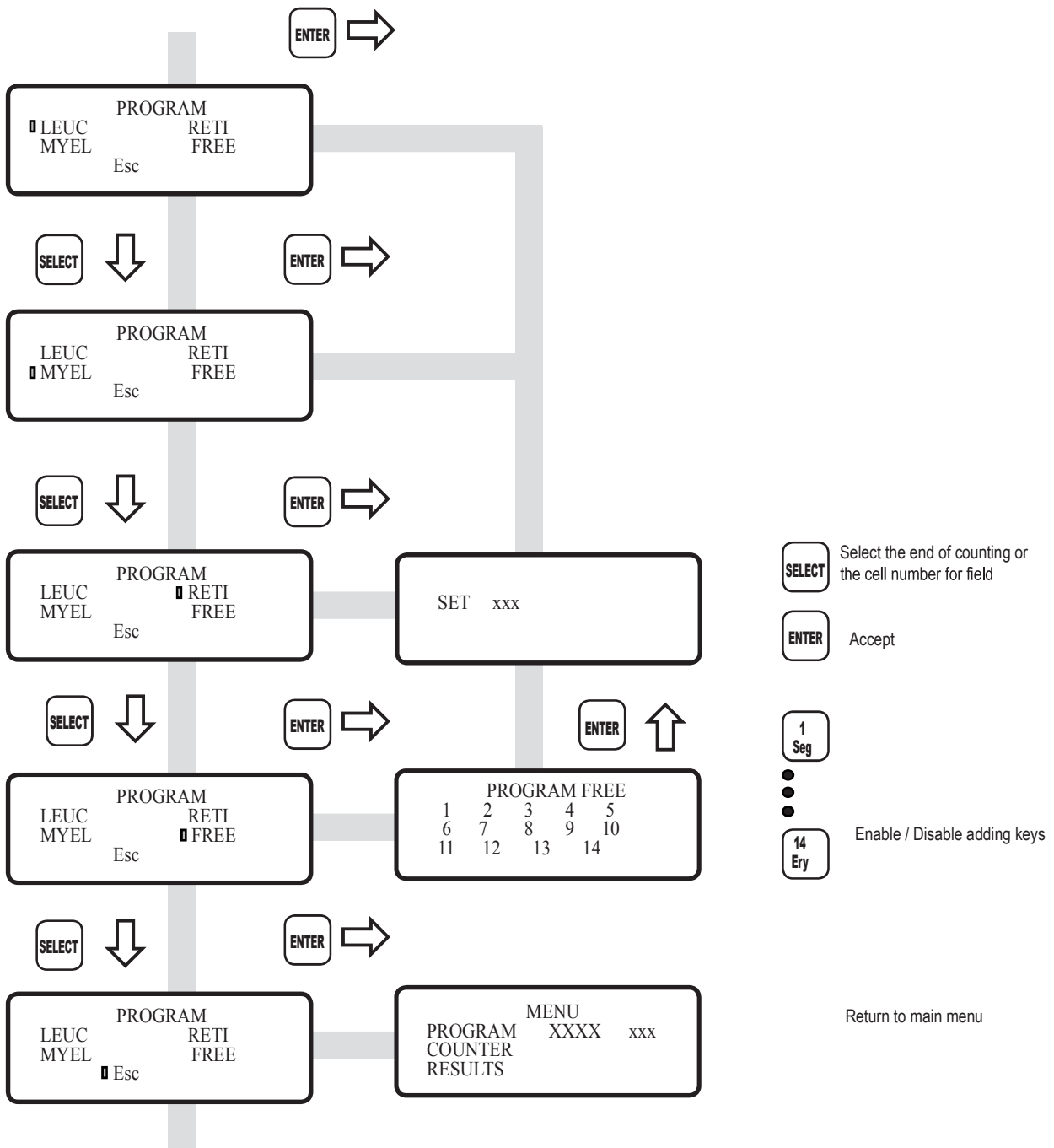
Select with **SELECT** and accept with **ENTER**

## 8 PROGRAMS MENU

In the programming screen, choose one of the four programs that includes the cells counter.

The following features are configurable:

- Total number of cells that we want to count in the programs LEUC, MYEL and FREE.
- Number of cells for field in the program RETI.
- Keys that increase total in the program FREE.

## 9 SCREEN COUNTER

In the counter function the sanguine cells counter is carried on. Each one of the keys has a name and a number associated.

Acoustic warnings:

Short beep: pressed key that it adds to the total.

Example: it programs the RETI key «14. Ery»

Long beep: Inactive key in the used program.

Example: it programs the RETI key «1. Seg»

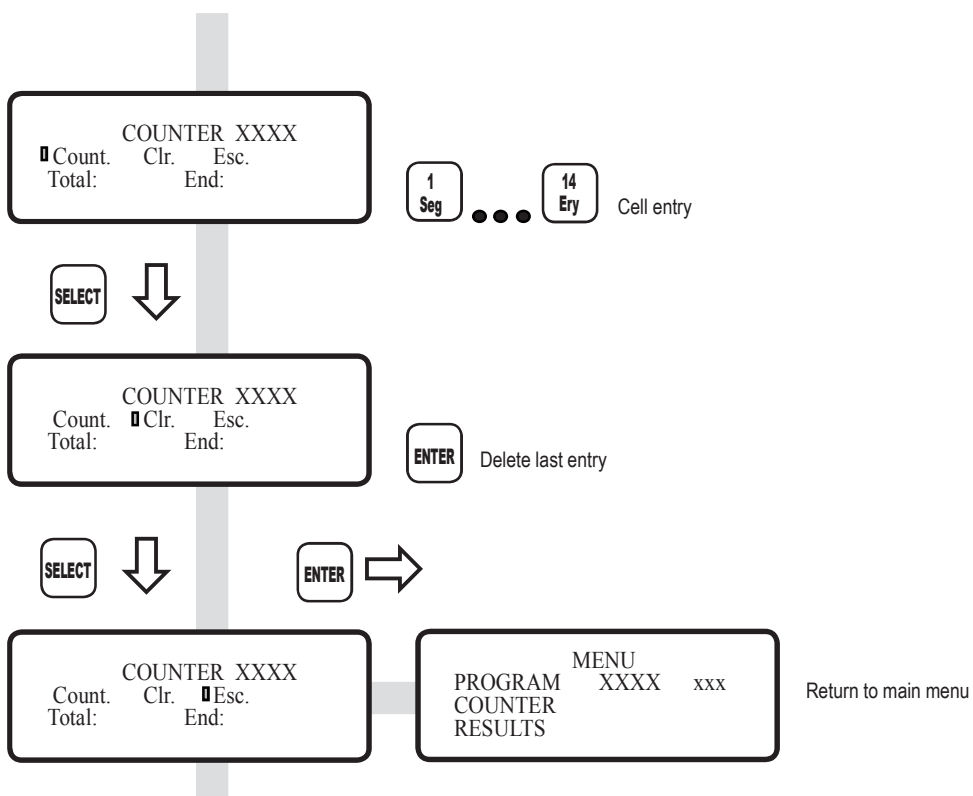
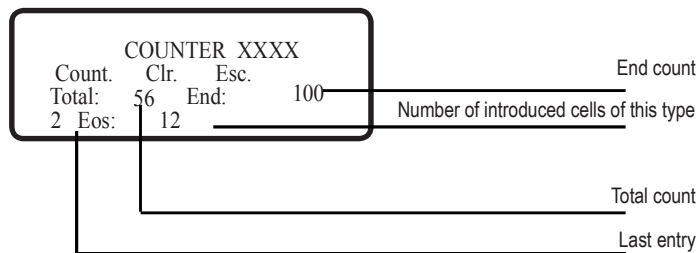
Two short beeps: Active key that doesn't add to the total and has a separate count.

Example: it programs the RETI key «13. RETI»

Triple short beep: end of count has been reached.

In the screen we can read the total number of counted cells up to the moment, the total number of cells to count, the last introduced cell and the quantity that has been introduced until the moment.

If pressing a wrong key, the program allows to correct the last introduced key.

## 10 RESULTS SCREEN

This function shows the results in screen and the current state of the count.

When pressing a key corresponding to a cell it will appear:

- The total number of cells of this type.
- The percentage of this cell on the total.

If «--,-» appears, it means that this cell does not affect the percentage.

The total amount of counted cells is always visible in this screen.