**REACTOR UNIVERSAL PARA LABORATORIO
UNIVERSAL REACTOR FOR LABORATORY**

RV-12

1001746

INDICE

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INDEX

1 General information	8
2 Features	9
3 Technical data	10
4 Parts and accessories description	11
5 Assembling	12

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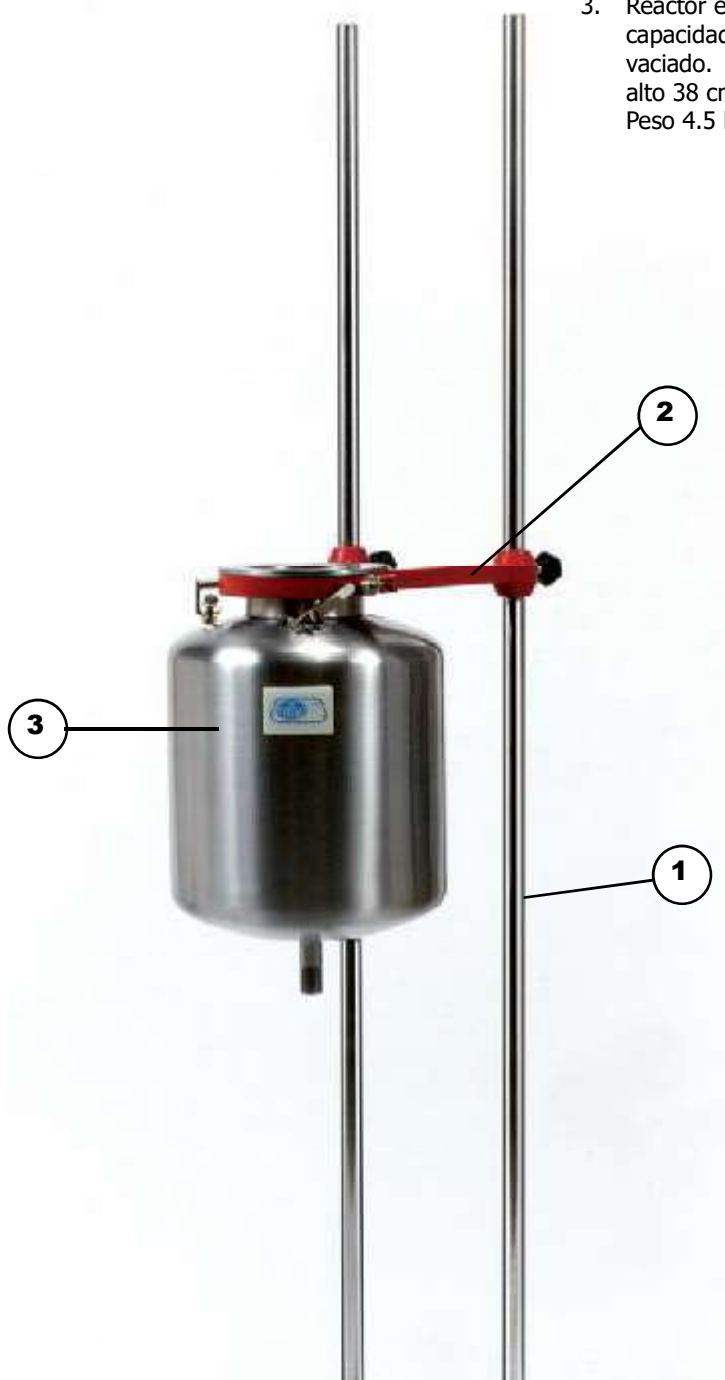
ENGLISH VERSION

1. General information

- Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.
- Do not install or use the equipment without reading this handbook before. This manual makes part of the equipment and must remain available to the users
- If you have any doubt inquiry to your J.P. Selecta products supplier.
- Do not use the equipment in explosive or corrosive atmospheres or with liquids which can produce explosive mixtures..

2. Features

This is a universal reactor of distillation or backflow-concentration suitable to laboratory and pilot plant operations. It has a standard entering necked able to lock of gless reactors DN100 type QVF or similar. Made in stainless steel AINSI 316 , 12 litres capacity and drainage manual valve able to be connected to a hose.

3. Technical data

Capacity	12 litres
Dimensions (Height x Diametre)	38 x 25 cm
Weigth	4.5 Kg
Inner diametre reactor necked	100 mm
External platform diametre	138 mm

4. Description of parts and accessories


Standard equipment

1. Standard support.
2. Clamp with support ring.
3. Reactor made in stainless steel AISI 316, capacity 12 litres with draining tap.
External dimensions:
height 38 cm, diameter 25 cm.
weight 4.5 kg.


Cod. 1001748

Cod. 1001747

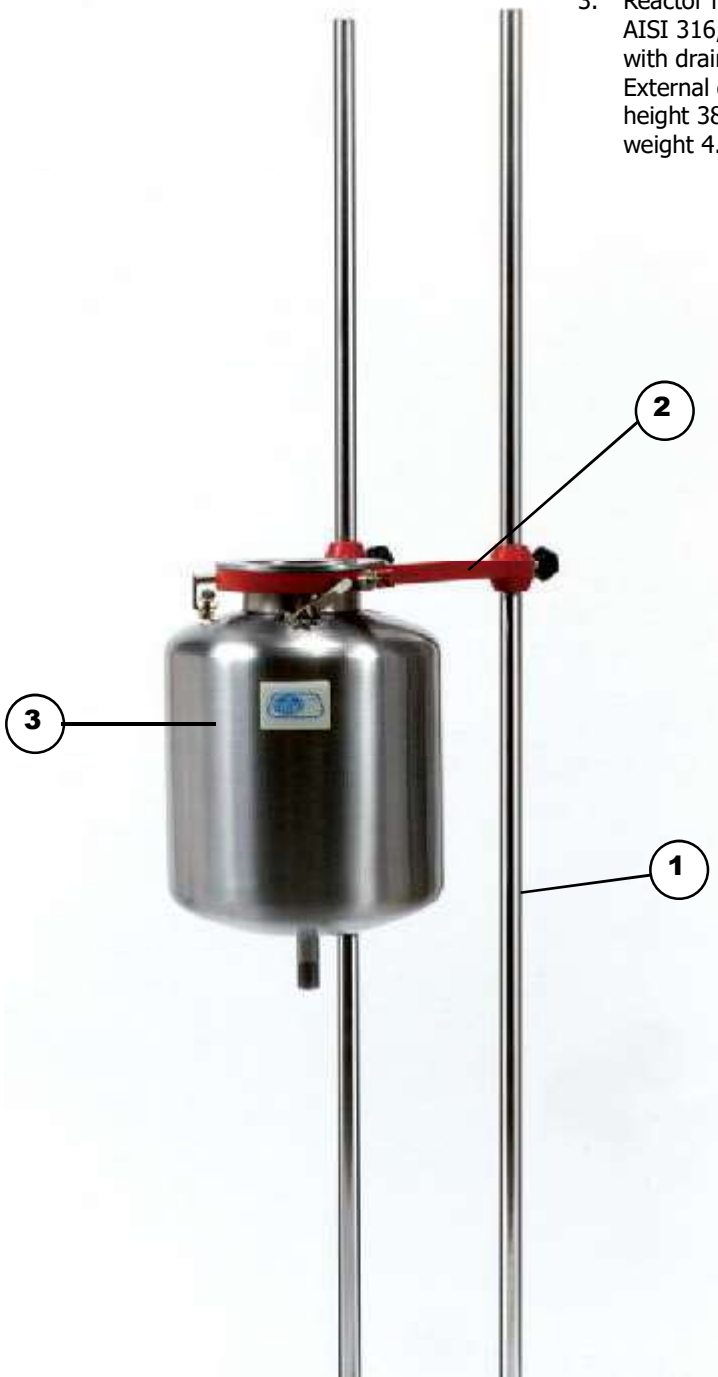Cod. 1001746

Accessories

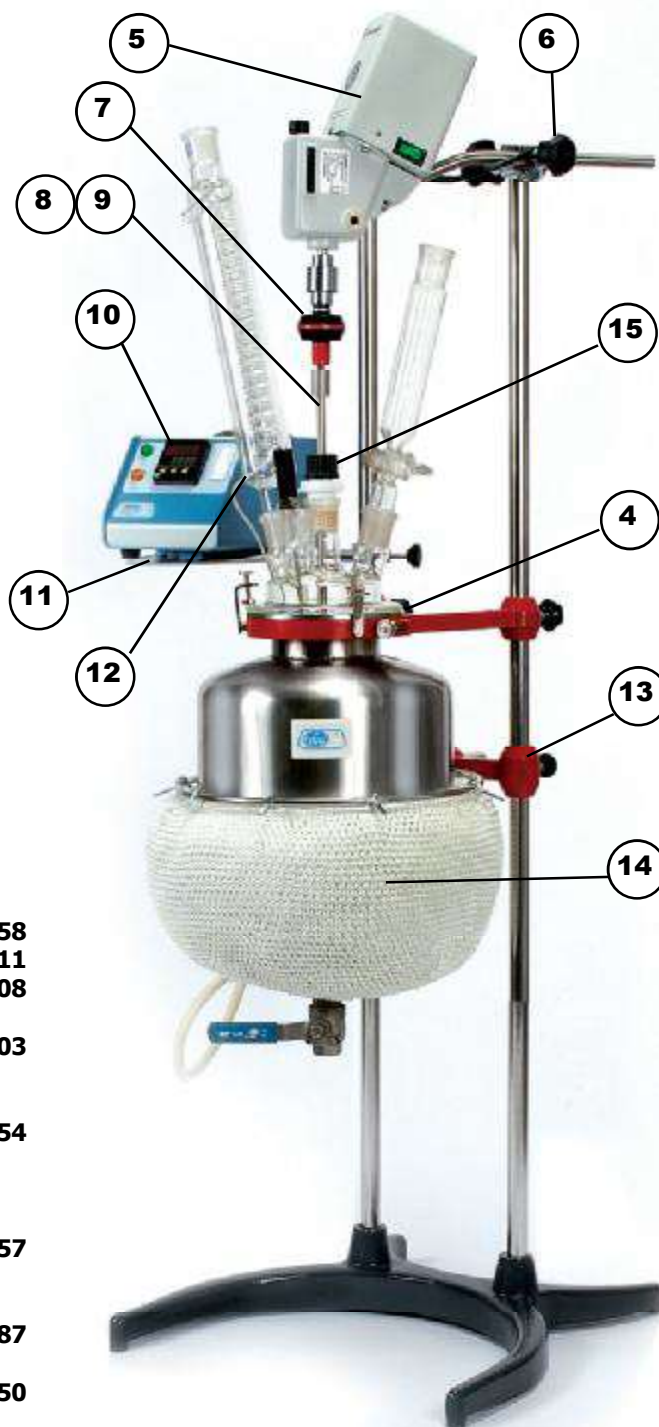- | | | |
|-----|---|---------------------|
| 4. | Clasp for fastening normalized flat gaskets for reactor. | Cod. 1001758 |
| 5. | Stirrer RZR-1. sdjustable speed | Cod. 5050111 |
| 6. | Double bosshead Maxim. | Cod. 7000408 |
| 7. | Flexible shaft holder for 6 to 10Ø stirring shafts | Cod. 5050903 |
| 8. | Half-moon shape stirring bar stainless steel AISI 304. Length 90 x shovel width 12 cm | Cod. 1001754 |
| 9. | Helix shape stirring shaft 8Ø in stainless steel AISI 304 and PTFE. Length 90 x shovel width 4 cm | Cod. 1001757 |
| 10. | Electronic controller Electemp width connector for inner temperature control. | Cod. 3000887 |
| 11. | Support plate with bosshead to regulator Electemp | Cod. 1001750 |
| 12. | Pt100 temperature probe 200 mm length x 4Ø mm with connection cord to the Electemp. | Cod. 1001749 |
| 13. | Ring holder with double bosshead for heating mantle | Cod. 1001751 |
| 14. | Heating mantle with double heating elements for temperatures up to 400°C. Power: 1400W | Cod. 3001752 |
| 15. | Guide NS29/32 adjustable to 8Ø stainless steel shafts made in PTFE/iton gasket | Cod. 5050909 |

5. Assembling

- 1 Place support (1)
- 2 Assemble support (13)
- 3 Place heating mantle (14)
- 4 Assemble clamp (2)
- 5 Place reactor (3)
- 6 Adjust the height of the reactor until making contact it to the heating mantle to manta (14) and level the support (2)
- 7 Assemble the stirring shaft (8)/(9)
- 8 Assemble the borosilicate lid 100Ø QVF type or similar
- 9 Adjust the system with the clamp (4) of three points
- 10 Assemble the guide NS29/32 to PTFE with adjustable gaskets (cod. 5050909) and adjust it.
- 11 Assemble the flexible shaft holder (7) cod. 5050907 or the hight viscosity one cod. 5050904
- 12 Place the clamp (6)
- 13 Asemble the stirrer (5) and adjust the shaft holder (7)
- 14 Assemble the support of the temperature regulator (11)
- 15 Place the temperature regulator (10)
- 16 Insert the temperature probe (12)

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